



Explorer Plus™ Balances

Instruction Manual



Version History

| Date | Version | Description |
|----------|---------|---|
| 2025/6/1 | A | <ul style="list-style-type: none"><li data-bbox="667 342 868 371">• Initial release |
| | | |
| | | |

TABLE OF CONTENTS

| | | |
|----------|---|-----------|
| 1 | INTRODUCTION | 8 |
| 1.1 | Description..... | 8 |
| 1.2 | Features | 8 |
| 1.2.1 | Display | 8 |
| 1.2.2 | Modular Design..... | 8 |
| 1.2.3 | Digital Loadcell with Advanced Environmental Sensors | 8 |
| 1.2.4 | Removable Draftshield Design | 8 |
| 1.2.5 | Automatic Draftshield Doors..... | 8 |
| 1.2.6 | Status Lights and Sample Spotlights..... | 8 |
| 1.2.7 | HID Data Transfer Function..... | 8 |
| 1.2.8 | Time Synchronization NTP (Network Time Protocol)..... | 8 |
| 1.2.9 | Fingerprint login System (Accessory)..... | 8 |
| 2 | Safety Precautions | 9 |
| 2.1 | Intended Use..... | 9 |
| 2.2 | Definition of Signal Warnings and Symbols | 9 |
| 2.3 | Safety Notes | 9 |
| 3 | Installation | 11 |
| 3.1 | Unpacking..... | 11 |
| 3.2 | Installing Components | 11 |
| 3.2.1 | For 0.01mg, 0.1mg, and 1mg Models..... | 11 |
| 3.2.2 | For 0.01mg, 0.01g, and 0.1g Models..... | 11 |
| 3.2.3 | For High-Capacity Models..... | 12 |
| 3.2.4 | Selecting a Location | 12 |
| 3.2.5 | Level the Balance Before Use | 13 |
| 3.3 | Connecting Power | 13 |
| 3.4 | Turn on the balance..... | 14 |
| 3.5 | Adjust the viewing angles | 14 |
| 3.6 | Disconnect the terminal from the weighing base | 14 |
| 3.7 | Remove Draftshield from the Base | 15 |
| 3.8 | Connecting the Interface | 15 |
| 3.9 | Overview of Parts and Features –Draft Shield Models..... | 16 |
| 3.9.1 | Side View..... | 16 |
| 3.9.2 | Rear View..... | 16 |
| 3.10 | Overview of Parts and Features –Non-Draft Shield Models | 17 |
| 3.10.1 | Side View..... | 17 |
| 3.10.2 | Rear View | 17 |
| 3.11 | Overview of High-Capacity Models | 17 |

| | | |
|----------|---|-----------|
| 3.11.1 | Side View..... | 17 |
| 3.11.2 | Rear View..... | 18 |
| 3.12 | Overview of Main Screen and Connectivity..... | 18 |
| 3.13 | Main Screen Features..... | 19 |
| 3.14 | Warning and Notifications..... | 19 |
| 3.14.1 | Calibration Failure..... | 19 |
| 3.14.2 | Leveling Required..... | 20 |
| 3.14.3 | Password Expiration..... | 20 |
| 3.15 | Hotkeys..... | 20 |
| 3.15.1 | Standard Hotkeys..... | 21 |
| 3.15.2 | Additional Hotkeys (Add+)..... | 22 |
| 3.16 | Shortcut Buttons..... | 22 |
| 3.17 | HID Connections..... | 22 |
| 3.17.1 | English Keyboard..... | 22 |
| 3.17.2 | French Keyboard..... | 23 |
| 4 | Weighing Application Modes..... | 24 |
| 4.1 | Introduction..... | 24 |
| 4.2 | Basic Weighing..... | 25 |
| 4.2.1 | Weighing Result..... | 25 |
| 4.2.2 | Reference Fields – Basic Weighing..... | 25 |
| 4.2.3 | Item Setting..... | 26 |
| 4.2.4 | Minimum Weight..... | 26 |
| 4.2.5 | Set minimum weight value..... | 26 |
| 4.2.6 | Auto Tare..... | 27 |
| 4.2.7 | Sample Library..... | 27 |
| 4.2.8 | Printing Settings..... | 28 |
| 4.2.9 | Shortcut Buttons..... | 28 |
| 4.3 | Parts Counting..... | 29 |
| 4.3.1 | Application buttons..... | 29 |
| 4.3.2 | Begin Parts Counting..... | 30 |
| 4.3.3 | Accumulate the data for Parts Counting..... | 30 |
| 4.3.4 | Sample Library – Parts Counting..... | 31 |
| 4.3.5 | Printing Settings..... | 32 |
| 4.4 | Check Counting..... | 32 |
| 4.4.1 | Application buttons and Reference Field..... | 33 |
| 4.4.2 | Begin Check Counting..... | 33 |
| 4.4.3 | Accumulate the data for Check Counting..... | 35 |
| 4.4.4 | Sample Library – Check Counting..... | 35 |

| | | |
|--------|--|-------------------------------------|
| 4.4.5 | Printing Settings | 36 |
| 4.5 | Percent Weighing | 36 |
| 4.5.1 | Application buttons and Reference Field | 37 |
| 4.5.2 | Begin Percent Weighing..... | 37 |
| 4.5.3 | Accumulate the data for Percent Weighing | 38 |
| 4.5.4 | Sample Library – Check Counting..... | 39 |
| 4.5.5 | Printing Settings | 39 |
| 4.6 | Check Weighing | 40 |
| 4.6.1 | Reference Field..... | 40 |
| 4.6.2 | Begin Check Weighing..... | 41 |
| 4.6.3 | Sample Library – Check Weighing | 42 |
| 4.6.4 | Printing Settings | 43 |
| 4.7 | Dynamic Weighing..... | 43 |
| 4.7.1 | Application buttons | 44 |
| 4.7.2 | Begin Check Weighing..... | 44 |
| 4.7.3 | Sample Library – Dynamic Weighing | 45 |
| 4.7.4 | Printing Settings | 45 |
| 4.8 | Totalization | 46 |
| 4.8.1 | Application buttons | 46 |
| 4.8.2 | Begin Totalization..... | 46 |
| 4.8.3 | Accumulate the data for Totalization | 47 |
| 4.8.4 | Sample Library – Totalization | 47 |
| 4.8.5 | Printing Settings..... | 48 |
| 4.9 | Formulation..... | 48 |
| 4.9.1 | Application buttons | 49 |
| 4.9.2 | Begin Free Formulation..... | 50 |
| 4.9.3 | Begin Recipe Based Formulation..... | 51 |
| 4.9.4 | Begin Recipe Based Formulation..... | Error! Bookmark not defined. |
| 4.9.5 | Begin Free Formulation..... | Error! Bookmark not defined. |
| 4.9.6 | Printing Settings..... | Error! Bookmark not defined. |
| 4.10 | Differential | 52 |
| 4.10.1 | Application Buttons..... | 53 |
| 4.10.2 | Begin Differential | Error! Bookmark not defined. |
| 4.10.3 | Item Settings – Differential | Error! Bookmark not defined. |
| 4.10.4 | Begin Differential | 54 |
| 4.10.5 | Printing Settings..... | 54 |
| 4.11 | Density Determination..... | 55 |
| 4.11.1 | ITEM SETTINGS | 56 |

| | | |
|---------|---|----|
| 4.11.2 | Reference Fields | 57 |
| 4.11.3 | Measuring the Density of a Solid Using Water | 57 |
| 4.11.4 | Measuring the Density of a Buoyant Material Using Water | 59 |
| 4.11.5 | Measuring the Density of a Solid Using an Auxiliary Liquid | 60 |
| 4.11.6 | Measuring the Density of Porous Material Using Oil | 61 |
| 4.11.7 | Measuring the Density of Porous Material Using Oil and Auxiliary Liquid | 62 |
| 4.11.8 | Measuring the Density of a liquid using a Calibrated Sinkers | 64 |
| 4.11.9 | Sample Library – Density Determination | 65 |
| 4.11.10 | Printing Settings | 67 |
| 4.12 | Peak Hold | 67 |
| 4.12.1 | ITEM SETTINGS | 67 |
| 4.12.2 | Reference Fields | 70 |
| 4.12.3 | Begin with Peak Hold | 70 |
| 4.12.4 | Begin with Display Hold | 71 |
| 4.12.5 | Sample Library – Totalization | 71 |
| 4.13 | Pipette Adjustment | 73 |
| 4.13.1 | ITEM SETTINGS | 74 |
| 4.13.2 | Application Buttons | 76 |
| 4.13.3 | Reference Fields | 76 |
| 4.13.4 | Test Method | 76 |
| 4.13.5 | Begin Pipette Adjustment in Manual Mode | 79 |
| 4.13.6 | Begin Pipette Adjustment in Auto Mode | 81 |
| 4.13.7 | Printing Settings | 83 |
| 4.14 | SQC | 83 |
| 4.14.1 | Reference Fields | 83 |
| 4.14.2 | Begin SQC | 83 |
| 4.14.3 | Acceptance Criterion | 88 |
| 4.14.4 | Determination of -T1, -T2, +T1, and +T2 | 89 |
| 4.14.5 | Printing Settings | 89 |
| 4.15 | Fill Weight Variation | 90 |
| 4.15.1 | Application buttons | 90 |
| 4.15.2 | Reference Fields | 90 |
| 4.15.3 | Item Settings – Fill Weight Variation | 91 |
| 4.15.4 | Begin Tests with USP <2091> | 91 |
| 4.15.5 | Begin Tests with Customized +/- Limits | 93 |
| 4.15.6 | Printing Settings | 95 |
| 4.16 | Flow Rate Control | 95 |
| 4.16.1 | Item Settings – Check Weighing | 96 |

| | | |
|----------|--|------------|
| 4.16.2 | Begin Flow Rate Control | 99 |
| 4.16.3 | Begin with Target Rate Mode | 101 |
| 4.16.4 | Sample Library – Flow Rate Control..... | 103 |
| 5 | Menu Settings..... | 106 |
| 5.1 | Menu Navigation | 106 |
| 5.2 | Menu Structure..... | 106 |
| 5.3 | Quick Setup | 107 |
| 5.3.1 | Language..... | 107 |
| 5.3.2 | Time Synchronization/ Network Server | 107 |
| 5.3.3 | Date and Time..... | 108 |
| 5.3.4 | Automatic Calibration..... | 108 |
| 5.3.5 | User Management | 108 |
| 5.3.6 | System Log..... | 108 |
| 5.3.7 | Balance Information..... | 109 |
| 5.3.8 | Motorized Leveling | 109 |
| 5.3.9 | Repeatability Test..... | 110 |
| 5.4 | Calibration | 111 |
| 5.4.1 | Calibration Settings | 111 |
| 5.4.2 | Internal Calibration | 112 |
| 5.4.3 | Span Calibration..... | 113 |
| 5.4.4 | Calibration History | 114 |
| 5.5 | Balance Setup..... | 114 |
| 5.5.1 | Language..... | 115 |
| 5.5.2 | Time Synchronization/ Network Server | 115 |
| 5.5.3 | Date and Time..... | 115 |
| 5.5.4 | Balance Name..... | 115 |
| 5.5.5 | Change Password | 116 |
| 5.5.6 | Fingerprint/ Fingerprint Setting | 116 |
| 5.5.7 | Filter Level | 118 |
| 5.5.8 | Stability Indicator Range | 118 |
| 5.5.9 | Auto Zero Tracking | 118 |
| 5.5.10 | Gross Indicator..... | 118 |
| 5.5.11 | Graduation | 118 |
| 5.5.12 | Ionizer | 119 |
| 5.5.13 | Approved Mode..... | 119 |
| 5.6 | Auto Door..... | 120 |
| 5.7 | Sensor..... | 120 |
| 5.8 | System Log..... | 121 |

| | | |
|------------|--|------------|
| 5.9 | ECO | 121 |
| 5.9.1 | Power Saving | 122 |
| 5.9.2 | Brightness | 122 |
| 5.9.3 | Volume | 122 |
| 5.9.4 | Status lights | 122 |
| 5.10 | User Management | 123 |
| 5.10.1 | Create, Edit Delete a User | 123 |
| 5.10.2 | Preset User Role Accessibility | 123 |
| 5.10.3 | Group User Permissions | 124 |
| 5.10.4 | Password Policy | 124 |
| 5.11 | Application Modes | 125 |
| 5.12 | Weighing Units | 125 |
| 5.13 | Communication | 127 |
| 5.13.1 | RS232 | 127 |
| 5.13.2 | RS232 (DB9) Pin Connections | 129 |
| 5.13.3 | USB | 129 |
| 5.13.4 | Ethernet | 129 |
| 5.13.5 | WI-FI & Bluetooth | 130 |
| 6 | Print Setting | 131 |
| 6.1 | Print Content | 131 |
| 6.2 | Connect to Printer | 132 |
| 6.3 | Connect to PC | 132 |
| 6.4 | Data to Excel | 133 |
| 6.5 | Save to USB | 133 |
| 6.5.1 | Application Print Out Template | 134 |
| 6.6 | Printout Examples | 139 |
| 6.6.1 | Calibration Report Template | 139 |
| 6.6.2 | Application Print Out Template | 140 |
| 6.6.3 | Batch Printing Output | 145 |
| 7 | Library | 146 |
| 7.1 | Library Data | 146 |
| 7.2 | Import and Export Library | 147 |
| 8 | Maintenance | 148 |
| 8.1 | Maintenance Menu | 148 |
| 8.2 | Software Upgrade | 149 |
| 8.2.1 | Software Upgrade Process | 149 |
| 8.2.2 | Balance Information | 150 |
| 8.3 | Service Menu | 150 |

| | | |
|-----------|---|------------|
| 8.4 | Service Log File | 150 |
| 8.5 | Factory Reset..... | 150 |
| 8.6 | Log Off | 150 |
| 8.7 | Power Off..... | 150 |
| 9 | Maintenance | 151 |
| 9.1 | Legal for Trade Setting | 151 |
| 9.2 | Settings Change When Approved Mode is ON | 152 |
| 9.3 | Verification..... | 152 |
| 9.4 | Sealing | 153 |
| 9.5 | Output Format..... | 153 |
| 10 | MAINTENANCE | 154 |
| 10.1 | Calibration | 154 |
| 10.2 | Cleaning | 154 |
| 10.3 | Battery Power | 154 |
| 10.4 | Troubleshooting | 154 |
| 10.5 | End of Life Instruction | 154 |
| 10.5.1 | Material Composition of Draftshield Models..... | 155 |
| 10.6 | Service Information..... | 155 |
| 11 | TECHNICAL DATA | 156 |
| 11.1 | Service Information..... | 156 |
| 11.2 | Model Specifications..... | 157 |
| 11.3 | Accessory Specifications | 164 |
| 11.4 | Drawings and Dimensions..... | 165 |
| 11.5 | Accessories | 166 |
| 11.6 | Interface Commands | 168 |
| 12 | COMPLIANCE | 170 |
| 13 | LIMITED WARRANTY | 172 |

1 INTRODUCTION

1.1 Description

The Explorer Plus series balance offers a premium weighing performance, compliance support and data traceability. The advanced features take the EXP to the next level by optimizing the user experience. This Flagship balance is crafted to meet the most demanding requirements of modern scientific research.

The Explorer Balance is a highly accurate weighing balance, with proper care, it can offer over a decade of reliable service. It comes in various capacities, ranging from 120 grams to 65 kilograms.

1.2 Features

1.2.1 Display

The 7-inch display features vibrant colors and a glass panel that is easy to clean and protect chemical spills.

1.2.2 Modular Design

The balances are composed of two interconnected modules: a Terminal and a Base. Detachable Terminal and Base design ready to use for small lab cabinet or fume hood.

1.2.3 Digital Loadcell with Advanced Environmental Sensors

The balances feature the digital load cell platform with three ambient sensors, which are loadcell temperature, humidity, and air pressure.

1.2.4 Removable Draftshield Design

Removable Draftshield for hassle-free cleaning and long-term care

1.2.5 Motorized Leveling System

The balances feature a Motorized leveling system equipped with a digital-level bubble.

1.2.6 Automatic Draftshield Doors

The balances feature an anti-pinch, half open, and full open auto door function for safe and easy sample access.

1.2.7 Status Lights and Sample Spotlights

Status lights for weighing visually indicate overload or underload behavior during the weighing process.

Sample spotlight for enhanced visibility and make a Clearview environment

1.2.8 HID Data Transfer Function

Support for HID (Human Interface Device) connection to a computer without the need for drivers

1.2.9 Time Synchronization NTP (Network Time Protocol)

The balance features Synchronization with Network (NTP function) which can use consistent time and data from the local network.

1.2.10 Fingerprint login System (Accessory)

The fingerprint login system offers a convenient and secure way to log in, eliminating the risk of forgetting passwords or unauthorized password leaks. These measures safeguard the weighing data and the defined weighing process configurations, preventing them from being altered or deleted.

2 Safety Precautions

2.1 Intended Use

This instrument is intended for use in laboratories, pharmacies, schools, businesses and light industry. It must only be used for measuring the parameters described in these operating instructions. Any other type of use and operation beyond the limits of technical specifications, without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use. If the instrument is not used according to these operating instructions, the intended protection provided by the instrument may be impaired.

2.2 Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

| | |
|-----------|--|
| WARNING | For a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided. |
| CAUTION | For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries if not avoided. |
| ATTENTION | For important information about the product. May lead to equipment damage if not avoided. |
| NOTE | For useful information about the product. |

Warning Symbols



General hazard



Explosion hazard



Electrical shock hazard

2.3 Safety Notes



CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Before connecting power, verify that the AC adapter's input voltage range and plug type are compatible with the local AC mains power supply.
- Only connect the power cord to a compatible grounded electrical outlet. (For Explorer Plus High-Capacity models only)
- Do not position the equipment such that it is difficult to reach the power connection.
- Make sure that the power cord does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in these instructions.
- This equipment is for indoor use only.
- Do not operate the equipment in wet, hazardous or unstable environments.
- Do not allow liquids to enter the equipment.
- Do not load the equipment above its rated capacity.
- Do not drop loads on the platform.
- Do not place the equipment upside down on the platform.
- Use only approved accessories and peripherals.

- Disconnect the equipment from the power supply when cleaning.
- Service should only be performed by authorized personnel.
- When shipping or transporting this product, follow the applicable regulations for equipment containing lithium-ion batteries.
-



WARNING: Never work in an environment subject to explosion hazards! The housing of the instrument is not gas tight. (Explosion hazard due to spark formation, corrosion caused by the ingress of gases).



WARNING: Electrical shock hazards exist within the housing! The housing should only be opened by authorized and qualified personnel. Remove all power connections to the unit before opening.



CAUTION: The protection provided by the unit may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.



WARNING: ELECTRICAL SHOCK HAZARD

Avoid pressing two carbon brushes while the lonizer is turned on.

3 Installation

3.1 Unpacking

Carefully unpack the Explorer Plus balance and its components. The components included may vary based on the balance model. Retain the packaging for safe storage and transportation.

- Balance
- Quick Start Guide
- Weighing Pan
- Power Adapter and Local Electronic Plug (except High-Capacity Models)

3.2 Installing Components

The Explorer Plus balance is designed for quick, no-assembly installation, allowing for immediate use. Simply follow the steps to assemble your Explorer Plus balance within seconds. Assembly of all components is required before using the balance.

3.2.1 For 0.01mg, 0.1mg, and 1mg Models

Unpack the balance and install the weighing pan.





3.2.2 For 0.01mg, 0.01g, and 0.1g Models

Unpack the balance and install the weighing pan.




3.2.3 For High-Capacity Models

Unpack the balance and install the weighing pan.

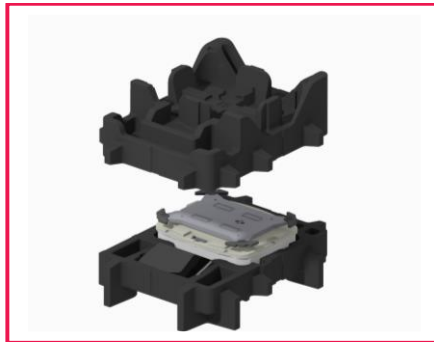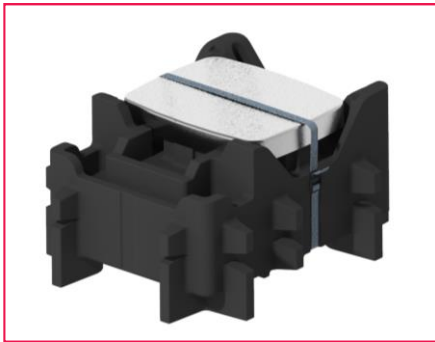

3.2.4 Selecting a Location

External factors such as temperature fluctuations, air currents, electromagnetic interference, and vibrations can all have an impact on the accuracy of the analytical balance. It is important to keep the analytical balance in a controlled environment to minimize these factors and ensure accurate readings.


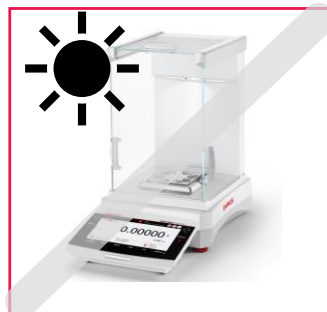


3.2.5 Level the Balance Before Use


The Explorer Plus analytical and precision models are equipped with a motorized leveling system and a digital level bubble. To level your Explorer Plus balance, simply follow the provided steps. The digital level indicator will notify you if the weighing table is not level. The Explorer Plus high-capacity model also features a built-in digital level bubble, but it has slightly different leveling steps.

The balance uses a built-in Motorized Leveling program to level the balance automatically.

- Go to main screen to find the Motorized Levelling button at the left corner, press it to start leveling the balance.
- The process is fully automated and do not need any user operation during the levelling.
- If the environment is unstable or the weighing table is tilting too much, the leveling process may fail. In such cases, you need to move the balance to a stable, level surface that can absorb vibrations.

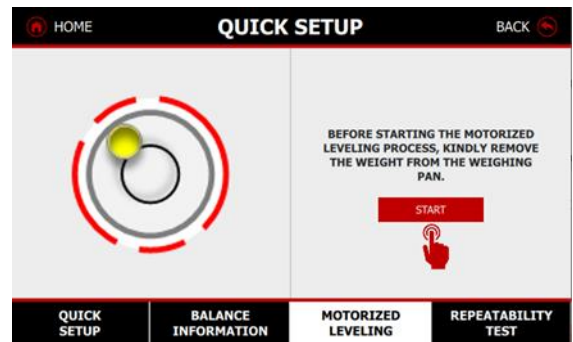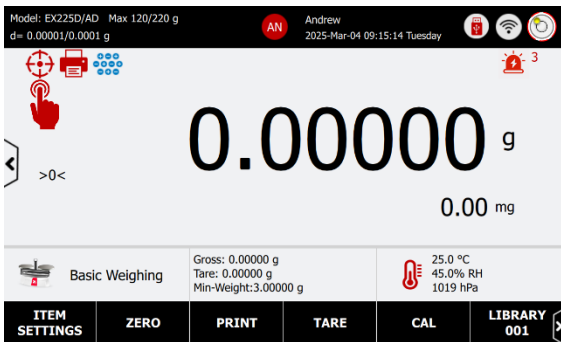

- For Explorer high-capacity Model, press the leveling button, and adjust the leveling feet according to the position of the digital bubble until the bubble is centered.

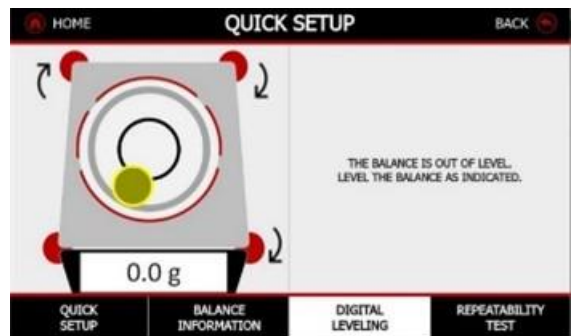

3.3 Connecting Power

- For Analytical and Precision models supplied with an AC adapter, connect the DC output connector to the power receptacle on the rear of the base. Then connect the AC power cord to a suitable electrical outlet.
- For High-Capacity models, you need to connect the power cord to a suitable electrical grounded outlet.
- When the balance run in sleep or standby or mode, press the **LABMAN** icon to start weighing.



Caution:

- For use with CSA certified (or equivalent approved) power source, which must have a limited current output.
- Allow equipment to warm up for 60 minutes for optimal weighing performance.

3.4 Turn on the balance

Long press the power button to turn on the balance.



3.5 Adjust the viewing angles

To adjust the display's angle, hold the base of the terminal and move it upward or downward. The internal connection cable has a maximum length of 1 meter.




3.6 Disconnect the terminal from the weighing base

- To remove the terminal from the base, simply lift the terminal straight up.
- To reattach the terminal, align the connection rods with the slots on the base frame and insert them.
- If you need a longer operating distance, the terminal extension cable and tower kit are available as optional accessories. Refer to the accessory list in Table xxxx.



3.7 Remove Draftshield from the Base

- Remove the weighing pan and unlock the draft shield fixed ring.

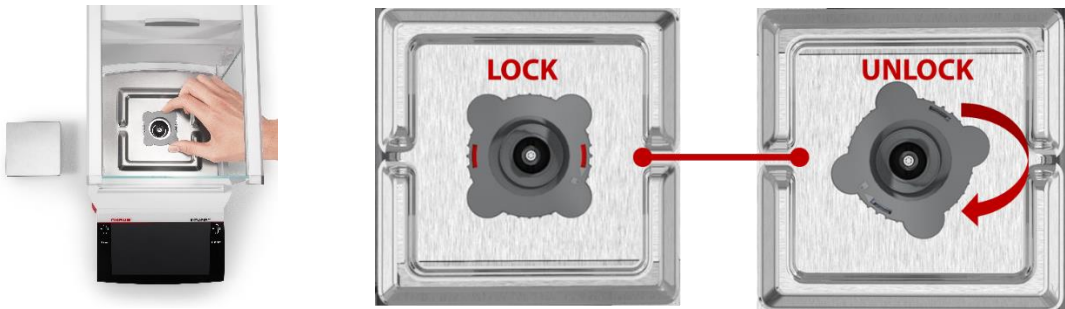

- Remove the EMC plate and take out the 4 screws beneath it. Raise the draft shield vertically.



3.8 Connecting the Interface

Enhanced communication options include 2 USB Host (Type A) x 2, USB Device (Type B) x 1, USB Device (Type C) x 1, Ethernet Lan port (RJ45) x 1, RS232 x 1; Optional Wi-Fi, Bluetooth Dongle.

USBA and C ports, Ethernet Lan port,
BT and WiFi connectivity




3.9 Overview of Parts and Features –Draft Shield Models

3.9.1 Side View



3.9.2 Rear View

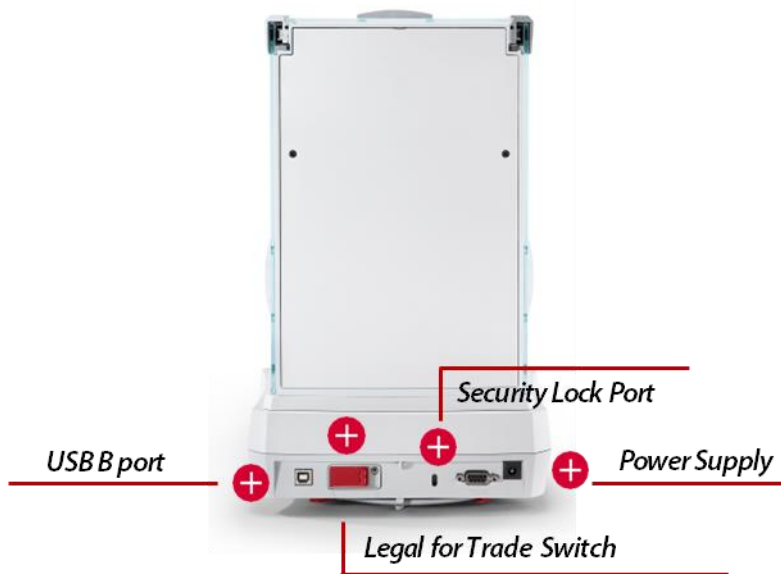

3.10 Overview of Parts and Features –Non-Draft Shield Models

3.10.1 Side View



3.10.2 Rear View



3.11 Overview of High-Capacity Models



3.11.1 Side View



3.11.2 Rear View




3.12 Overview of Main Screen and Connectivity

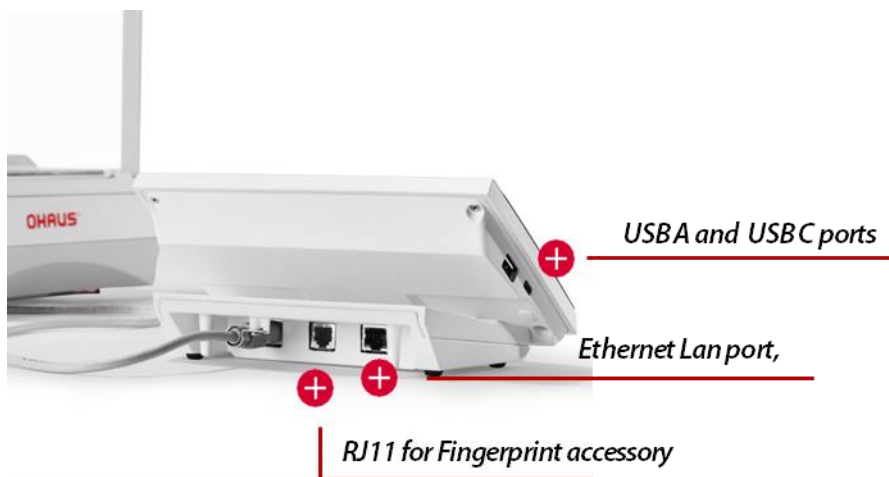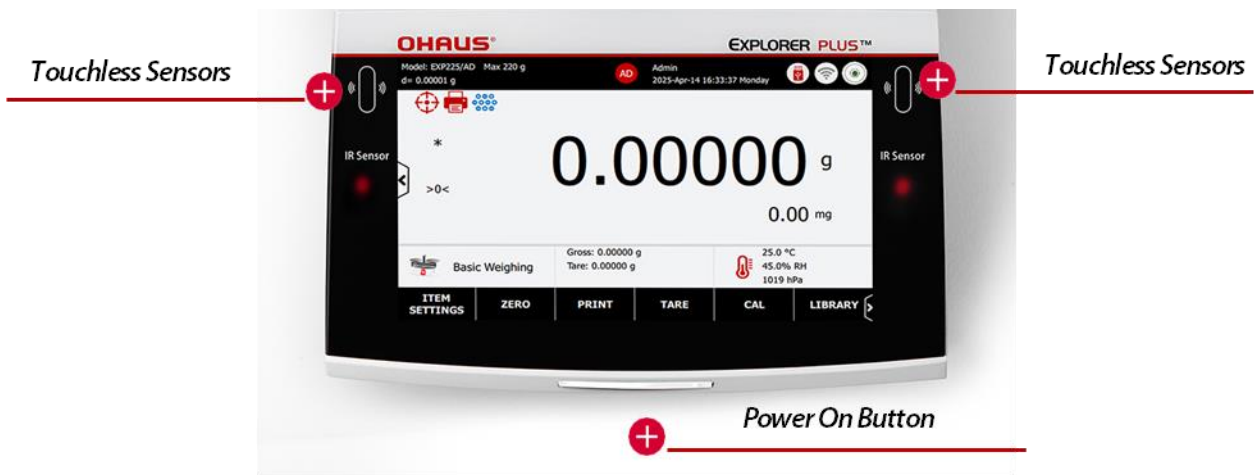

3.13 Main Screen Features

This Explorer Plus balance has built a 7-inch display features vibrant colors and a glass panel that is easy to clean and protect chemical spills.
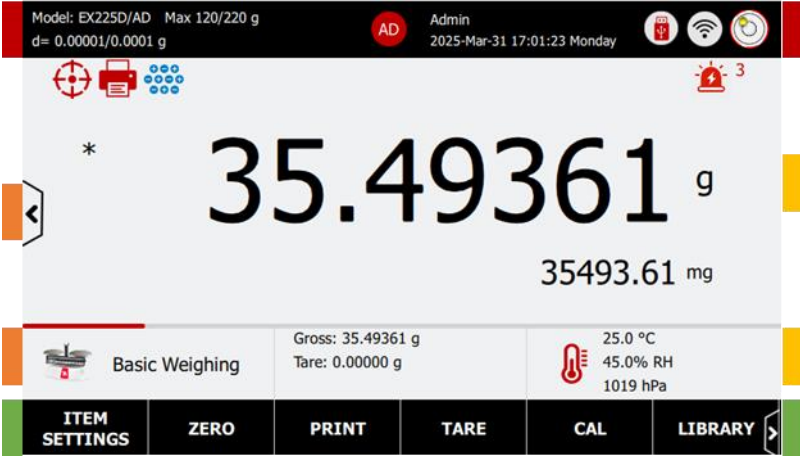
| | | |
|---|---|--|
| <p>Upper Left (Shortcut buttons)</p> <ul style="list-style-type: none"> ✓ Motorized Leveling ✓ Batch Print ✓ Ionizer | <p>Top Line</p> <ul style="list-style-type: none"> ✓ Display the essential balance information, such as model name, the max capacity, d value and e value. ✓ User Name and Date/Time | <p>Upper Right</p> <ul style="list-style-type: none"> ✓ Peripherals Icons connected to balance ✓ Real time leveling bubble ✓ Warning Sign |
| <p>Press the trapezoidal key to switch between main menus.</p> <ul style="list-style-type: none"> ✓ Quick Setup ✓ Calibration ✓ Balance Setup ✓ User Management ✓ Application Modes ✓ Weighing Units ✓ Communication ✓ Library ✓ Maintenance ✓ Factory Reset ✓ Log Off ✓ Power Off |  | <p>Main Display Area</p> <ul style="list-style-type: none"> ✓ Click “g” to switch an alternative unit ✓ Display the parameter for the current application on the reference field ✓ Digital loadcell Environmental Sensor |
| <p>Press the “Basic Weighing” key to switch weighing application modes</p> | <p>Hotkeys</p> <ul style="list-style-type: none"> ✓ Item Settings, Zero, Print, Tare, Calibration, Library/Method, Sensors, and press “ADD+” key to add customized hotkeys | <p>ADD+</p> <ul style="list-style-type: none"> ✓ 1d/10, Open Left Door, Open Right Door Open, Motorized Leveling, Repeatability Test |


3.14 Warning and Notifications

The Explorer Plus balance features an internal warning and notification system that alerts users to take action in the following circumstances:

- Calibration failure: The balance has not calibrated successfully.
- Leveling required: The balance needs to be leveled when the digital level bubble is not centered.
- Password expiration: The password has expired in accordance with the password policy.

3.14.1 Calibration Failure


- Press on the warning sign, and the notification will show up on the screen, press the message, you will find “Initial calibration” key.

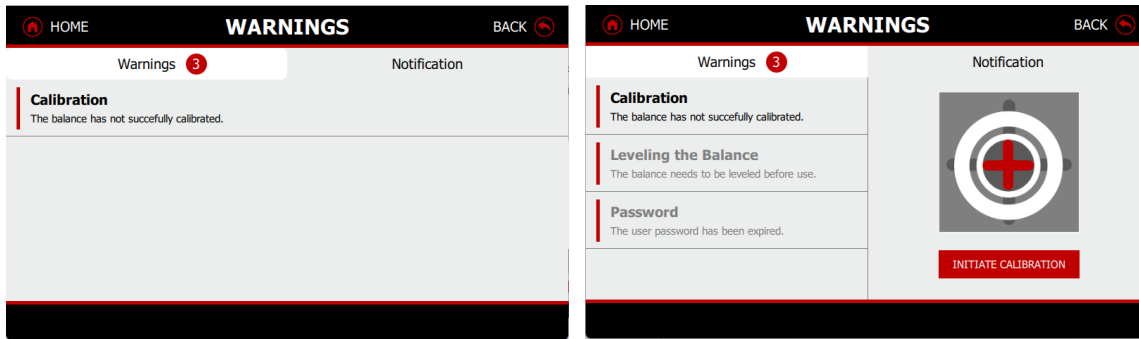

- Perform internal calibration or span calibration. For details, please refer to [5.4 Calibration](#).


3.14.2 Leveling Required

- Press on the warning sign, and the notification will show up on the screen, press the message, you will find “Leveling” key.

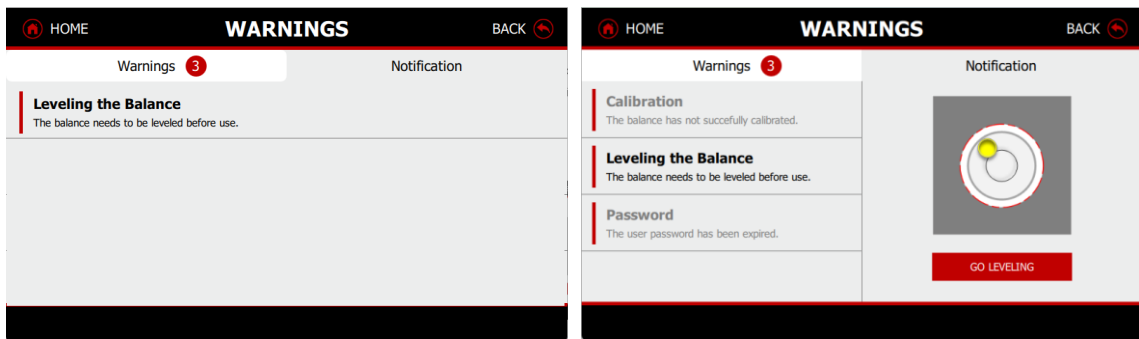

3.14.3 Password Expiration

Press on the warning sign, and the notification will show up on the screen, press the message, you will find “Reset Password” key.

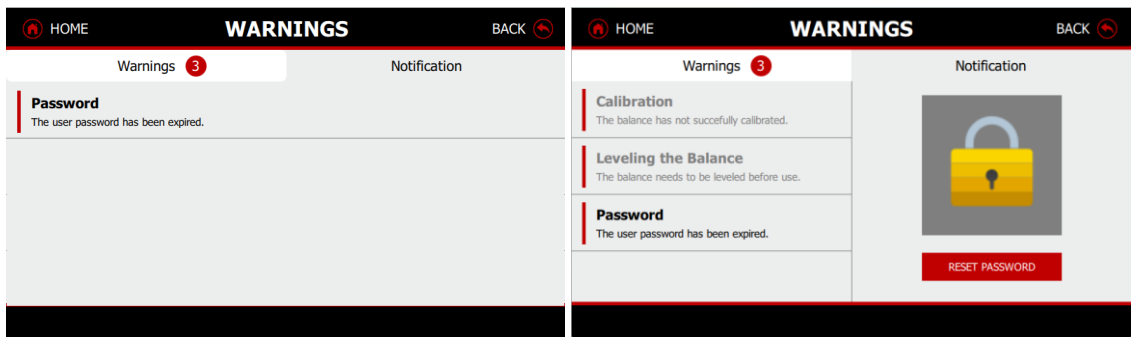



Note: If you forget your password, please contact Ohaus or your local Ohaus dealer for assistance.

3.15 Hotkeys

At the bottom of the main screen, there are eight hotkeys for basic weighing applications. Users can customize an additional five hotkeys by using the ADD+ feature.

3.15.1 Standard Hotkeys
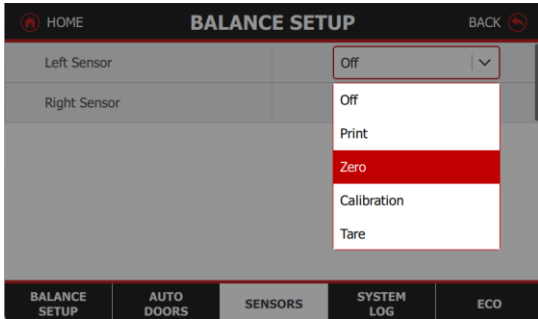
| Hotkeys | Description |
|----------------------|---|
| Item Settings | Press to configure current application mode settings. |
| Zero | <p>Remove the load from the pan and press Zero to set the display to zero. When weighing pan is empty, the >0< indicator turns on when the measurement is within $\pm \frac{1}{4}$ division (d) of the zero setting.</p> <p>Note: The balance also features Auto Zero Tracking (AZT) that automatically maintains a Center of Zero within the tolerances set in the AZT menu (see Balance Setup).</p> |
| Print | <p>Press Print to send the displayed value to a printer or computer via the active COM port.</p> <p>Make sure that the balance has connected with printer or PC and interface parameters are set up correctly.</p> |
| Tare | Tare key is to tare the weight on the container to zero, and the net weight will show on the reference field |
| CAL | Press CAL to calibrate the balance or to set the calibration parameters. |
| Library | Press "Library" key to either create a new library or retrieve an existing one for the current application. |
| Sensors | <p>The balances have two pressless sensors that can be assigned a unique function when activated. The Sensor setting are: Off, Print, Zero, Calibration, Tare, Auto Doors, Open Left Door, Open Right Door, Ionizer, Ionizer + Tare, and Draft Shield Light.</p>  <p>Once the sensor is set up for a specific function, waving your hand over it will trigger that function and cause the sensor light to turn green. If the sensor fails to activate, verify the settings. The sensor lights will remain red.</p> |
| Add+ | <p>Additional five hotkey to be customized.</p> <p>Click "Right Arrow" key to move the available hotkey to selected area and press "Save" to exit.</p> <p>The new hotkey will appear on the right side of the "Sensor" section. Swipe along the bottom line, you will find the new hotkey.</p> |

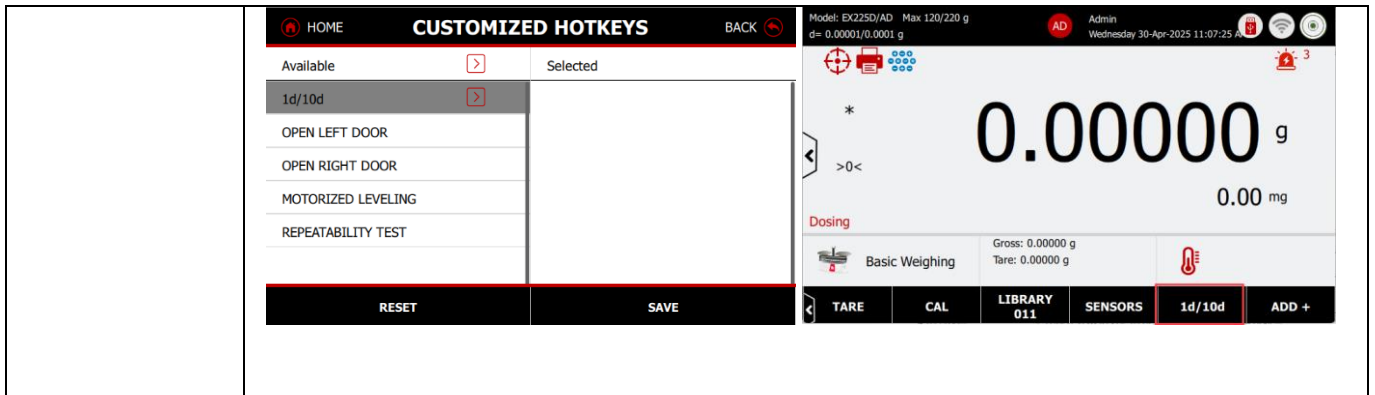

3.15.2 Additional Hotkeys (Add+)

| | |
|--------------------|---|
| 1d /10d | Press 1/10 to change 1d or 10d diviation for weighing result |
| Motorized Leveling | Press to Motorized Leveling the balance |
| Repeatability Test | Press Repeatability Test to perform repeatability test |
| Left Door Open | Open the draftshield left door |
| Right Door Open | Open the draftshield right door |

3.16 Shortcut Buttons


The shortcut buttons are available only in Basic Weighing application. For details, please refer to shortcut button in xxx.

3.17 HID Connections

The Explorer Plus balance supports HID (Human Interface Device) connection to a computer without needing drivers. This allows users to use a mouse or keyboard to enter information such as User Name, USER ID, sample name, batch name, and other input details. The input field supports both English and French characters, as well as numbers, and symbols.


3.17.1 English Keyboard

- English keyboard is default setting, click “Upward arrow” key, you will find the lower-case letter and additional symbol such as @, %, -, ^ and etc.
- Repeat three times, and you can find all the characters you need to input.
- The currently selected language is displayed on the SPACE key.

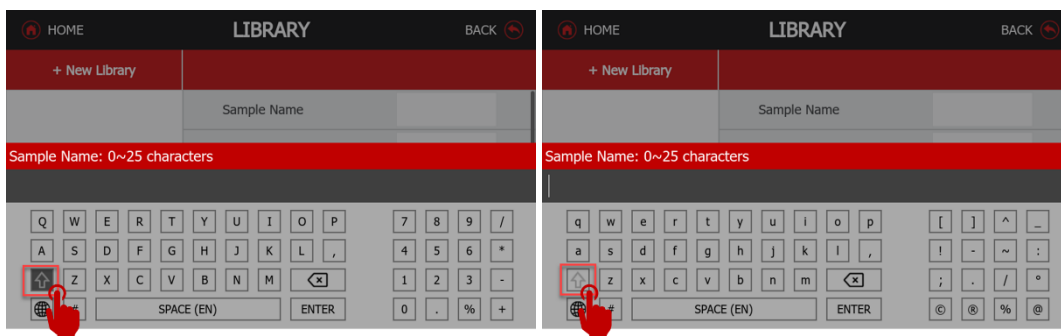

Uppercase Letter/ Numeric Input


Lowercase Letter/ Symbol Input

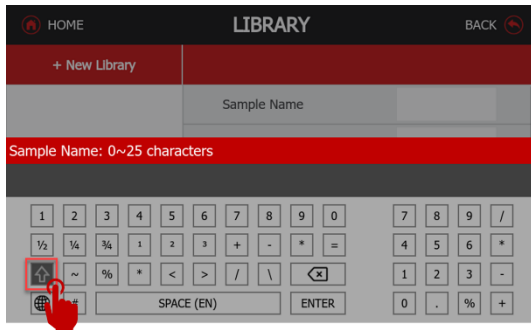


Numeric/ Symbol Input

3.17.2 French Keyboard


- Click “Earth” key, the keyboard changes to French language.
- Press the Upward arrow to access French characters, as well as symbols like %, 1/2, 1/4, 3/4, √ 2, ∅, ml, and more.
- Repeat three times, and you can find all the characters you need to input.
- The currently selected language is displayed on the SPACE key.

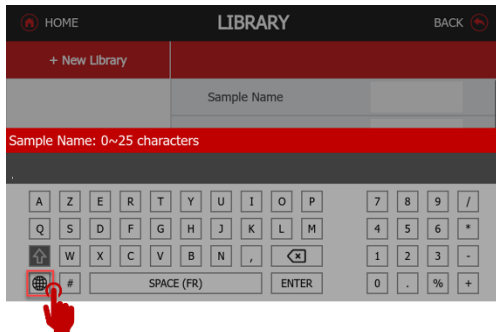

Uppercase Letter/ Numeric Input

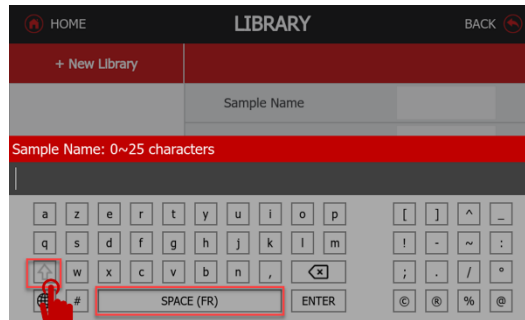

Lowercase Letter/ Symbol Input

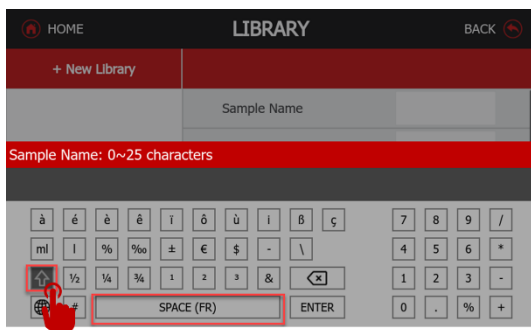


French character/ Number Input

4 Weighing Application Modes

4.1 Introduction

The Explorer Plus balance has 15 weighing application modes. All applications are by default enabled. You can disable the weighing application modes in Menu Menu/Application Modes. Before using any weighing application, ensure that the balance has been leveled and calibrated successfully.

The application mode switch button is located in the lower left corner. Press the “Basic Weighing” key to switch weighing application modes.

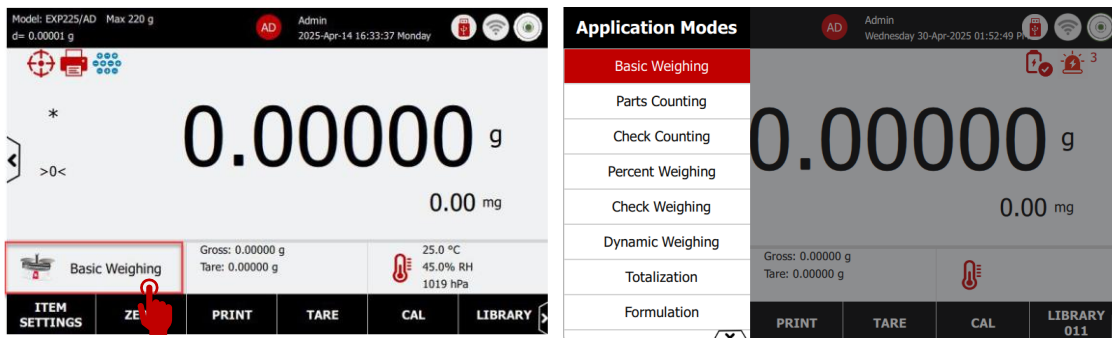


- **The Basic Weighing** application is used to determine the weight of items in the selected unit of measure.
- **Parts Counting** displays the number of pieces or parts based on an average piece weight. Checks if the current sample pieces are within tolerances (e.g. an over and under limit).
- **Check Counting** checks if the current sample pieces are within tolerances (e.g. an over and under limit).
- **Percent Weighing** displays the current weight as a percentage of a reference weight.
- **Check Weighing** checks if the current weight is within tolerances (e.g. an over and under limit).
- **Dynamic Weighing** is set to weigh objects that are not stable, such as animals.
- **Totalization** is used to sum the weights of multiple samples and report the statistical data for the series of samples.
- **Formulation** is used to combine various elements in proportionate amounts.
- **Differential** is used to calculate the difference in weights of multiple samples taken at different times.
- **Density Determination** is used to determine the density of a solid or a liquid.
- **Peak Hold** captures the maximum weight in a series of weighing.
- **Pipette Adjustment** calculates the inaccuracy and imprecision of pipettes, used to check if a pipette's dosage is within tolerances.
- **SQC** is used to determine the homogeneity of items in a batch and batches over time.
- **Fill Weight Variation** Checks the homogeneity of a series of items.
- **Flow Rate Control** aims at managing the volume or mass of fluid (liquid or gas) that passes through a pump system over a specific period.

4.2 Basic Weighing

- Basic weighing application modes are displayed by default on the main screen.
- Press **Tare** or **Zero** to start weighing.
- Place sample on the pan to display the weight. When stable, the * appears.
- The resulting value is displayed in the main Weighing Line in the active unit of measure.

4.2.1 Weighing Result

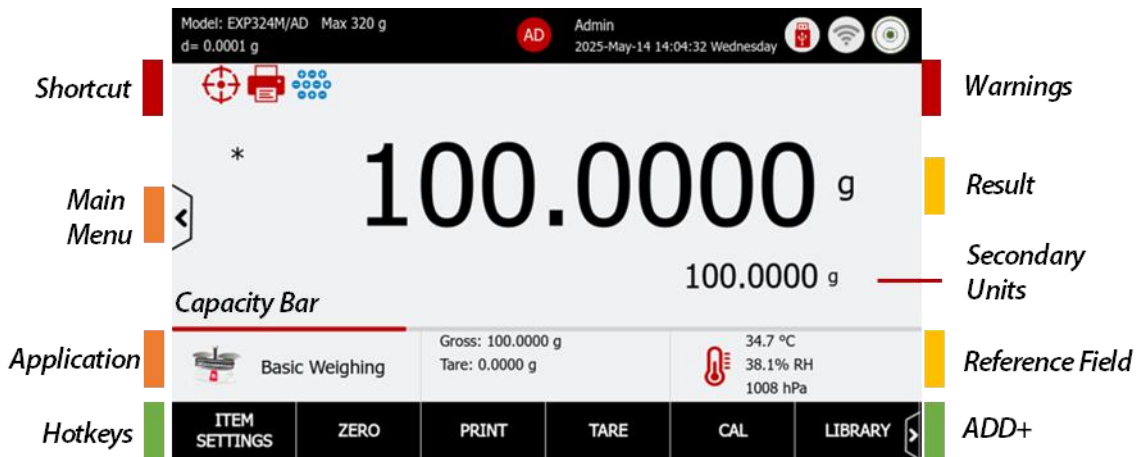


4.2.2 Reference Fields – Basic Weighing

| Reference Fields | Description |
|--------------------------|--|
| Gross/ Tare | Display the Gross weight and tare value when the container is tared |
| Min Weight (Conditional) | When the Minimum Weight is set to On or Below Minimum Weight. Display the minimum weight value. |
| Environment Sensor | Load cell temperature, humidity and air pressure. Default setting is on. Click off to hide this field. |
| Capacity Bar | The red line represents the current load as a proportion of the balance maximum capacity. |
| Weighing Mode | Standard and Sample Dosing modes are two weighing modes in the balance filter settings. <ul style="list-style-type: none"> • Sample Dosing Mode is specifically designed for applications involving powder samples." |
| Item Settings | <ul style="list-style-type: none"> • Primary Weighing Unit: The default unit is grams. The operator can switch to alternative weighing units and two custom units. • Secondary Weighing Unit: able to alternative weighing units and 2 custom units • Weighing Mode: Standard, Sample Dosing Mode <ul style="list-style-type: none"> ■ Sample Dosing Mode is design for power sample, or any sample are sensitive to be stable • Minimum Weighing • Auto Tare: Automatic tare the container value |

4.2.3 Item Setting

Press the ITEM SETTINGS button to change the application settings.

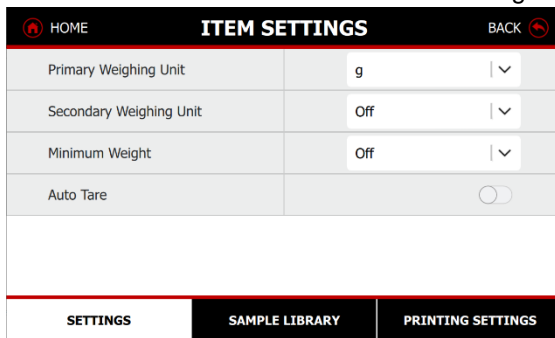


Primary Weighing Unit/ Secondary Weighing Unit

- Click “g” to switch an alternative primary weighing unit. The default unit is gram.
- Select a secondary unit to display below the primary line.

Switch to alternative Unit

Click “g” to switch an alternative unit

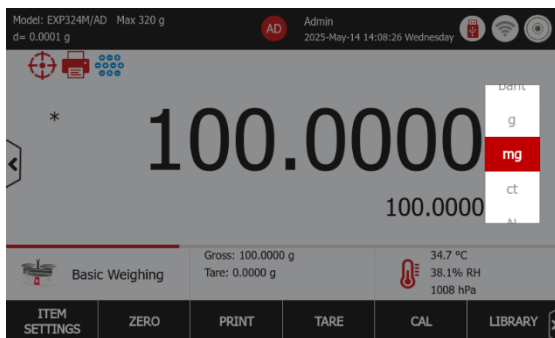

4.2.4 Minimum Weight


Determining the minimum weight in laboratories is crucial for ensuring accurate and reliable weighing results. The Explorer balance has the min-weight function that the user has already get the result of the manual weight from repeatability test. In the basic weighing, user can input the known min-weight in this balance. When the result is lower than min-weight, the screen will show “Below Minimum Weight”, and the operator has to add more samples.

The Explorer Plus balance offers three options of minimum weight settings: Off, On, and Below Minimum Weight.

- **Off:** Disable function
- **On:** When the result is lower than min-weight, the screen will show “Below Minimum Weight”.
 - The operator had to add more samples. The result will not be printed in paper printouts.
- **Below Minimum Weight:** When the result is lower than min-weight, the screen will show “Below Minimum Weight”. The operator had to add more samples. The result will be printed in paper printouts.


4.2.5 Set minimum weight value

- Press to the Sample Library tab to input the min-weight value when you turn on the min-weight in settings.
- Input the min-weight value, and the range of minimum weight is from 0g to max capacity. If the number is out of the range, it will display an error message “invalid entry”.

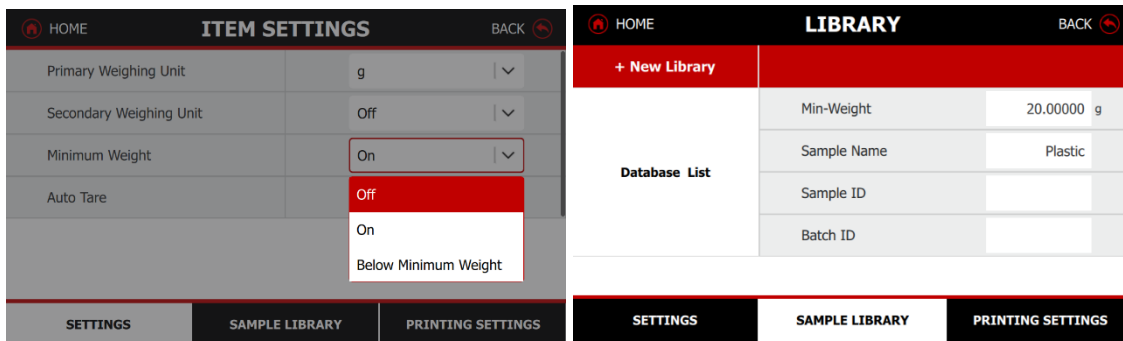

4.2.6 Auto Tare

- Auto tare allows users to tare the container weight
- When Auto Tare is enabled, the screen would display the message “Place container on the pan.”
- After a container is placed on the weighing pan, its weight will be stored in the balance. The net weight and gross weight values will then be displayed in the reference field.

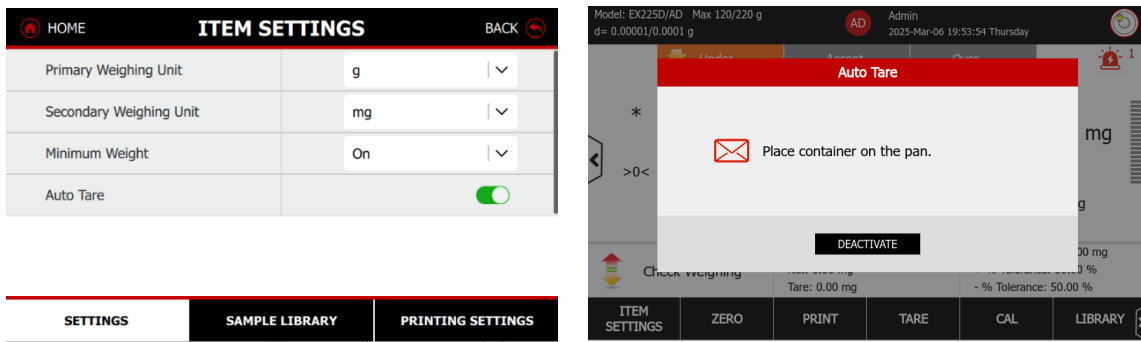


4.2.7 Sample Library

The Explorer Plus balance features a built-in library for managing multiple sample profiles. Up to 3,000 library records can be saved and recalled for use. The total library capacity and records are accessible across various application modes and through the library menu. When the library usage exceeds 80%, a pop-up message will alert the user to export the library or delete all items. The entire library menu, please refer to the section of Library.

Create, Active and Delete a Library

- Press “+New Library” button to create a new Library.
- When a new library is created, it remains inactive. The operator has to activate the library after ticking the box.
- Press on the RUBISH BIN button to delete a library.
- After the library item is activated, the main screen will show the activated library number.
- Other sample information can be configured in the library, including sample name, sample ID, batch ID, lot ID, project ID, and the creation of 10 more customized ID.

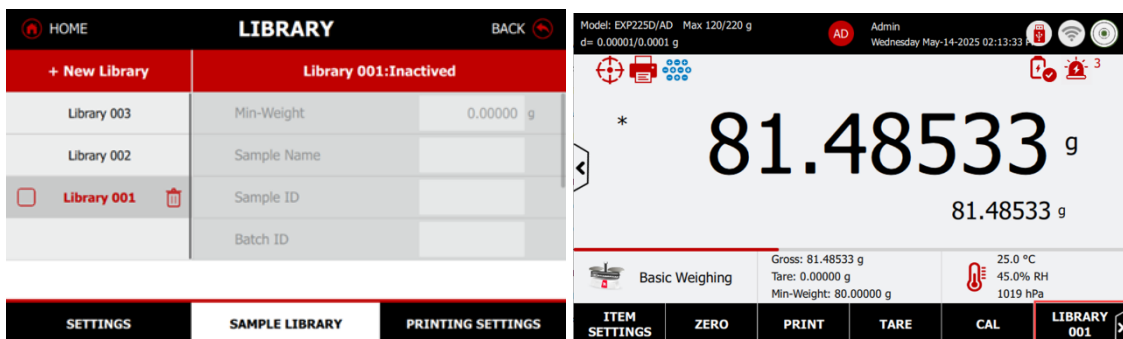


4.2.8 Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to 6.0 Printer section

4.2.9 Shortcut Buttons

The Explorer Plus balance offers three shortcut buttons for easy leveling, print batch sample result, and turn on lonizer function.

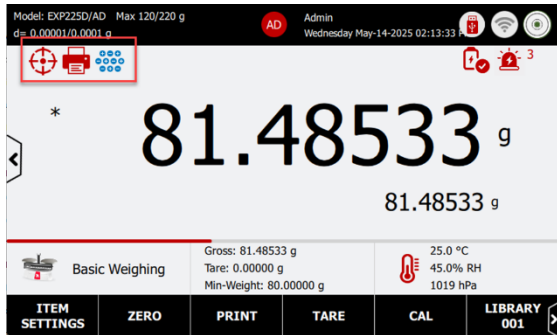


Motorized Leveling

Press the Motorized Leveling icon to initiate the automatic leveling of the balance. The initial digital level bubble will display the real-time status at the upper right corner. See detail in section of Motorized leveling in Balance setup xxx

Batch Printing

Batch printing allows users to record and print a list of sample weight. It is available only in the Basic Weight application.

- Press the button to start measure batch samples. The batch sample information will show on the left side of screen. The right side of screen shows the current sample result.
- Place a sample onto the weighing pan, press accept when the weight is stable.
- Press “End Batch” button to complete the sample batch weighing, then the operator can print all the batch weighing data in once.
- Before printing, user can go back to edit when the result has errors.
- User can export the data to PDF or print to PC/ Printer. After printing, operator can review the history result
- Note: If users would like to continue with batch printing, tap "BACK TO EDIT" to exit the result screen. The result will neither be printed nor saved to the history.

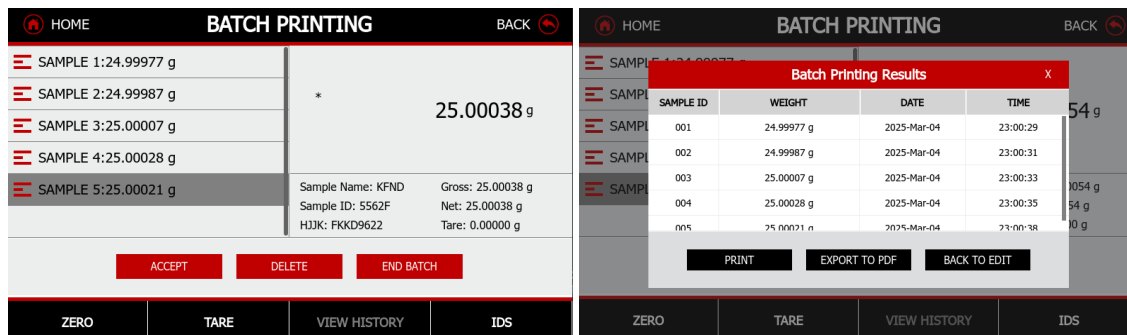


Ionizer

Press the Ionizer icon to use Electrode brush to neutralize the weighing chamber. See details in the section of ionizer in Balance setup XXX.

4.3 Parts Counting

- Use this application to count samples of uniform weight.
- In the lower portion of the home screen, select Parts Counting.
- Press **Tare** or **Zero** to start weighing.
- Place objects on the pan to display the number of pieces. The default (or last) Average Piece Weight (APW) is displayed.
- Setup the SAMPLE SIZE, Average Piece Weight (APW) before starting Parts Counting application.

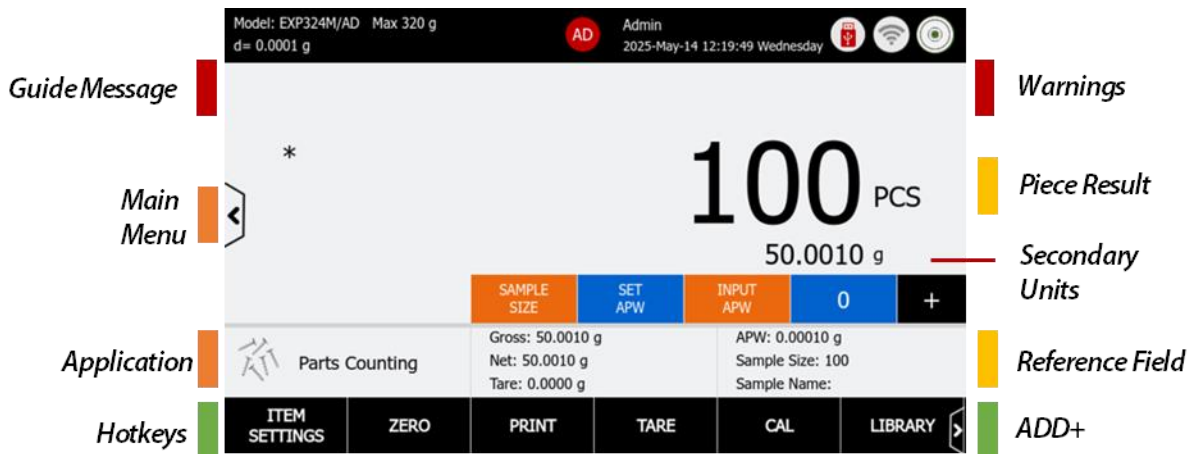

4.3.1 Application buttons

| Application Button | Description |
|--------------------|--|
| Sample Size | Input sample size |
| Set APW | Place the sample on the pan, the balance will calculate Average Piece Weight (APW) |
| Input APW | Input Average Piece Weight (APW) |
| 0 | The statistic key is to recall the number of parts counting, and press this button, operator can review the statics results. |
| + | Press this button to manually accumulate weight. In Automatic Statistics mode, this button will be inactive. |
| Item Settings | <ul style="list-style-type: none"> • Auto Optimization: able to automatically optimized APW Auto Optimization improves counting accuracy. • Statistics: Automatic, Manual • Secondary Weighing Units: able to use 15 weighing units and 2 customer units • Auto Tare: Automatic tare the container value |


4.3.2 Begin Parts Counting

Step 1: Establish an Average Piece Weight (APW)

Each time a new type of part is counted, the nominal weight of one piece (Average Piece Weight or APW) must be established using a small quantity of pieces.


Set a Known Average Piece Weight (APW)

- Press the INPUT APW button and enter the desired APW value.
- Tap ENTER to save the value

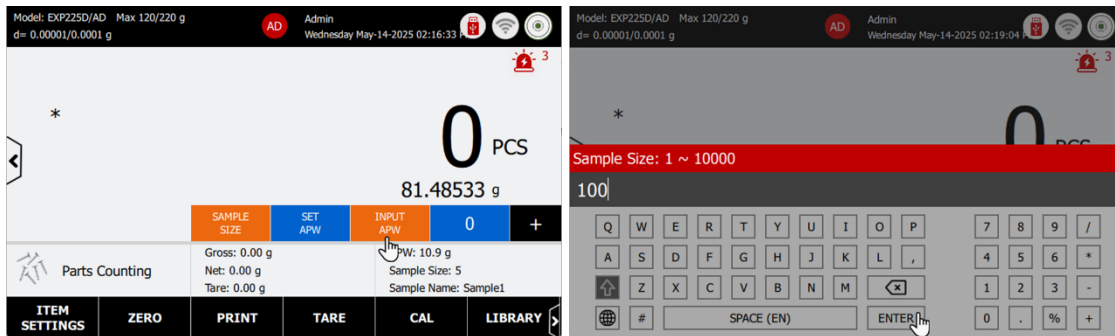

Set a new Average Piece Weight (APW) – Derived

- Press the Sample Size button and enter the desired Sample Size.
- Tap ENTER to save the value
- Place reference weight on the pan, and press Accept to establish a new APW.
- Example: The home screen displays 2.50006 g at the new APW.

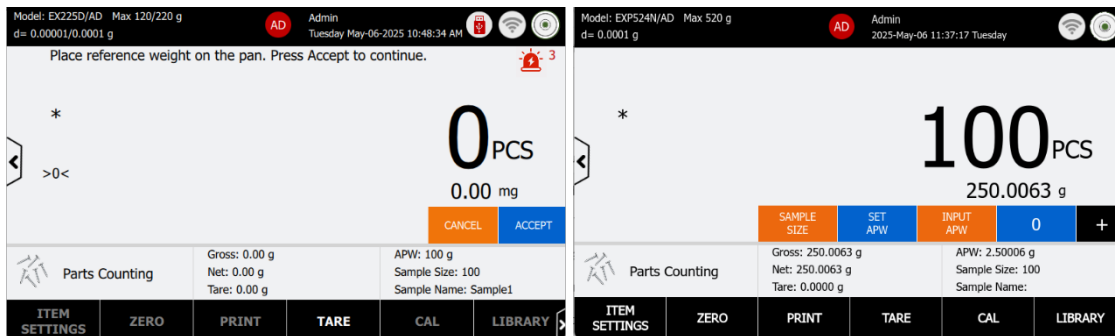

- The sample size can be 1 to 10 000 pieces. The default sample size is 10.
- Once a sample size is changed, the balance will immediately recalculate APW screen, expecting to establish a new APW. The home screen shows 100 pieces at the new APW

Step 2: Counting the Samples

Place samples on the pan, the balance will display the number of pieces.

4.3.3 Accumulate the data for Parts Counting


Utilize the Statistics "+" function to aggregate samples based on counting numbers and generate statistical reports.

The balance can store up to 999 accumulation records.

Step 1: Set the Statistic Mode in Parts Counting Mode


- Statistics can be determined Manual mode or automatically
- Automatic Mode: Weights are automatically recorded when stable
- To use the statistics function, enable it first in the Item Settings.
- Add items and press the "+ Statistic button to accumulate the statistical data

- Tap the statistics number to view the statistical results.

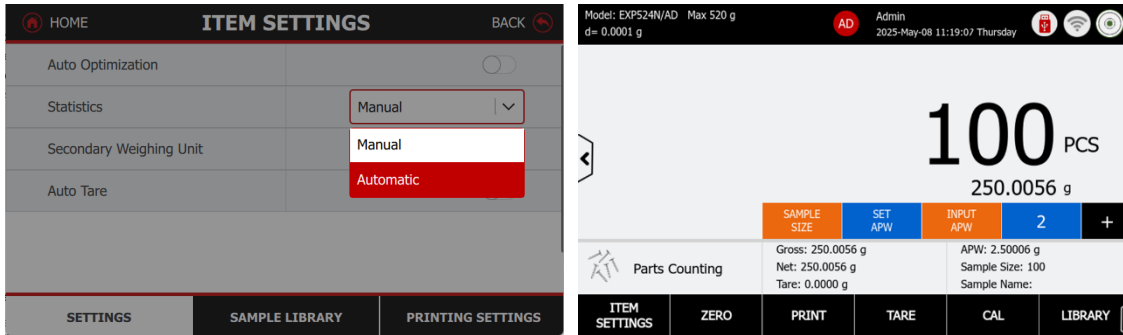

Step 2: Statistics Report View

- The report content includes Start Time, Number of Samples, Total, Average, Maximum, Minimum, Range, Standard Deviation, Relative Deviation % and individual item weighing values.

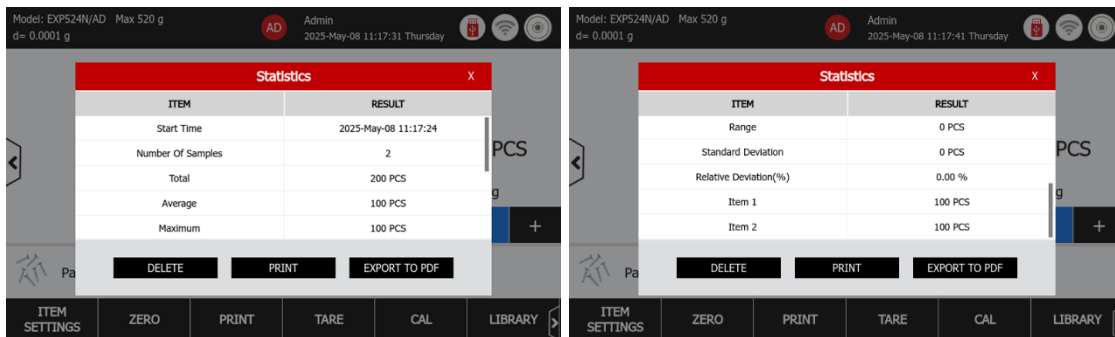

Step 3: Print Report

After reviewing the report, users can choose from several actions:


- Delete the statistics and restart the percent weighing process
- Print the report to a printer or PC based on the Print Settings
- Export the report as a PDF file to a USB flash drive

4.3.4 Sample Library – Parts Counting


The Explorer Plus balance features a built-in library for managing multiple sample profiles. Up to 3,000 library records can be saved and recalled for use. The total library capacity and records are accessible across various application modes and through the library menu. When the library usage exceeds 80%, a pop-up message will alert the user to export the library or delete all items. The entire library menu, please refer to 6.5.1 Library.

Create, Active and Delete a Library

- Press “+New Library” button to create a new Library.
- When a new library is created, it remains inactive. The operator has to activate the library after ticking the box.
- Press on the RUBISH BIN button to delete a library.

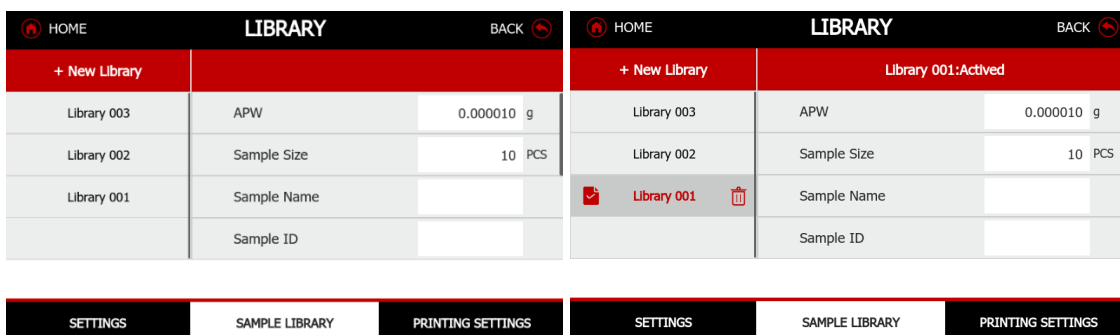

- After the library item is activated, the main screen will show the activated library number.

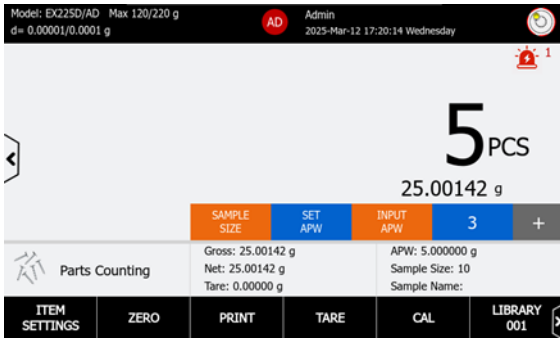

- Other sample information can be configured in the library, including sample name, sample ID, batch ID, lot ID, project ID, and the creation of 10 more customized ID.


4.3.5 Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.


For details of Printing Setting, please refer to 6.0 Printing.

4.4 Check Counting

- Use this application to checks if the current sample pieces are within tolerances (e.g. an over and under limit).
- In the lower portion of the home screen, select Check Counting.
- Press **Tare** or **Zero** to start weighing.
- Place objects on the pan to check if the number of pieces within the tolerances. The default over and under limit pieces is displayed.
- Setup the SAMPLE SIZE, Average Piece Weight (APW) before starting Check Counting application.

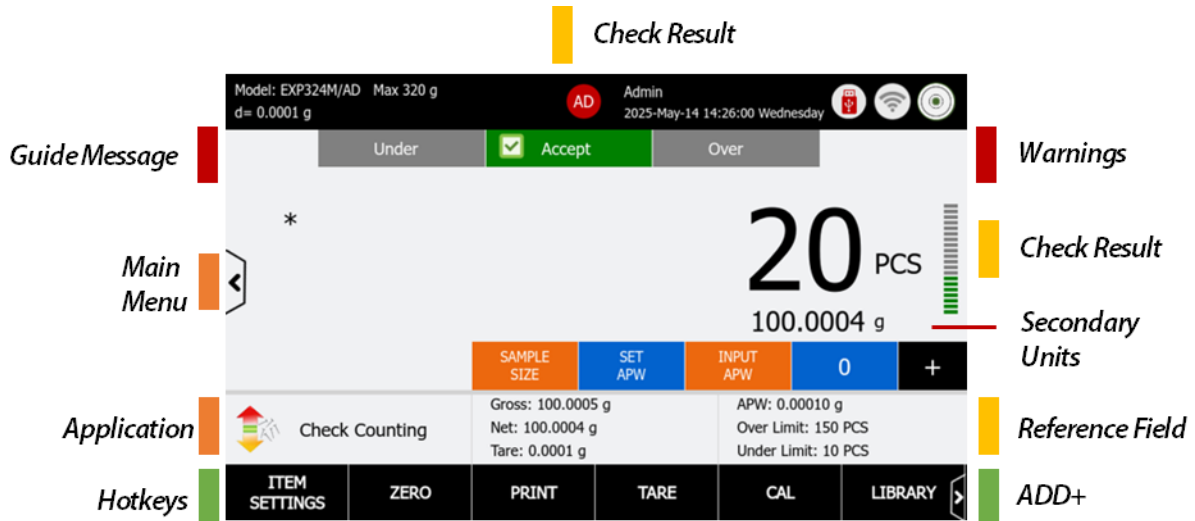

4.4.1 Application buttons and Reference Field

| Application buttons/ Reference Field | Description |
|--|---|
| Sample Size | Input sample size |
| Set APW | Place the sample on the pan, the balance will calculate Average Piece Weight (APW) |
| Input APW | Input Average Piece Weight (APW) |
| 0 | The statistic key is to recall the number of check counting, and press this button, operator can review the statics results. |
| + | Press this button to manually accumulate weight. In Automatic Statistics mode, this button will be inactive. |
| Item Settings | <ul style="list-style-type: none"> • Auto Optimization: able to automatically optimized APW Auto Optimization improves counting accuracy. • Statistics: Automatic, Manual • Secondary Weighing Units: able to use 15 weighing units and 2 customer units • Audible Signal: The balance beeper will sound to alert the user of the check status. <ul style="list-style-type: none"> ■ Available Setting: Off, Under, Accept, Over, Under and Over • Auto Tare: Automatic tare the container value |
| Check Statuses  | <ul style="list-style-type: none"> • Under-Orange indicator • Accept-Green Indicator • Over-Red Indicator • The check statuses are divided into 20 segments to display the proportion of the current load relative to the total capacity. Each segment corresponds to 5% of the total capacity. |
| Statuses Lights | <ul style="list-style-type: none"> • If the check status is either Under or Over, the lights will turn red to indicate this. |
| Over Limit | <ul style="list-style-type: none"> • The piece value exceeds the maximum allowed threshold. |
| Under Limit | <ul style="list-style-type: none"> • The piece value exceeds the maximum allowed threshold. |

4.4.2 Begin Check Counting


Step 1: Establish an Average Piece Weight (APW)

Each time a new type of part is counted, the nominal weight of one piece (Average Piece Weight or APW) must be established using a small quantity of pieces.

Set a Known Average Piece Weight (APW)


- Press the INPUT APW button and enter the desired APW value.

- Tap ENTER to save the value

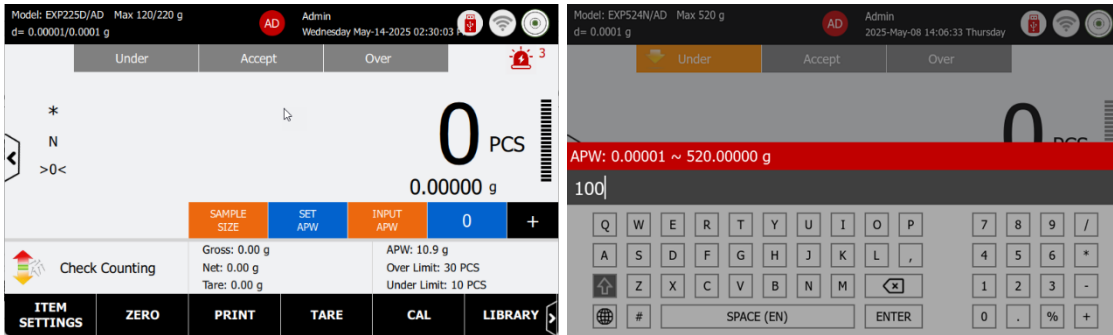


Set a new Average Piece Weight (APW) – Derived

- Press the Sample Size button and enter the desired Sample Size.
- Tap ENTER to save the value
- Place reference weight on the pan, and press Accept to establish a new APW.
- The sample size can be 1 to 10 000 pieces. The default sample size is 10.
- Once a sample size is changed, the balance will immediately recalculate APW screen, expecting to establish a new APW. The home screen shows 100 pieces at the new APW
- Example: The home screen displays 10.00000 g at the new APW.

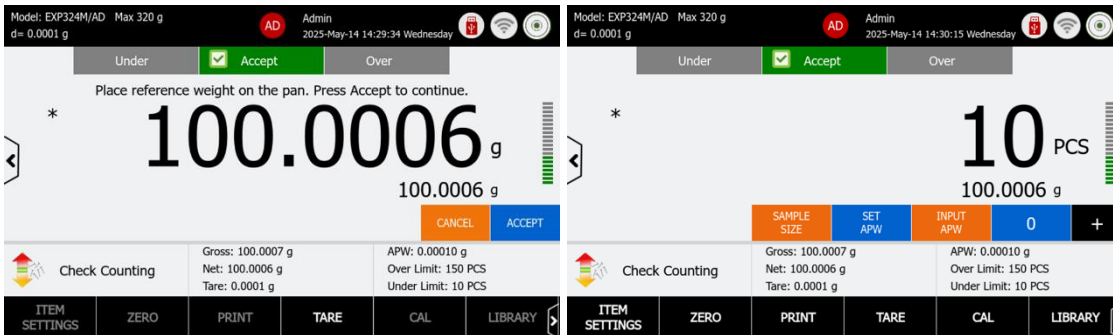

Step 2: Establish Over Limit and Under Limit and check the result

- The Over Limit and Under Limit are set in Library.
- The main screen will display the check status of the current sample in the library.
 - If the check status is either Under or Over, the lights will turn red to indicate this.
 - When the piece value exceeds the over limits, the screen will display Over.
 - When the piece value exceeds the under limits, the screen will display Under.
 - When the piece value is in the accept range, the screen will display Accept.
- Example: The home screen shows 30 pieces at the **Accept Range**.

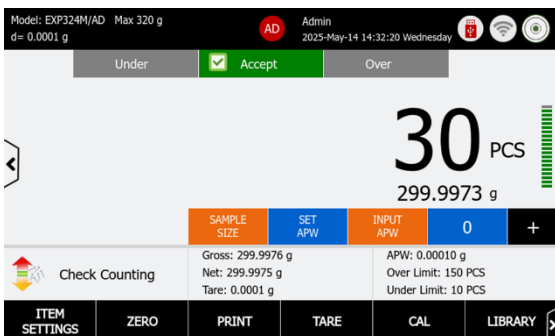


4.4.3 Accumulate the data for Check Counting

Utilize the Statistics "+" function to aggregate samples based on counting numbers and generate statistical reports.

The balance can store up to 999 accumulation records.


Step 1: Set the Statistic Mode in Parts Counting Mode

- Statistics can be determined Manual mode or automatically
- Automatic Mode: Weights are automatically recorded when stable

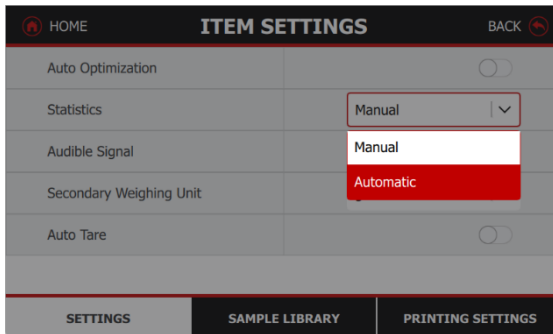

Step 2: Statistics Report View

- To use the statistics function, enable it first in the Item Settings.
- Add items and press the "+ Statistic button to accumulate the statistical data
- Tap the statistics number to view the statistical results.
- Statistics Report View
 - The report content includes Start Time, Number of Samples, Total, Average, Maximum, Minimum, Range, Standard Deviation, Relative Deviation % and individual item weighing values.

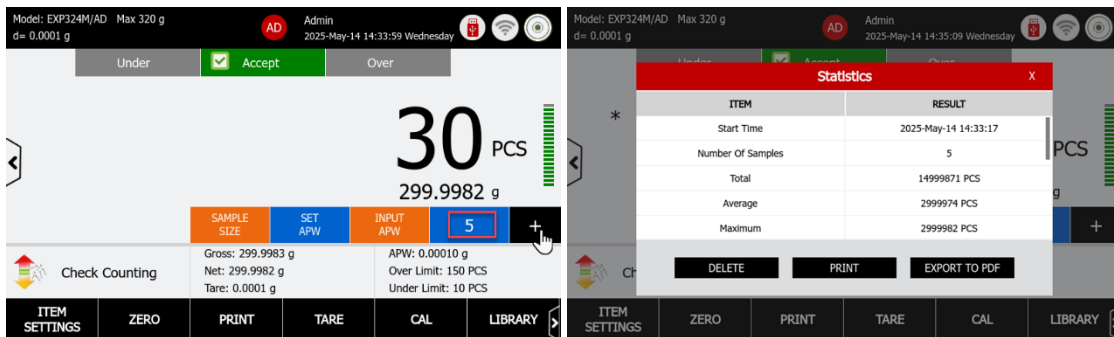

Step 3: Print Report

After reviewing the report, users can choose from several actions:


- Delete the statistics and restart the percent weighing process
- Print the report to a printer or PC based on the Print Settings
- Export the report as a PDF file to a USB flash drive

4.4.4 Sample Library – Check Counting

The Explorer Plus balance features a built-in library for managing multiple sample profiles. Up to 3,000 library records can be saved and recalled for use. The total library capacity and records are accessible across various application modes and through the library menu. When the library usage exceeds 80%, a pop-up message will alert the user to export the library or delete all items. The entire library menu, please refer to 6.5.1

Create, Active and Delete a Library

- Press “+New Library” button to create a new Library.
- When a new library is created, it remains inactive. The operator has to activate the library after ticking the box.
- Press on the RUBISH BIN button to delete a library.
- After the library item is activated, the main screen will show the activated library number.
- Other sample information can be configured in the library, including sample name, sample ID, batch ID, lot ID, project ID, and the creation of 10 more customized ID.

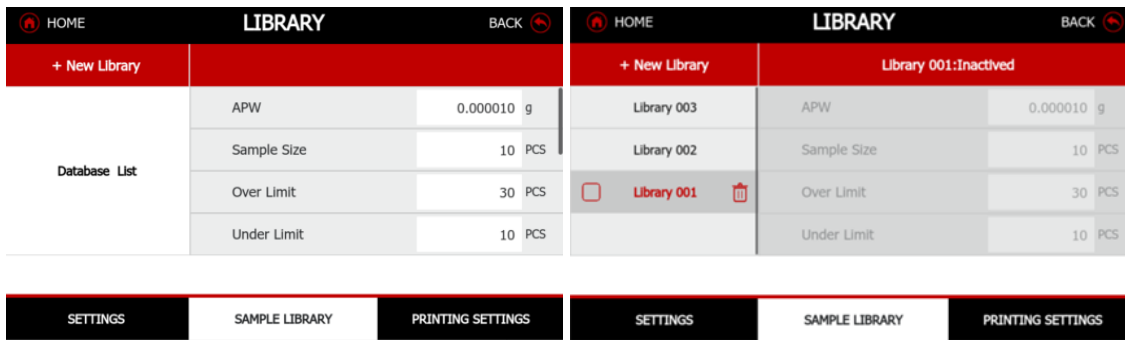

4.4.5 Printing Settings


The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to 6.1 Print Setting

4.5 Percent Weighing

Use this application to displays the current weight as a percentage of a reference weight.

- In the lower portion of the home screen, select Percent Weighing.
- Place objects on the pan to display the percentage of a reference weight.
- Press **Tare** or **Zero** to start weighing.
- The default Reference weight is displayed.
- Setup the Reference Weight, Reference Factor or Set Reference Weight before starting Percent Weighing application.

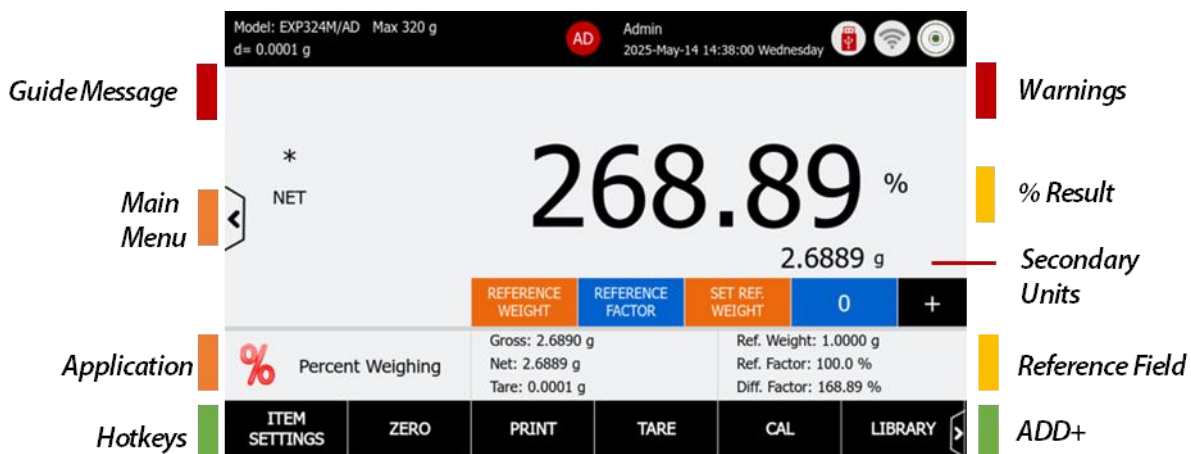


4.5.1 Application buttons and Reference Field

| Application Button/ Reference Field | Description |
|--|--|
| Reference Weight | In manufacturing or quality control processes, a reference weight can be used to verify that products meet specified weight standards. The reference weight can be set within the range of 100d to the maximum capacity. |
| Reference Factor | a reference factor is a known value used to adjust the Reference Weight |
| Set Ref. Weight | Place the sample on the pan, the balance will store the reference weight. |
| 0 | The statistic key is to recall the number of percent weighing, and press this button, operator can review the statics results. |
| + | Press this button to manually accumulate weight. In Automatic Statistics mode, this button will be inactive. |
| Item Settings | <ul style="list-style-type: none"> • Statistics: Automatic, Manual • Secondary Weighing Units: able to use 15 weighing units and 2 customer units • Auto Tare |
| Ref. Factor | <ul style="list-style-type: none"> • It displays the percentage of the current reference weight, with the unit in %, accurate to one decimal place. |
| Diff. Factor | <ul style="list-style-type: none"> • It displays the percentage difference between the sample weight and the predetermined reference weight in weighing calculations. • The unit is %, accurate to two decimal places. |

4.5.2 Begin Percent Weighing


Set a Known Reference Weight

- Press the REFERENCE WEIGHT button and enter the desired Reference Weight value.
- Tap ENTER to save the value
- Example: The home screen displays 30.00000 g as the Reference Weight.

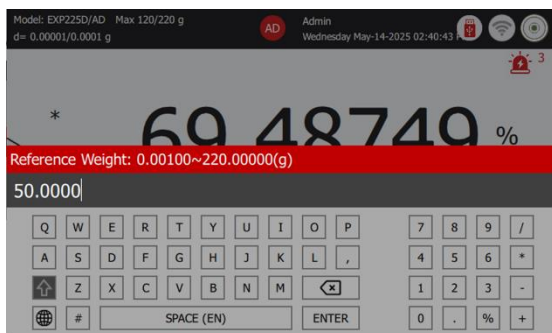

Set a new Reference Weight – Derived


- Place reference weight on the pan, and press Accept to establish a new Reference Weight
- The home screen displays 50.0013 g at the new Reference Weight

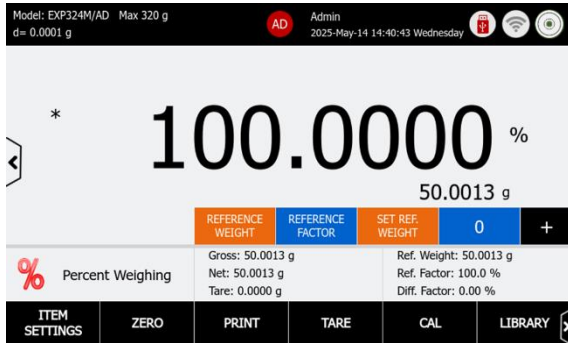

Step 2: Establish a Reference Factor

Adjust the reference in a desired factor.

- Press the REFERENCE FACTOR button and set the factor percentage.
- Tap ENTER to save the value
- If a library is activated, the new Reference Weight will overwrite the existing value in the library.
- Example: The home screen displays an 80% Ref. Factor for adjusting the Reference Weight.

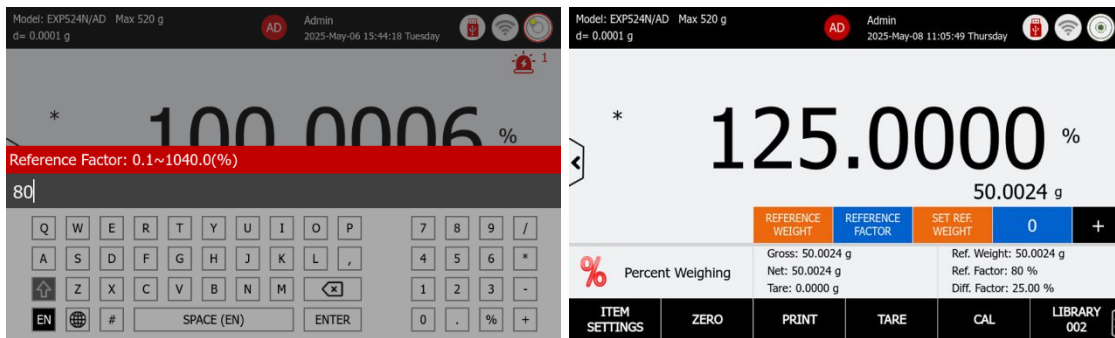

Step 3: Weighing the Samples

Place an object on the pan. The difference between the sample and the Reference Weight is displayed as both a weight and a percentage.

Place samples on the pan, the balance will display the percentage of reference weight. The reference weight value, the set of ref factor and difference factor will display in the reference field.

4.5.3 Accumulate the data for Percent Weighing


Utilize the Statistics "+" function to aggregate samples based on the numbers of samples and generate statistical reports. The balance can store up to 999 accumulation records.

Step 1: Set the Statistic Mode in Percent Weighing Mode


- Statistics can be determined Manual mode or automatically
- Automatic Mode: Weights are automatically recorded when stable

Step 2: Statistics Report View

- To use the statistics function, enable it first in the Item Settings.
- Add items and press the "+ Statistic button to accumulate the statistical data

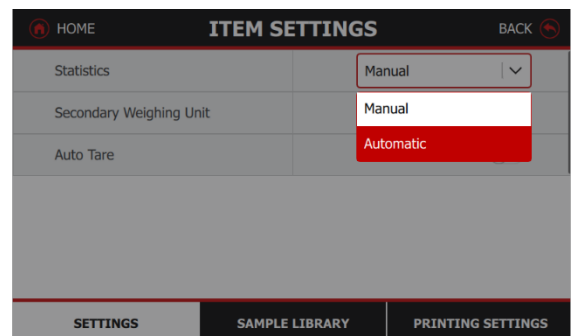

- Tap the statistics number to view the statistical results.
- Statistics Report View
 - The report content includes Start Time, Number of Samples, Total, Average, Maximum, Minimum, Range, Standard Deviation, Relative Deviation % and individual item percent weighing values.

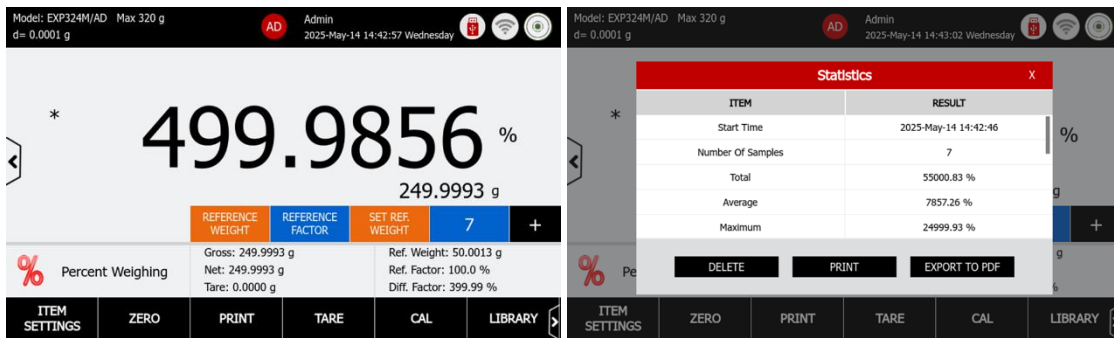

Step 3: Print Report

After reviewing the report, users can choose from several actions:


- Delete the statistics and restart the percent weighing process
- Print the report to a printer or PC based on the Print Settings
- Export the report as a PDF file to a USB flash drive

4.5.4 Sample Library – Check Counting

The Explorer Plus balance features a built-in library for managing multiple sample profiles. Up to 3,000 library records can be saved and recalled for use. The total library capacity and records are accessible across various application modes and through the library menu. When the library usage exceeds 80%, a pop-up message will alert the user to export the library or delete all items. The entire library menu, please refer to 6.5.1

Create a Library Record

- Press “+New Library” button to create a new Library.
- When a new library is created, it remains inactive. The operator has to activate the library after ticking the box.
- Press on the RUBISH BIN button to delete a library.
- After the library item is activated, the main screen will show the activated library number.

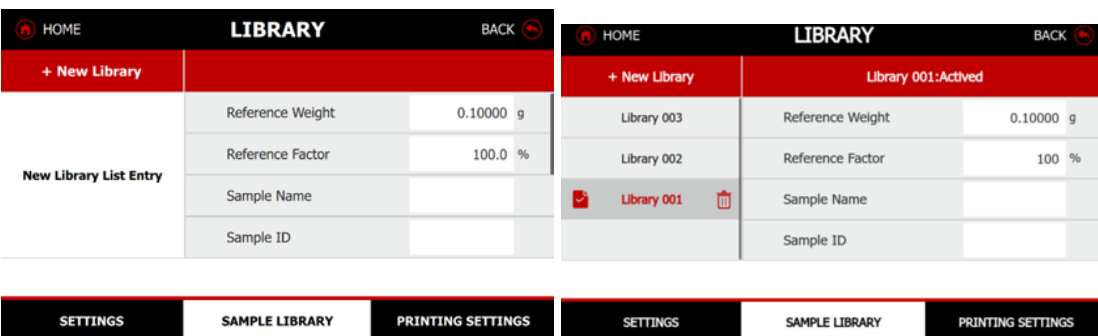

- Other sample information can be configured in the library, including sample name, sample ID, batch ID, lot ID, project ID, and the creation of 10 more customized ID.


4.5.5 Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to 6.1 Print Setting


4.6 Check Weighing

- Use this application to checks if the current sample pieces are within tolerances (e.g. an over and under limit, nominal weight vs. \pm tolerance).
 - This function serves to alert the operator of the check status and overload lights rather than requiring them to read the weighing value from the working distance
- In the lower portion of the home screen, select Check Weighing.
- Press **Tare** or **Zero** to start weighing.
- Place objects on the pan to check if the number of pieces within the tolerances. The default over and under limit pieces is displayed.
- Setup the Over and Under Limits, and or Nominal weight & Tolerance before starting Check Weighing application.

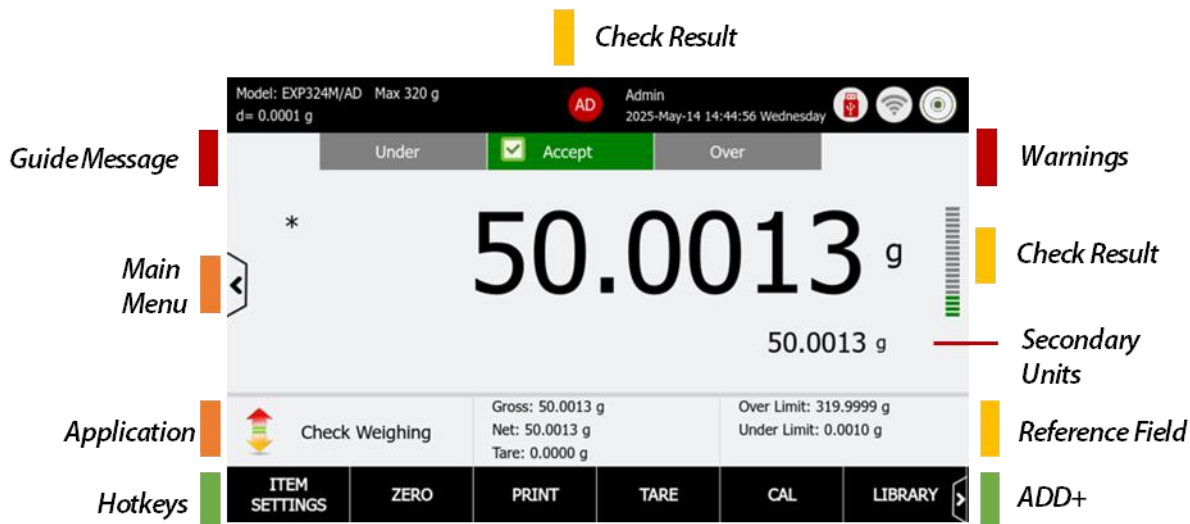

4.6.1 Reference Field

| Reference Field | Description |
|-----------------|--|
| Over Limit | The piece value exceeds the maximum allowed threshold. |
| Under Limit | The piece value exceeds the maximum allowed threshold. |
| Item Settings | <ul style="list-style-type: none"> • Limits Setting Mode: check result with different limits setting. <ul style="list-style-type: none"> ■ Under and Over ■ Nominal weight and Tolerance ■ Nominal Weight and Percent Tolerance • Display: Weight or Check Statues • Audible Signal: The balance beeper will sound to alert the user of the check status. <ul style="list-style-type: none"> ■ Available Setting: Off, Under, Accept, Over, Under and Over • Primary Weighing Unit: The default unit is grams. The operator can switch to alternative weighing units and two custom units. |

| | |
|---|--|
| | <ul style="list-style-type: none"> • Secondary Weighing Units: able to use 15 weighing units and 2 custom units |
|  | <ul style="list-style-type: none"> • The check statuses are divided into 20 segments to display the proportion of the current load relative to the total capacity. Each segment corresponds to 5% of the total capacity. • Under-Orange indicator • Accept-Green Indicator • Over-Red Indicator |


4.6.2 Begin Check Weighing

Step 1: Establish an Over Limits and Under Limits

- The Over Limit and Under Limit are set in Library.
- The main screen will display the check status of the current sample in the library.
- If the check status is either Under or Over, the status lights will turn red to indicate this.
- When the piece value exceeds the maximum, the screen will display Over.
 - When the piece value exceeds the maximum, the screen will display Under.
 - When the piece value is in the accept range, the screen will display Accept.


In ITEM SETTINGS, select the Mode as Over and Under, Nominal Weight Tolerance, or Nominal Percent Tolerance.




Example: The home screen displays 319.9999 gram as Over Limited Value, and 50.0013 g is in the range of Accept.

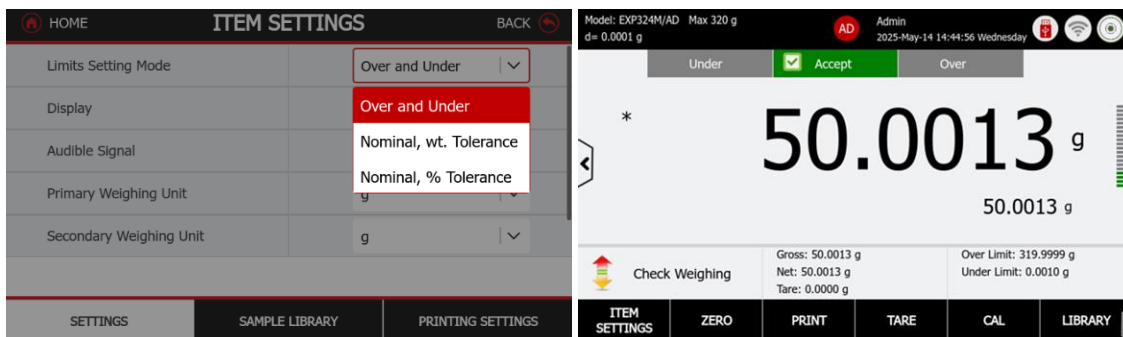

Step 2: Set the check result display as Under, Accept and Over


This function serves to alert the operator of the check status and overload lights rather than requiring them to read the weighing value from the working distance

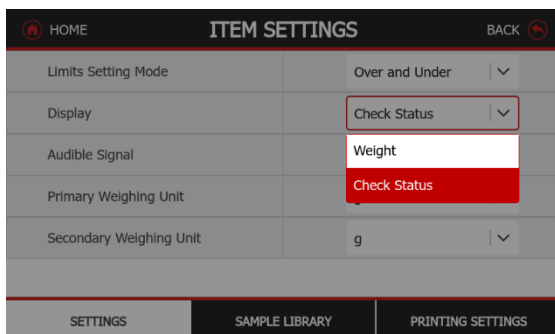

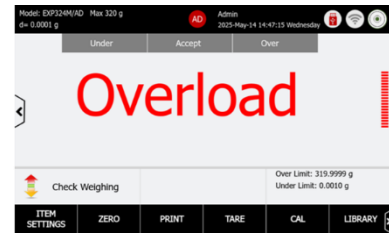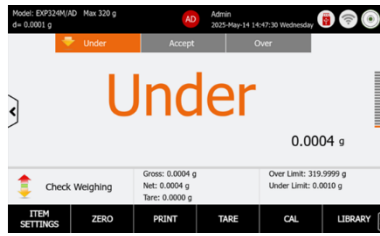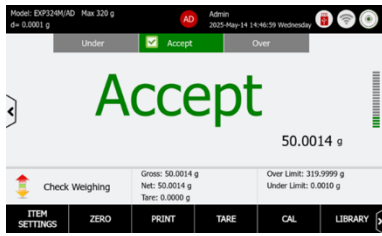

Step 3: Set up the Audible Signal

The operator can configure check sounds for Under, Accept, and Over in the Item Settings menu. Available Setting: Off, Under, Accept, Over, Under and Over

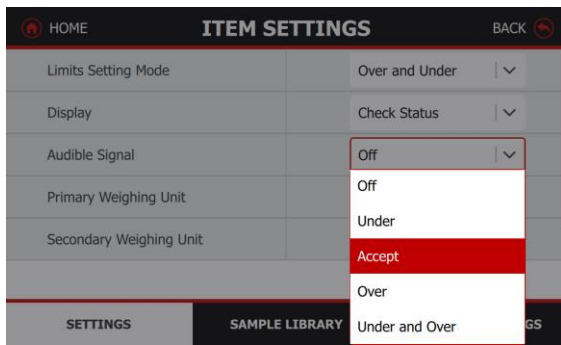


4.6.3 Sample Library – Check Weighing


The Explorer Plus balance features a built-in library for managing multiple sample profiles. Up to 3,000 library records can be saved and recalled for use. The total library capacity and records are accessible across various application modes and through the library menu. When the library usage exceeds 80%, a pop-up message will alert the user to export the library or delete all items. The entire library menu, please refer to 6.5.1

Create, Active and Delete a Library


- Press “+New Library” button to create a new Library.
- When a new library is created, it remains inactive. The operator has to activate the library after ticking the box.
- Press on the RUBISH BIN button to delete a library.
- After the library item is activated, the main screen will show the activated library number.
- Example: Limits setting mode is Under and Over

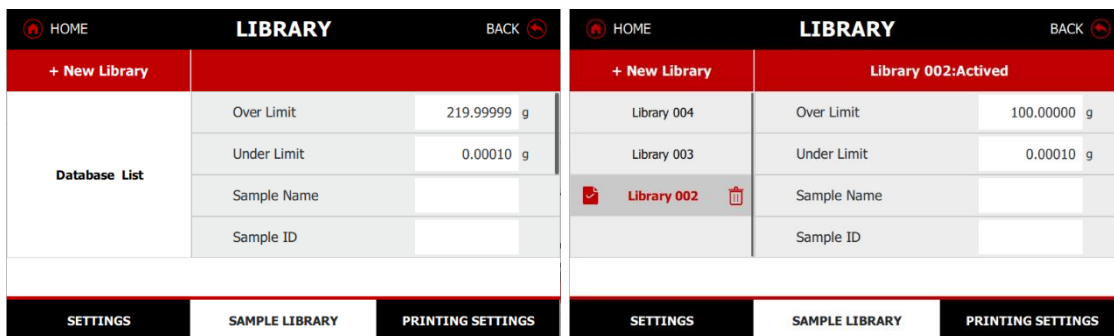

Example: Nominal weight and Tolerance

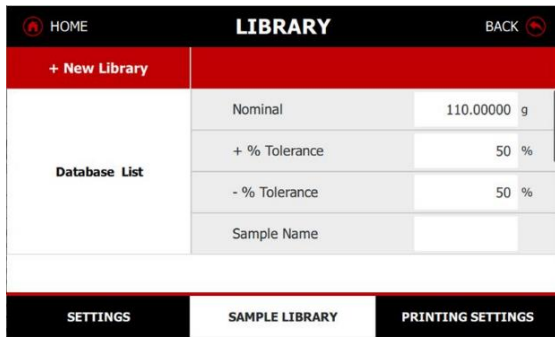

Example: Nominal Weight and Percent Tolerance

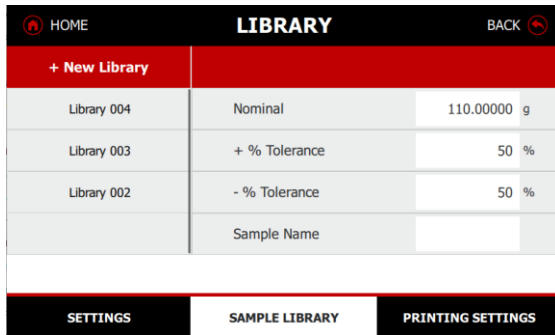

Other sample information can be configured in the library, including sample name, sample ID, batch ID, lot ID, project ID, and the creation of 10 more customized ID.

4.6.4 Printing Settings


The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to 6.1 Printing Setting.

4.7 Dynamic Weighing

This function serves to is set to weigh objects that are not stable, such as animals.

- In the lower portion of the home screen, select Dynamic Weighing.
- Press **Tare** or **Zero** to start weighing.
- Place the moving sample on the pan to achieve an averaged weight reading within seconds. The default weighing average time is displayed.
- Setup the Start Mode, Weighing Average Time, Auto Print Result before starting Check Weighing application.

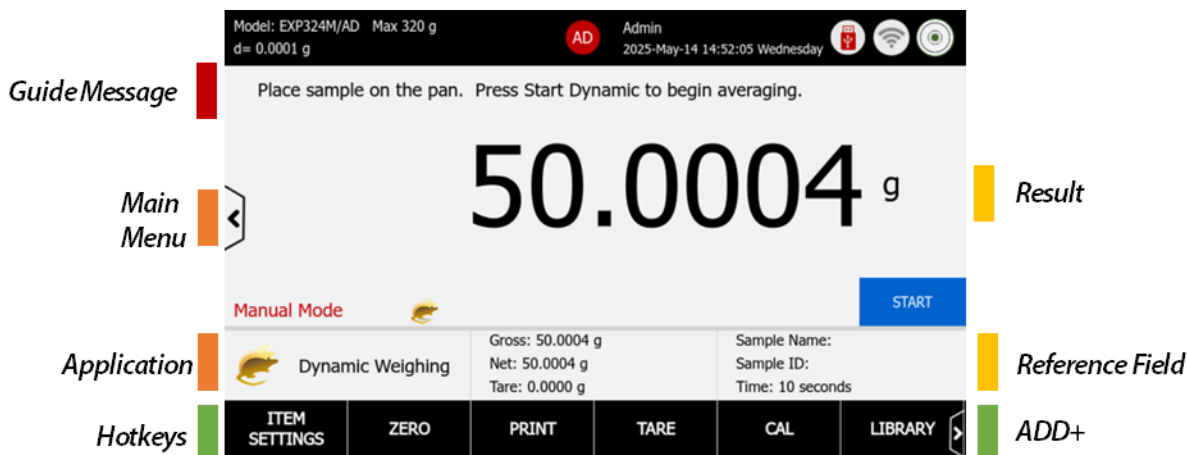

4.7.1 Application buttons


| Application Button/ Reference Field | Description |
|--|--|
| Start | Press “Start” button to begin averaging. |
| Time | It displays the averaging time. |
| Item Settings | <ul style="list-style-type: none"> Start Mode: Automatic and Manual Automatic start without press Start button after second dynamic weighing Auto Print Result: The averaged result will be printed immediately, without press “Print” button in the main screen. Main Unit: The default setting is gram. It is able to change to other weighing units Auto Tare: Automatic tare the container value |

4.7.2 Begin Check Weighing

Step 1: Setup the weighing averaging time in the Sample Library

- The weighing averaging time are set in Library.
- The main screen will display the averaging weight of the current sample.


Example: The screen displays 10 seconds as the averaging time.

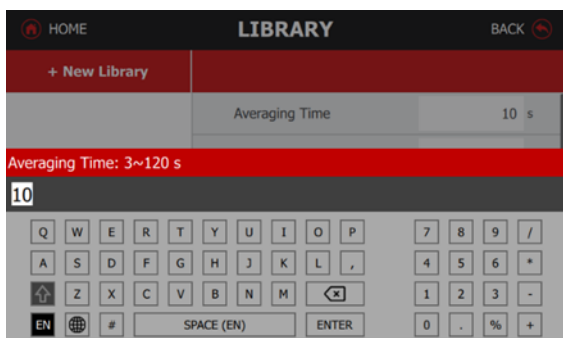

Step 2: Weigh movement sample on the weighing pan

- Place a sample on the weighing pan, and the balance will immediately calculate the average weight.

- The process will then commence a countdown based on the pre-set time configuration.
- The moving mouse icon will be stop at the end of the process, the result will be highlighted.

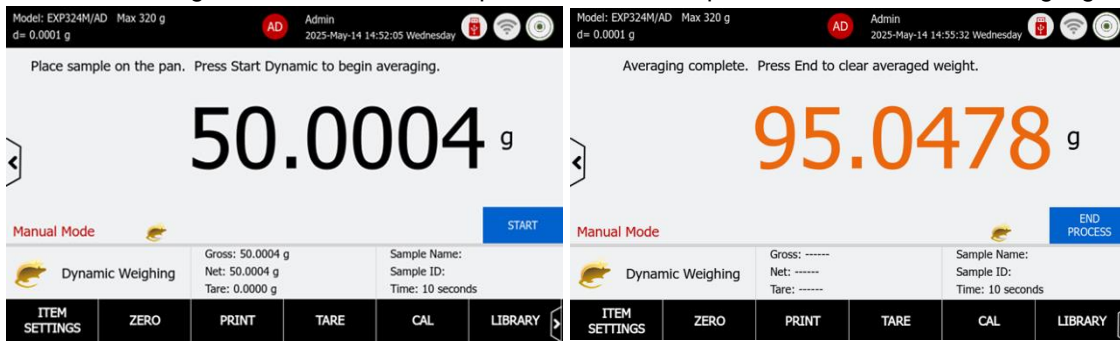

4.7.3 Sample Library – Dynamic Weighing

The Explorer Plus balance features a built-in library for managing multiple sample profiles. Up to 3,000 library records can be saved and recalled for use. The total library capacity and records are accessible across various application modes and through the library menu. When the library usage exceeds 80%, a pop-up message will alert the user to export the library or delete all items. The entire library menu, please refer to 6.5.1 Application Print Out Template

| Basic Weighing | | Parts Counting | | Check Counting | |
|-----------------------|----------------|-----------------------|----------------|-----------------------|----------------|
| 12-24-2024 | 09:31:46 | 12-24-2024 | 10:13:34 | 12-24-2024 | 10:17:42 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN058 | Project Name: | PN070 | Project Name: | PN081 |
| ----- | | ----- | | ----- | |
| Application Name: | Basic Weighing | Application Name: | Parts Counting | Application Name: | Check Counting |
| Sample Name: | SN058 | Sample Name: | SN070 | Sample Name: | SN081 |
| Sample ID: | S058 | Sample ID: | S070 | Sample ID: | S081 |
| Batch ID: | B058 | Batch ID: | B070 | Batch ID: | B081 |
| Lot ID: | L058 | Lot ID: | L070 | Lot ID: | L081 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | Customer ID2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 3.5275 | Result: | 25 PCS | Result: | 74 PCS |
| oz N | | Gross: | 78.48 g | Status: | Accept |
| Gross: | 6.2960 | Net: | 78.48 g N | Gross: | 178.49 g |
| oz | | Tare: | 0.00 g T | Net: | 178.49 g N |
| Net: | 3.5275 oz N | | | Tare: | 0.00 g T |
| Tare: | 2.7690 oz T | Library: | Library 070 | | |
| | | APW: | 3.124 g | Library: | Library 081 |
| Library: | Library 058 | Samples: | 10 PCS | APW: | 2.412 g |
| | | | | Samples: | 81 PCS |
| Signature: _____ | | Signature: _____ | | Over: | 254 PCS |
| | | | | Under: | 51 PCS |
| Verify By: _____ | | Verify By: _____ | | | |
| | | | | Signature: _____ | |
| | | | | Verify By: _____ | |

| Percent Weighing | | Check Weighing | | Dynamic | |
|-----------------------|------------------|-----------------------|----------------|-----------------------|------------------|
| 12-26-2024 | 13:30:04 | 12-26-2024 | 13:33:25 | 12-26-2024 | 13:36:21 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID : | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN010 | Project Name: | PN028 | Project Name: | PN055 |
| ----- | | ----- | | ----- | |
| Application Name: | Percent Weighing | Application Name: | Check Weighing | Application Name: | Dynamic Weighing |
| Sample Name: | SN010 | Sample Name: | SN028 | Sample Name: | SN055 |
| Sample ID: | S010 | Sample ID: | S028 | Sample ID: | S055 |
| Batch ID: | B010 | Batch ID: | B028 | Batch ID: | B055 |
| Lot ID: | L010 | Lot ID: | L028 | Lot ID: | L055 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | C2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 361.69 % | Result: | Accept | Result: | 367.00 g |
| Gross: | 198.93 g | Gross: | 198.92 g | Gross: | 198.93 g |
| Net: | 198.93 g N | Net: | 198.92 g N | Net: | 198.93 g N |
| Tare: | 0.00 g T | Tare: | 0.00 g T | Tare: | 0.00 g T |
| Library: | Library 010 | Library: | Library 028 | Library: | Library 055 |
| Reference Weight: | 550.00 g | Over Limit: | 214.25 g | Averaging Time: | 3 seconds |
| Reference Factor: | 10.0 % | Under Limit: | 12.54 g | Signature: | _____ |
| Difference Weight: | 143.93 g | Signature: | _____ | Verify By: | _____ |
| Difference Factor: | 261.69 % | Verify By: | _____ | | |
| Signature: | _____ | | | | |
| Verify By: | _____ | | | | |

| Totalization | | Formulation | | Differential | |
|---------------------------|---------------------|-----------------------|---------------------|-----------------------|----------------------|
| 12-26-2024 | 15:11:42 | 12-26-2024 | 16:30:04 | 2025-Apr-08 | 19:58:46 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN072 | Project Name: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Name: | Totalization | Application Name: | Formulation | Application Mode: | Differential |
| Sample Name: | SN072 | Result: | 199.09 g N | Result: | 3.89[3] g N |
| Sample ID: | S072 | Gross: | 299.09 g | Gross: | 18.42[0] g |
| Batch ID: | B072 | Net: | 199.09 g N | Net: | 3.89[3] g N |
| Lot ID: | L072 | Tare: | 100.00 g T | Tare: | 14.52[7] g T |
| Customer ID1: | C1 | Recipe: | Recipe 001 | Start Time: | 2025-Apr-08 19:57:07 |
| Customer ID2: | C2 | Item Name | Item Weight | End Time: | 2025-Apr-08 19:58:41 |
| Customer ID3: | C3 | Item 1 | 100.00 g | Reciprocal Proportion | On |
| Customer ID4: | C4 | Item 2 | 199.08 g | Absolute Value | On |
| Customer ID5: | C5 | Start Time: | 12-26-2024 16:29:44 | Item Name: | ITEM1 |
| Customer ID6: | C6 | End Time: | 12-26-2024 16:30:01 | Initial Weight: | 0.766 g |
| Customer ID7: | C7 | -----Sample Data----- | | Final Weight: | 0.760 g |
| Customer ID8: | C8 | Item Name: | Item 1 | Difference Weight: | 0.006 g |
| Customer ID9: | C9 | Target Weight: | 100.00 g | Different Percent: | 0.8 % |
| Customer ID10: | C10 | Actual Weight: | 100.00 g | Item Name: | ITEM2 |
| Result: | 278.33 g | Difference: | 0.00 % | Initial Weight: | 1.528 g |
| Gross: | 0.00 g | Item Name: | Item 2 | Final Weight: | 1.534 g |
| Net: | 0.00 g N | Target Weight: | 199.08 g | Difference Weight: | 0.006 g |
| Tare: | 0.00 g T | Actual Weight: | 199.08 g | Different Percent: | 0.4 % |
| Library: | Library 072 | Difference: | 0.00 % | Item Name: | ITEM3 |
| Start Time: | 12-26-2024 15:08:29 | Total Weight: | 299.08 g | Initial Weight: | 2.292 g |
| End Time: | 12-26-2024 15:11:42 | Signature: _____ | | Final Weight: | 2.325 g |
| -----Sample Data (g)----- | | Verify By: _____ | | Difference Weight: | 0.033 g |
| '1: | 198.93 | | | Different Percent: | 1.4 % |
| '2: | 19.85 | | | Item Name: | ITEM4 |
| '3: | 19.86 | | | Initial Weight: | 3.094 g |
| '4: | 19.87 | | | Final Weight: | 3.110 g |
| '5: | 19.88 | | | Difference Weight: | 0.016 g |
| Samples: | 5 | | | Different Percent: | 0.5 % |
| Average: | 55.67 g | | | Item Name: | ITEM5 |
| Maximum: | 198.93 g | | | Initial Weight: | 3.871 g |
| Minimum: | 19.85 g | | | Final Weight: | 3.893 g |
| Range: | 179.08 g | | | Difference Weight: | 0.022 g |
| Standard Deviation: | 71.63 g | | | Different Percent: | 0.6 % |
| Relative Deviation(std%): | 128.68 % | | | Signature: _____ | |
| Signature: _____ | | | | Verified By: _____ | |
| Verify By: _____ | | | | | |

| Density Determination | | Peak Hold | | Pipette Adjustment | |
|-----------------------|---------------------------|-----------------------|-----------|---------------------------------------|---------------------------|
| 2025-Apr-08 | 20:01:25 | 2025-Apr-08 | 20:25:41 | 025-Apr-08 | 20:11:58 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| ----- | | ----- | | ----- | |
| Application Mode: | Density Determination | Application Mode: | Peak Hold | Application Mode: | Pipette Adjustment |
| Result: | 7.435 g/cm ³ | Result: | 355.512 g | Result: | Pass |
| Gross: | 19.44[1] g | Gross: | 0.000 g | Gross: | 24.76[4] g ? |
| Net: | 4.91[4] g N | Net: | 0.000 g N | Net: | 1.01[1] g ? N |
| Tare: | 14.52[7] g T | Tare: | 0.000 g T | Tare: | 23.75[3] g T |
| Weight in air.: | 5.67[4] g | On Stability: | No | Start Time: | 2025-Apr-08 20:09:03 |
| Weight in liquid.: | 4.91[3] g | Signature: _____ | | End Time: | 2025-Apr-08 20:11:52 |
| Auxiliary liquid.: | Water | Verified By: _____ | | Liquid Density: | 0.99823 g/cm ³ |
| Liquid Density: | 0.99823 g/cm ³ | | | Pressure: | 1.0 ATM |
| Water Temp.: | 20.0 °C | | | Pipette Test Method: | Method 001 |
| Porous Material: | Off | | | Nominal Inaccuracy Imprecision Number | 1000 ul 5.00 % 5.00 % 3 |
| Signature: _____ | | | | Pipette Test Results | |
| Verified By: _____ | | | | ----- | |
| | | | | Pipette Name: | pip001 |
| | | | | Pipette Number: | |
| | | | | Nominal Volume: | 1000 ul Pass |
| | | | | -----Sample Data----- | |
| | | | | 1 | 1023 ul 1.021 g |
| | | | | 2 | 1032 ul 1.030 g |
| | | | | 3 | 1016 ul 1.014 g |
| | | | | Inaccuracy | |
| | | | | E%: | 2.64 % |
| | | | | E% Limit: | 5.00 % |
| | | | | Average: | 1026 ul |
| | | | | Imprecision | |
| | | | | CV%: | 0.97 % |
| | | | | CV% Limit: | 5.00 % |
| | | | | Standard Deviation: | 10 ul |
| | | | | > +2S: | 0.00 % , 0 |
| | | | | > +1S: | 0.00 % , 0 |
| | | | | +1S > Mean > -1S: | 0.00 % , 0 |
| | | | | < -1S: | 0.00 % , 0 |
| | | | | < -2S: | 100.00 % , 10 |
| | | | | Signature: _____ | |
| | | | | Verified By: _____ | |

| SQC | | Fill Weight Variation | | Flow Rate Control | |
|------------------------------------|-----------------|-----------------------|-----------------------|------------------------|-------------------|
| 2025-Apr-08 | 20:13:55 | 2025-Apr-08 | 20:17:46 | 2025-Apr-08 | 19:49:45 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | | Project ID: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Mode: | SQC | Application Mode: | Fill Weight Variation | Application Mode: | Flow Rate Control |
| Result: | Accept | Sample Name: | | Result: | |
| Average: | 100.010 g | Sample ID: | | Max Flow Rate: | 0.543 ml/min |
| Minimum: | 100.010 g | Batch ID: | | Min Flow Rate: | 0.526 ml/min |
| Maximum: | 100.011 g | Lot ID: | | Average Flow Rate: | 0.533 ml/min |
| Standard Deviation: | 0.000 g | Samples: | 5 | Pump Cycle: | 3 times |
| Test Result: | | Unit1: | 100.01[1] g | Test Result: | |
| Sample ID Weight T1/T2 | | Unit2: | 100.01[1] g | 1: | 0.529 ml/min |
| Sample 001 100.011 g >=-T1 & <=+T1 | | Unit3: | 100.01[1] g | 2: | 0.526 ml/min |
| Sample 002 100.010 g >=-T1 & <=+T1 | | Unit4: | 100.01[1] g | 3: | 0.543 ml/min |
| Sample 003 100.010 g >=-T1 & <=+T1 | | Unit5: | 100.01[1] g | Pump Speed: | 0.500 ml/min |
| Batch: | Batch 01 | Difference: | | Liquid Density: | 0.99823 g/ml |
| Number in Batch: | 100% Inspection | 1: 0.00[3] g 0.00 % | | Target Volume: | 10.00 ml |
| Inspection Quantity: | 3 | 2: 0.00[3] g 0.00 % | | Number of Pump Cycles: | 3 times |
| Tare Mode: | Off | 3: 0.00[3] g 0.00 % | | Mode: | Flow Rate Control |
| Pretare: | 0.000 g | 4: 0.00[3] g 0.00 % | | Signature: _____ | |
| Sample Type: | Solid | 5: 0.00[3] g 0.00 % | | Verified By: _____ | |
| EU Criteria: | Off | Result: | | | |
| Nominal Weight: | 100.000 g | Accept: | 5 Units | | |
| +T1: | 0.300 g | Fail: | 0 Units | | |
| -T1: | -0.300 g | Acceptance Criteria: | Off | | |
| +T2: | 0.600 g | Average Weight: | 100.00[8] g | | |
| -T2: | -0.600 g | +/- Limit %: | 7.50 % | | |
| Signature: _____ | | Signature: _____ | | | |
| Verified By: _____ | | Verified By: _____ | | | |

4.8 Printout Examples

Here are the examples for each application are displayed with all items enabled in the Print Content menu. The default values for Header lines 1-5 are also indicated.

Notes:

- When the user management and electronic signature are active, the User ID will display in the Signature Line. The electronic signature is disabled by default.
- When a library is activated, "Library Name:" will appear below "Application" in the printout.

4.8.1 Calibration Report Template

| Internal Calibration | Span Calibration | Repeatability Test | | | | | | | | | | | | | | | | | | | | | |
|--|--|---|--------|-----------|-----------|---|-----------|------------|---|------------|------------|---|------------|------------|---|------------|------------|---|-----------|------------|---|-----------|------------|
| Calibration Report -----Internal Calibration----- Date and Time: 2025-May-12 10:34:02 Balance ID: Balance Name: EXP224/AD Result: 0.0000g Difference Internal Calibration Successful Signature: User ID Verified By: _____ | Calibration Report -----Span Calibration----- Date and Time: 2025-May-12 10:34:02 Balance ID: Balance Name: EXP224/AD Result: 0.0000g Difference Span Calibration Successful Signature: User ID Verified By: _____ | Routine Test: Repeatability Test Date: 2025-Apr-11 Start Time: 19:13:45 End Time: 19:15:52 Test Weights ID: Test Weights Value: 25.00000 g Test Weights Class: Test Result: <table border="1"> <thead> <tr> <th>Number</th> <th>Zero Load</th> <th>Full Load</th> </tr> </thead> <tbody> <tr> <td>1</td> <td>0.00000 g</td> <td>25.00496 g</td> </tr> <tr> <td>2</td> <td>-0.00002 g</td> <td>25.00262 g</td> </tr> <tr> <td>3</td> <td>-0.00013 g</td> <td>25.00063 g</td> </tr> <tr> <td>4</td> <td>-0.00006 g</td> <td>25.01753 g</td> </tr> <tr> <td>5</td> <td>0.00001 g</td> <td>25.00375 g</td> </tr> <tr> <td>6</td> <td>0.00002 g</td> <td>25.00014 g</td> </tr> </tbody> </table> SD (Span): 0.006441 g Signature: _____ Verified By: _____ | Number | Zero Load | Full Load | 1 | 0.00000 g | 25.00496 g | 2 | -0.00002 g | 25.00262 g | 3 | -0.00013 g | 25.00063 g | 4 | -0.00006 g | 25.01753 g | 5 | 0.00001 g | 25.00375 g | 6 | 0.00002 g | 25.00014 g |
| Number | Zero Load | Full Load | | | | | | | | | | | | | | | | | | | | | |
| 1 | 0.00000 g | 25.00496 g | | | | | | | | | | | | | | | | | | | | | |
| 2 | -0.00002 g | 25.00262 g | | | | | | | | | | | | | | | | | | | | | |
| 3 | -0.00013 g | 25.00063 g | | | | | | | | | | | | | | | | | | | | | |
| 4 | -0.00006 g | 25.01753 g | | | | | | | | | | | | | | | | | | | | | |
| 5 | 0.00001 g | 25.00375 g | | | | | | | | | | | | | | | | | | | | | |
| 6 | 0.00002 g | 25.00014 g | | | | | | | | | | | | | | | | | | | | | |

4.8.2 Application Print Out Template

| Basic Weighing | | Parts Counting | | Check Counting | |
|-----------------------|----------------|-----------------------|----------------|-----------------------|----------------|
| 12-24-2024 | 09:31:46 | 12-24-2024 | 10:13:34 | 12-24-2024 | 10:17:42 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN058 | Project Name: | PN070 | Project Name: | PN081 |
| ----- | | ----- | | ----- | |
| Application Name: | Basic Weighing | Application Name: | Parts Counting | Application Name: | Check Counting |
| Sample Name: | SN058 | Sample Name: | SN070 | Sample Name: | SN081 |
| Sample ID: | S058 | Sample ID: | S070 | Sample ID: | S081 |
| Batch ID: | B058 | Batch ID: | B070 | Batch ID: | B081 |
| Lot ID: | L058 | Lot ID: | L070 | Lot ID: | L081 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | Customer ID2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 3.5275 | Result: | 25 PCS | Result: | 74 PCS |
| oz N | | Gross: | 78.48 g | Status: | Accept |
| Gross: | 6.2960 | Net: | 78.48 g N | Gross: | 178.49 g |
| oz | | Tare: | 0.00 g T | Net: | 178.49 g N |
| Net: | 3.5275 oz N | | | Tare: | 0.00 g T |
| Tare: | 2.7690 oz T | Library: | Library 070 | | |
| | | APW: | 3.124 g | Library: | Library 081 |
| Library: | Library 058 | Samples: | 10 PCS | APW: | 2.412 g |
| | | | | Samples: | 81 PCS |
| Signature: _____ | | Signature: _____ | | Over: | 254 PCS |
| | | | | Under: | 51 PCS |
| Verify By: _____ | | Verify By: _____ | | Signature: _____ | |
| | | | | Verify By: _____ | |

| Percent Weighing | | Check Weighing | | Dynamic | |
|-----------------------|------------------|-----------------------|----------------|-----------------------|------------------|
| 12-26-2024 | 13:30:04 | 12-26-2024 | 13:33:25 | 12-26-2024 | 13:36:21 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID : | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN010 | Project Name: | PN028 | Project Name: | PN055 |
| ----- | | ----- | | ----- | |
| Application Name: | Percent Weighing | Application Name: | Check Weighing | Application Name: | Dynamic Weighing |
| Sample Name: | SN010 | Sample Name: | SN028 | Sample Name: | SN055 |
| Sample ID: | S010 | Sample ID: | S028 | Sample ID: | S055 |
| Batch ID: | B010 | Batch ID: | B028 | Batch ID: | B055 |
| Lot ID: | L010 | Lot ID: | L028 | Lot ID: | L055 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | C2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 361.69 % | Result: | Accept | Result: | 367.00 g |
| Gross: | 198.93 g | Gross: | 198.92 g | Gross: | 198.93 g |
| Net: | 198.93 g N | Net: | 198.92 g N | Net: | 198.93 g N |
| Tare: | 0.00 g T | Tare: | 0.00 g T | Tare: | 0.00 g T |
| Library: | Library 010 | Library: | Library 028 | Library: | Library 055 |
| Reference Weight: | 550.00 g | Over Limit: | 214.25 g | Averaging Time: | 3 seconds |
| Reference Factor: | 10.0 % | Under Limit: | 12.54 g | Signature: | _____ |
| Difference Weight: | 143.93 g | Signature: | _____ | Verify By: | _____ |
| Difference Factor: | 261.69 % | Verify By: | _____ | | |
| Signature: | _____ | | | | |
| Verify By: | _____ | | | | |

| Totalization | | Formulation | | Differential | |
|---------------------------|---------------------|-----------------------|---------------------|-----------------------|----------------------|
| 12-26-2024 | 15:11:42 | 12-26-2024 | 16:30:04 | 2025-Apr-08 | 19:58:46 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN072 | Project Name: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Name: | Totalization | Application Name: | Formulation | Application Mode: | Differential |
| Sample Name: | SN072 | Result: | 199.09 g N | Result: | 3.89[3] g N |
| Sample ID: | S072 | Gross: | 299.09 g | Gross: | 18.42[0] g |
| Batch ID: | B072 | Net: | 199.09 g N | Net: | 3.89[3] g N |
| Lot ID: | L072 | Tare: | 100.00 g T | Tare: | 14.52[7] g T |
| Customer ID1: | C1 | | | | |
| Customer ID2: | C2 | Recipe: | Recipe 001 | Start Time: | 2025-Apr-08 19:57:07 |
| Customer ID3: | C3 | Item Name | Item Weight | End Time: | 2025-Apr-08 19:58:41 |
| Customer ID4: | C4 | Item 1 | 100.00 g | | |
| Customer ID5: | C5 | Item 2 | 199.08 g | Reciprocal Proportion | On |
| Customer ID6: | C6 | | | Absolute Value | On |
| Customer ID7: | C7 | Start Time: | 12-26-2024 16:29:44 | | |
| Customer ID8: | C8 | End Time: | 12-26-2024 16:30:01 | Item Name: | ITEM1 |
| Customer ID9: | C9 | -----Sample Data----- | | Initial Weight: | 0.766 g |
| Customer ID10: | C10 | Item Name: | Item 1 | Final Weight: | 0.760 g |
| Result: | 278.33 g | Target Weight: | 100.00 g | Difference Weight: | 0.006 g |
| Gross: | 0.00 g | Actual Weight: | 100.00 g | Different Percent: | 0.8 % |
| Net: | 0.00 g N | Difference: | 0.00 % | | |
| Tare: | 0.00 g T | | | Item Name: | ITEM2 |
| | | Item Name: | Item 2 | Initial Weight: | 1.528 g |
| Library: | Library 072 | Target Weight: | 199.08 g | Final Weight: | 1.534 g |
| Start Time: | 12-26-2024 15:08:29 | Actual Weight: | 199.08 g | Difference Weight: | 0.006 g |
| End Time: | 12-26-2024 15:11:42 | Difference: | 0.00 % | Different Percent: | 0.4 % |
| | | | | | |
| -----Sample Data (g)----- | | Total Weight: | 299.08 g | Item Name: | ITEM3 |
| '1: | 198.93 | | | Initial Weight: | 2.292 g |
| '2: | 19.85 | Signature: _____ | | Final Weight: | 2.325 g |
| '3: | 19.86 | | | Difference Weight: | 0.033 g |
| '4: | 19.87 | Verify By: _____ | | Different Percent: | 1.4 % |
| '5: | 19.88 | | | | |
| | | | | Item Name: | ITEM4 |
| Samples: | 5 | | | Initial Weight: | 3.094 g |
| Average: | 55.67 g | | | Final Weight: | 3.110 g |
| Maximum: | 198.93 g | | | Difference Weight: | 0.016 g |
| Minimum: | 19.85 g | | | Different Percent: | 0.5 % |
| Range: | 179.08 g | | | | |
| Standard Deviation: | 71.63 g | | | Item Name: | ITEM5 |
| Relative Deviation(std%): | 128.68 % | | | Initial Weight: | 3.871 g |
| | | | | Final Weight: | 3.893 g |
| Signature: _____ | | | | Difference Weight: | 0.022 g |
| | | | | Different Percent: | 0.6 % |
| Verify By: _____ | | | | | |
| | | | | Signature: _____ | |
| | | | | Verified By: _____ | |

| Density Determination | | Peak Hold | | Pipette Adjustment | |
|-----------------------|---------------------------|-----------------------|-----------|---------------------------------------|---------------------------|
| 2025-Apr-08 | 20:01:25 | 2025-Apr-08 | 20:25:41 | 025-Apr-08 | 20:11:58 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| ----- | | ----- | | ----- | |
| Application Mode: | Density Determination | Application Mode: | Peak Hold | Application Mode: | Pipette Adjustment |
| Result: | 7.435 g/cm ³ | Result: | 355.512 g | Result: | Pass |
| Gross: | 19.44[1] g | Gross: | 0.000 g | Gross: | 24.76[4] g ? |
| Net: | 4.91[4] g N | Net: | 0.000 g N | Net: | 1.01[1] g ? N |
| Tare: | 14.52[7] g T | Tare: | 0.000 g T | Tare: | 23.75[3] g T |
| Weight in air.: | 5.67[4] g | On Stability: | No | Start Time: | 2025-Apr-08 20:09:03 |
| Weight in liquid.: | 4.91[3] g | Signature: _____ | | End Time: | 2025-Apr-08 20:11:52 |
| Auxiliary liquid.: | Water | Verified By: _____ | | Liquid Density: | 0.99823 g/cm ³ |
| Liquid Density: | 0.99823 g/cm ³ | | | Pressure: | 1.0 ATM |
| Water Temp.: | 20.0 °C | | | Pipette Test Method: | Method 001 |
| Porous Material: | Off | | | Nominal Inaccuracy Imprecision Number | |
| Signature: _____ | | | | 1000 ul 5.00 % 5.00 % 3 | |
| Verified By: _____ | | | | Pipette Test Results | |
| | | | | ----- | |
| | | | | Pipette Name: | pip001 |
| | | | | Pipette Number: | |
| | | | | Nominal Volume: | 1000 ul Pass |
| | | | | -----Sample Data----- | |
| | | | | 1 1023 ul 1.021 g | |
| | | | | 2 1032 ul 1.030 g | |
| | | | | 3 1016 ul 1.014 g | |
| | | | | Inaccuracy | |
| | | | | E%: | 2.64 % |
| | | | | E% Limit: | 5.00 % |
| | | | | Average: | 1026 ul |
| | | | | Imprecision | |
| | | | | CV%: | 0.97 % |
| | | | | CV% Limit: | 5.00 % |
| | | | | Standard Deviation: | 10 ul |
| | | | | > +2S: | 0.00 % , 0 |
| | | | | > +1S: | 0.00 % , 0 |
| | | | | +1S > Mean > -1S: | 0.00 % , 0 |
| | | | | < -1S: | 0.00 % , 0 |
| | | | | < -2S: | 100.00 % , 10 |
| | | | | Signature: _____ | |
| | | | | Verified By: _____ | |


| SQC | | Fill Weight Variation | | Flow Rate Control | |
|------------------------------------|-----------------|-----------------------|-----------------------|------------------------|-------------------|
| 2025-Apr-08 | 20:13:55 | 2025-Apr-08 | 20:17:46 | 2025-Apr-08 | 19:49:45 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | | Project ID: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Mode: | SQC | Application Mode: | Fill Weight Variation | Application Mode: | Flow Rate Control |
| Result: | Accept | Sample Name: | | Result: | |
| Average: | 100.010 g | Sample ID: | | Max Flow Rate: | 0.543 ml/min |
| Minimum: | 100.010 g | Batch ID: | | Min Flow Rate: | 0.526 ml/min |
| Maximum: | 100.011 g | Lot ID: | | Average Flow Rate: | 0.533 ml/min |
| Standard Deviation: | 0.000 g | Samples: | 5 | Pump Cycle: | 3 times |
| Test Result: | | Unit1: | 100.01[1] g | Test Result: | |
| Sample ID Weight T1/T2 | | Unit2: | 100.01[1] g | 1: | 0.529 ml/min |
| Sample 001 100.011 g >=-T1 & <=+T1 | | Unit3: | 100.01[1] g | 2: | 0.526 ml/min |
| Sample 002 100.010 g >=-T1 & <=+T1 | | Unit4: | 100.01[1] g | 3: | 0.543 ml/min |
| Sample 003 100.010 g >=-T1 & <=+T1 | | Unit5: | 100.01[1] g | Pump Speed: | 0.500 ml/min |
| Batch: | Batch 01 | Difference: | | Liquid Density: | 0.99823 g/ml |
| Number in Batch: | 100% Inspection | 1: 0.00[3] g 0.00 % | | Target Volume: | 10.00 ml |
| Inspection Quantity: | 3 | 2: 0.00[3] g 0.00 % | | Number of Pump Cycles: | 3 times |
| Tare Mode: | Off | 3: 0.00[3] g 0.00 % | | Mode: | Flow Rate Control |
| Pretare: | 0.000 g | 4: 0.00[3] g 0.00 % | | Signature: _____ | |
| Sample Type: | Solid | 5: 0.00[3] g 0.00 % | | Verified By: _____ | |
| EU Criteria: | Off | Result: | | | |
| Nominal Weight: | 100.000 g | Accept: | 5 Units | | |
| +T1: | 0.300 g | Fail: | 0 Units | | |
| -T1: | -0.300 g | Acceptance Criteria: | Off | | |
| +T2: | 0.600 g | Average Weight: | 100.00[8] g | | |
| -T2: | -0.600 g | +/- Limit %: | 7.50 % | | |
| Signature: _____ | | Signature: _____ | | | |
| Verified By: _____ | | Verified By: _____ | | | |

4.8.3 Batch Printing Output

| Printout | Export to PDF |
|-----------------------------|---------------|
| User ID: Admin | |
| First Name/Last Name: Admin | |
| Sample Name: | |
| Sample ID: | |
| Customized ID: | |
| -----Sample Data----- | |
| 1 25.00120 g 23:19:39 | |
| 2 24.99835 g 23:19:46 | |
| 3 25.00070 g 23:19:47 | |
| 4 25.00087 g 23:19:48 | |
| 5 25.00106 g 23:19:49 | |
| Signature: _____ | |
| Verified By: _____ | |

LibraryCreate, Active and Delete a Library

- Press "+New Library" button to create a new Library.
- When a new library is created, it remains inactive. The operator has to activate the library after ticking the box.
- Press on the RUBISH BIN button to delete a library.
- After the library item is activated, the main screen will show the activated library number.
- Example: Limits setting mode is Under and Over

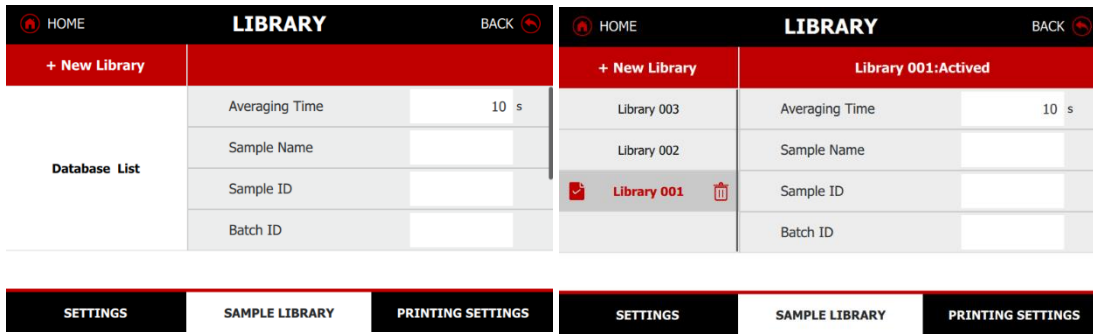

4.8.4 Printing Settings


The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to 6.1

4.9 Totalization

This function serves to is used to sum the weights of multiple samples and report the statistical data for the series of samples.

- In the lower portion of the home screen, select Totalization.
- Press **Tare** or **Zero** to start weighing.
- Place the series sample on the pan. Press the Accumulate button to add up the weights. The total weight of the samples will be displayed in the reference field.
- Setup the Start Mode before starting Check Weighing application.

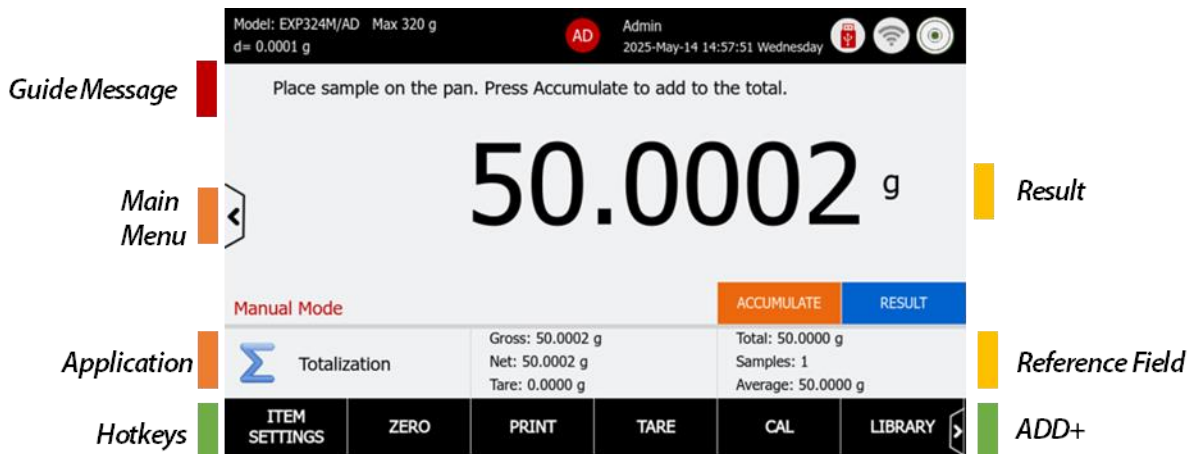

4.9.1 Application buttons


| Application Button/ Reference Field | Description |
|--|--|
| Accumulate | Press the “Accumulate” button to add up the sample weight. |
| Result | After the weight series samples, press “Result“ button to review the total and other statistic datas |

| | |
|----------------------|--|
| Item Settings | <ul style="list-style-type: none"> Start Mode: Automatic and Manual. Automatic start without press Accumulate button after second sample weighing Primary Weighing Unit: The default unit is grams. The operator can switch to alternative weighing units and two custom units. Auto Tare: Automatic tare the container value |
|----------------------|--|

4.9.2 Begin Totalization

Step 1: Setup the start mode in the Item Settings


- Start Mode: Automatic and Manual. Automatic start without press Accumulate button after second sample weighing e.

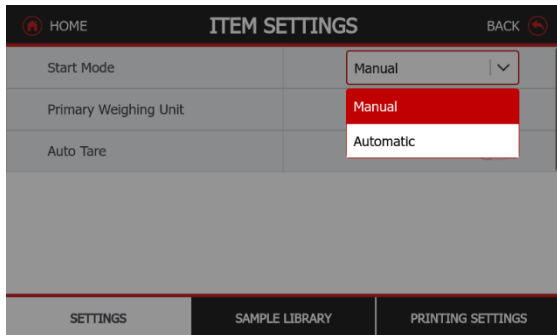

Step 2: Weigh the samples on the weighing pan

- Place a sample on the weighing pan,
- Press the Accumulate button to add up the weights.

Example: The home screen displays 50.0002 gram as the fourth weight value and press Result button to review the data

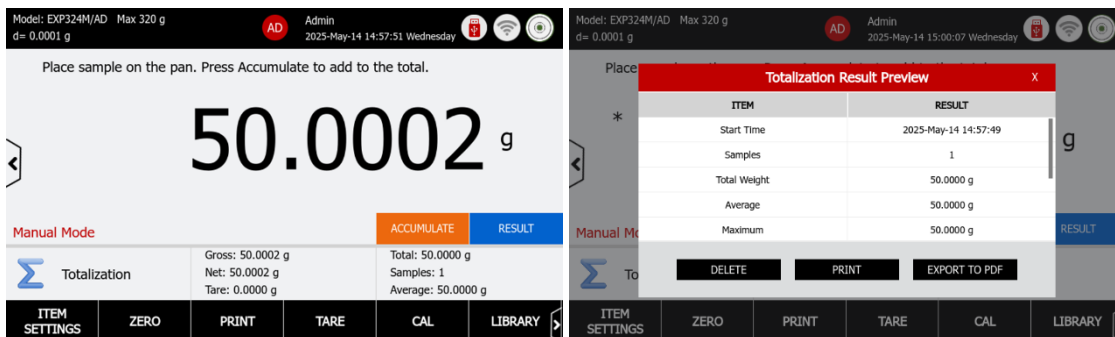


4.9.3 Accumulate the data for Totalization

Utilize the Accumulate function to aggregate samples and generate statistical reports.

The balance can store up to 999 accumulation records. Tap the Result to view the totalization and other statistical results.

The report content includes Start Time, Number of Samples, Total, Average, Maximum, Minimum, Range, Standard Deviation, Relative Deviation % and individual item weighing values.

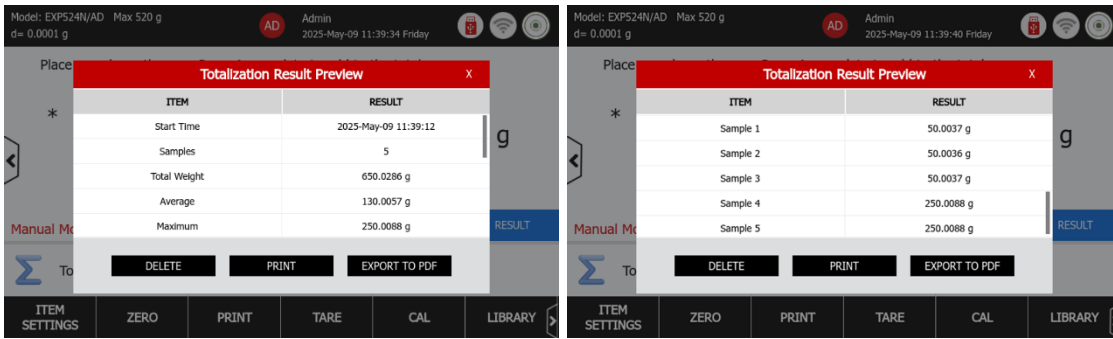


4.9.4 Sample Library – Totalization

The Explorer Plus balance features a built-in library for managing multiple sample profiles. Up to 3,000 library records can be saved and recalled for use. The total library capacity and records are accessible across various application modes and through the library menu. When the library usage exceeds 80%, a pop-up message will alert the user to export the library or delete all items. The entire library menu, please refer to 6.5.1

Create, Active and Delete a Library

- Press “+New Library” button to create a new Library.
- When a new library is created, it remains inactive. The operator has to activate the library after ticking the box.
- Press on the RUBISH BIN button to delete a library.
- After the library item is activated, the main screen will show the activated library number.
- Example: Limits setting mode is Under and Over

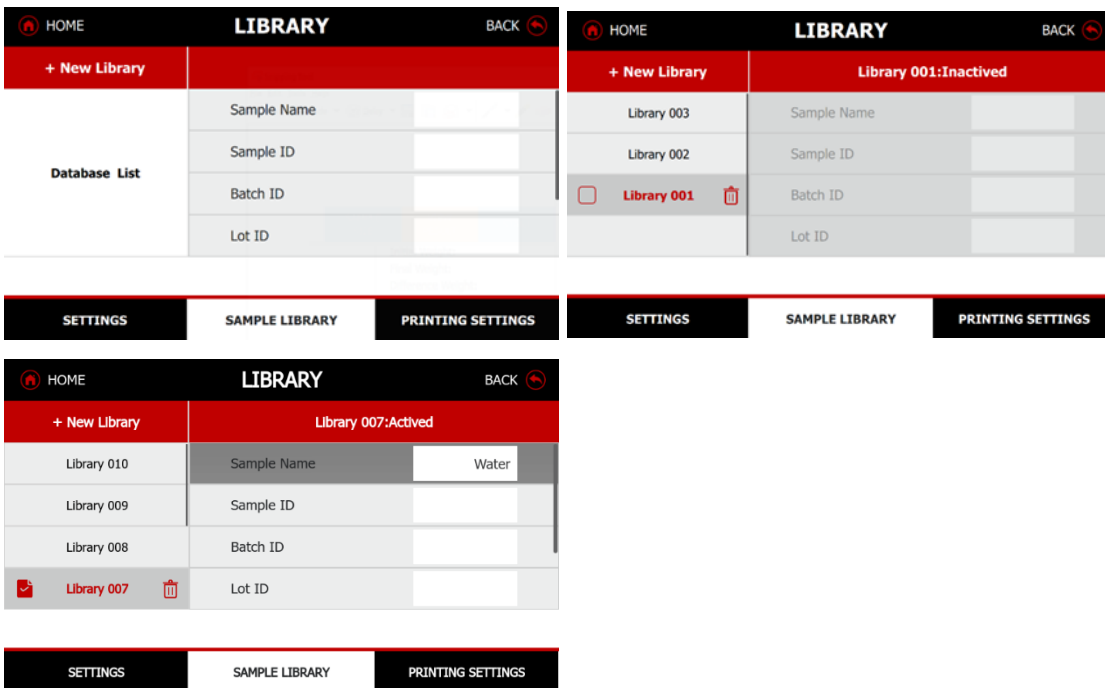

4.9.5 Printing Settings


The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to

4.10 Formulation

This function serves to is used to combine various elements in proportionate amounts. It has two formulation modes, free recipe and recipe-based formulation. Both modes allow filerto increase the volume of the formulation, enabling the dosage form to achieve the desired size and weight. This is particularly important for formulations containing small amounts of active pharmaceutical ingredients (APIs).

- In the lower portion of the home screen, select Formulation.
- Press **Tare** or **Zero** to start weighing.
- Please the ingredients on the pan to start the process. The total weight of the samples will be displayed in the reference field.
- Setup the Filler weight before starting Formulation application.

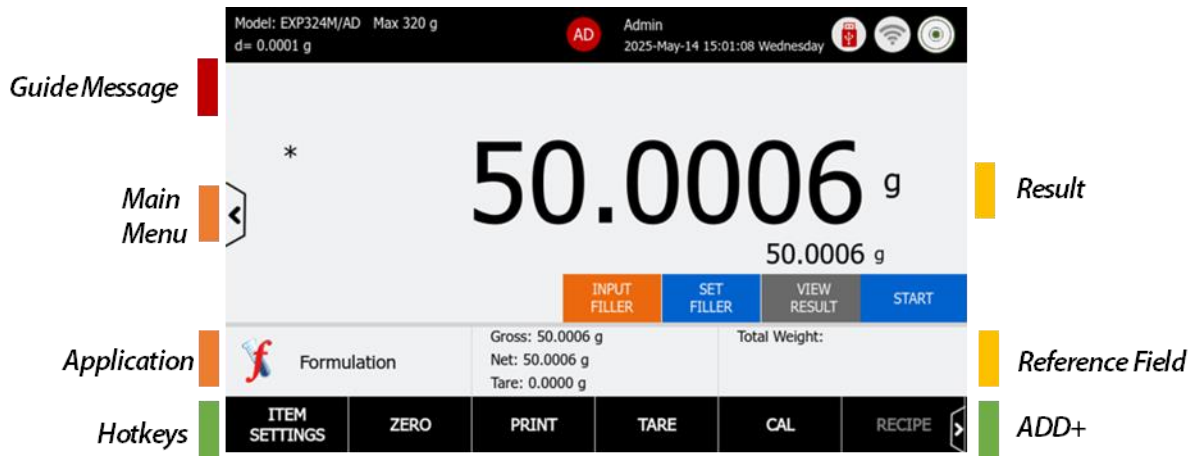


4.10.1 Application buttons

| Application Button/ Reference Field | Description |
|--|---|
| Input Filler | Input Filler value |
| Set Filler | Place Filler on the pan and accept the value |
| View Result | Press this button to view the formulation results once the process has finished. |
| Start | Press the button to start formulation. |
| Diff. (wt.): | Display the weight difference. |
| Diff. (%) | Display the difference in percentage. |
| Target | Display the target weight based on the recipe library. |
| Item Settings | <ul style="list-style-type: none"> • Formulation Mode: Recipe Based and Free Formulation |

| | |
|--|---|
| | <p>The system can store up to 99 ingredients in a recipe, and there are 25 recipes available for setup.</p> <ul style="list-style-type: none"> • Filler: Turn on and off • Primary Weighing Unit: The default unit is grams. The operator can switch to alternative weighing units and two custom units. • Secondary Weighing Unit: The operator can switch to alternative weighing units and two custom units. • Auto Tare: Automatic tare the container value |
|--|---|


4.10.2 Begin Free Formulation

Step 1: Setup the formulation mode in the Item Settings

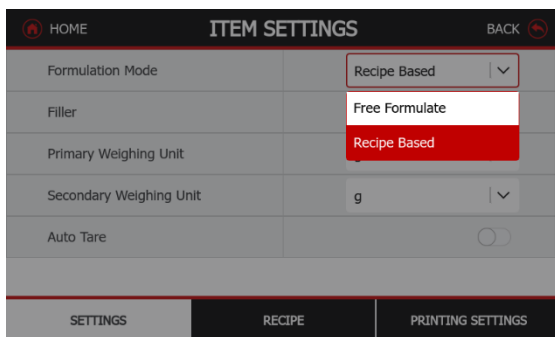

Step 2: Input Filler


- If the user has enabled the Filler option in the item settings, a filler value must be entered before creating a Free Recipe.
- The user can input the filler value by pressing the "Input Filler" button.
- Alternatively, the user can place the filler on the weighing pan and store the displayed value as the Filler weight.

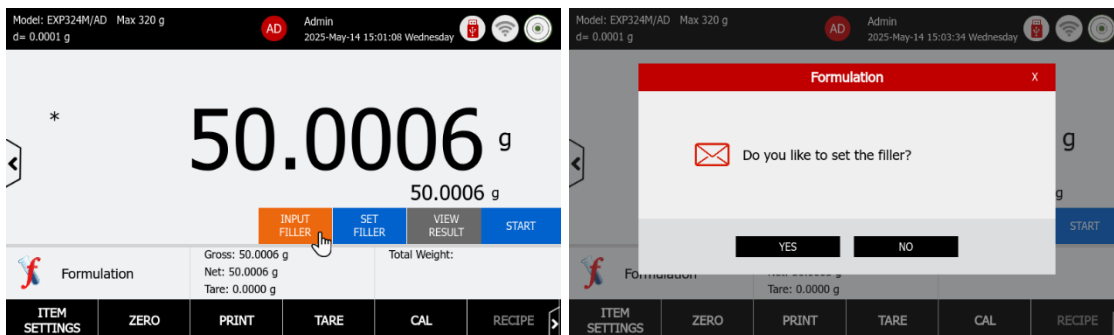

Weigh the samples on the weighing pan


- Place a sample on the weighing pan,
- Press the Accumulate button to add up the weights.
- The user can choose to Add this free formulation to Recipe, Print the Result or transfer the data to a PDF file using a USB flash drive.

Example: The home screen displays 199.9982 gram as the fourth weight value and press **View Result** button to review the data

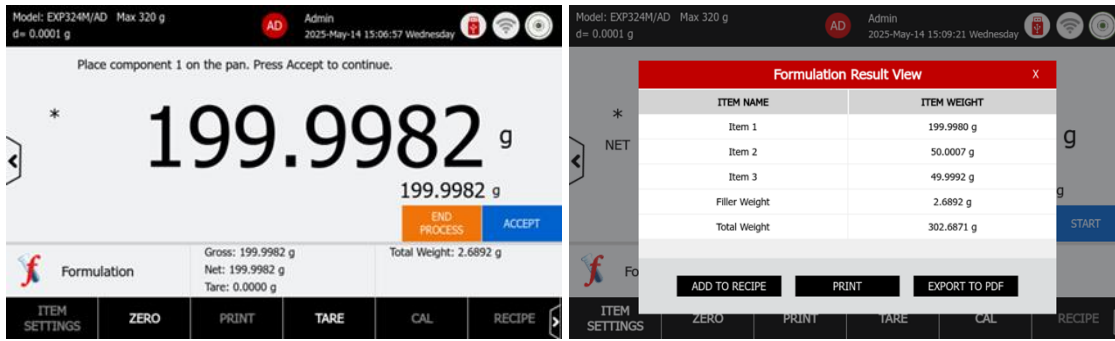


4.10.3 Begin Recipe Based Formulation

Step 1: Setup the formulation mode in the Item Settings

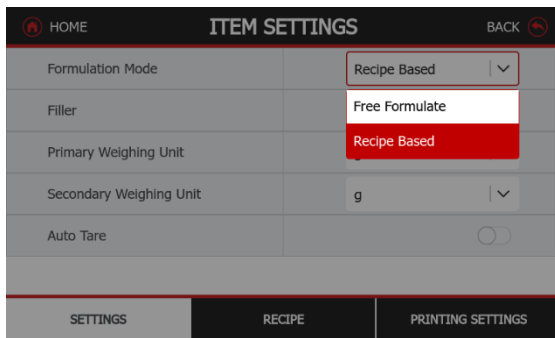

Step 2: Create Recipe ID and Name

- Press the **Create Recipe** button and create a New Recipe Name
- Click the tick box to activate Recipe 01
- Press the **Edit** button to go to next step

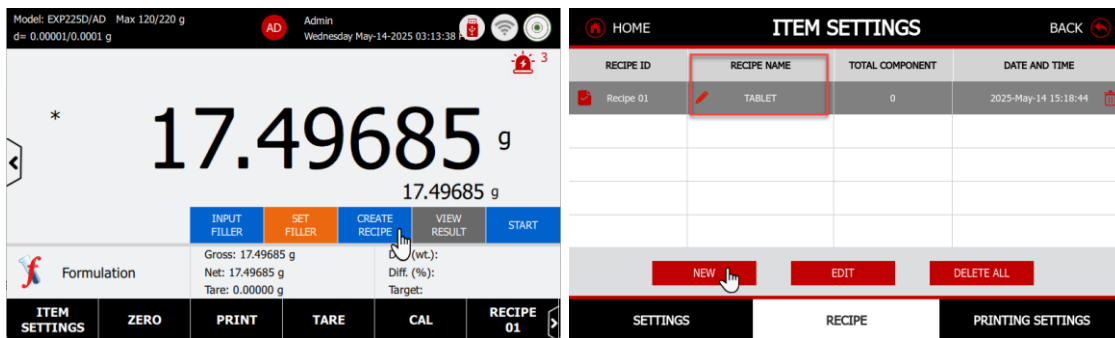

Step 2: Input Filler


- If the user has enabled the Filler option in the item settings, a filler value must be entered before creating a Free Recipe.
- The user can input the filler value by pressing the "Input Filler" button.

Alternatively, the user can place the filler on the weighing pan and store the displayed value as the Filler weight.

Step 3: Create Recipe Items


- Press **Add New** button to create ingredient names and their corresponding weight in gram.
- Use the **Up** and **Down** buttons to adjust the sequence order of the ingredients.

- Press the **Save** button before exiting the Recipe screen.
- Once the recipe has been saved, the user can press the **Start** button to begin the formulation process.

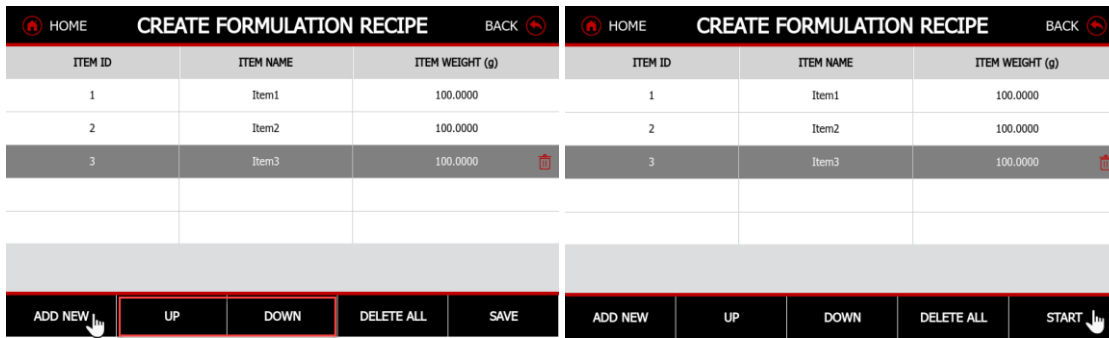

Step 4: Weigh the ingredients in the order by activated Recipe.

- Place the sample on the weighing pan and press the Accept button to confirm the weight.
- Follow the on-screen guide messages and repeat this process until all ingredients have been weighed.
- Once the process is completed, a result screen will appear with a button to review the data.
- The user can choose to Print the Result or transfer the data to a PDF file using a USB flash drive.

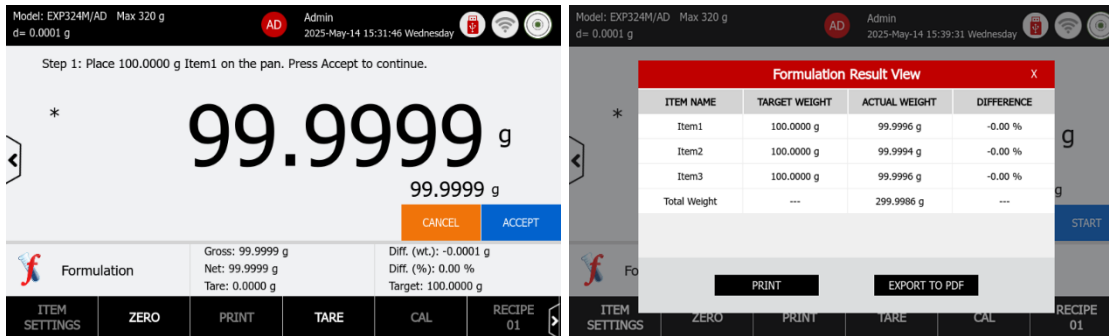

4.10.4 Printing Settings


The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to xxxx

4.11 Differential

This function serves to is used to calculate the difference in weights of multiple samples taken at different times.

- In the lower portion of the home screen, select Differential.
- Press **Tare** or **Zero** to start weighing.
- Setup the Sample Groups number before starting Differential Weighing. The initial weight, final weight and difference weight will be displayed in the reference field.

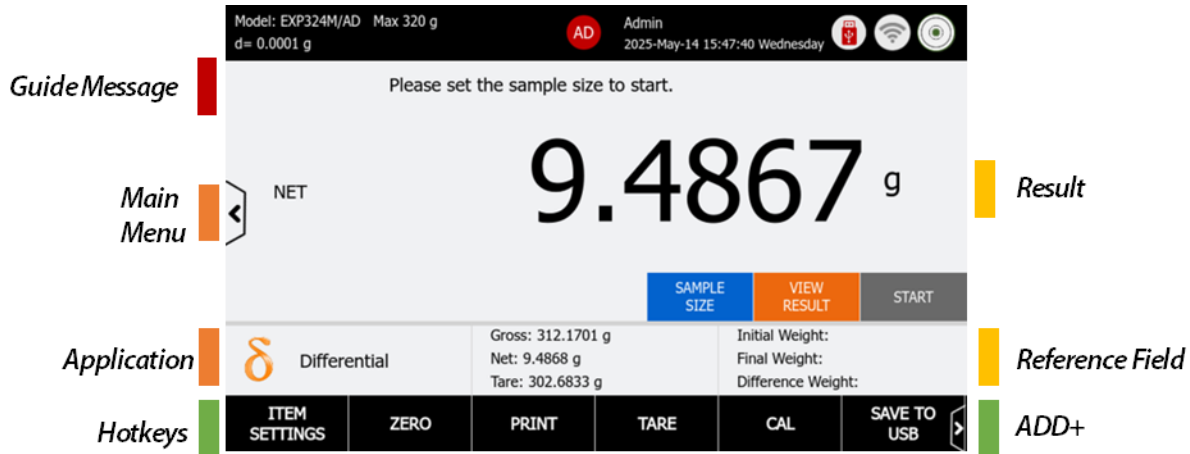


4.11.1 Application Buttons

| Application Button | Description |
|--------------------|--|
| Sample Groups | The number of sample groups requiring both initial and final weight measurements. |
| View Result | Press this button to view the Differential results once the process has finished. |
| Start | Press Start button to start the process |
| Accept | Accept the weight on the pan. |
| End Process | Press this button to complete the process |
| Initial Weight | Display the sample initial weight value. |
| Final Weight | Display the sample final weight value. |
| Difference Weight | Display the difference weight between initial and final measurement. |
| Item Settings | Sample Sequence: Off, A set of the initial and final weight Reciprocal Proportion: ON/OFF Absolute Value Auto Tare: |

Sample Sequence mode

- Off: After weight the series samples initial weight, and then weight the final weight. e.g. Weigh the initial weight of items 1, 2, 3, 4, 5, and then weigh the final weight of items 1, 2, 3, 4, 5.
- The set of sample measurement is design to place sample initial and final weight in a set
 - Example: Weigh the initial weight of items 1 and then final weight, and then weigh the initial weight of items 2 and then the final weight of items 2

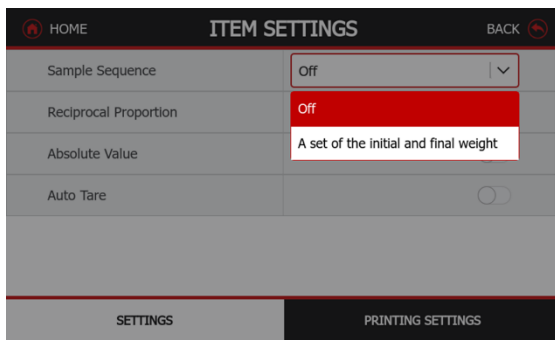

Reciprocal Mode

- Set on off:
 - $Difference\ Weight = Final\ Weight - Initial\ Weight$
- Set it on: The difference result=Final Weight -Initial Weight.
 - $Difference\ Weight = Initial\ Weight - Final\ Weight$

Absolute Value


The difference result will be positive no matter initial weight is heavier than final weight

- Set it on:
 - $Difference\ Weight = |Final\ Weight - Initial\ Weight|$

4.11.2 Begin Differential Weighing


Step 1: Setup Group Sample: Multiple samples can be tested in the differential weighing application.

- Input group sample number

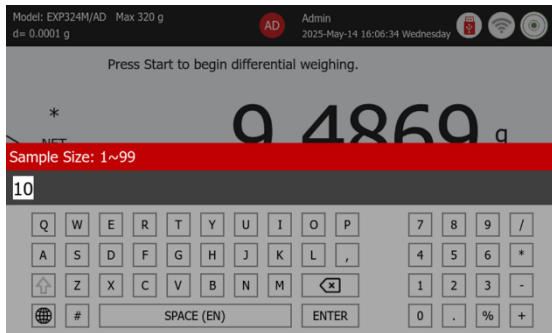

Step 2: Establish the initial weight of each sample

- Press START button and place the sample on the pan
- Press Accept to confirm the initial weight
- Follow the on-screen guide messages and repeat this process until all samples initial weight have been weighed.
- After that process, weigh the samples final weight in order, and the result screen will pop up at the end.

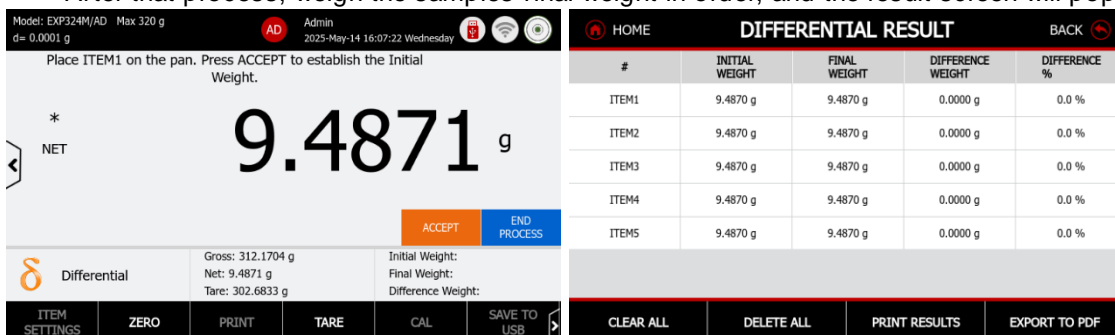

- The results include the initial weight, the final weight, the difference in weight, and the percentage difference.
- Users can clear the information, print the results to a PC or printer, and transfer the data to a PDF file via a USB flash drive.

4.11.3 Printing Settings


The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to xxx

4.12 Density Determination

This function serves to is used to determines the density of a solid or a liquid..

- In the lower portion of the home screen, select Density Determination.
- Press Tare or Zero to start weighing.
- User Help is available for review before setting up the process.

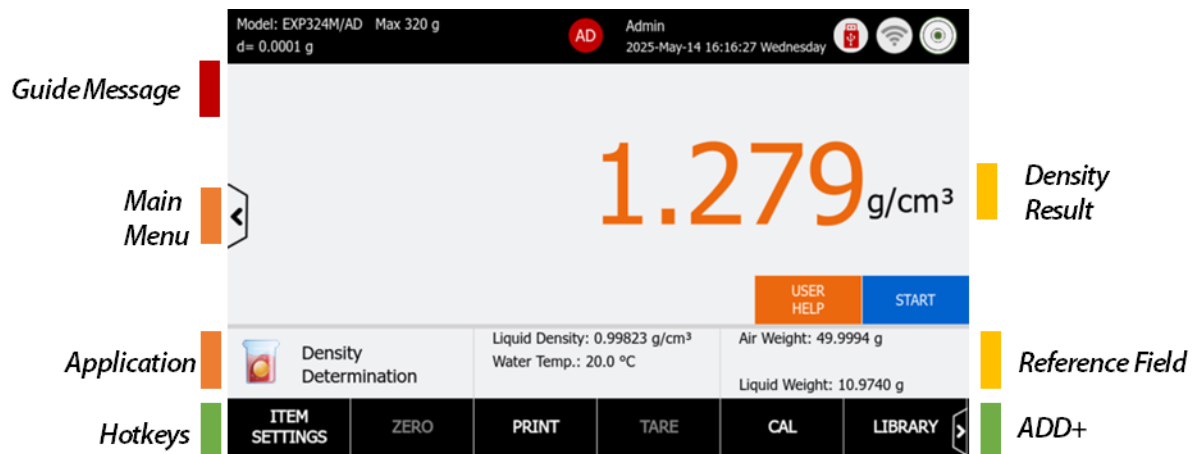


4.12.1 Application Buttons/ Reference Field

| Application Button | Description |
|--------------------|-------------|
| User Help | |
| Start | |
| Liquid Density | |
| Water Temp. | |
| Air Weight | |
| Liquid Weight | |

| | |
|---------------|--|
| Sinker | |
|---------------|--|

4.12.2 ITEM SETTINGS


Press the ITEM SETTINGS button to enter the application settings.

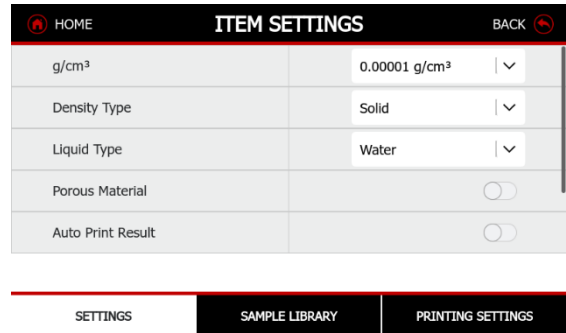

g/cm³


Select the displayed digits in g/cm³.

- 0.1 g/cm³
- 0.01 g/cm³
- 0.001 g/cm³
- 0.0001 g/cm³
- 0.00001 g/cm³

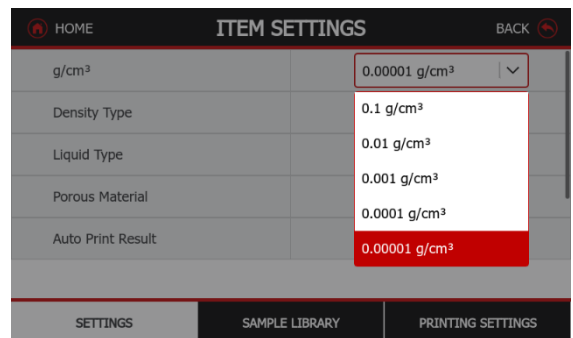


Density Type

Select the material type to be tested.

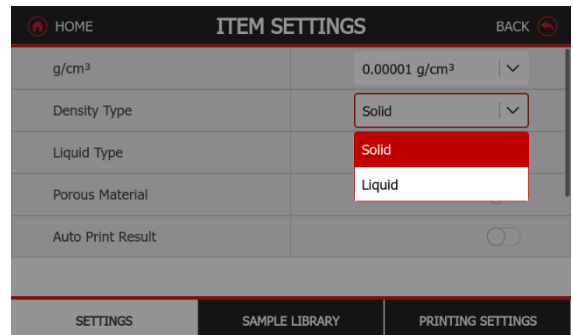


Liquid Type [Conditional]

This field will be visible when Density Type is selected to Solid.
 Liquid type can be selected to Water or Other.
 When it is selected to Other, users can set up the density of this liquid in Library.

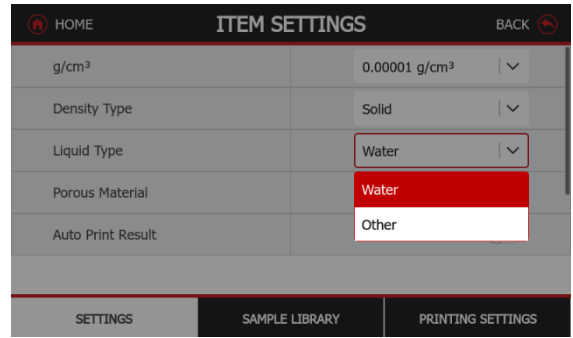


Porous Material [Conditional]

This field will be visible when Density Type is selected to Solid.
 User can select if the solid is porous material.
 When it is selected to Other, users can set up the oil density in Library.

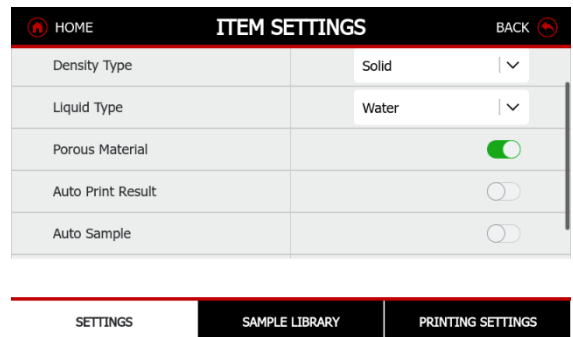

Auto Sample

Auto Sample ON: Density can be determined with a button press after each step.
 Auto Sample Off: Weights are automatically recorded in each step.

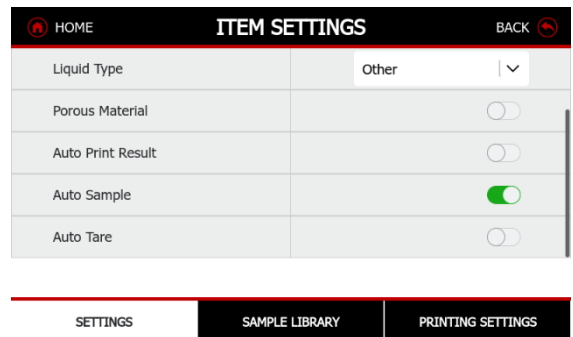

4.12.3 Reference Fields

The reference fields are different depending on Item Settings.

| Density Type | Liquid Type | Porous Material | Reference Values | Weighing Values |
|--------------|-------------|-----------------|---|---|
| Solid | Water | Off | Liquid Density Water Temp | Air Weight Liquid Weight |
| Solid | Water | On | Liquid Density Water Temp Oil Density | Dry Weight Oiled Weight Liquid Weight |
| Solid | Other | Off | Liquid Density | Air Weight Liquid Weight |

| | | | | |
|--------|-------|----|-------------------------------|---|
| Solid | Other | On | Liquid Density Oil Density | Dry Weight Oiled Weight Liquid Weight |
| Liquid | \ | \ | Sinker Volume | Air Weight Liquid Weight |

4.12.4 Measuring the Density of a Solid Using Water

Prerequisites

The density kit is installed. For details, please refer to [Density Determination Kitt Instruction Manual](#).


Step 1: Item Settings

| Density Type | Liquid Type | Porous Material |
|--------------|-------------|-----------------|
| Solid | Water | Off |

Step 2: Edit in Library


Edit the water temperature value in the Library.

Note: For more details about Library operations, please refer to [4.12.10 Sample Library – Density Determination](#).

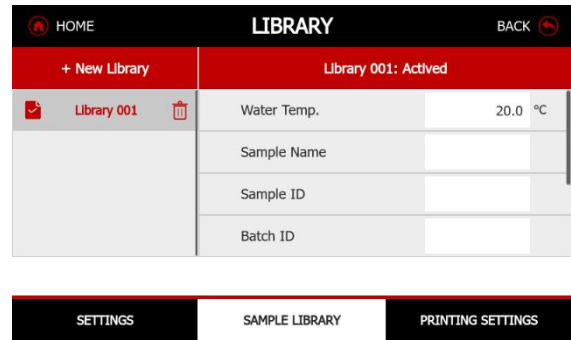


Step 3: Weigh the Sample

1. Weigh sample in air.

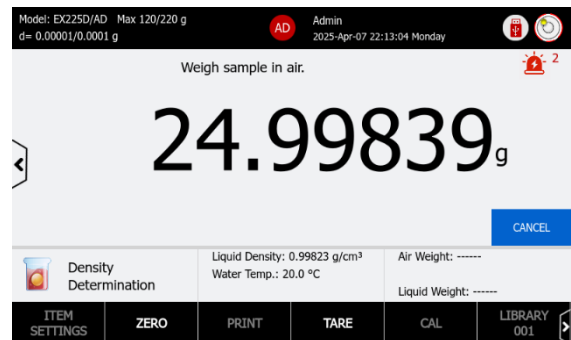


2. When finished, clear the pan.

3. Weigh sample in liquid.

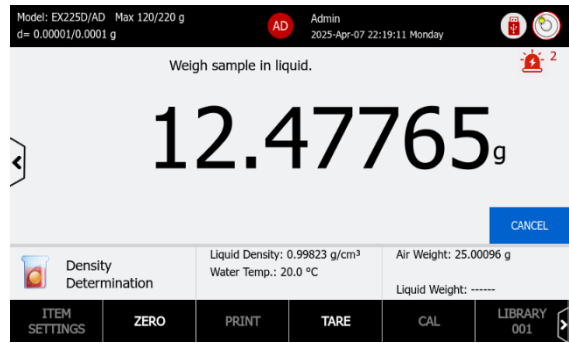

4. The result will be displayed in orange font color.

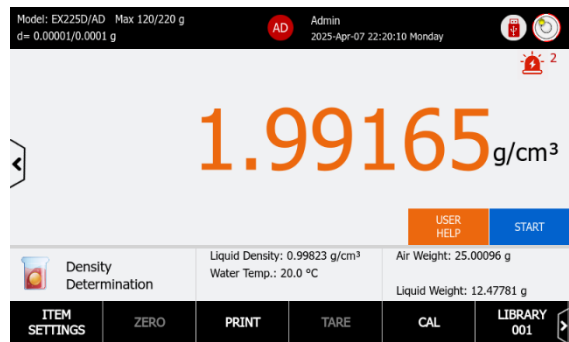

4.12.5 Measuring the Density of a Buoyant Material Using Water

Prerequisites

The density kit is installed. For details, please refer to [Density Determination Kitt Instruction Manual](#).


Step 1: Item Settings

| Density Type | Liquid Type | Porous Material |
|--------------|-------------|-----------------|
| Solid | Water | Off |

Step 2: Edit in Library


Edit the water temperature value in the Library.

Note: For more details about Library operations, please refer to [4.12.10 Sample Library – Density Determination](#).

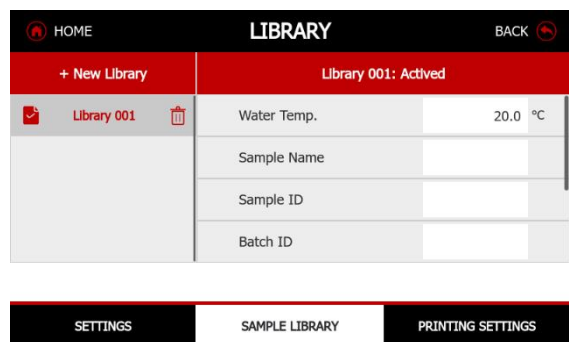


Step 3: Weigh the Sample


1. Weigh sample in air.

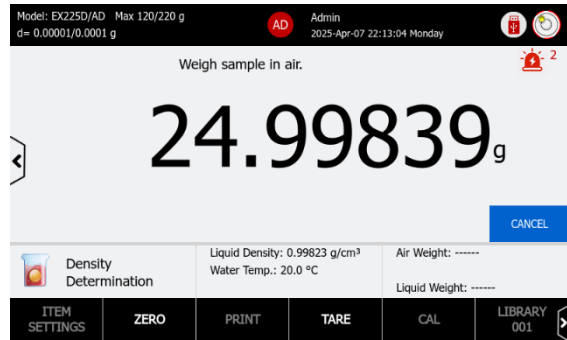

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2. When finished, clear the pan.
3. Weigh sample in liquid. Push the sample down into the liquid until it is fully submerged.



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4. The result will be displayed in orange font color.



4.12.6 Measuring the Density of a Solid Using an Auxiliary Liquid

Prerequisites

The density kit is installed. For details, please refer to [Density Determination Kitt Instruction Manual](#).


Step 1: Item Settings

| Density Type | Liquid Type | Porous Material |
|--------------|-------------|-----------------|
| Solid | Other | Off |

Step 2: Edit the Library


Edit the liquid density value in the Library.

Note: For more details about Library operations, please refer to [4.12.10 Sample Library – Density Determination](#).

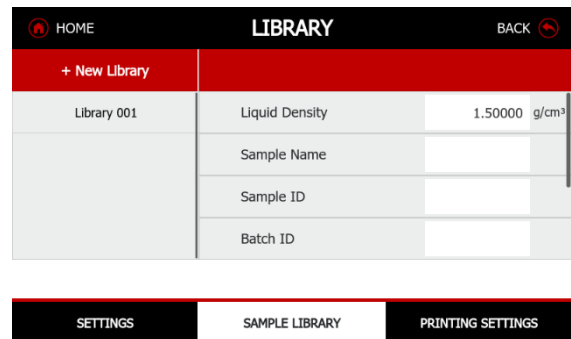


Step 3: Weigh the Sample

1. Weigh sample in air.

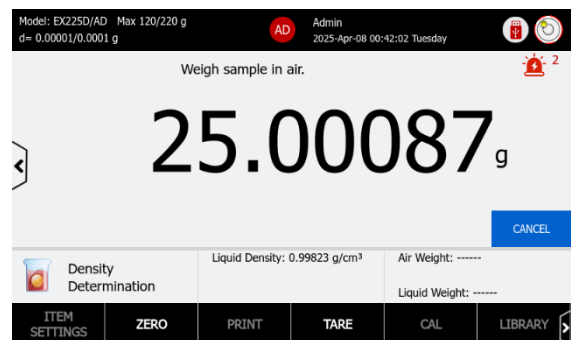


2. When finished, clear the pan.

3. Weigh sample in liquid.

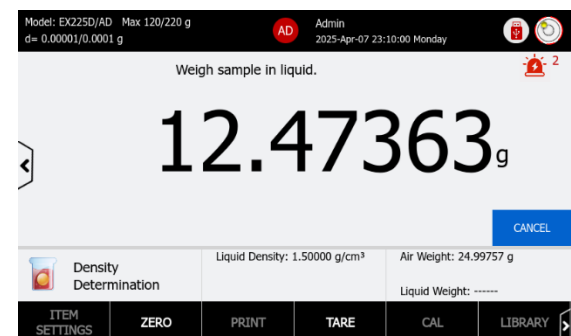

4. The result will be displayed in orange font color.

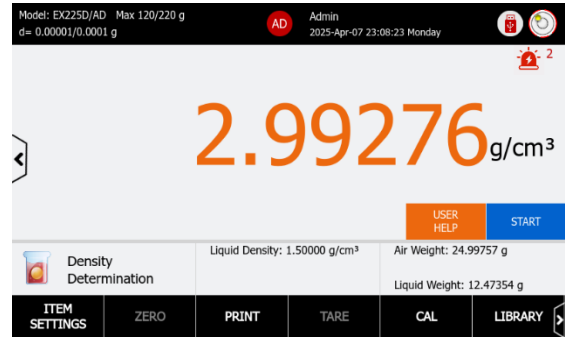

4.12.7 Measuring the Density of Porous Material Using Oil

Prerequisites

The density kit is installed. For details, please refer to [Density Determination Kitt Instruction Manual](#).


Step 1: Item Settings

| Density Type | Liquid Type | Porous Material |
|--------------|-------------|-----------------|
| Solid | Water | On |

Step 2: Edit in Library


Edit the water temperature and oil density values in the Library.

Note: For more details about Library operations, please refer to [4.12.10 Sample Library – Density Determination](#).

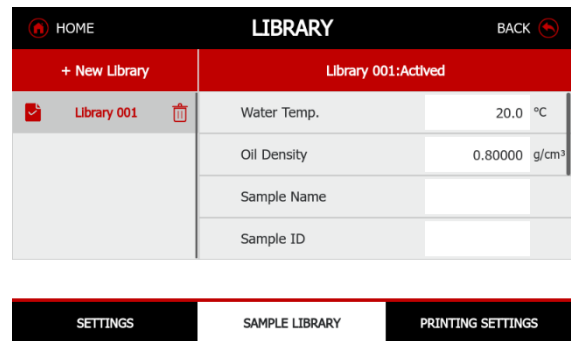


Step 3: Weigh the Sample

1. Weigh the un-oiled sample in air.

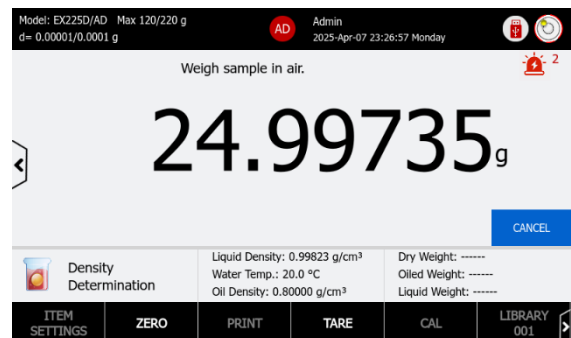


2. Weigh the oiled sample in air.

3. Weigh oiled sample in liquid.

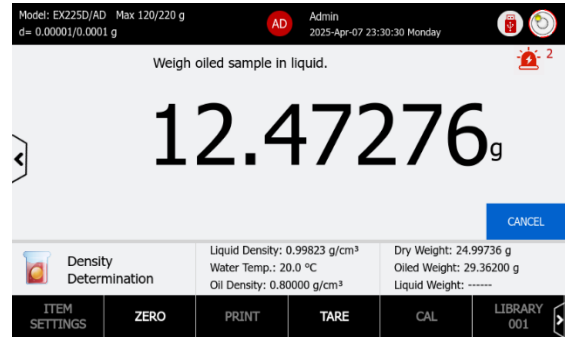

4. The result will be displayed in orange font color.

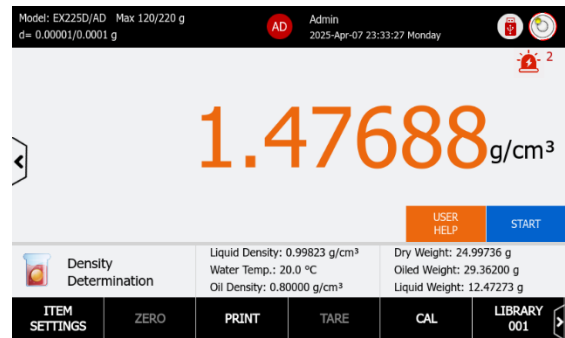

4.12.8 Measuring the Density of Porous Material Using Oil and Auxiliary Liquid

Prerequisites

The density kit is installed. For details, please refer to [Density Determination Kitt Instruction Manual](#).


Step 1: Item Settings

| Density Type | Liquid Type | Porous Material |
|--------------|-------------|-----------------|
| Solid | Other | On |

Step 2: Edit in Library


Edit the liquid density and oil density value in the Library.

Note: For more details about Library operations, please refer to [4.12.10 Sample Library – Density Determination](#).

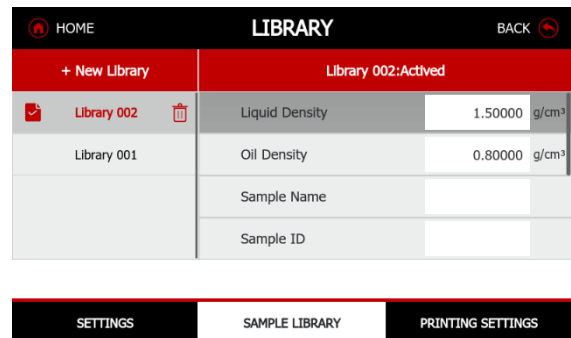


Step 3: Weigh the Sample


1. Weigh the un-oiled sample in air.

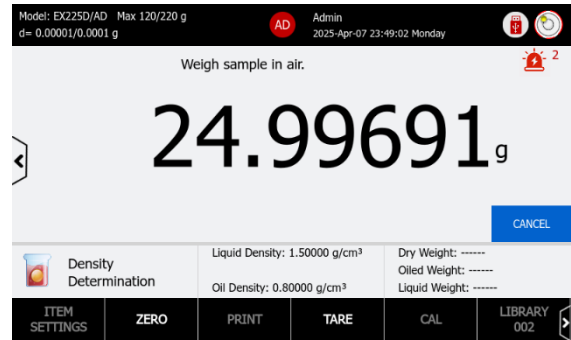


2. Weigh the oiled sample in air.

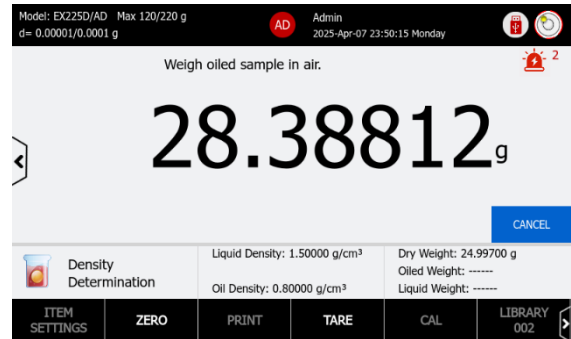

3. Weigh oiled sample in liquid.

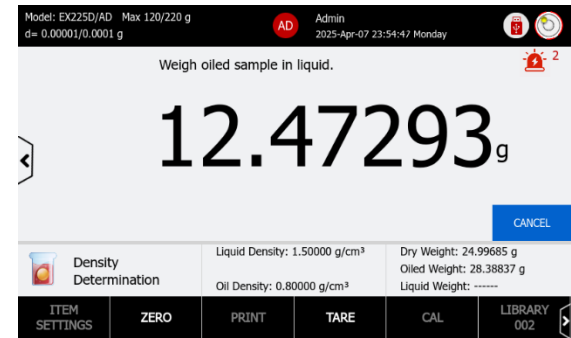

4. The result will be displayed in orange font color.

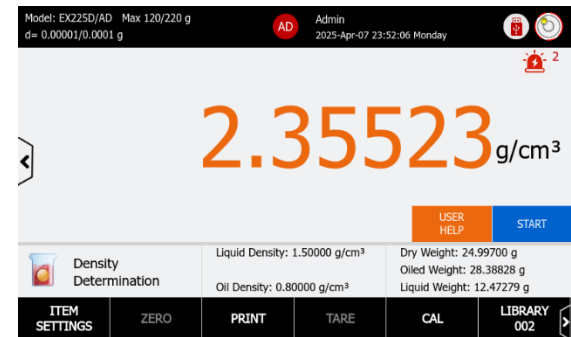

4.12.9 Measuring the Density of a liquid using a Calibrated Sinkers

Prerequisites

The density kit is installed. For details, please refer to [Density Determination Kitt Instruction Manual](#).


Step 1: Item Settings

| Density Type | Liquid Type | Porous Material |
|--------------|-------------|-----------------|
| Liquid | \ | \ |

Step 2: Edit in Library


Edit the sinker volume in the Library.

Note: For more details about Library operations, please refer to [4.12.10 Sample Library – Density Determination](#).

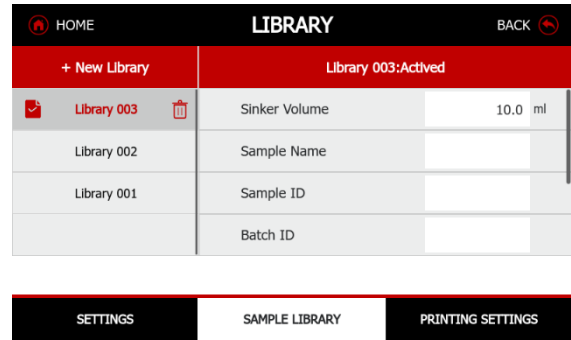


Step 3: Weigh the Sample


1. Weigh the sinker in air.

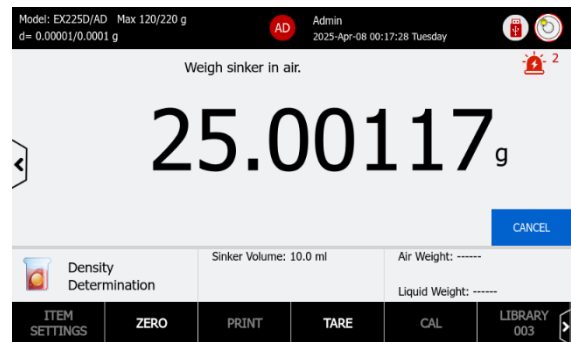

2. Weigh sinker in liquid.

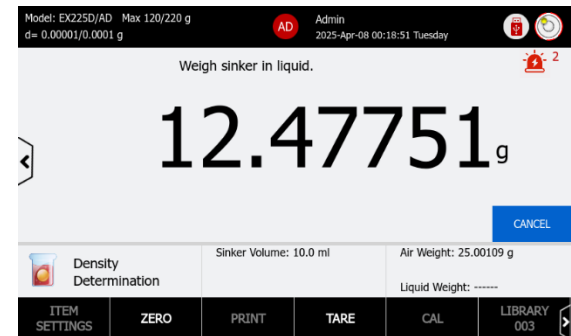

3. The result will be displayed in orange font color.

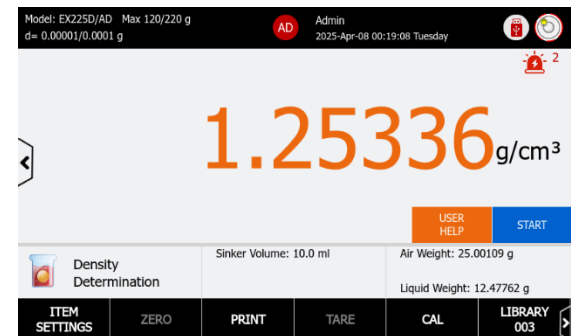

4.12.10 Sample Library – Density Determination



Explorer Plus balance built in library simplifies working with multiple samples, allowing application settings to be saved and recalled for reuse.

The Sample Library can store up to 3000 records, without limitation on individual applications. To check the status of library usage, navigate to Main Menu > Library.

Note: In Main Menu > Library, the balance would notify users to delete or export libraries when the storage is used 80%.


Create a Library Record

1. Press the +New Library button. When a new library is created, it is inactivated and not editable.

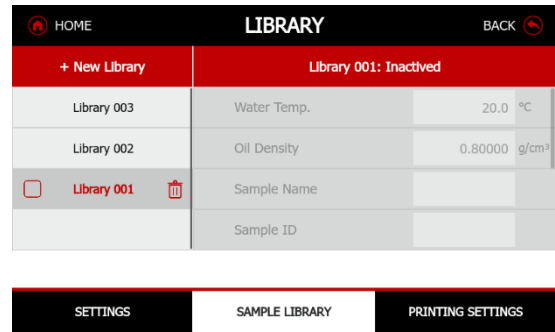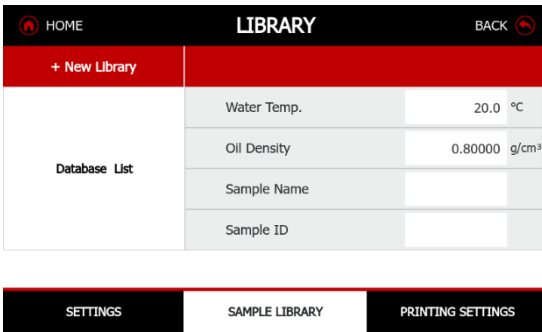


2. Activate the new library by check the box in front of the library id.

The library status will be Activated and editable.


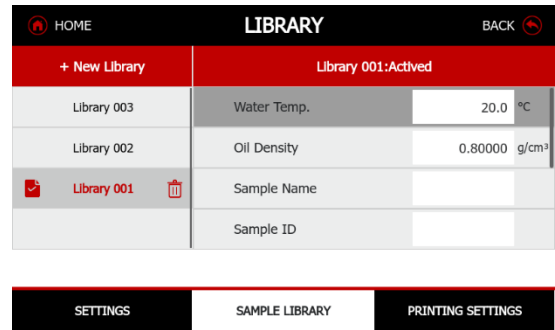


3. Edit the data fields using the keyboard. Press **ENTER** to confirm editing.

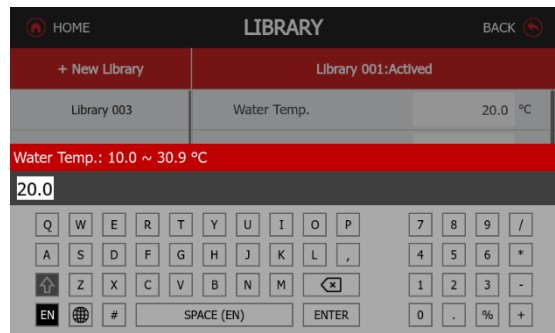

Activate a Library


To activate a library item, check the box in front of the library item.

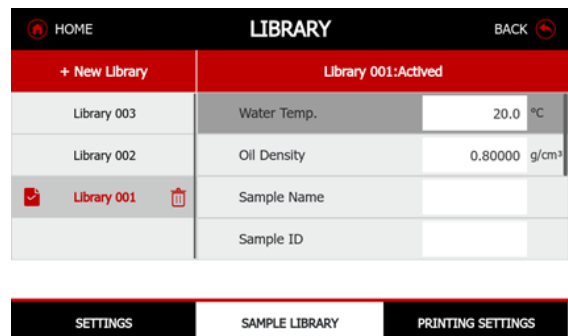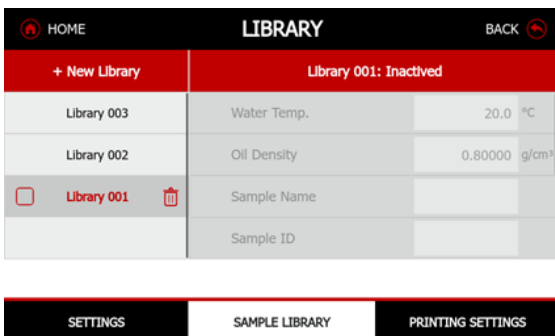

After the library item is loaded,

- On the screen, it would notify that the library is activated.

- On the application main screen, the library button will become the activated library id.

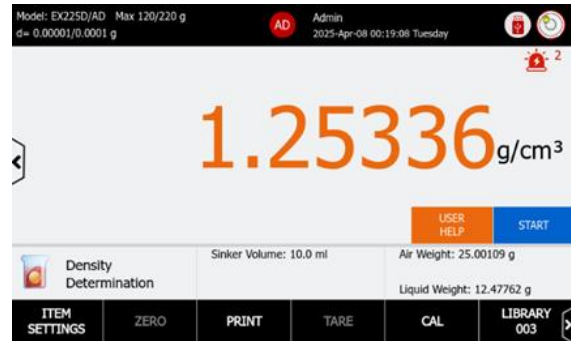

Delete a Library of Current Application

Press on the RUBISH BIN button behind the library ID to delete a library.

Manage Libraries

The balance allows users to review, delete, export or import libraries in Main Menu > Library.

For details, please refer to [6.5.1](#)

4.12.11 Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.


For details of Printing Setting, please refer to [Error! Reference source not found. Error! Reference source not found.](#)

4.13 Peak Hold

Peak Hold captures the maximum weight in a series of weighings.



Application Icon

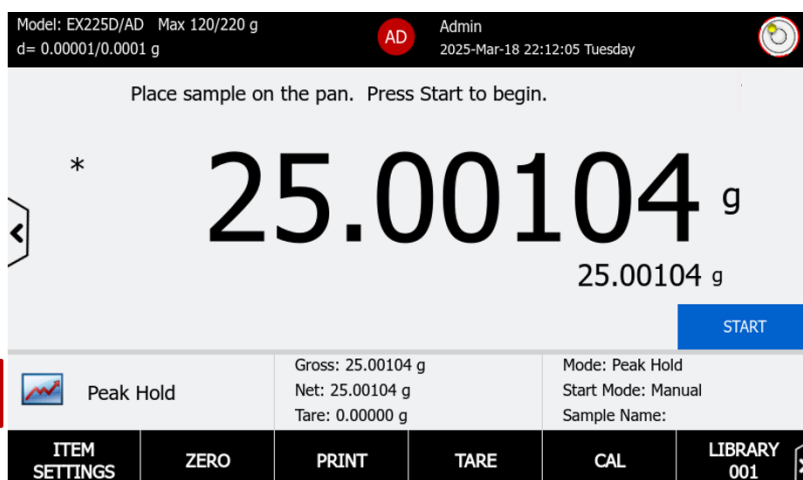

Application Guide Message

Main Display Line

Secondary Unit

Application Buttons


Reference Fields

Hotkeys

Application Name


4.13.1 ITEM SETTINGS

Press the ITEM SETTINGS button to enter the application settings.

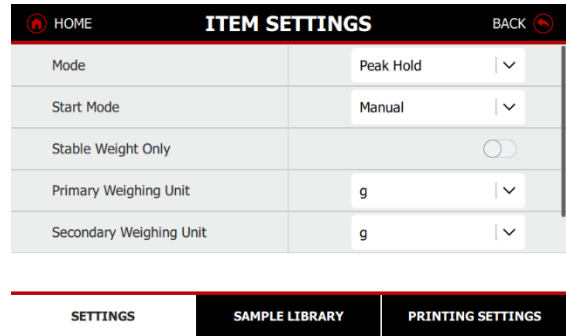

Mode

Select Start Mode as Peak Hold or Display Hold.

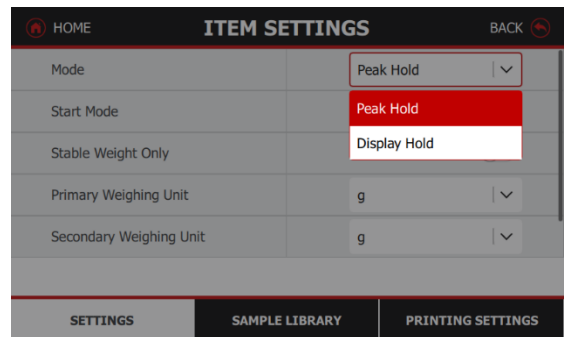


Peak Hold

The Peak Hold mode allows the user to capture and store the highest stable or unstable weight in a series of weight measurements.

Start Mode

Three different Start Modes can be selected in Peak Hold:

- **Manual** (start and stop the capture via button press).
- **Semi-Automatic** (start automatically/reset manually).
- **Automatic** (start and stop the capture automatically).

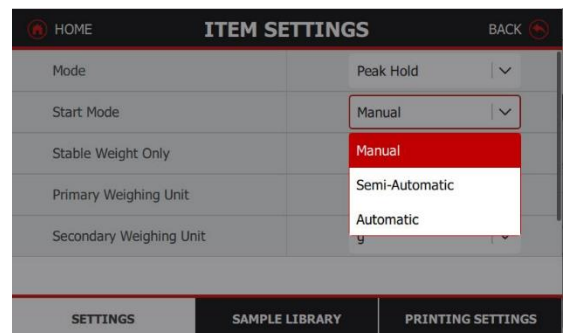

Manual

Application Guide Message

Main Display Line

Secondary Unit

Application Buttons

Application Name

Reference Fields

Hotkeys

Semi-Automatic

Application Guide Message

Main Display Line

Secondary Unit

Application Buttons

Application Name

Reference Fields

Hotkeys

Automatic

Application Guide Message

Main Display Line

Secondary Unit


Application Name

Reference Fields

Hotkeys

Display Hold

The Display Hold mode allows the user to capture and store a stable or unstable weight in a series of weight measurements by pressing the HOLD button.

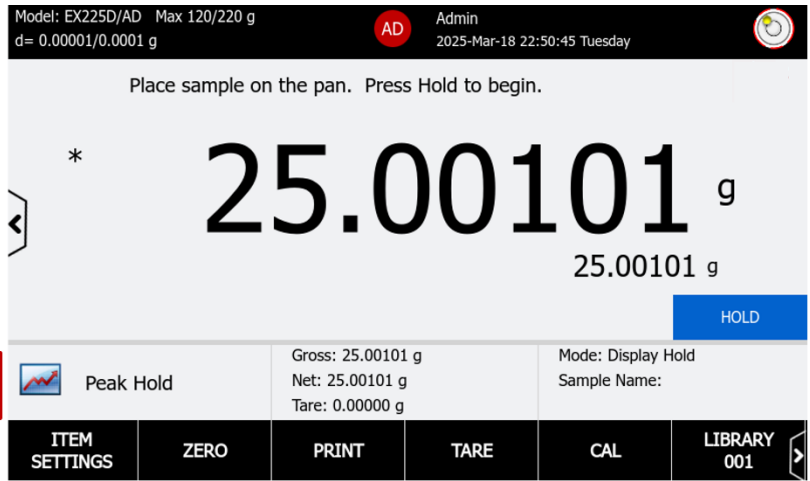

Application Guide Message
 Main Display Line
 Secondary Unit
 Application Buttons
 Reference Fields
 Hotkeys

Application Name


4.13.2 Reference Fields

| Reference Fields | Description |
|------------------|---|
| Mode | Display the selected mode. |
| Start Mode | The Start Mode field is visible when Mode is selected to Peak Hold. |


4.13.3 Begin with Peak Hold

Weigh Samples in Manual Mode

1. Place samples onto the weighing pan.

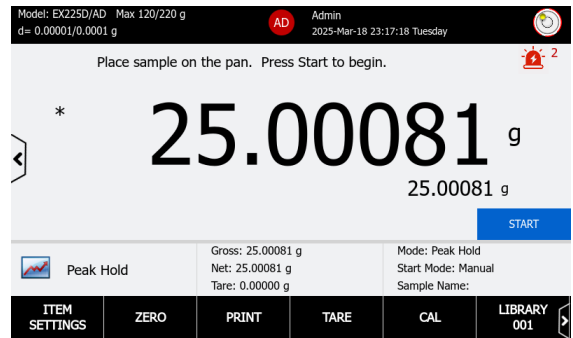


1. Press START to begin the capture process.
 The highest stable weight reading is held indefinitely.

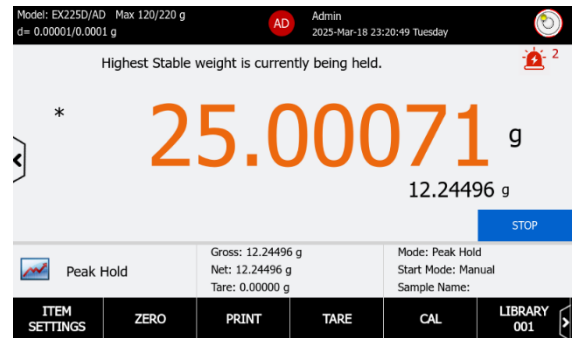

2. Press STOP to end and reset the process.

Weigh Samples in Semi Automatic Mode


1. Place samples onto the weighing pan. The balance will start the capture process automatically.
The highest stable weight reading is held indefinitely.

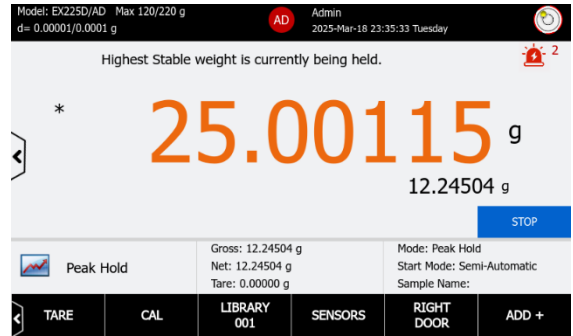

2. Press STOP to end and reset the process.

Weigh Samples in Automatic Mode



1. Place samples onto the weighing pan. The balance will start the capture process automatically.
The highest stable weight reading is held indefinitely.

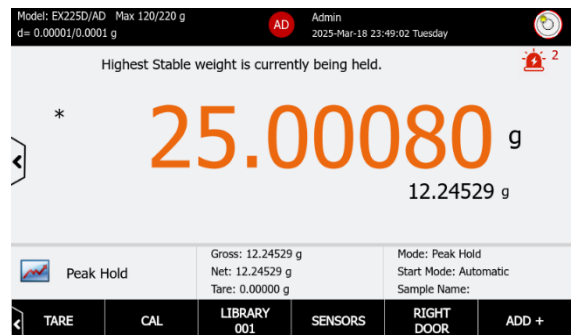

2. Remove all items from the pan to end and reset the process.

4.13.4 Begin with Display Hold

1. Place samples onto the weighing pan.
2. Press HOLD to capture and hold the current weight reading.

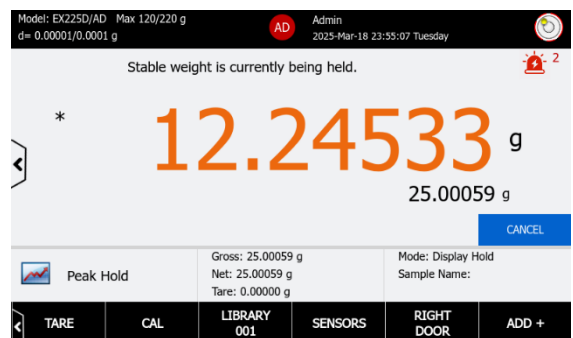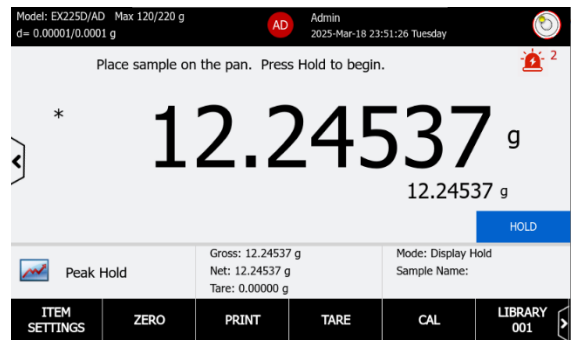

3. Press CANCEL to end and reset the process.

4.13.5 Sample Library – Totalization



Explorer Plus balance built in library simplifies working with multiple samples, allowing application settings to be saved and recalled for reuse.

The Sample Library can store up to 3000 records, without limitation on individual applications. To check the status of library usage, navigate to Main Menu > Library.

Note: In Main Menu > Library, the balance would notify users to delete or export libraries when the storage is used 80%.


Create a Library Record

1. Press the +New Library button. When a new library is created, it is inactivated and not editable.

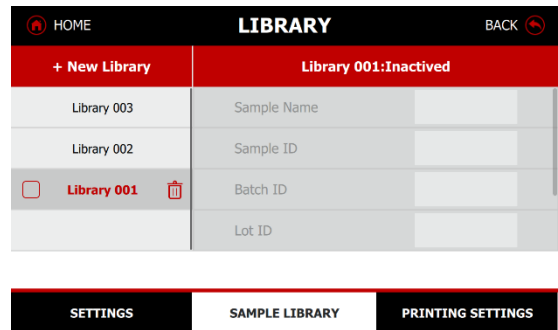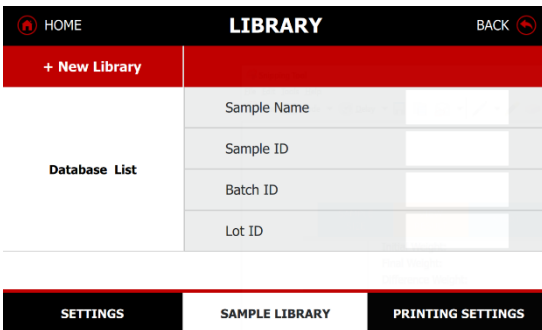


2. Activate the new library by check the box in front of the library id.

The library status will be Activated and editable.


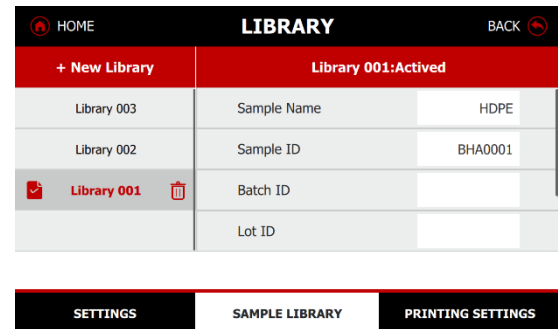


3. Edit the data fields using the keyboard. Press **ENTER** to confirm editing.

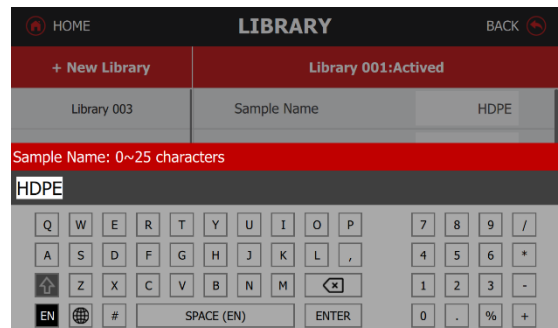


Activate a Library

To activate a library item, check the box in front of the library item.

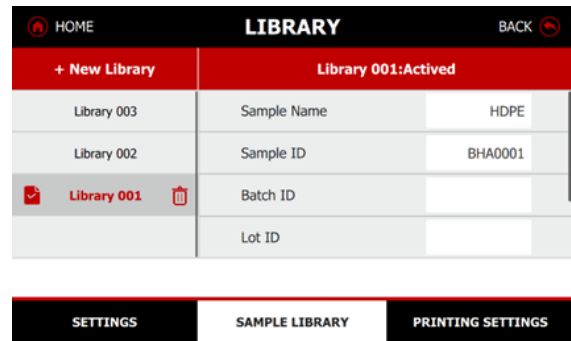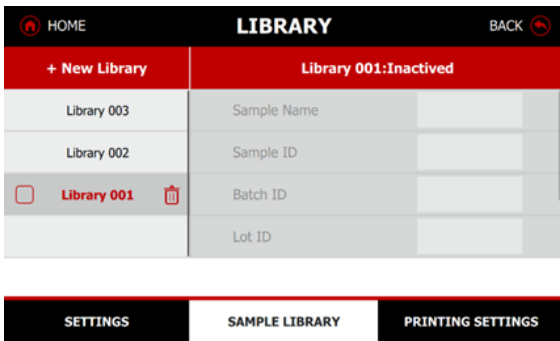

After the library item is loaded,

- On the screen, it would notify that the library is activated.
- On the application main screen, the library button will become the activated library id.

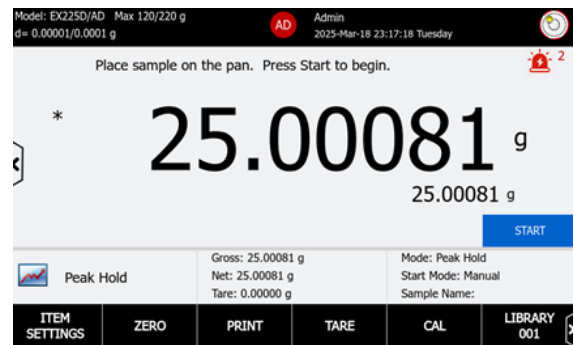

Delete a Library of Current Application

Press on the RUBISH BIN button behind the library ID to delete a library.

Manage Libraries

The balance allows users to review, delete, export or import libraries in Main Menu > Library.


For details, please refer to [6.5.1](#) Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to

4.14 Pipette Adjustment

Pipette Adjustment calculates the inaccuracy and imprecision of pipettes. The application is used to check if a pipette's dosage is within tolerances.

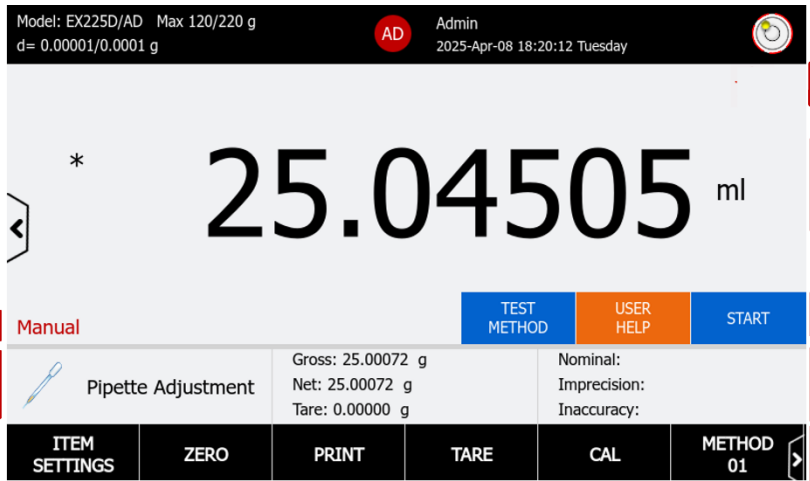

Application Guide Message

Main Display Line

Process Mode **Manual**

Application Buttons


Application Name

Reference Fields

Hotkeys


4.14.1 ITEM SETTINGS

Press the ITEM SETTINGS button to enter the application settings.

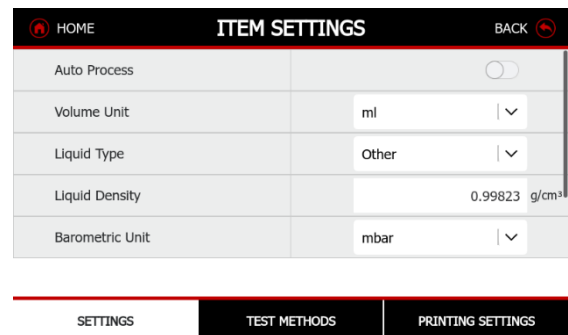


Auto Process

When Auto Process is enabled, the weight value will be recorded automatically when the reading is stable.

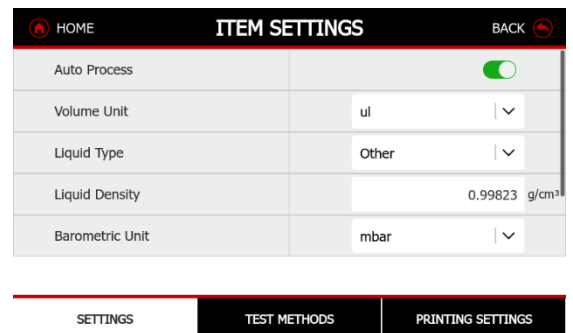


Volume Unit

The volume unit can be selected as ul or ml.

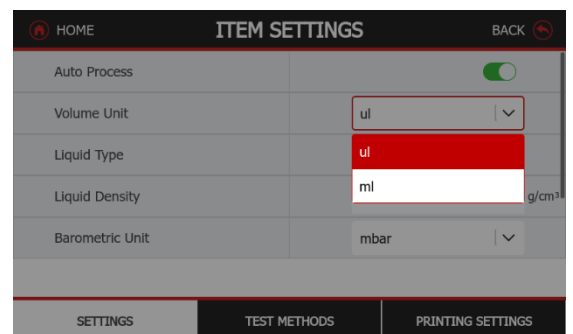


Liquid Type

Select the liquid used during measurements. The liquid can be selected as Water or Other.

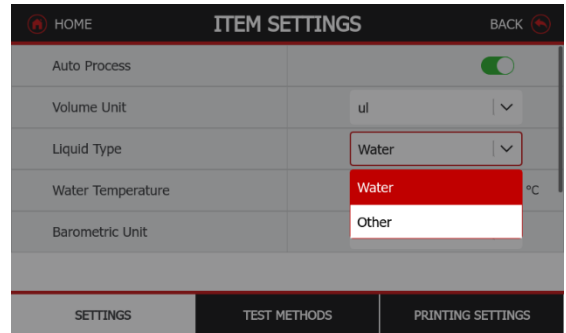


Water Temperature [Conditional]

When the liquid type is selected as Water, users can input the water temperature.

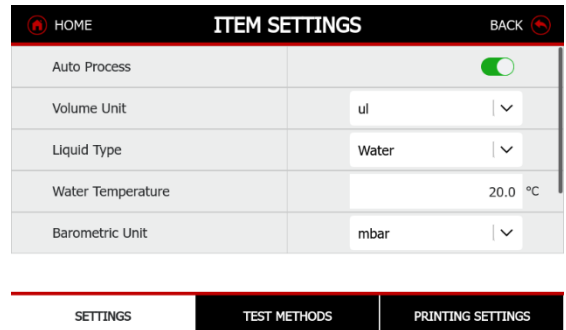


Liquid Density [Conditional]

When the liquid type is selected as Other, users can input the liquid density.

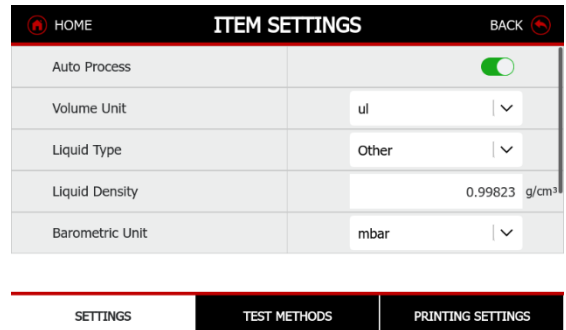

Barometric Unit

To select Atmospheric Pressure unit as ATM, inHg, KPa, mbar, mmHg, or PSIA

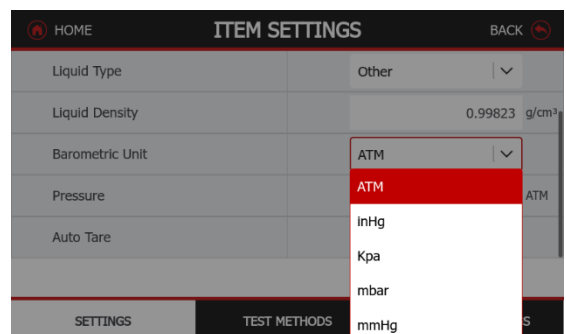

Pressure

User can set up the air pressure value in the selected barometric unit.

The screenshot shows the 'ITEM SETTINGS' screen with the following configuration:

| Setting | Value |
|-----------------|---------------------------|
| Liquid Type | Other |
| Liquid Density | 0.99823 g/cm ³ |
| Barometric Unit | ATM |
| Pressure | 1.0 ATM |
| Auto Tare | <input type="checkbox"/> |

Navigation buttons at the bottom: SETTINGS, TEST METHODS, PRINTING SETTINGS.

4.14.2 Application Buttons

| Application Button | Icon | Description |
|-------------------------|------|---|
| TEST METHOD | | Set up Test Methods. A Test Method must be selected to start the test. |
| USER HELP | | View user help instructions. |
| START | | Press the button to start the test. Note: If no method is selected, the START button will be greyed and inactivated. |
| ACCEPT (Manual Only) | | When a test is started, press the button to accept and record the weight value. |
| VIEW RESULT | | After a test is completed, press the button to view the test result. |
| CANCEL | | In a test, press the button to cancel the test. |

4.14.3 Reference Fields

| Reference Fields | Description |
|------------------|---|
| Nominal | Display the nominal volume of the test. |
| Imprecision | Display the imprecision rate. |
| Inaccuracy | Display the inaccuracy rate. |

4.14.4 Test Method


The balance allows users to set up maximum 25 methods for pipette adjustment. Each method can include 5 tests and include the following parameters:

- Nominal Volume
- Pipette Name

- Pipette Number
- Inaccuracy (%)
- Imprecision (%)
- Number of Samples in a method


Create a Method

1. Press the TEST METHOD button or the METHOD hotkey to enter the Method screen.

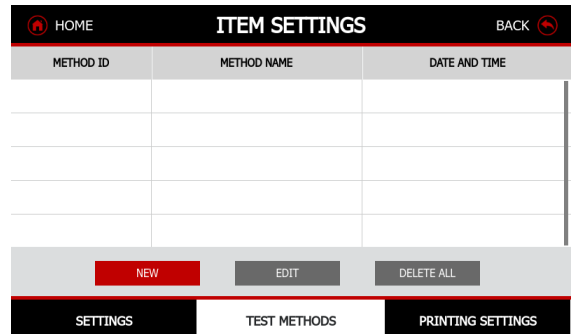


2. Press the NEW button to create a method.

Users can create maximum 25 methods for pipette adjustment.

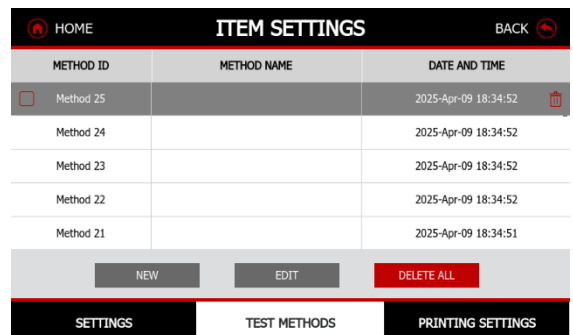


3. To rename the method,

- Check the box in front of the method ID
- Press the METHOD NAME field, and the EDIT icon would be visible.
- Press the EDIT icon to rename the method.

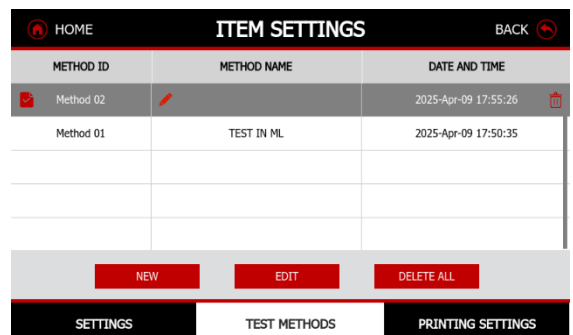


Edit a Method

1. Select the method by check the box in front of the METHOD ID.

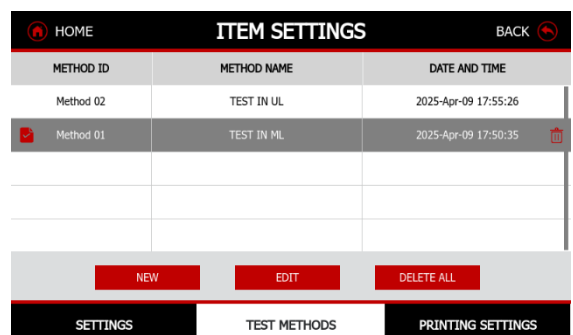


2. Press the EDIT button to enter the editing screen.

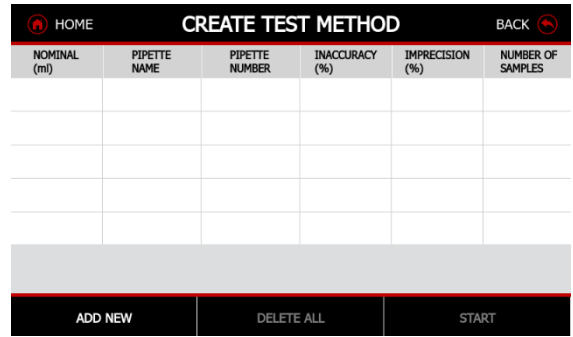


3. Press the ADD NEW to create a new test. A test includes the following parameters:


- Nominal (in the unit selected in ITEM SETTINGS): Set up the nominal capacity of the pipette.
- Pipette Name
- Pipette Number
- Inaccuracy (%)
- Imprecision (%)
- Number of Samples

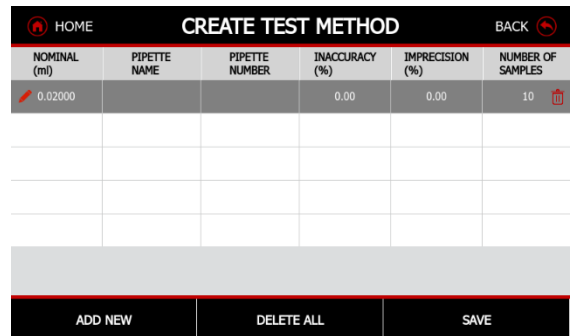

4. Press the field and the EDIT icon will be visible. Press the icon to edit the parameters of a test.

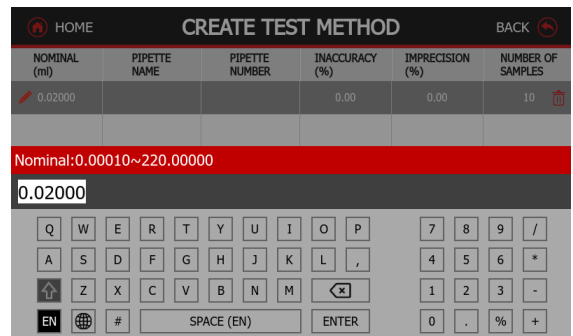

5. Continue to press ADD NEW button to add more tests. Users can create maximum 5 tests in a Method.

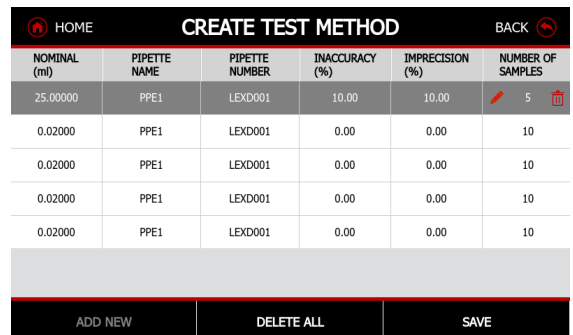


6. To delete a test, press the RUBISH BIN icon at the end of the row.

7. Press the SAVE button to save the changes.

Activate a Method

1. To activate a method in the adjustment, check the box in front of the Method ID.
2. This method will be activated for pipette adjustment. In the main screen, the METHOD hotkey will display the method ID.

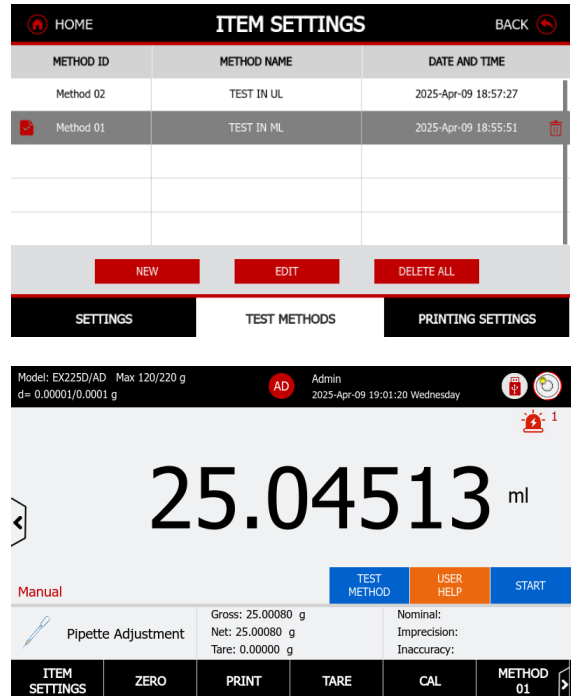


Delete a Method

To delete a method, press the RUBBISH BIN icon at the end of the row. Users can also delete all methods by pressing DELETE ALL button.

4.14.5 Begin Pipette Adjustment in Manual Mode

Step 1: Item Settings

1. Turn On the Auto Process toggle.

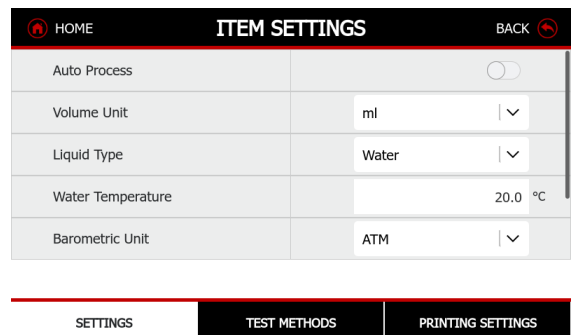

2. Set up Volume Unit, Liquid Type, Water Temperature, Liquid Density (if Liquid type is Other), Barometric Unit, and Pressure in ITEM SETTINGS.

Step 2: Set up a Method


Set up and activate a method for testing. For details, please refer to [4.14.4 Test Method](#).

This instruction uses a method as the right image displayed.


| HOME | | CREATE TEST METHOD | | | | BACK |
|--------------|--------------|--------------------|----------------|-----------------|-------------------|------|
| NOMINAL (ml) | PIPETTE NAME | PIPETTE NUMBER | INACCURACY (%) | IMPRECISION (%) | NUMBER OF SAMPLES | |
| 0.50000 | PPE1 | LDUA001 | 10.00 | 10.00 | 3 | |
| 2.00000 | PPE2 | LDUA002 | 10.00 | 10.00 | 3 | |
| | | | | | | |
| | | | | | | |
| | | | | | | |
| | | | | | | |
| ADD NEW | | DELETE ALL | | START | | |

Step 3: Start Pipette Adjustment


1. Press the START button at this page or at the Main Screen.
2. Place a container on the pan, then press TARE
3. Dispense sample 1 into the container. Then press Accept to record the value.

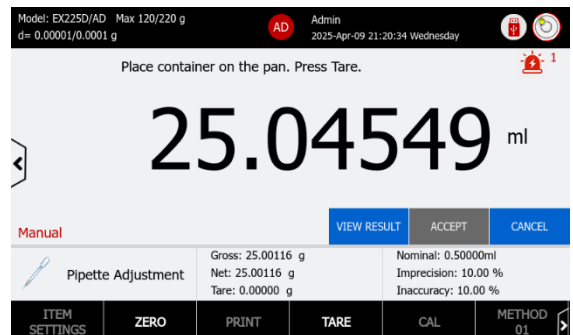

4. Follow the instruction to dispense the next sample. Continue the step 3 until all samples are dispensed.
5. When the first test is completed, the balance will ask if you want to start the next test.

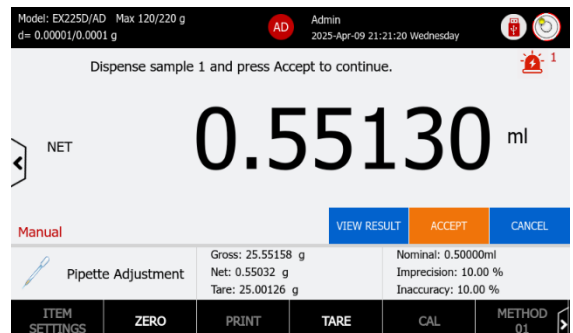

6. Press Yes to continue the next test. Follow the step 3 and 4 until all tests are completed. Press No to finish the test and to view the result.

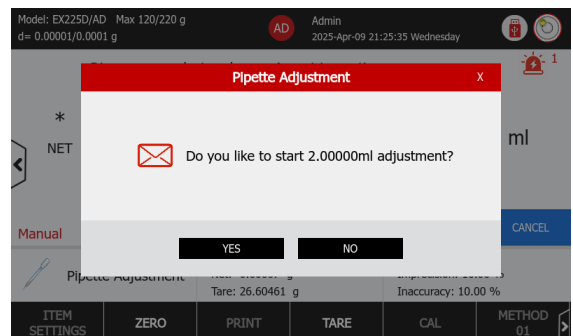

Step 4: View the Results

| HOME | | PIPETTE ADJUSTMENT RESULT | | | BACK | |
|---|--------------|---------------------------|---|---|--------|--|
| NOMINAL (ml) | PIPETTE NAME | PIPETTE NUMBER | INACCURACY (%) | IMPRECISION (%) | STATUS | |
| 0.50000 | PPE1 | LDUA001 | E%: 7.11 % E% Limit: 10.00 % Average:0.53556 ml | CV%: 4.90 % CV% Limit: 10.00 % Std:0.02624 ml | Pass | |
| 2.00000 | PPE2 | LDUA002 | E%: 6.73 % E% Limit: 10.00 % Average:2.13460 ml | CV%: 2.46 % CV% Limit: 10.00 % Std:0.05258 ml | Pass | |
| <div style="display: flex; justify-content: space-around;"> VIEW RESULT PRINT RESULTS DELETE ALL EXPORT TO PDF </div> | | | | | | |

4.14.6 Begin Pipette Adjustment in Auto Mode

Step 1: Item Settings

1. Turn off the Auto Process toggle.

| HOME | | ITEM SETTINGS | | BACK | |
|---|--|---------------|-------------------------------------|------|--|
| Auto Process | | | <input checked="" type="checkbox"/> | | |
| Volume Unit | | ml | v | | |
| Liquid Type | | Water | v | | |
| Water Temperature | | 20.0 | °C | | |
| Barometric Unit | | ATM | v | | |
| <div style="display: flex; justify-content: space-around;"> SETTINGS TEST METHODS PRINTING SETTINGS </div> | | | | | |

2. Set up Volume Unit, Liquid Type, Water Temperature, Liquid Density (if Liquid type is Other), Barometric Unit, and Pressure in ITEM SETTINGS.

Step 2: Set up a Method


Set up and activate a method for testing. For details, please refer to [4.14.4 Test Method](#).

This instruction uses a method as the right image displayed.


| HOME | | CREATE TEST METHOD | | | BACK | |
|--|--------------|--------------------|----------------|-----------------|-------------------|--|
| NOMINAL (ml) | PIPETTE NAME | PIPETTE NUMBER | INACCURACY (%) | IMPRECISION (%) | NUMBER OF SAMPLES | |
| 0.50000 | PPE1 | LDUA001 | 10.00 | 10.00 | 3 | |
| 2.00000 | PPE2 | LDUA002 | 10.00 | 10.00 | 3 | |
| <div style="display: flex; justify-content: space-around;"> ADD NEW DELETE ALL START </div> | | | | | | |

Step 3: Start Pipette Adjustment


1. Press the START button at this page or at the Main Screen.
2. Place a container on the pan. The balance will perform tare automatically.

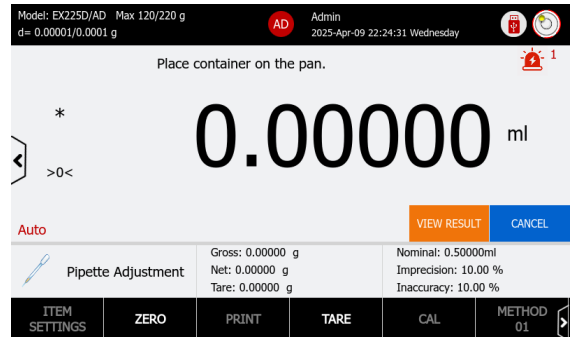

3. Dispense sample 1. The balance will accept the stable weight automatically.

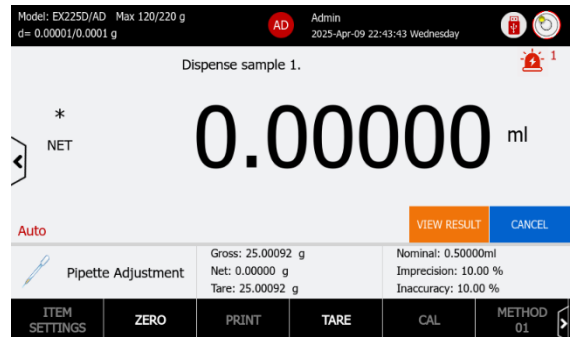

4. Follow the instruction to dispense the next sample. Continue the step 3 until all samples are dispensed.
5. When the first test is completed, the balance will ask if you want to start the next test.

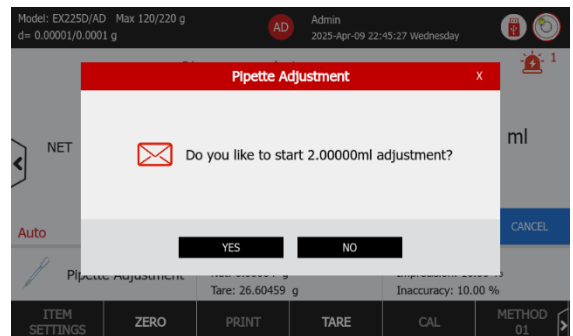

6. Press Yes to continue the next test. Follow the step 3 and 4 until all tests are completed. Press No to finish the test and to view the result.

Step 4: View the Results

| HOME | | PIPETTE ADJUSTMENT RESULT | | | | BACK |
|--------------|--------------|---------------------------|--|--|--------|---------------|
| NOMINAL (ml) | PIPETTE NAME | PIPETTE NUMBER | INACCURACY (%) | IMPRECISION (%) | STATUS | |
| 0.50000 | PPE1 | LDUA001 | E%: 7.11 % E% Limit: 10.00 % Average: 0.53556 ml | CV%: 4.90 % CV% Limit: 10.00 % Std: 0.02623 ml | Pass | |
| 2.00000 | PPE2 | LDUA002 | E%: 6.73 % E% Limit: 10.00 % Average: 2.13463 ml | CV%: 2.47 % CV% Limit: 10.00 % Std: 0.05263 ml | Pass | |
| | | | | | | |
| | | | | | | |
| VIEW RESULT | | PRINT RESULTS | | DELETE ALL | | EXPORT TO PDF |

4.14.7 Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to [Error! Reference source not found. Error! Reference source not found.](#)

4.15 SQC

The SQC application is used to determine the homogeneity of items in a batch and batches over time.



Application Icon

The screenshot shows the SQC application interface with the following components labeled:

- Application Guide Message:** A black header bar at the top containing model information (EX225D/AD, Max 120/220 g), user info (Admin, 2025-Apr-08 18:21:19 Tuesday), and a red 'AD' indicator.
- Main Display Line:** A large central display showing a weight of 25.00080 g with a '*' symbol above it.
- Application Buttons:** A row of buttons including 'VIEW RESULT' (grey) and 'START' (blue).
- Reference Fields:** A section showing 'Sample Status' as 'Recorded Sample: 0/30' and 'Application Name' as 'SQC'. It also displays weight statistics: Gross: 25.00080 g, Net: 25.00080 g, Tare: 0.00000 g, and Samples: 0. Below these are two rows of statistics: '+T1>N>-T1: 0, 0.00%' and '+T2>N>-T2: 0, 0.00%'.
- Hotkeys:** A bottom navigation bar with buttons for 'ITEM SETTINGS', 'ZERO', 'PRINT', 'TARE', 'CAL', and 'BATCH BATCH 25'.

4.15.1 Reference Fields

| Reference Fields | Description |
|---------------------|--|
| Samples | Count of tested samples |
| +T1>N>-T1: 0, 0.00% | Statistics of samples that lies between +T1 and -T1. |
| +T2>N>-T2: 0, 0.00% | Statistics of samples that lies between +T2 and -T2. |

4.15.2 Begin SQC

Step 1: Set up SQC Batch.

SQC Batch Library allows users to set up sample number of the batch, sample type, EU Criteria, Nominal Weight, Tolerances (when EU Criteria is off), Tare Type, Sample Name, Sample ID, Batch ID, Lot ID, Project Name, and Customized IDs.


The SQC Batch Library can store up to 25 records.

Users must load a batch to start SQC tests.


Create a Batch Library Record

1. Press the +New Batch button. When a new library is created, it is inactivated and not editable.
2. Activate the new batch by check the box in front of the batch id.

The batch status will be Activated and editable.


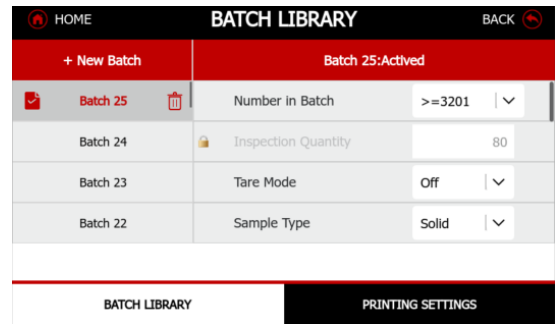


3. Edit the data fields

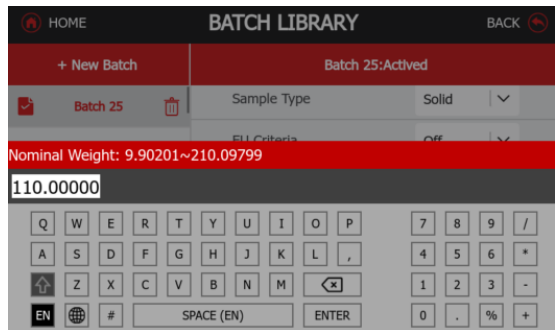


Activate a Batch

To activate a library item, check the box in front of the library item.

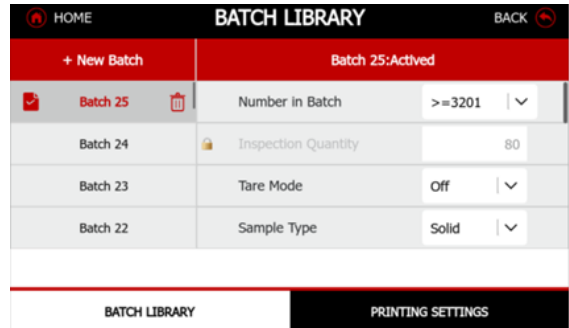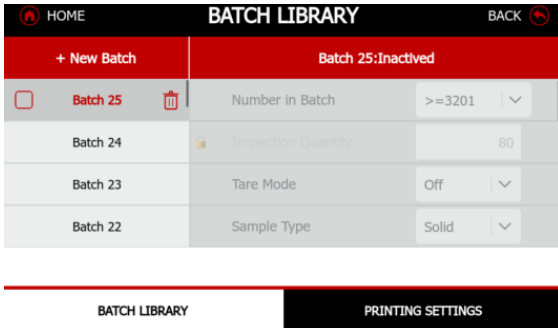

After the batch is loaded,

- On the screen, it would notify that the library is activated.
- On the application main screen, the library button will become the activated library id.

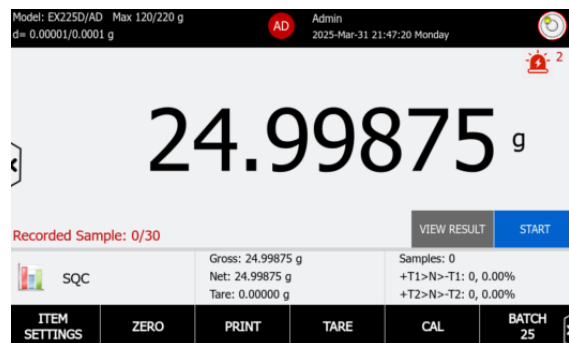

Delete a Library of Current Application





Press on the RUBISH BIN button to delete a library.

SQC Batch Library Settings

| Library Item | Options | Description |
|-----------------------------------|--|---|
| Number in Batch | <ul style="list-style-type: none"> From 100 ~ 500 From 501 ~ 3200 >= 3200 100% Inspection | Set up the total item number in the batch. |
| Inspection Quantity [Conditional] | 3~500 | Set up the inspection quantity when Number in Batch is set to 100% Inspection. |
| Tare Mode | <ul style="list-style-type: none"> Off Pretare | When tare mode is off, users can tare the container by pressing Tare button When set to Preset, Users can set up a fixed tare value. |
| Pretare [Conditional] | / | Visible when Tare Mode is set to Preset. Set up a fixed tare value. |
| Sample Type | <ul style="list-style-type: none"> Solid Liquid | Set up the sample type. |
| Density [Conditional] | Input liquid density in g/ml | Visible when Sample Type is set to Liquid. User can set up the density value to auto convert the sample weigh to ml. |
| EU Criteria | <ul style="list-style-type: none"> On Off | On: Evaluate samples based on tolerances stated in the legal directive (refer to -T1 / +T1 (main tolerance) based on the target.) Off: Evaluate samples based on the tolerance set up by users. |
| Nominal Weight | Input nominal target weight in g or ml. | Set the value of nominal target weight in g or ml. <ul style="list-style-type: none"> To weigh in g: Set the Sample Type to Solid. To weigh in ml: Set the Sample type to Liquid. |
| +T2 | Legal: The field is locked Free: Set Target Over 2 value | In Legal Mode, +T2 (Target Over 2) is twice of the +T1 value. In Free mode, +T2 can be defined freely. |
| +T1 | Legal: The field is locked Free: Set Target Over 1 value | In Legal Mode, +T1 (Target Over 1) is determined based on the target value according to -T1 / +T1 (main tolerance) based on the target In Free mode, +T1 can be defined freely. |
| -T1 | Legal: The field is locked Free: Set Target Under 2 value | In Legal Mode, -T1 (Target Under 1) is determined based on the target value according to -T1 / +T1 (main tolerance) based on the target In Free mode, -T1 can be defined freely. |
| -T2 | Legal: The field is locked Free: Set Target Under 1 value | In Legal Mode, -T2 (Target Under 2) is twice of the -T1 value. In Free mode, -T2 can be defined freely. |

Step 2: Weigh Individual Samples


1. Press the Start button to begin.
2. If there's an item on the weighing pan, remove the item from the pan.
3. Follow the screen instruction to weigh a sample.
4. Remove the sample from the pan
7. Follow the step 3 and 4 until all samples are weighed.

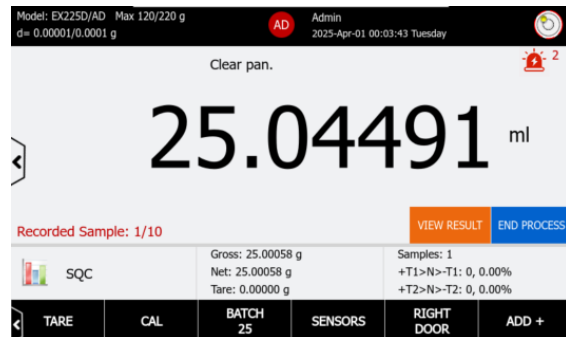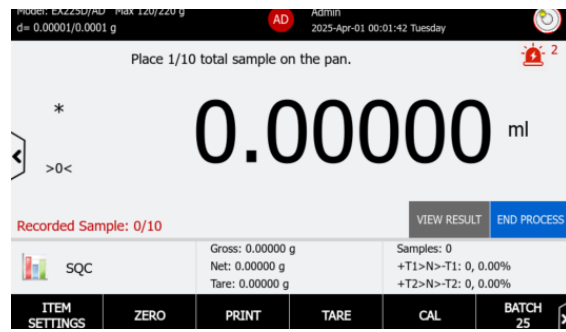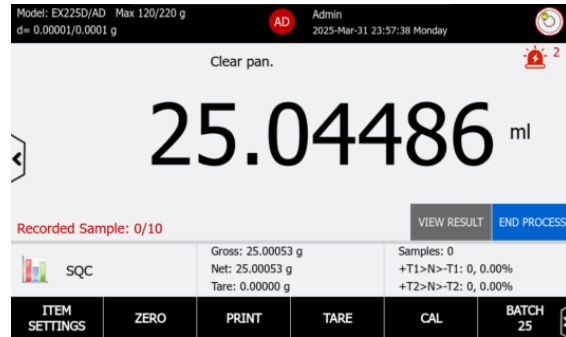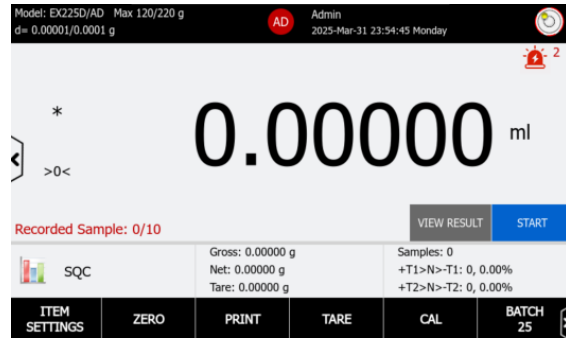


Step 3: Lot Accepted or Refused.

The scale would auto determine the lot to be accepted or refused based on [4.15.3 Acceptance Criterion](#).

If a 2nd samples need to be provided, SQC would update the required samples number directly.

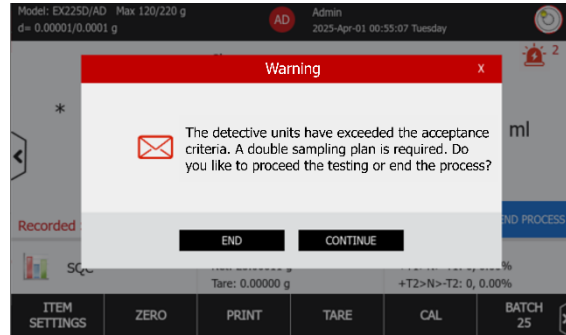

The balance would inform users if the test was already failed before all samples are weighed.

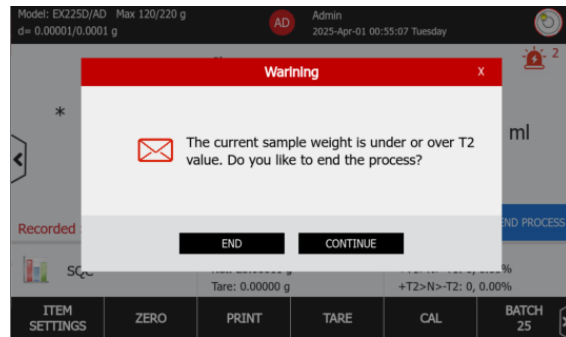


Step 4: View and Print Statical Data.

The result screen will be displayed when the test is ended. The balance presents the data in both list and graphic views.

List View

| HOME | | VIEW RESULT | | | BACK |
|----------------|--------------|---------------|------------|---------------|------|
| # | BATCH NUMBER | SAMPLE ID | WEIGHT (g) | T1/T2 | |
| 001 | Batch 03 | Sample 001 | 249.99 | >+T2 | |
| 002 | Batch 03 | Sample 002 | 199.99 | >=-T1 & <=+T1 | |
| 003 | Batch 03 | Sample 003 | 199.99 | >=-T1 & <=+T1 | |
| 004 | Batch 03 | Sample 004 | 200.00 | >=-T1 & <=+T1 | |
| Result: Reject | | | | | |
| VIEW GRAPHICS | | PRINT RESULTS | | EXPORT TO PDF | |

Graphic View

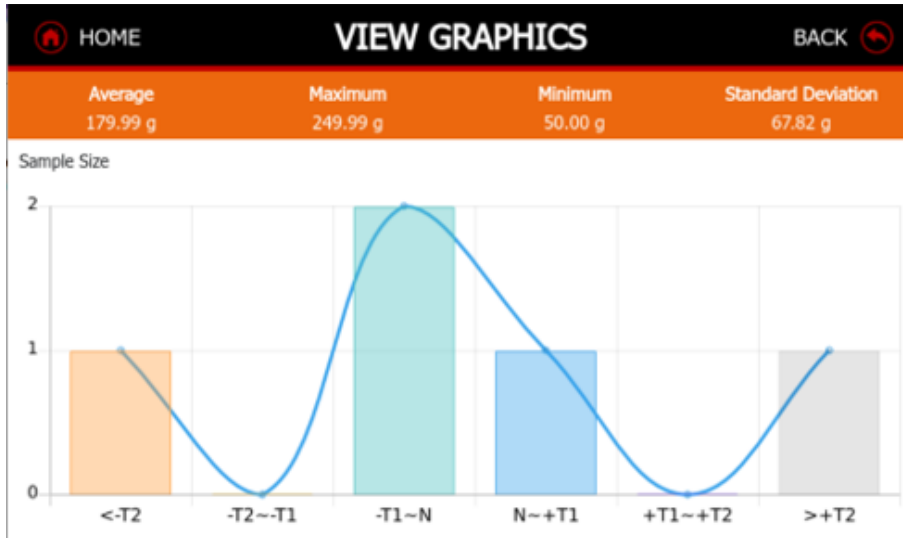

4.15.3 Acceptance Criterion

| SQC Result | Description |
|--------------|--|
| Lot Accepted | <p>The lot will be considered accepted when all conditions are met:</p> <ul style="list-style-type: none"> The average weight of samples must be greater than the nominal weight. The defective sample¹ number in the 1st samples is less than or equal to the 1st acceptance criterion. <p>If the defective samples number lies between the 1st Acceptance and Rejection criterion, a 2nd samples should be tested.</p> <p>The lot will be considered accepted when all conditions are met:</p> <ul style="list-style-type: none"> The average weight of aggregated samples must be greater than the nominal weight. The aggregated defective samples number is less than or equal to the 2nd acceptance criterion. |
| Lot Rejected | <p>The lot will be rejected when any of the condition is met:</p> <ul style="list-style-type: none"> if any of the sample weight is not greater than the -T2 limits; Defective sample number is equal to or greater than Rejection criterion. The average weight of samples is less than the nominal weight |

Note: 1 Defective sample are samples with a weight that lies between -T1 and -T2.

Sample Number and Criterion Table

| Lot Piece Number | Samples | | | Number of defective samples | |
|------------------|-----------------|--------|-------------------|-----------------------------|---------------------|
| | Order | Number | Aggregated Number | Acceptance criterion | Rejection criterion |
| 100 to 500 | 1 st | 30 | / | 1 | 3 |
| | 2 nd | 30 | 60 | 4 | 5 |
| 501 to 3200 | 1 st | 50 | / | 2 | 5 |

| | | | | | |
|---------------|-----------------|----|-----|---|---|
| | 2 nd | 50 | 100 | 6 | 7 |
| 3201 and over | 1 st | 80 | / | 3 | 7 |
| | 2 nd | 80 | 160 | 8 | 9 |

4.15.4 Determination of -T1, -T2, +T1, and +T2

In EU Criteria On mode, -T1 and +T1 are determined based on the target according to the below table.

-T2 and +T2 are the value twice of -T1 and +T1.

In Free mode, -T1, -T2, +T1, and +T2 values can be customized by the users.

Samples with a weight that lies between -T1 and -T2 are determined as defective.

-T1 / +T1 (main tolerance) based on the target

| TARGET IN GRAMS OR IN MILLILITRES | -T1 / +T1 TOLERANCE ON THE TARGET | |
|--------------------------------------|-----------------------------------|------------|
| | in % | in g or ml |
| From 5 to 50 | 9 | - |
| From 50 to 100 | - | 4,5 |
| From 100 to 200 | 4,5 | - |
| From 200 to 300 | - | 9 |
| From 300 to 500 | 3 | - |
| From 500 to 1000 | - | 15 |
| From 1000 to 10000 | 1,5 | - |
| From 10000 to 15000 | - | 150 |
| Beyond 15000 | 1 | - |

4.15.5 Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to [Error! Reference source not found. Error! Reference source not found.](#)

4.16 Fill Weight Variation

Fill Weight Variation checks the homogeneity of a series of items.



Application Icon

Model: EX225D/AD Max 120/220 g
d= 0.00001/0.0001 g

Andrew
2025-Mar-19 10:42:21 Wednesday

* 3.99928 g

SAMPLES AVERAGE +/- LIMIT % VIEW RESULT START

Application Name: Fill Weight Variation

Reference Fields:
Gross: 10.0 g
Net: 10.0 g
Tare: 0.0 g
Samples: 20
Average: ---
+/- Limit %: 7.50 %

Hotkeys: ITEM SETTINGS ZERO PRINT TARE CAL SENSORS

4.16.1 Application buttons


| Application Button | Icon | Description |
|--------------------|------|---|
| SAMPLES AVERAGE | | Press to set sample quantity and establish sample average weight. |
| +/- LIMIT % | | Press to set the +/- Limit %. Note: When Acceptance Criteria is set to USP <2091> mode, the button will be greyed and inactivated. |
| VIEW RESULT | | Press to the button to establish the reference weight derived from a sample weight. |
| START | | The icon directly shows the number of data that have been accumulated for statistics, Press the button to view statistics report. |

4.16.2 Reference Fields

| Reference Fields | Description |
|------------------|--|
| Samples | Display the sample quantity for setting the sample average weight. |
| Average | Display the sample average weight. |
| +/- Limit % | Display the +/- Limit % value. |

4.16.3 Item Settings – Fill Weight Variation

Press the Item Settings button to enter the application settings.

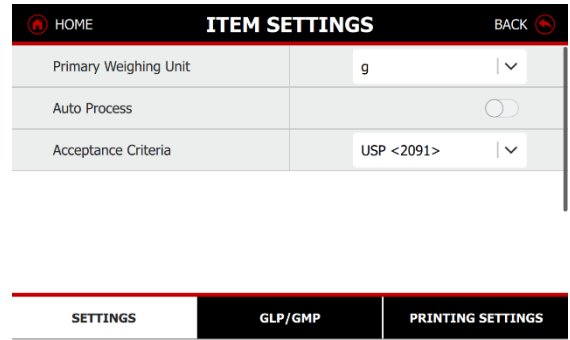

| Item | Description |
|---------------------|---|
| Auto Process | <ul style="list-style-type: none"> On: The balance will accept the sample weight automatically. Then it will tare and progress to weigh the next sample. Off: The sample weight will be recorded by pressing the ACCEPT button. |
| Acceptance Criteria | Off: User can customize the +/- Limit % value. USP <2091>: +/- Limit % value is auto calculated according to the USP guidelines. |

4.16.4 Begin Tests with USP <2091>


Step 1: Set up Acceptance Criteria

1. Press the Item Settings button to enter the application settings.
2. Set Acceptance Criteria to Off.


Step 2: Set up Samples Average

1. Press the Item Settings button to enter the application settings.
Set Acceptance Criteria to USP <2091>.
2. Press the SAMPLES AVERAGE button.

3. Set up the sample quantity.

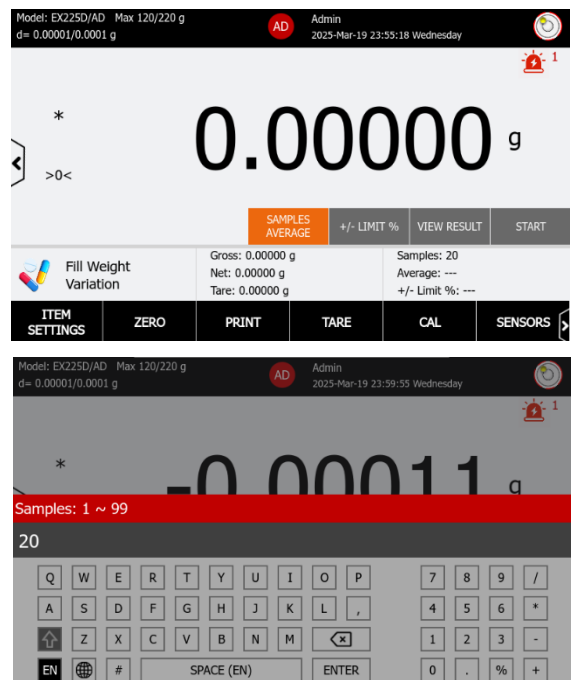


4. Place the samples onto the weighing pan for averaging.

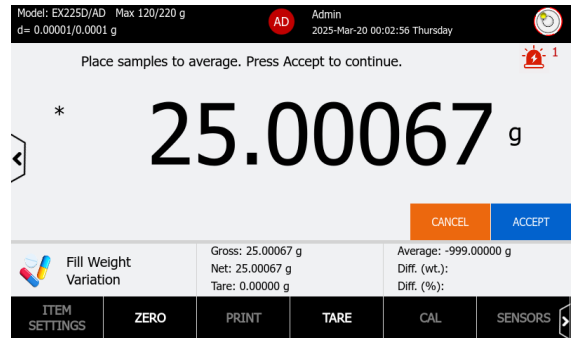


5. Press the ACCEPT button to confirm.

The value of sample quantity, average weight and +/- limit will be calculated and displayed in the reference fields. The START button will be activated when samples average is set up.

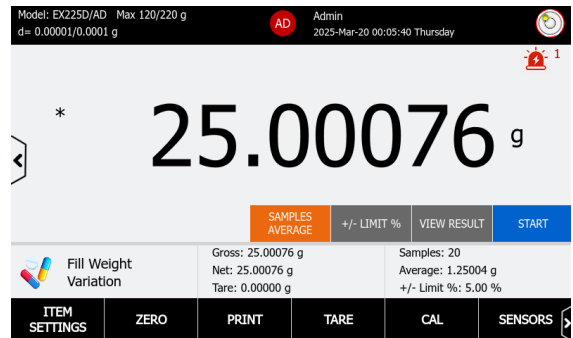


Step 3: Weigh the Samples

1. Press the START button.

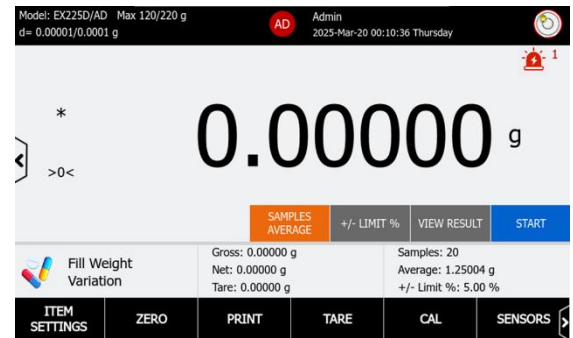

2. Please a sample onto the weighing pan.

- Auto Process Off: Press the ACCEPT button to record the weight value, and progress to the next sample.
- Auto Process On: Stable weight will be recorded automatically. The balance will tare and progress to the next sample.

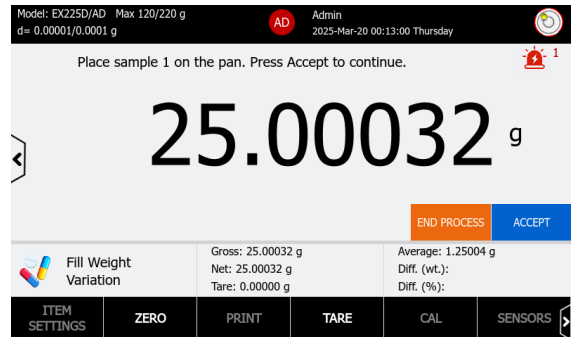


3. Repeat the step 2 until all samples are weighed.

4. When all samples are weighed, the process will end and the screen will display the test results.

Step 4: Check the Results

The result screen will be displayed when a text is finished. User can also press the VIEW RESULT button to view the result of the latest test.

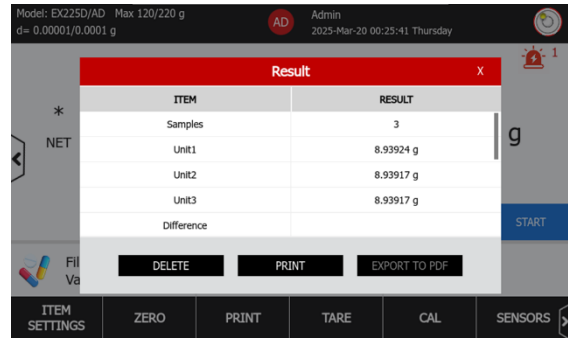

The test result would include the following information:

- Sample quantity
- Sample x weight
- Difference in weight and percentage
- Result (Accept and Fail numbers)
- Average Weight
- +/- Limit %


4.16.5 Begin Tests with Customized +/- Limits

Step 1: Set up Acceptance Criteria


5. Press the Item Settings button to enter the application settings.
6. Set Acceptance Criteria to Off.

Step 2: Set up Samples Average


1. Press the SAMPLES AVERAGE button.

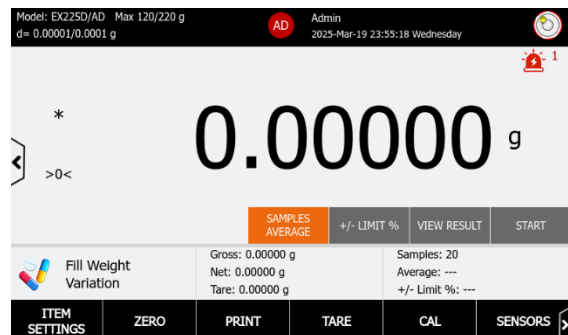


2. Set up the sample quantity.

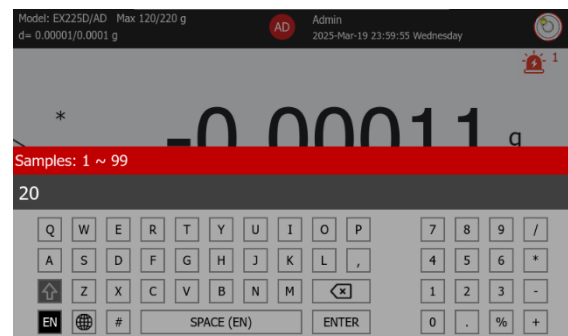

3. Place the samples onto the weighing pan for averaging.

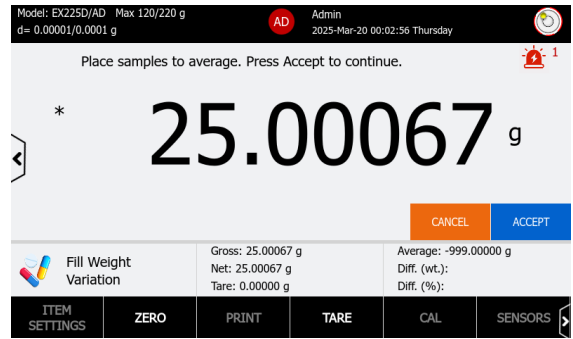


4. Press the ACCEPT button to confirm.
 The value of sample quantity and average weight will be calculated and displayed in the reference fields.
 The START button will be activated when samples average is set up.

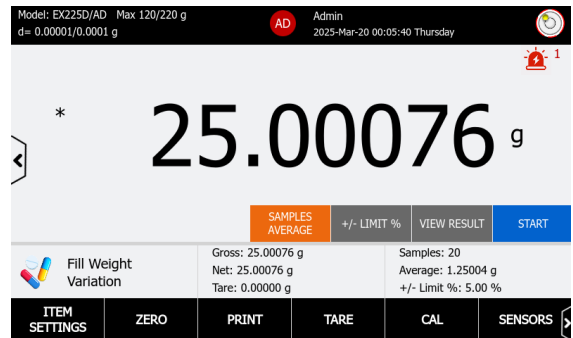


Step 3: Set up +/- Limits %

1. Press the +/- LIMIT % button.

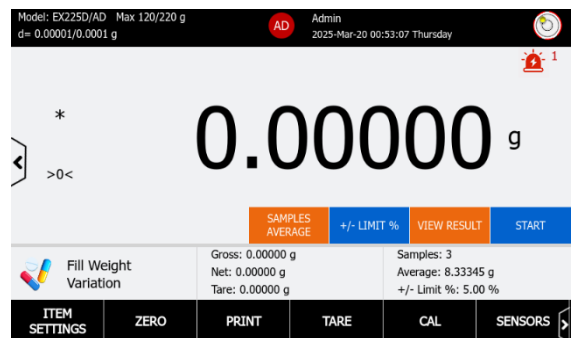


2. Set the +/- LIMIT value in %.

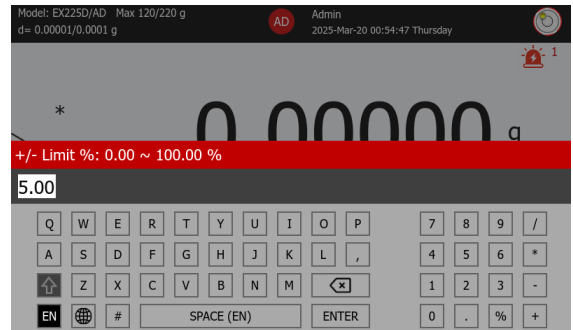


Step 4: Weigh the Samples

1. Press the START button.

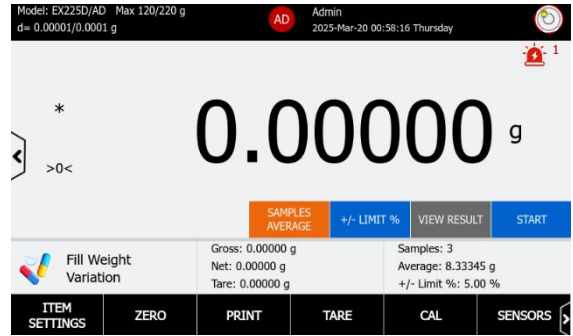

2. Please a sample onto the weighing pan.

- Auto Process Off: Press the ACCEPT button to record the weight value, and progress to the next sample.
- Auto Process On: The balance will record the stable weight value and progress to the next sample automatically.

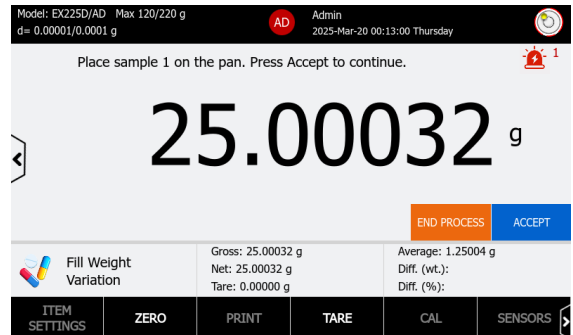

3. Repeat the step 2 until all samples are weighed.

4. When all samples are weighed, the process will end and the screen will display the test results.

Step 5: Check the Results

The result screen will be displayed when a text is finished. User can also press the VIEW RESULT button to view the result of the latest test.

4.16.6 Printing Settings

The Explorer Plus balance offers advanced print settings. Users can customize the output format, and content and direct it to Excel or save it in a USB-compatible format. Additionally, users can review the print content to ensure all items are correctly selected before printing on paper or PC.

For details of Printing Setting, please refer to 6.9.5 Printing Settings.

4.17 Flow Rate Control

Flow Rate Control aims at managing the volume or mass of fluid (liquid or gas) that passes through a pump system over a specific period.



Application Icon

The screenshot shows the main display of the Explorer Plus™ Balances. At the top, it displays the model (EX225D/AD), maximum capacity (Max 120/220 g), and density (d= 0.00001/0.0001 g). The user is logged in as Admin, and the date and time are 2025-Mar-24 18:34:37 Monday. A message at the top says "Press the Start button to initiate the flow rate measurement." The main display shows a large weight of 1.10067 g and a secondary unit of 1.10067 ml. Below the display are four buttons: TARGET VOLUME (orange), PUMP SPEED (blue), VIEW RESULT (grey), and START (blue). The mode is set to Target Control. The application name is Flow Rate Control. Reference fields show Gross: 1.10067 g, Net: 1.10067 g, Tare: 0.00000 g, Target Volume: 10.00 ml, Pump Speed: 25.00 ml/min, and Number of Pump Cycles: 6. At the bottom are hotkeys: ITEM SETTINGS, ZERO, PRINT, TARE, CAL, and LIBRARY.

Annotations on the right side of the screenshot:

- Application Guide Message
- Main Display Line
- Secondary Unit
- Application Buttons
- Reference Fields
- Hotkeys

4.17.1 Item Settings – Check Weighing

Press the ITEM SETTINGS button to enter the application settings.


The screenshot shows the ITEM SETTINGS screen. It has a HOME button on the left and a BACK button on the right. The screen is divided into three sections: Mode (Flow Rate Control), Pump Speed Unit (ml/min), and Audible Signal (off). At the bottom, there are three buttons: SETTINGS, SAMPLE LIBRARY, and PRINTING SETTINGS.

Mode

Users can select the Mode as Flow Rate Control or Target Control.

This screenshot shows the ITEM SETTINGS screen with the Mode dropdown menu open. The dropdown menu has two options: Flow Rate Control (highlighted in red) and Target Control. The other settings and buttons remain the same as in the previous screenshot.

Flow Rate Control

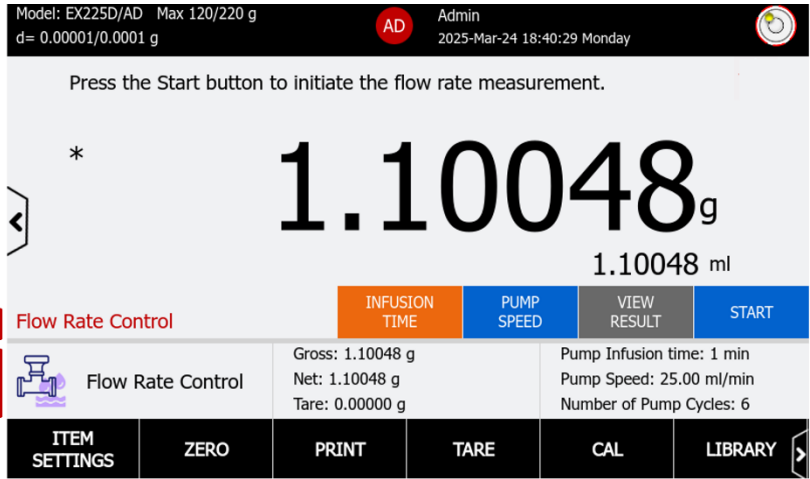

Application Guide Message
 Main Display Line
 Secondary Unit
 Application Buttons
 Reference Fields
 Hotkeys


Application buttons

| Application Button | Icon | Description |
|--------------------|------|---|
| INFUSION TIME | | Press to set up a testing time of each cycle. |
| PUMP SPEED | | Press to set up the pump speed. |
| VIEW RESULT | | Press to view the result of the latest test. |
| START | | Press the button to begin testing. |

Reference Fields

| Reference Fields | Description |
|-----------------------|--|
| Pump Infusion time | Display the pump infusion time. |
| Pump Speed | Display the theoretical pump speed. |
| Number of Pump Cycles | Display the pump cycles value configured in Library. |

Target Control

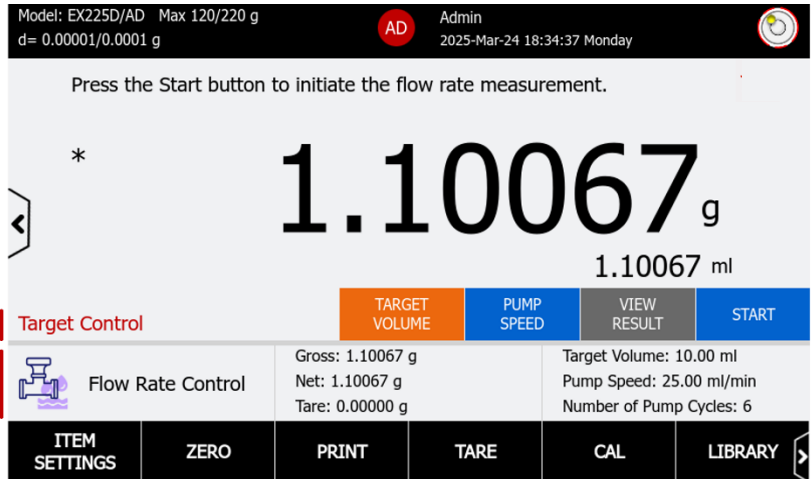

Application Guide Message

Main Display Line

Secondary Unit

Application Buttons

Reference Fields

Hotkeys

Application buttons

| Application Button | Icon | Description |
|--------------------|------|--|
| Target VOLUME | | Press to set up a target volume. |
| PUMP SPEED | | Press to set up the pump speed. |
| VIEW RESULT | | Press to view the result of the latest test. |
| START | | Press the button to begin testing. |

Reference Fields

| Reference Fields | Description |
|-----------------------|---|
| Target Volume | Display the target volume. |
| Pump Speed | Display the theoretical pump speed. |
| Number of Pump Cycles | Display the pump cycles number configured in Library. |

Pump Speed Unit

Users can set up the pump speed unit in g/s, g/min, ml/s, ml/min, ul/s, or ul/min.

Volume Unit

Users can set up the volume unit in ul, ml, or l.

The field is visible in Target Control mode.


4.17.2 Begin Flow Rate Control

Flow Rate Calculation


$$FlowRate = \frac{EndWeight - StartWeight}{LiquidDensity \cdot InfusionTime}$$

Step 1: Set Infusion Time

1. Press the INFUSION TIME button.

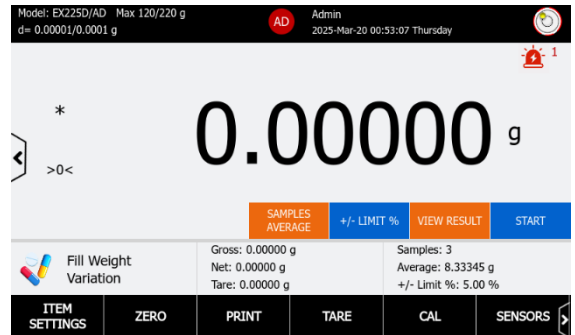

2. Set up the infusion time in minute.

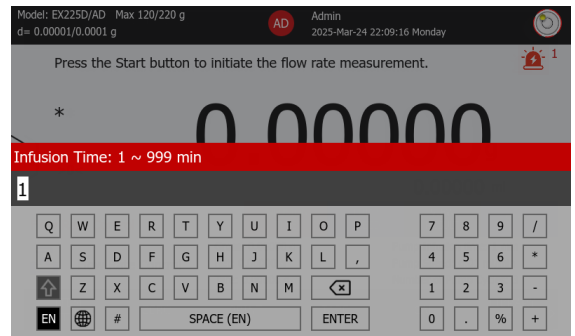


Note: Pump Time can also be set up in Library.


Step 2: Set up Number of Pump Cycles

Number of pump cycles can be setup in Library.

1. Press the Library button.

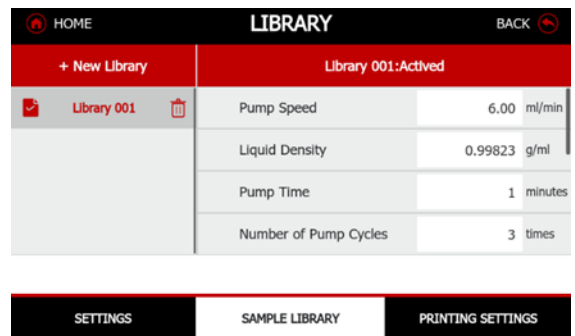


- Set the number of pump cycles.

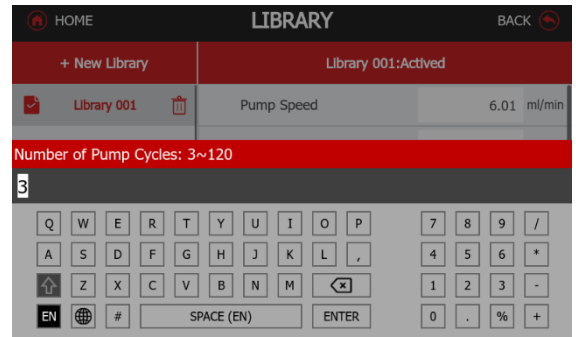


Step 3: Weigh the Samples

- Press the START button to initiate the flow rate measurement. The current weight will be recorded as Start Weight.

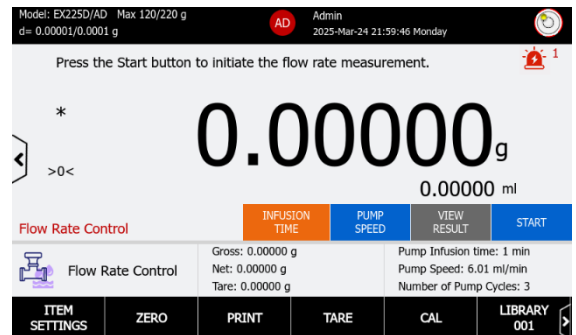

- Begin flow rate calculation. When reaching the time, the current weight will be recorded as the End Weight.
- The flow rate is calculated and displayed in orange font color.

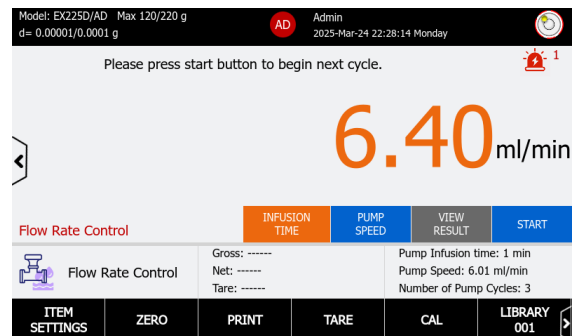


- Press the START button to begin the next cycle. Repeat the step 2 and 3 until all cycles are completed.
- When all cycles are finished, the result will be displayed.

Step 3: Check Flow Rate Result

The result pages displays information include Pump Speed, Max Flow Rate, Min Flow Rate, Average Flow Rate, and Flow Rate values presented graphically.

Each page will display six pump cycle results.

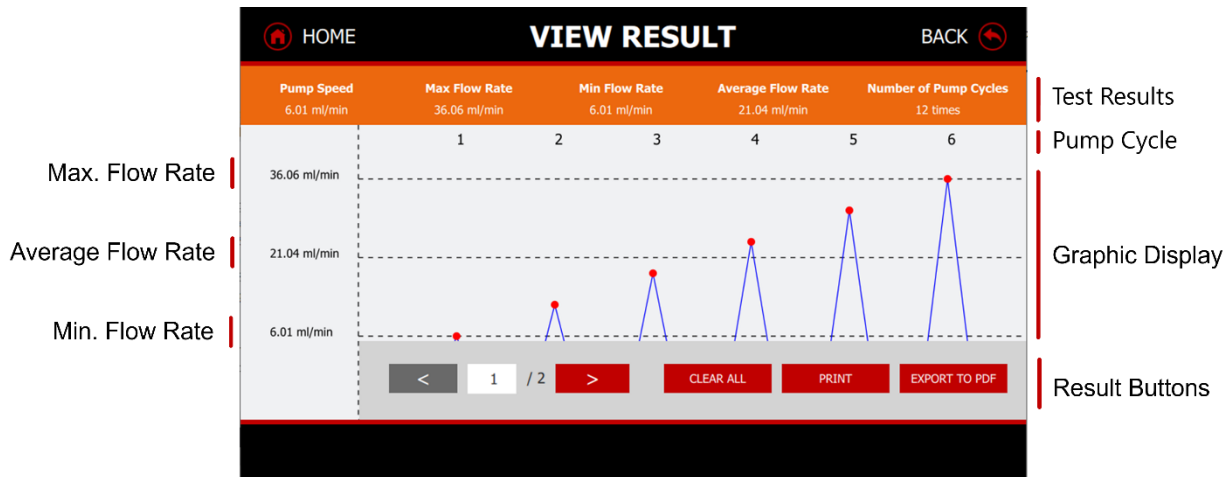


4.17.3 Begin with Target Rate Mode

Flow Rate Calculation in Target Mode


$$FlowRate = \frac{EndWeight - StartWeight}{LiquidDensity} / EndTime$$

Step 1: Set Infusion Time

1. Press the Target Volume button.

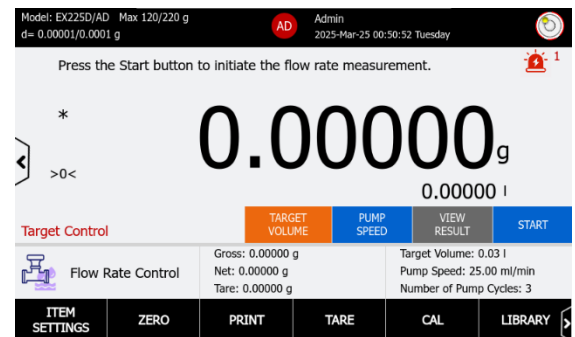

2. Set up the target volume.

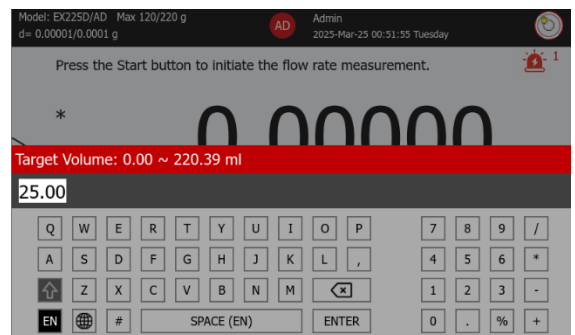


Note: Target volume can also be set up in Library.


Step 2: Set up Number of Pump Cycles

Number of pump cycles can be setup in Library.

1. Press the Library button.

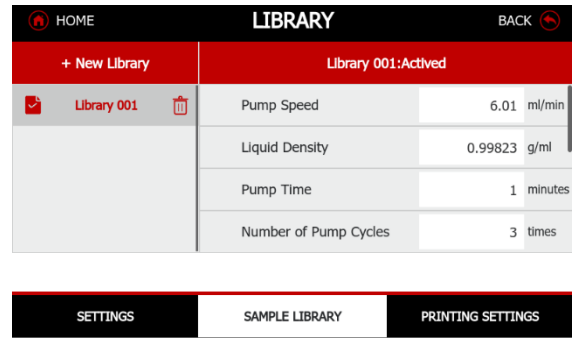


2. Set the number of pump cycles.

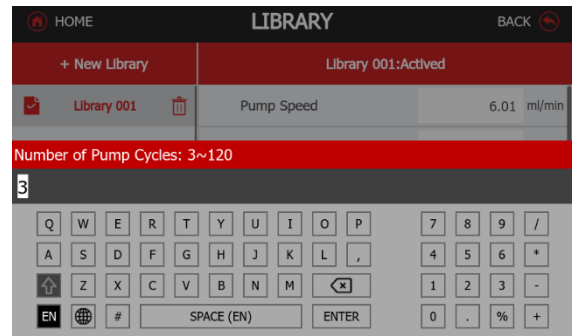


Step 3: Weigh the Samples

1. Press the START button to initiate the flow rate measurement.
The current weight will be recorded as Start Weight.

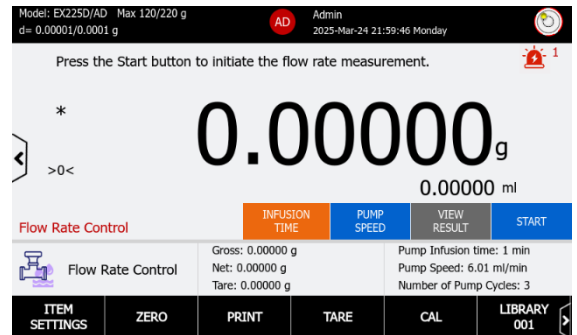

2. Begin flow rate calculation.
When reaching the time, the current weight will be recorded as the End Weight.

3. The flow rate is calculated and displayed in orange font color.

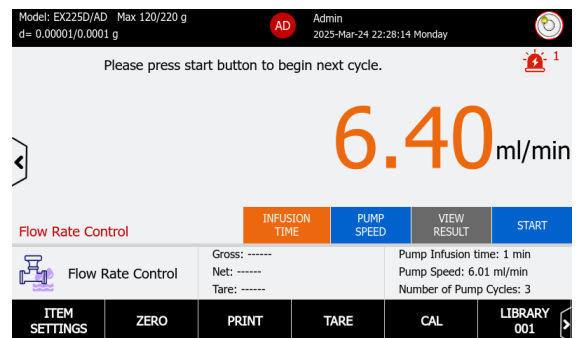


4. Press the START button to begin the next cycle.
Repeat the step 2 and 3 until all cycles are completed.

5. When all cycles are finished, the result will be displayed.

Step 3: Check Flow Rate Result

The result pages displays information include Pump Speed, Max Flow Rate, Min Flow Rate, Average Flow Rate, and Flow Rate values presented graphically.

Each page will display six pump cycle results.

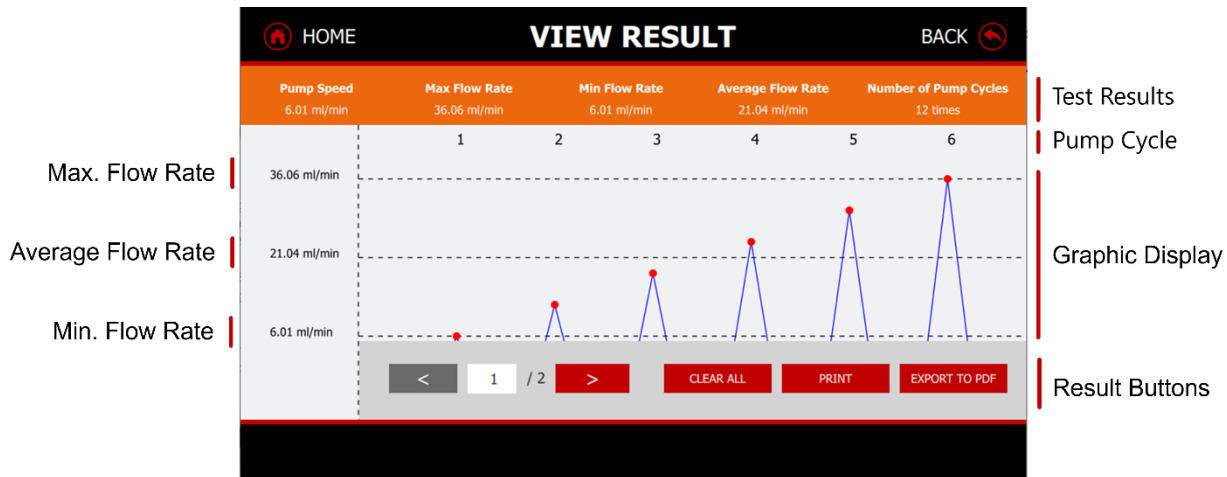

4.17.4 Sample Library – Flow Rate Control



Explorer Plus balance built in library simplifies working with multiple samples, allowing application settings to be saved and recalled for reuse.

The Sample Library can store up to 3000 records, without limitation on individual applications. To check the status of library usage, navigate to Main Menu > Library.

Note: In Main Menu > Library, the balance would notify users to delete or export libraries when the storage is used 80%.


Create a Library Record

1. Press the +New Library button. When a new library is created, it is inactivated and not editable.

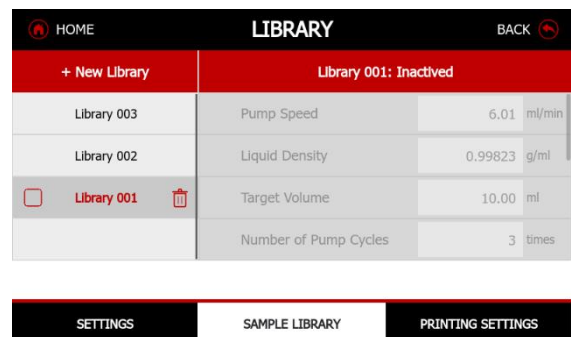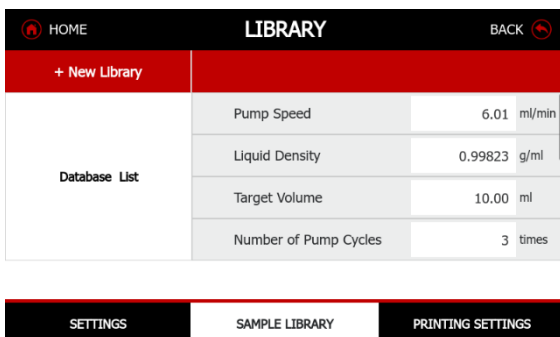


2. Activate the new library by check the box in front of the library id.

The library status will be Activated and editable.


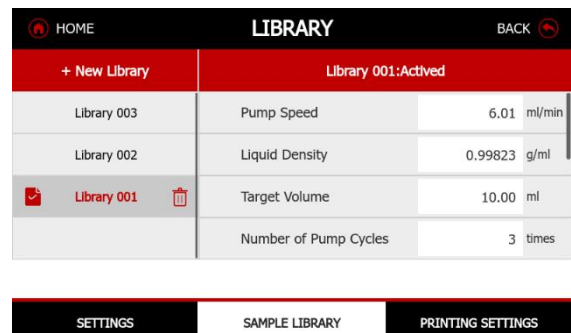


3. Edit the data fields using the keyboard. Press **ENTER** to confirm editing.

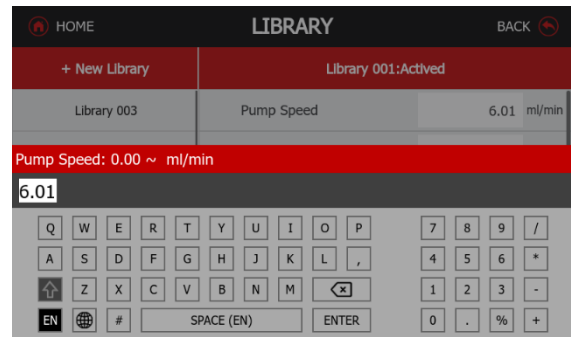


Activate a Library

To activate a library item, check the box in front of the library item.

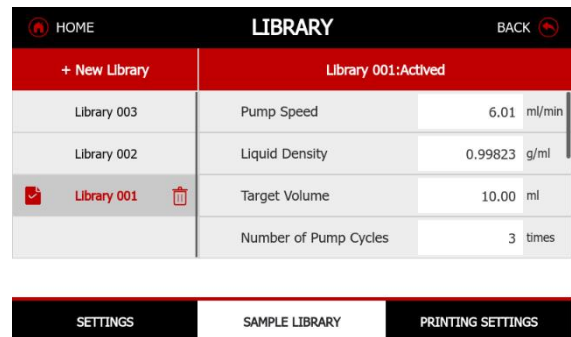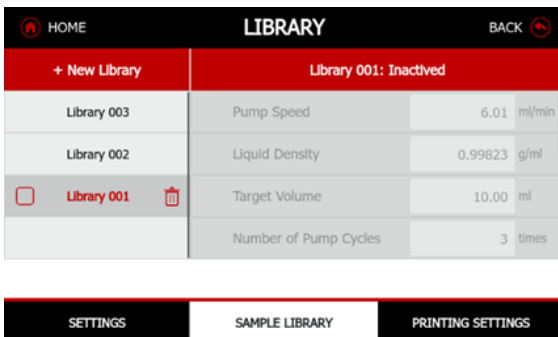

After the library item is loaded,

- On the screen, it would notify that the library is activated.
- On the application main screen, the library button will become the activated library id.

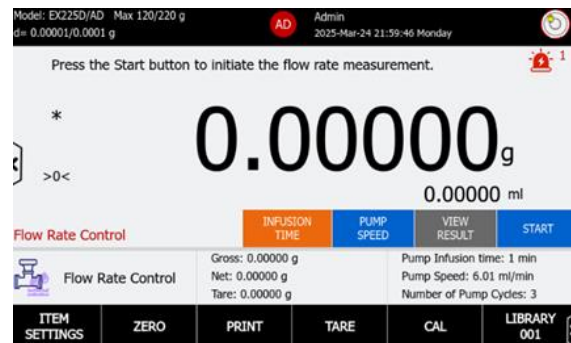

Delete a Library of Current Application

Press on the RUBISH BIN button to delete a library.


Manage Libraries

The balance allows users to review, delete, export or import libraries in Main Menu > Library.

For details, please refer to [6.5.1](#) Printing Settings

Users can set the Printing Settings for the Totalization application.

For details of Printing Setting, please refer to

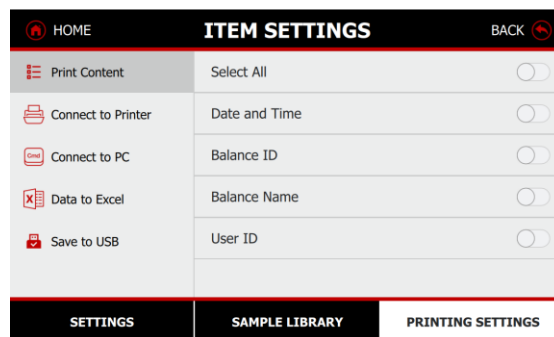


5.3 Quick Setup

The Quick Setup menu is designed for first-time users. By quickly navigating through this menu, you can easily configure the most desired functions.

5.3.1 Language


Set the language for menu display and pop-up messages. The default setting is configured based on the country of sale. The 13 available languages are English, German, French, Spanish, Italian, Polish, Czech, Hungarian, Portuguese, Chinese, Japanese, Korean, and Turkish.

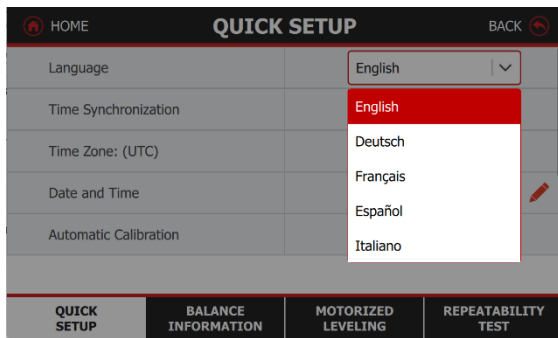

5.3.2 Time Synchronization/ Network Server


The new balance features Synchronization with Network (NTP function) which can use consistent time and data from the local network. The NTP system supports DNS domain name resolution.

- When you input the address of a public domain server, it will automatically convert it into an IP address. Here are some examples of NTP public time server address
 - “0.europe.pool.ntp.org” when the location is in Europe
 - “cn.ntp.org.cn” for China
 - “0.us.pool.ntp.org” for the US

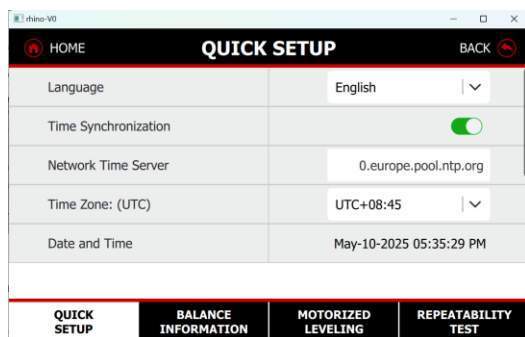

5.3.3 Date and Time

If you are offline use the balance, the new data/time setting fulfill your documentation needs with UTC time zone and customized the time format MMM-DD-YY HH-MM-SS and working weekdays setting.

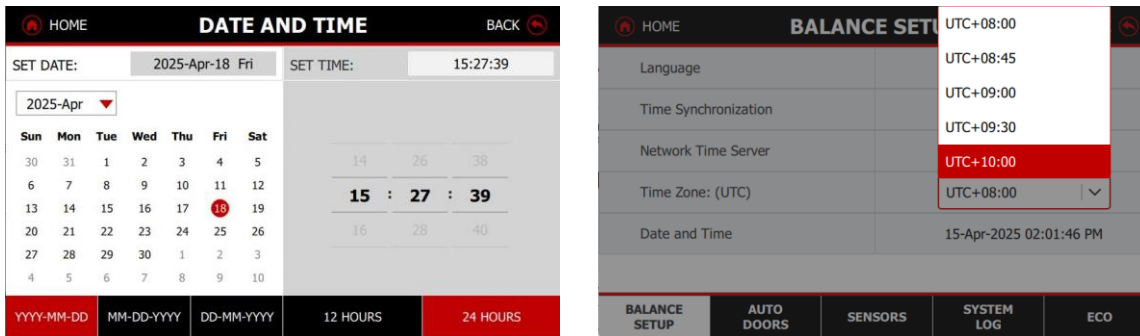

5.3.4 Automatic Calibration

When Automatic Calibration is set ON, the balance performs a self-calibration: AutoCal will automatically calibrate the Balance (using the internal mass) each time there is a change in temperature significant enough to affect accuracy or every 11 hours.

An information screen will appear when an Automatic Calibration is about to start. Three option buttons will be displayed:

- Now – Press to perform the calibration directly.
- After 5 min – Press to perform the calibration after 5 minutes.
- Deactivate – Press to deactivate the Automatic Calibration function

5.3.5 User Management

When **User Management** is set ON, the balance will be able to create up to 200 user accounts, featuring four predefined roles. 4-level user management with an underdefined user group with access rights to the balance.

- Administrator (1)
- Supervisor (maximum of 20)
- 179 allocated to other roles like operator, log viewer, and group users

See detail setting in User Management section 5.6

5.3.6 System Log


When the **System Log** is enabled, the balance can store up to 100,000 log entries. It has consist with System Log, Calibration Log, Printed Data Log, and Failure log. These electronic records include changes made on balance setting will keep records in system log file. e.g. every printed data, date/time change, balance setting changes, perform calibration action, user log in/log out, user account create/edit/delete and etc.

See detail setting in Maintenance section 5.6

5.3.7 Balance Information

The balance information will display the Information displayed includes:


- Balance Information: Balance ID, Balance Model, Balance Software Version, Digital Loadcell Type, and Digital Loadcell Software Version
 - The balance max capacity, readability, d and (e) value are displayed in the Upper left corner of the main screen.
- Environmental Parameters: Loadcell Temperature, Humidity, and Atmospheric Pressure
 - The loadcell temperature, humidity, and air pressure are displayed in the reference field.
 - The user can click on the display to turn it off.

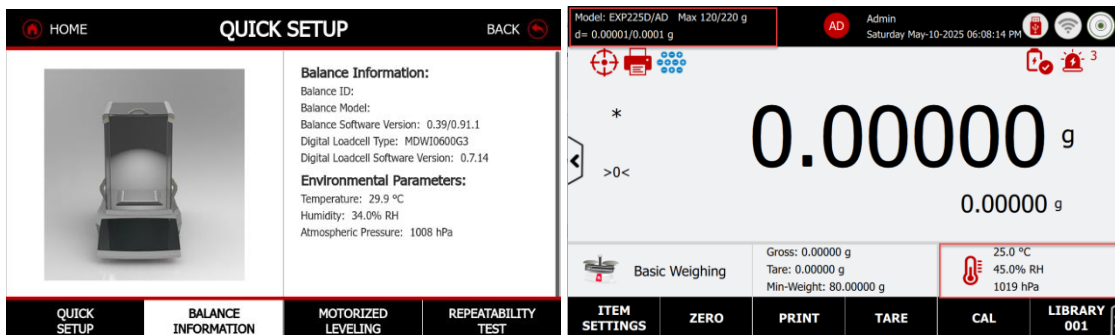


5.3.8 Motorized Leveling

The Motorized Leveling System is a simple and effective balance feature for the first setup user and the technician dealers. Press Start button to level the balance automatically. The initial digital level bubble will indicate the bubble real time statuses.

- Before starting the motorized leveling process, kindly remove the weight from the weighing pan.

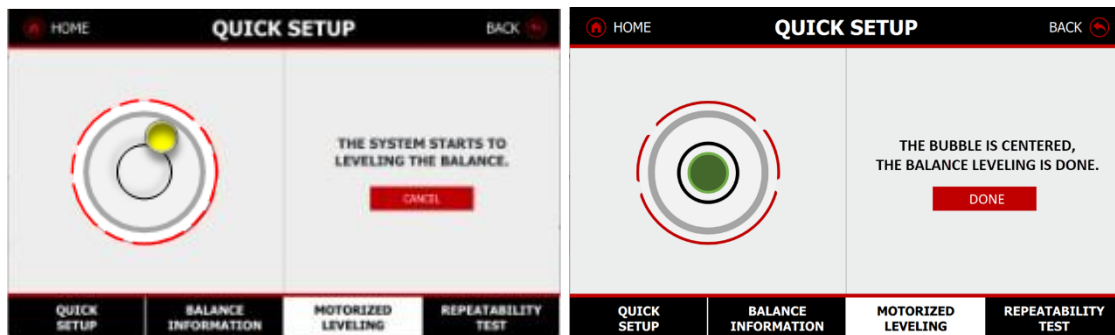

- After the level the balance, the digital bubble will display green at the upper right corner in the main screen
- If the leveling system fail to level the balance successfully, move the balance to a relatively flat surface. The warning notification will pop up

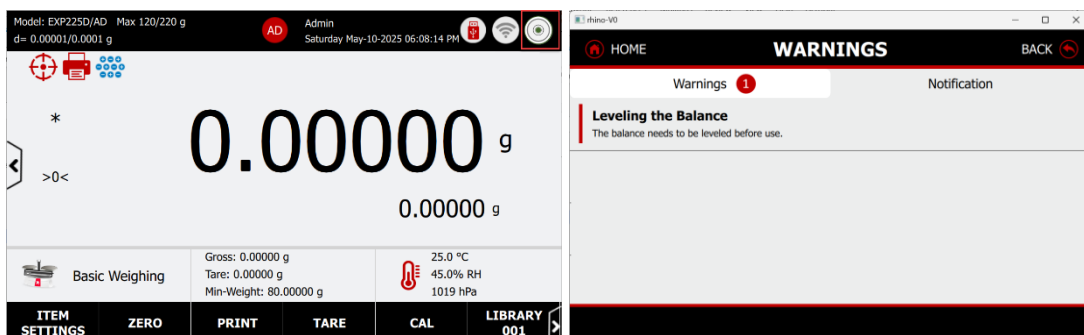

5.3.9 Repeatability Test

This function is designed to conduct the daily repeatability test of the balance. Users can perform a 10-time repeatability test using weights that are close to their daily working point.


Note: To determine the minimum weight in accordance with <USP41>, test environment must be equivalent to standard lab <USP1205> and the test weights must be certified and their values checked periodically.

- Press Repeatability Test tab and then follow the onscreen instruction.

Step 1: Set Number of Repeatability Tests, Test Weight ID, Test Weight Value and Test Weight Class in this menu.

The Number of Tests, Test Weight ID, Test Weight Value, and Test Weight Class can be set in accordance with the workshop's standard procedure.

- The default setting is 10 time in Number of Test

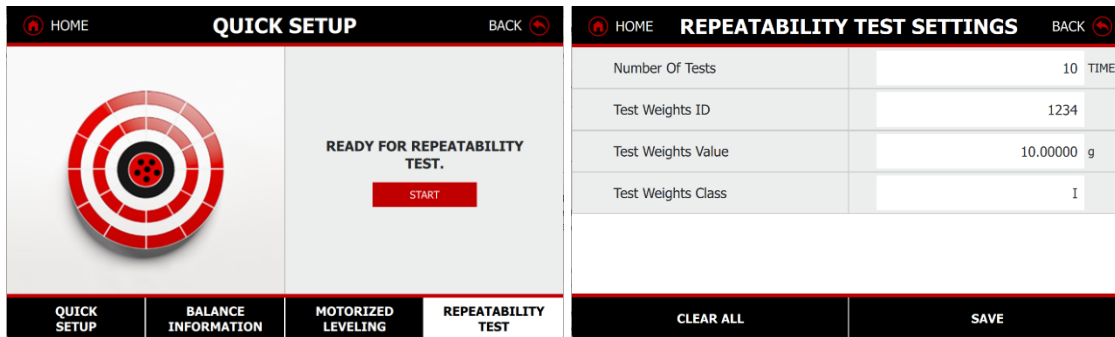


- Press Save button to store all the settings.

Step 2: Perform repeatability Test

- Place the balance in a stable environment and press Zero or Tare before starting.
- Press the Start button to place weights on the pan one by one.
- During the testing, do not tare or zero the pan, as the load will calculate all factors to produce accurate results.
- Example: After 10 tests with 10-gram weights, the balance will automatically display the repeatability test results.

Step 3: Review report

- Press the View History button to review the last repeatability test result.
- Press the Print button to print the repeatability test report and or export it to a PDF file when a USB flash drive is inserted.
- Press Min-weight to display the reference minimum weight.

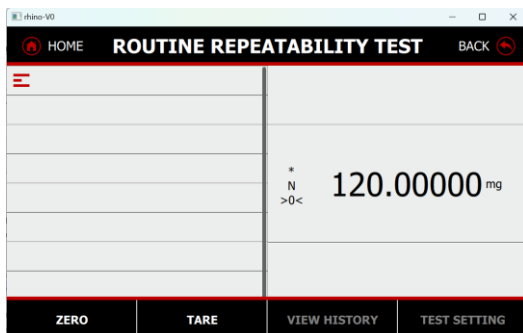


5.4 Calibration

Explorer Plus Balances build Automatic Calibration System. offer a choice of three calibration methods: Calibration Method: Automatic Calibration, Internal Calibration (User triggered), Span Calibration, and 3000 x Calibration History/ Log.

- Calibration is accomplished with the internal calibration mass. Internal calibration can be performed at any time, provided the balance has warmed up to operating temperature and is level.
- In the main screen, user can press CAL button to perform Internal Calibration.


Note: Do not disturb the balance during any calibration.

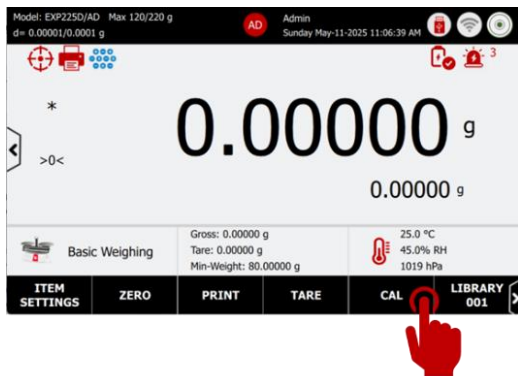

5.4.1 Calibration Settings

The Calibration Setting is used to configure the calibration method and the calibration points for Span Calibration.

- Internal Calibration
 - On/Off Internal Calibration function
 - Default setting: ON
- Automatic Calibration
 - When Internal Calibration is set ON, the balance performs a self calibration:
 - The environment temperature change of 1.5°C
 - Every 3 hours for Semi-Micro model, Every 11 hours for rest of the models
 - Default setting: ON
 - AutoCal Interval Hours: The automatic calibration is triggered by a time-based schedule.
 - Enable the Automatic Calibration function, user can defined the time.
 - The value can be set from 3-11 hours.
- Print Automatic Calibration
 - Enable the auto-print function. When the AutoCal is completed, the balance will send the calibration report to the PC or printer if connected.
- Span Calibration Points
 - The Span calibration uses two calibration points, one at zero load and the other at specified full load (span). For detailed calibration mass information please see the specification tables in section 9.
 - Refer to Table xxx for Span values and factory default span calibration points in the table xxx.
 - The user can defined the Span Calibration weights value by preset value and alternative weights.
 - ◆ Alternative Span Calibration Points: The weight value between 20% and 100% of the maximum balance capacity.

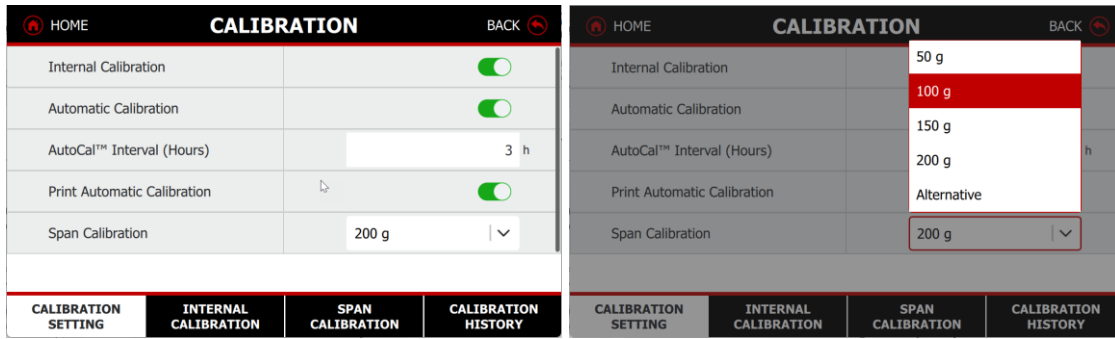


5.4.2 Internal Calibration

Internal calibration is performed using the built-in calibration weight. This process can be initiated at any time, as long as the balance has reached its operating temperature and balance is properly leveled.

Step 1: Start Calibration Process

- Press **Internal Calibration tab**, and press **Start**, the Balance will begin to calibrate. To cancel at any time, press **Cancel**.

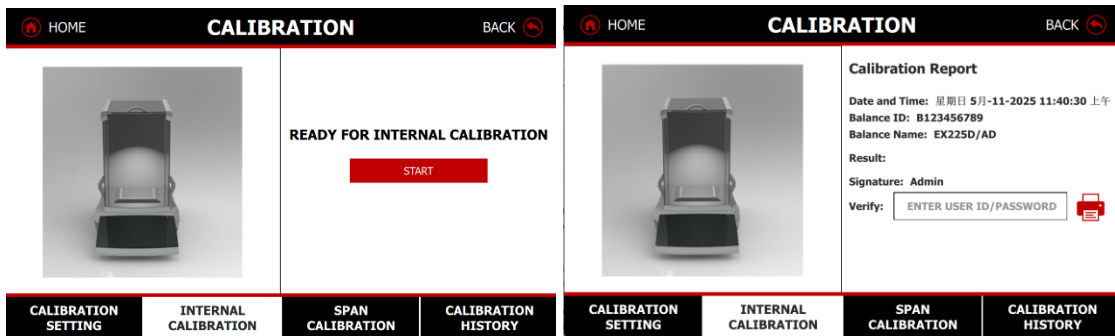

Step 2: Verify the calibration result

- After calibration, place the test mass on the pan and verify that the mass value now matches the displayed value. If not, repeat the procedure until the reading agrees with the test mass.

Step 3: Calibration Report

- When the calibration process is completed, the balance will generate a calibration report that includes the following information:
 - Date/Time: Week, Date and Time
 - Balance ID: Balance series number
 - Balance Name: Balance Model
 - Result: Successful or Failed
 - Signature: User ID (when the user management is set on)
 - Verified by: User can input the User ID/ Password or use the Fingerprint accessory to input the User information automatically. Fingerprint accessory order number is in the table XXXX
- Press the Print button to print the calibration report, and the results will be saved in the Calibration History for future reference.

5.4.3 Span Calibration

Step 1: Select Span Calibration weights value in the Calibration Setting

- Different capacity balances have different Span Calibration points. Users can choose the best Span Calibration points by using either the full load or half load of the balance's capacity.
- If the weights are certified with a tolerance value, users can input an **Alternative Weight Value** before starting the Span Calibration process.


Calibration Masses Table

| Model | Span Calibration Points | Weight Class | |
|---|--|------------------------------|--------------------|
| EXP125D; EXP125DM | 50g, 100g | ASTM Class 1 | OIML E2 |
| EXP125; EXP125M EXP225D; EXP225DM | 50g, 100g 100g, 150g, 200g | ASTM Class 1 ASTM Class 1 | OIML E2 OIML E2 |
| EXP225D/AD; EXP225DM/AD | 100g, 150g, 200g | ASTM Class 1 | OIML E2 |
| EXP225/AD; EXP225M/AD | 100g, 150g, 200g | ASTM Class 1 | OIML E2 |
| EXP124 | 50g, 100g | ASTM Class 1 | OIML E2 |
| EXP224 | 100g, 150g, 200g | ASTM Class 1 | OIML E2 |
| EXP324, M, N | 150g, 200g, 300g | ASTM Class 1 | OIML E2 |
| EXP223 | 100g, 150g, 200g | ASTM Class 1 | OIML E2 |
| EXP423 | 200g, 300g, 400g | ASTM Class 1 | OIML E2 |
| EXP623 | 300g, 400g, 500g , 600g | ASTM Class 1 | OIML E2 |
| EXP1203, M, N | 500g, 1000g | ASTM Class 1 | OIML E2 |
| EXP2202 | 1000g, 2000g | ASTM Class 1 | OIML E2 |
| EXP4202 | 2000g, 3000g, 4000g | ASTM Class 1 | OIML E2 |
| EXP6202 | 3000g, 4000g, 5000g , 6000g | ASTM Class 1 | OIML E2 |
| EXP10202, M, N | 5000g, 10,000g | ASTM Class 1 | OIML E2 |
| EXP6201 | 3000g, 4000g, 5000g , 6000g | ASTM Class 2 | OIML F1 |
| EXP10201 | 5000g, 10,000g | ASTM Class 1 | OIML E2 |
| EXP24001 | 10000g, 15000g, 20000g, 24000g | ASTM Class 1 | OIML F1 |
| EXP35001 | 10000g, 20000g, 30000g, 35000g | ASTM Class 1 | OIML F1 |
| EXP65001 | 10000g, 20000g, 30000g, 35000g | ASTM Class 1 | OIML F1 |

Step 1: Start Calibration Process

- Press **Span Calibration tab**, and press **Start**, the Balance will begin to calibrate. To cancel at any time, press **Cancel**.
- Follow the guide message to place the weights on the pan and wait until the balance is stable.
- **Step 2: Verify the calibration result**

After calibration, place the test mass on the pan and verify that the mass value now matches the displayed value. If not, repeat the procedure until the reading agrees with the test mass.

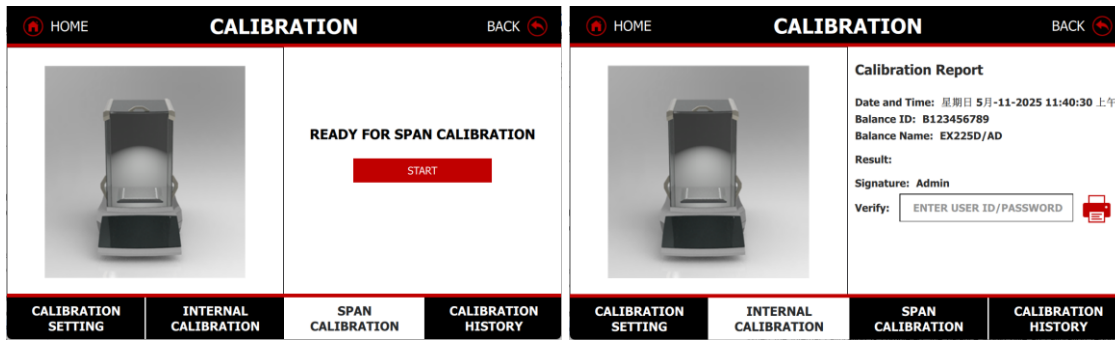

Step 3: Calibration Report


- When the calibration process is completed, the balance will generate a calibration report that includes the following information:
 - Date/Time: Week, Date and Time
 - Balance ID: Balance series number
 - Balance Name: Balance Model
 - Result: Successful or Failed
 - Signature: User ID (when the user management is set on)
 - Verified by: User can input the User ID/ Password or use the Fingerprint accessory to input the User information automatically. Fingerprint accessory order number is in the table XXXX
- Press the Print button to print the calibration report, and the results will be saved in the Calibration History for future reference.

5.4.4 Calibration History

The balance can store up to 3000 x Calibration History/ Log whenever the calibration is performed.

The total memory is displayed as xx/xx% in the upper right corner. At any time, users can export the calibration records to a PDF file when a USB flash drive is inserted.

Note: When the user activates the **System Log** function, the calibration history will be recorded as **Calibration Log** in the Service Menu.

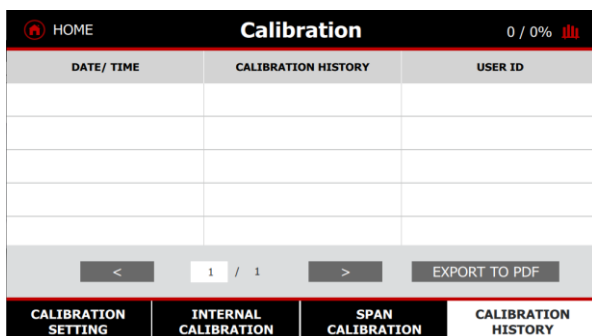


5.5 Balance Setup

Enter this sub-menu to customize Balance functionality

5.5.1 Language


Set the language for menu display and pop-up messages. The default setting is configured based on the country of sale. The 13 available languages are English, German, French, Spanish, Italian, Polish, Czech, Hungarian, Portuguese, Chinese, Japanese, Korean, and Turkish.

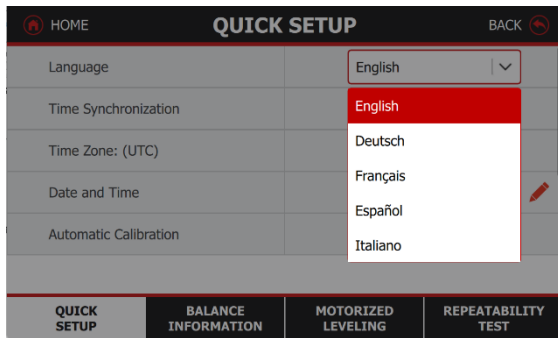

5.5.2 Time Synchronization/ Network Server


The new balance features Synchronization with Network (NTP function) which can use consistent time and data from the local network. The NTP system supports DNS domain name resolution.

- When you input the address of a public domain server, it will automatically convert it into an IP address. Here are some examples of NTP public time server address
 - “0.europe.pool.ntp.org” when the location is in Europe
 - “cn.ntp.org.cn” for China
 - “0.us.pool.ntp.org” for the US

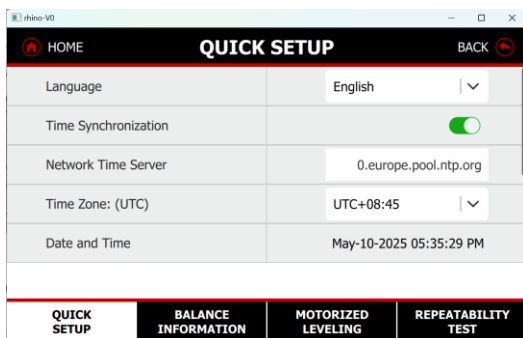

5.5.3 Date and Time

If you are offline use the balance, the new data/time setting fulfill your documentation needs with UTC time zone and customized the time format MMM-DD-YY HH-MM-SS and working weekdays setting.

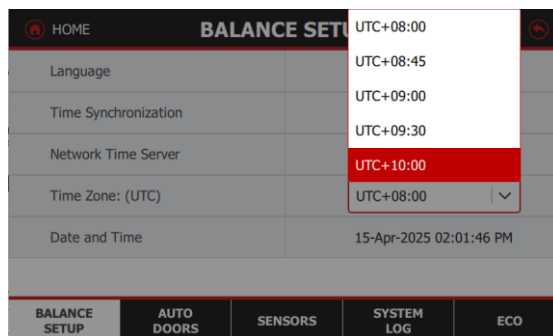

5.5.4 Balance Name

Set the balance identification. Alphanumeric settings up to 25 characters are available. The default setting is the model name of the balance

5.5.5 Change Password

- Change the password for the currently logged-in user. The Password Policy is defined in User Management:
- Option 1: Alphanumeric password containing 8 to 10 characters, combining letters and numbers.
- Option 2: Numeric password containing up to 25 characters.

See detail in section of User Management

Note: If you forget your password, please contact Ohaus or your local Ohaus dealer for assistance.

5.5.6 Fingerprint/ Fingerprint Setting

The fingerprint data is stored within the module itself. The OHAUS Balance cannot decrypt the original biometric data. A legal statement will be displayed on the terminal. The fingerprint data cannot be extracted from the module. This is because the module supplier defines the identification protocol for the fingerprint data. The OHAUS balance only manages the fingerprint ID derived from the original fingerprint data. The balance cannot transfer or copy the fingerprint ID from other balances.

2 Fingerprints per user, and total 100 x Fingerprint Identifications are available for user management control. Refer to the accessory list in Table xxxx.


Step 1: Connect terminal with the fingerprint accessory


Connect the cable to the terminal RJ11 port, and the fingerprint accessory will be ready for use.



Step 2: Enroll the first fingerprint


- Place your finger on the sensor and follow the prompts to add your fingerprint.
- Repeat this process five times. After five successful readings, the balance will enroll your fingerprint data.
- As a final step, place your finger on the sensor again to verify that the stored data matches your fingerprint.

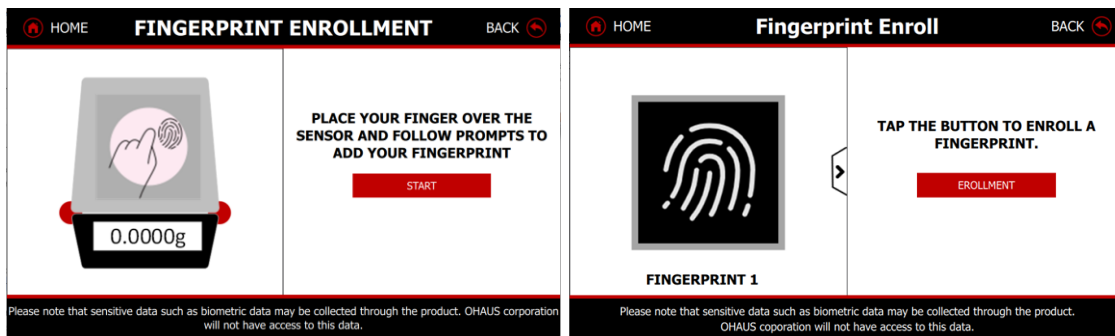

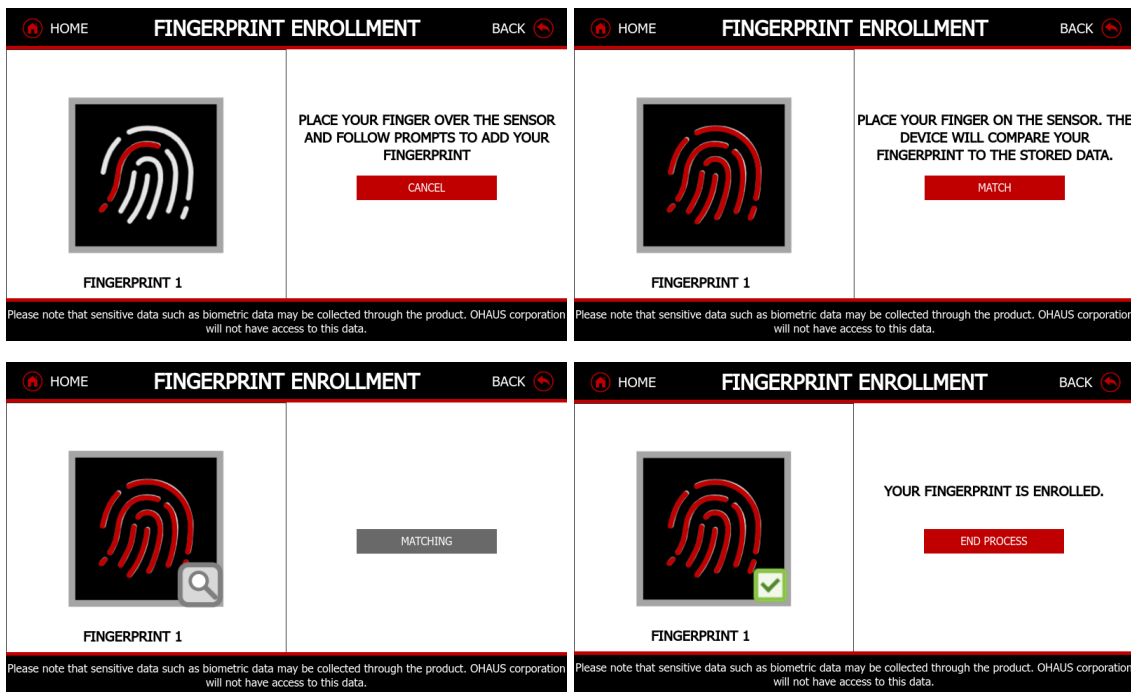

Step 2: Enroll the second fingerprint

- Tap the right arrow button to add new Fingerprint.
- The screen will show the FINGERPRINT 2.

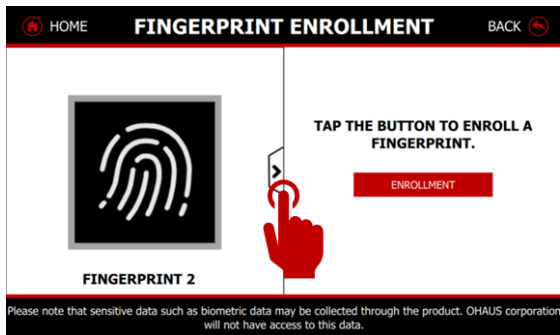


- Place your finger on the sensor and follow the prompts to add your fingerprint.
- Repeat this process five times. After five successful readings, the balance will enroll your fingerprint data.
- As a final step, place your finger on the sensor again to verify that the stored data matches your fingerprint.

Deactivated the Fingerprint Function

In user management, Administrator will turn on/off on user's fingerprint. See detail in the section of User Management xxx

Delete Fingerprint

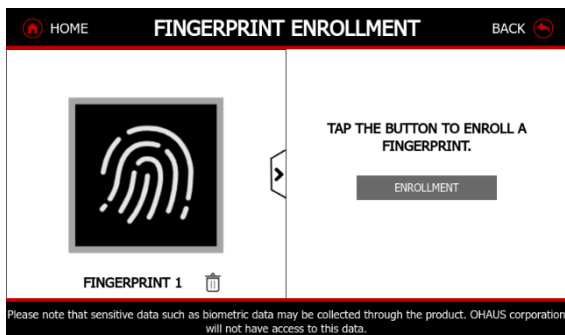

Legal Information

Legal Information will stay at the setting screen to comply with biometric authority regulation

Notes: The sensitive data such as biometric data may be collected through the product. OHAUS corporation will not have the access to this data.

5.5.7 Filter Level

Set the amount of signal filtering: 1, 2, 3, 4, **5** (default), 6, 7, 8, 9. The higher the number means slower the stabilization time with more stability.

Note: The setting is locked at the current setting when the Legal for Trade switch is set to the ON position.

5.5.8 Stability Indicator Range

Set the amount the reading can vary while the stability symbol remains on.

- 0.5 D = 0.5 graduations
- 1 D = 1 graduation
- 2D = 2 graduations
- 3D = 3 graduations
- D = 5 graduations

5.5.9 Auto Zero Tracking

Set the automatic zero tracking functionality.

- Off = disabled.
- 0.5 D = display maintains zero up to a drift of 0.5 graduation per second
- 1 D = display maintains zero up to a drift of 1 graduation per second.
- 2D = display maintains zero up to a drift of 3 graduations per second.

Note: When Legal for Trade is set to ON, the AZT setting is forced to 0.5 D.

The OFF setting is still available. The setting is locked at the current setting when the Legal for Trade switch is set to the ON position.

5.5.10 Gross Indicator

Set the symbol displayed for gross weights.

- Off = Disabled
- GROSS = the G symbol is displayed.
- BRUTTO = the B symbol is displayed.

The setting is locked at the current setting when the Legal for Trade switch is set to the ON position.

5.5.11 Graduation

Set the displayed readability of the balance.

- 1d = standard readability.
- 10d = readability is increased by a factor of 10.

For example, if the standard readability is 0.01g, selecting 10 Divisions will result in a displayed reading of 0.1g.

The setting is locked at the current setting when the Legal for Trade switch is set to the ON position.

5.5.12 Ionizer

Electrostatic charges on weighing containers and samples can result in false readings or instability as they create attractive and repulsive forces. The Explorer Plus Balance has built-in ionizers (Electrode brush) to neutralize the weighing chamber.

Set the ionizer's working time to either 5 or 10 seconds. Typically, 5 seconds is sufficient to eliminate static electricity from a sample. However, in cases where the air is extremely dry or the sample volume is large, a longer working time of 10 seconds may be necessary.

Note: For optimal results, it is recommended to place the lab spoon at the center of the ionizer electrodes before weighing.



5.5.13 Approved Mode

Use this menu to set the Legal for Trade mode.

- OFF = standard operation.
- ON = operation complies with Weights and Measures regulations.

Note:

- The security switch must be in the locked position to set Legal for Trade to ON.
- When Legal for Trade is set to ON, the menu settings are affected as follows:

- **Calibration Menu:**





- ◆ Automatic Calibration is set to ON and the menu is hidden. Internal Calibration is available, while all other functions are concealed.
- ◆ For EXP...N... models: Automatic Calibration will be locked at its current setting.

If you set Internal Calibration to be ON before you turn on Approved Mode, Internal Calibration menu will still be available. If you set Internal Calibration to be OFF before you turn on Approved Mode, Internal Calibration menu will be locked.

- **Balance Setup Menu:**

- ◆ Filter Level is locked at the current setting.
- ◆ Stability Indicator Range is locked at 1 Division.
- ◆ Auto Zero Tracking is limited to 0.5 Division and OFF.
- ◆ Auto Tare and Gross Indicator are locked at their current setting.
- ◆ Graduations is forced to 1 Division and the menu item is hidden.
- ◆ For EX...N...models, graduations will be locked at its current setting.


- Weighing Units menu: all units are locked at their current settings.
- **Communication Menu:**
 - ◆ Stable Weight Only is locked ON.
 - ◆ For EX...N...models, auto print mode selections are limited to OFF, On Stability, and Interval. Continuous is not available. Numeric Value Only is locked OFF.
- **Legal for Trade Switch:**

| | Explorer Plus Analytical and Precision Balance | Explorer Plus High-Capacity Balance |
|-------------|--|---|
| Position |  |  |
| Lock/Unlock |  |  |

5.6 Auto Door

The balances feature an anti-pinch, half open, and full open auto door function for safe and easy sample access.


- - If the user accidentally places their hands or other objects in the Draftshield doors while they are automatically opening, the door will stop to prevent hitting the object.
- - If the user accidentally places their hands or other objects at the back of the doors while they are automatically opening, the door will retract to prevent hitting the object.
- To achieve a faster stabilization time when weighing powder samples, the door can be set to either half-open or fully open.

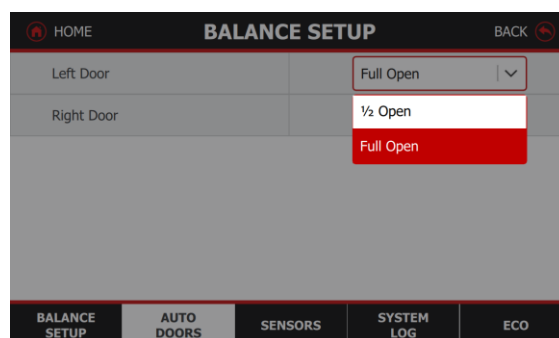

5.7 Sensor


There are 2 touchless sensors on the top of balance terminal.

When the user waves their hand over the sensor, the function will be activated, and a green light will flash for a second. The available settings are as follows: Print, Zero, Tare, Calibrate, Ionizer, Automatic Draft Shield Door, or Draft Shield Light. If the sensor is not activated, the sensor light will be red.

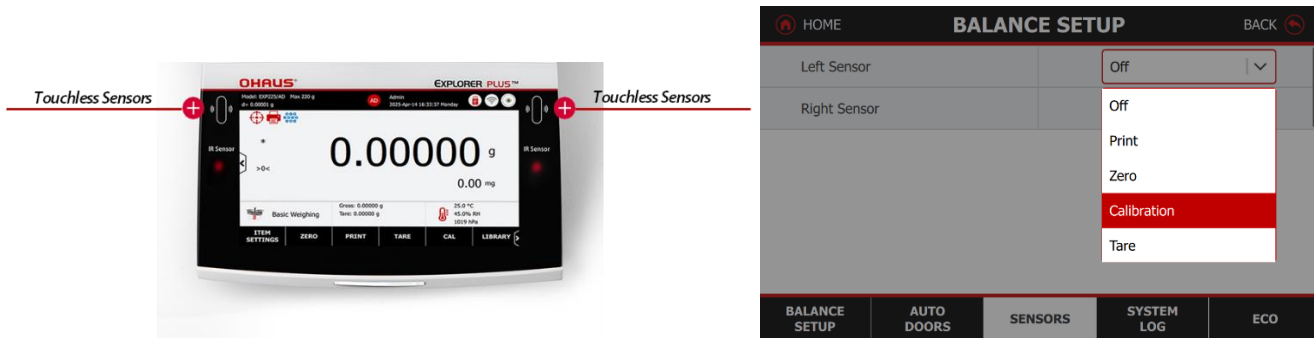

5.8 System Log

- All the changes made on balance setting will keep records in system log file. e.g. every printed data, date/time change, balance setting changes, perform calibration action, user log in/log out, user account create/edit/delete and etc. These system log files can be reviewed and exported as a PDF to a USB flash drive.
- Internal system log capacity of 100,000 entries records can be stored in balance memories. When the memories are full, a message will pop up for exporting the log files to a USB flash diver.
- System Log is saved in the Service Menu. See the detail function in section of Maintenance/ System Log.

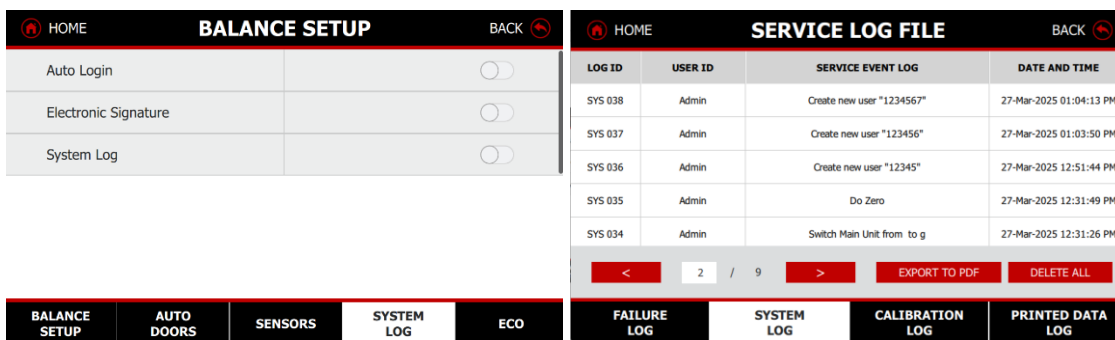

Auto Login

- Auto login is beneficial for Admin users who have not set a password, as no login is required after pressing the standby button.
- Other users cannot use this function when user management is enabled.

Electronic Signature

When the Electronic Signature feature is enabled, the User ID will be used in the printout field when the User ID content option is selected. For details, refer to the printout template in the Printing section.

System Log

Users can enable or disable the System Log as needed. When it is deactivated, changes to the balance will not be recorded in the balance memory.

5.9 ECO


OHAUS places energy-saving design as one of its top priorities, evident in all of its laboratory weighing products which feature power-saving capabilities. As a leading laboratory manufacturer, this small step forward demonstrates our commitment to these important causes.

To be specific, OHAUS electronic balance has designed a three-tiered energy-saving strategy for laboratory use.

- **Off** means the touchscreen display will operate with default settings.
- **Auto Dim** means the display will shift to a darker shade when there is inactivity.
- **Auto Standby** indicates the machine will log itself off automatically.
- **Sleep Mode** indicates that the display will be turned off at a predetermined time, while only retaining the base operation power.

5.9.1 Power Saving

- When the user selects Auto Dim and Auto Standby, the ECO time will automatically be set to 20 minutes. Additionally, the brightness, volume, and light functions will be locked to conserve power.
- When the user selects Sleep mode, the ECO time will automatically be set to 0 minutes. Additionally, the brightness, volume, and light functions will be locked to conserve power.

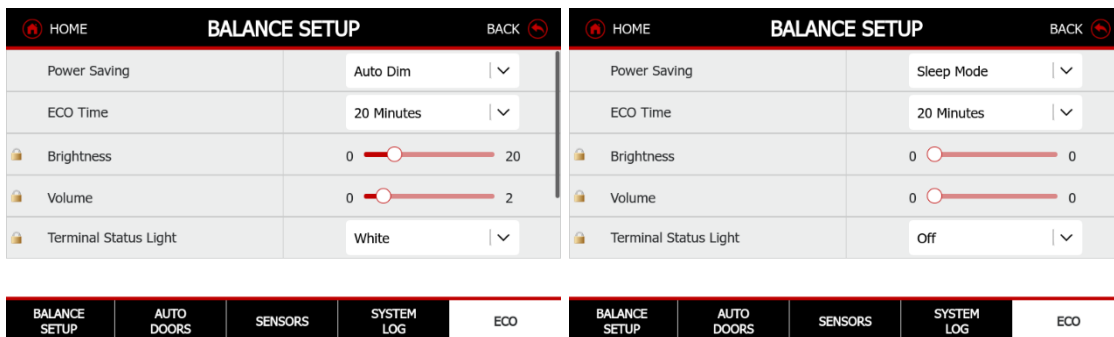

5.9.2 Brightness

Screen Brightness: Adjust the brightness from 0 to 100. Default value= 90

5.9.3 Volume

Adjust the volume from 0 to 10. Default value= 2

The Explorer Plus balance offers three sound options: a sound for click sound, successful actions, and an error sound.

5.9.4 Status lights

The lighting system can be customized into rainbow colors.

Terminal Status Lights: Users can change the light to different colors, such as Red, Pink, Yellow, Green, Cyan, Blue, and White.

Rainbow Draftshield Lights: Users can change the light to different colors, such as Red, Pink, Yellow, Green, Cyan, Blue, and White.

- **Note:** Overload, Accept, Underload color applies to application mode: check weighing and check counting



Sample spotlight for enhanced visibility and make a “ClearView” environment. Users can turn on and off this light. Spotlight light color: White.

5.10 User Management


The enhanced user management system enables administrators to create up to 200 user accounts, featuring four predefined roles. 4-level user management with an underdefined user group with access rights to the balance.

- Administrator (1)
- Supervisor (maximum of 20)
- 179 allocated to other roles like operator, log viewer, and group users

User Management: The User Name is generated from First Name and Last Name entries.

User ID is unique, can't change once its confirmed once you turn on Electronic Signature.

Group User: It's easy to set the group user with the same access to the balance setting

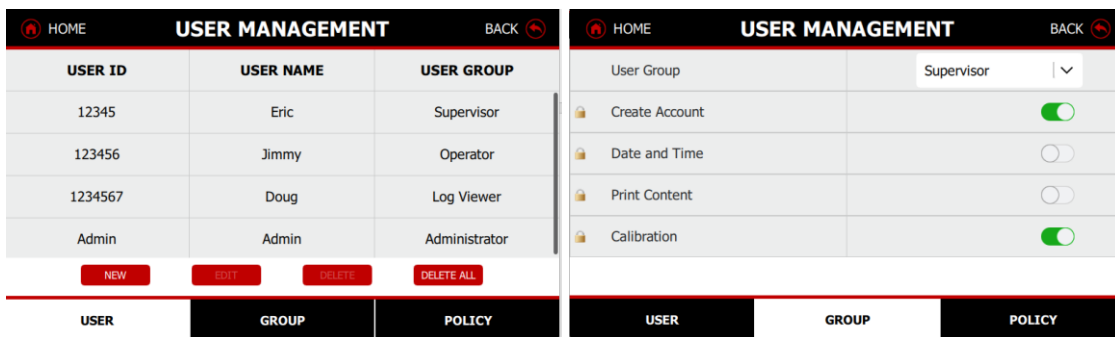

5.10.1 Create, Edit Delete a User

Admin can create, edit or delete Supervisor, Operator and Log Viewer, while Supervisor can create, edit or delete Operator. Operator and Log Viewer cannot access User management menu.

To create new users or edit or delete the current users

- Press to New button to add new user account, press Edit button to modify the user profiles
- Press delete or delete all user account by Administrator or Supervisor.

Note: Admin user cannot be deleted.

5.10.2 Preset User Role Accessibility


| Settings | Administrator | Supervisor | Operator | Log Viewer (Auditor) | Group 1 User |
|-------------------------------|---------------|----------------|------------------|----------------------|----------------|
| Number of Users | 1 person | 20 people | Total 179 people | | |
| Create Account/Delete Account | √ | Create Account | x | x | Create Account |
| Date and Time | | x | x | x | x |
| Print Content | | √ | x | x | x |
| Calibration Function | | √ | x | x | √ |
| Weighing Mode | | √ | √ | x | √ |
| Weighing Units | | √ | √ | x | √ |
| Communication Print Setting | | √ | √ | x | √ |

| | | | | | |
|--------------------|--------------------|--------------------|---|-----------|--------------------|
| Library | | √ | Allow to load existing library | x | √ |
| Factory Reset | | √ | x | x | √ |
| Routine Test | | √ | Allow to operate, not allow to change the Weights setting | x | √ |
| Balance Name | | x | x | x | x |
| IDs | | √ | v | x | √ |
| Setup a Min-Weight | | √ | x | x | √ |
| Service Mode | Print/Save/Read | Print/Save/Read | x | x | Print/Save/Read |
| System Log | Read/Export/Delete | Read/Export/Delete | x | Read only | Read/Export/Delete |
| Customized Hot Key | √ | √ | √ | x | √ |
| Graduation 1d/10d | √ | √ | x | x | x |

5.10.3 Group User Permissions

The Group User Permission function allows for the streamlined assignment of identical access permissions to multiple users. Four predefined groups are available: Administrator, Supervisor, Operator, and Log Viewer. In addition, a customizable Group 1 can be configured with unique access permissions.

All access options, except for Date and Time, Print Content, and Balance Name, are pre-configured and ready for use.

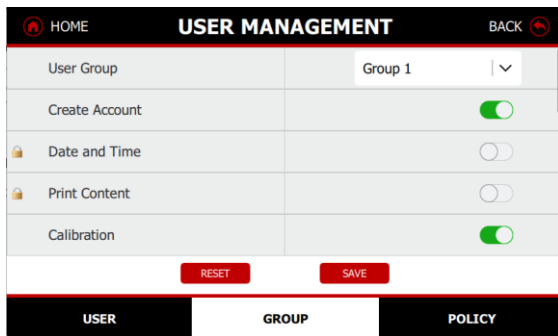

5.10.4 Password Policy

The system not only features alphanumeric password protection, but also incorporates a fingerprint login system. The fingerprint login system offers a convenient and secure way to log in, eliminating the risk of forgetting passwords or unauthorized password leaks.


These measures safeguard the weighing data and the defined weighing process configurations, preventing them from being altered or deleted.

Expiration Days:

To set when password will expire. There are three selections, 30 days, 60 days and 90 days.

Password Requirement:

- Option 1: Alphanumeric password containing 8 to 10 characters, combining letters and numbers.
- Option 2: Numeric password containing up to 25 characters.


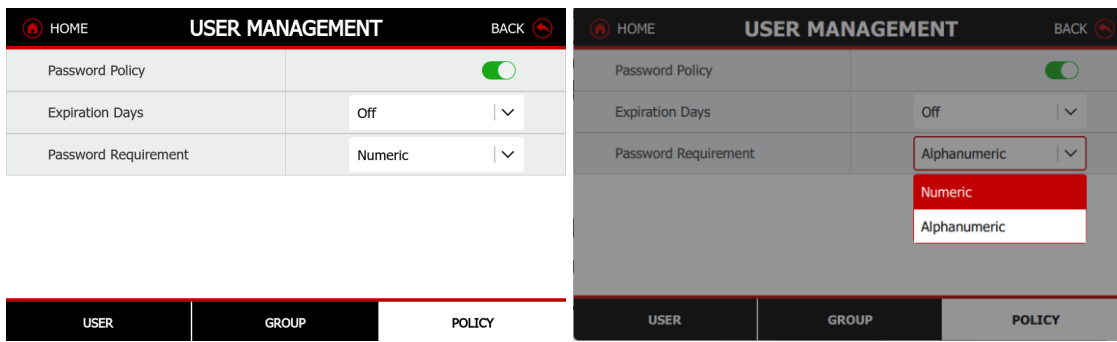


Note: If you forget your password, please contact Ohaus or your local Ohaus dealer for assistance.

5.11 Application Modes


The balance can be configured to operate in various Application modes.

Note: If an application mode does not appear in the list of available Applications, it must be enabled in the **Application Menu** touch Menu, then select Application Modes. The full list of modes appears. Touch the one you want to enable.

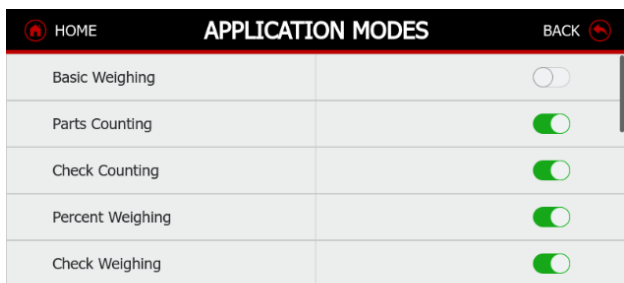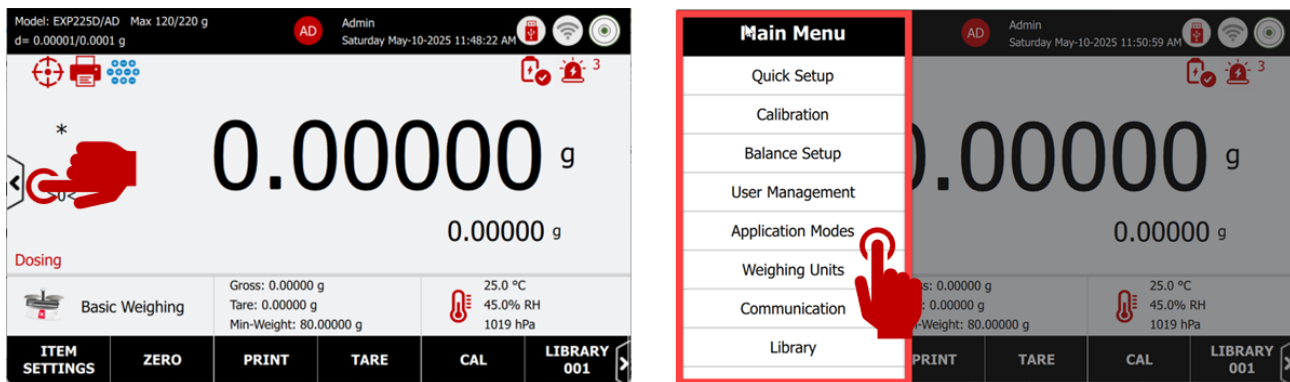

5.12 Weighing Units

The Explorer Plus balance can be configured to measure in a variety of weighing units, including three (2) custom units. In the main screen, user can click “g” to switch an alternative weighing unit. The default unit is gram. In this menu, you can turn on/off the weighing unit according to the sample application.

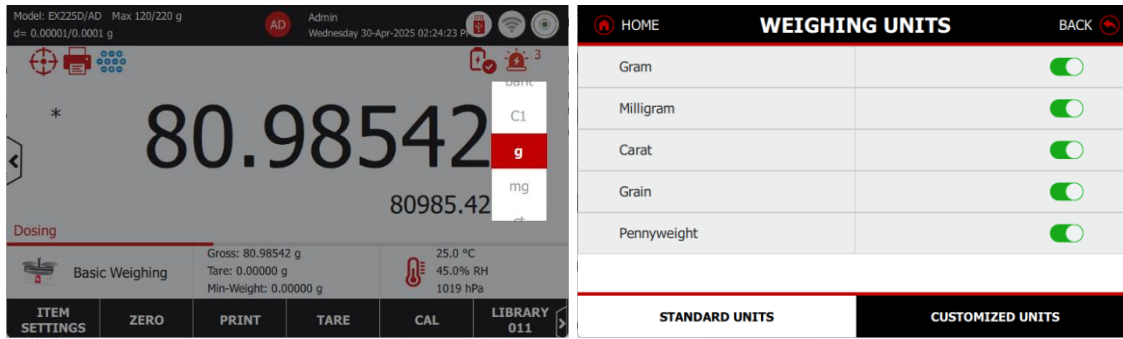


Customized Units

Use each Customized Unit to display weight in an alternative unit of measure. The custom unit is defined using a conversion factor, where the conversion factor is the number of custom units per kilogram expressed in scientific notation (Factor x 10^Exponent). For example: To display weight in troy ounces (32.15075 troy ounces per kilogram) enter a Factor of 0.321508 and an Exponent of 2. Set the status.

- OFF = disabled
- ON = enabled

When Customized Unit is set to ON, the Unit Name Factor, Exponent and Least Significant Digit must be set.

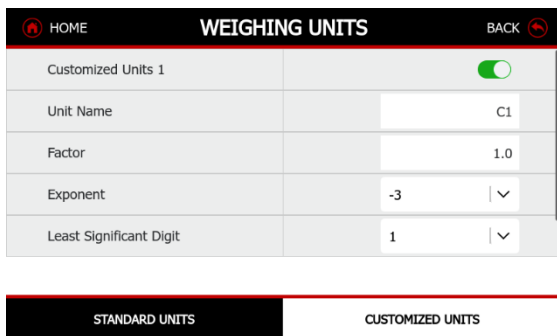

Factor: Set the conversion factor using the numeric keypad.

Settings of 0.00001 to 1.9999999 are available. The default setting is 1.000000

Exponent: Set the factor multiplier.

- 0 = multiply the Factor by 1 (1x10⁰)
- 1 = multiply the Factor by 10 (1x10¹)
- = multiply the Factor by 100 (1x10²)
- = multiply the Factor by 1000 (1x10³)
- - 3 = divide the Factor by 1000 (1x10⁻³)
- - 2 = divide the Factor by 100 (1x10⁻²)
- - 1 = divide the Factor by 10 (1x10⁻¹)


Least Significant Digit: Set the graduation.

Settings of 0.00001, 0.00002, 0.00005, 0.0001, 0.0002, 0.0005, 0.001, 0.002, 0.005, 0.01, 0.02, 0.05, 0.1, 0.2, 0.5, 1, 2, 5, 10, 20, 50, 100, 200, 500 and 1000 are available.

Note: Least Significant Digit setting selections are dependent on the Factor and Exponent settings. Selections are limited.

5.13 Communication

The enhanced communication options include 2 USB Type-A ports, 1 USB Type-C port, 1 Ethernet LAN, 1 RS232, 1 RJ11, and optional Bluetooth and WiFi functionalities. Additionally, the Explorer Plus balance supports HID (Human Interface Device) connection to a computer without requiring drivers. This enables users to use a mouse or keyboard to enter information such as User Name, USER ID, sample name, batch name, and other input details. The input field supports English and French characters, as well as numbers and symbols.

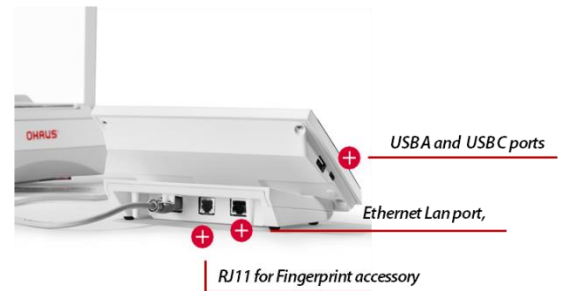

Enter this menu to define external communication methods and to set printing parameters. Data may be output to either a Printer, PC and or Fingerprint accessory.

•

| | |
|-------------------|---|
| USB Type A | USB Type A – Connect to USB flash driver, RFID Reader, Barcode Scanner and Wi-Fi & Bluetooth Dongle |
| USB Type B | Connect Balance to PC |
| USB Type C – | Connect Balance to PC |
| Ethernet Lan port | Connect Balance to PC |
| RS232 | Connect Balance to Printer or PC |
| RJ11 | Connect Balance to Fingerprint accessory |

5.13.1 RS232

The RS232 port can connect to various peripheral devices such as printers, PCs, and label printers, each with different output formats. Ensure you select the appropriate peripheral device before transferring data.

• The connection settings for Printer and PC are as follows. Factory default settings are indicated in bold.

- **Baud Rate** Set the baud rate (bits per second).
 - ◆ 1200 = 1200 bps, 2400 = 2400 bps, 4800 = 4800 bps
 - ◆ **9600 = 9600 bps**
 - ◆ 19200 = 19200 bps, 38400 = 38400 bps
- **Transmission** Set the data bits, stop bit, and parity.
 - ◆ 7 EVEN 1 = 7 data bits, even parity, 7 ODD 1 = 7 data bits, odd parity
 - ◆ 7 NONE 1 = 7 data bits, no parity
 - ◆ **8 NONE 1 = 8 data bits, no parity**

- ◆ 7 EVEN 2 = 7 data bits, even parity, 7 ODD 2 = 7 data bits, odd parity, 7 NONE 2 = 7 data bits, no parity, 8 NONE 2 = 8 data bits, no parity


■ **Handshake Set the flow control method.**

- ◆ NONE = no handshaking
- ◆ **XON-XOFF = XON/XOFF handshaking**
- ◆ HARDWARE = hardware handshaking

■ **12V Output**

- ◆ Enable the function to allow the RS232 port to output 5V, which supports the Bluetooth Adaptor.
- ◆ Example: When users connect an SF40A/BT printer to the balance, they need to enable this function.


Note: The purpose of this switch is to prevent damage to external devices from the 12V output.

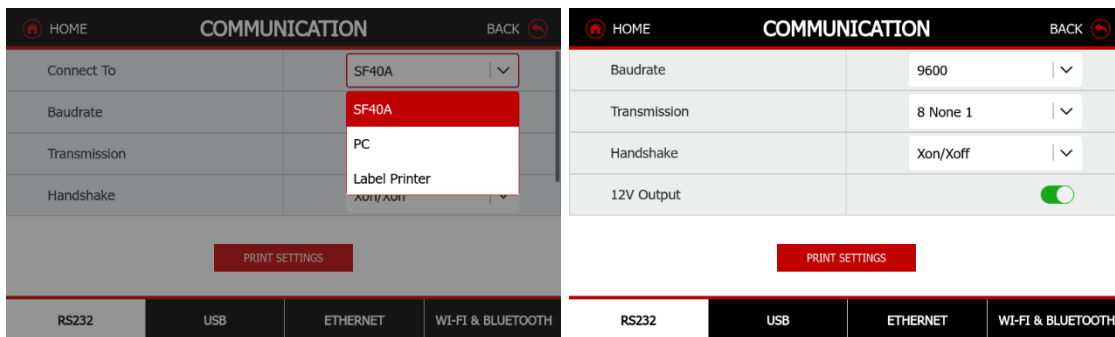

- The connection settings for Label Printer are as follows.


All templates including the default one can be edited via OHAUS Label Designer. Please go to the website below to download the software. For how to use the software, please contact an authorized dealer to obtain the software's instruction. Overall, 5 label templates can be stored in the balance.

- When user select connect to Label Printing, the Label Template menu appears with default selection, "100mm x 75mm".
 - ◆ This function is compatible with any label printer using the ZPLII programming language. We recommend Zebra GC420t label printer.
 - ◆ **#1-100mm x 75mm**, which is designed based on label size "100mm x 75mm". This template includes all basic information "Data/Time", "Sample ID", "Sample Name", "Batch ID", "Result", "Gross Weight", "Tare Weight", "Net Weight", and "User Name".
 - ◆ **#2-100mm x 10mm**, which is designed for small label size "100mm x 10mm" with weighing result only.
- To select another label template, touch the other numbers. Templates #3 - #5 are empty templates in Explorer until they are edited by OHAUS label Designer and written to the balance.

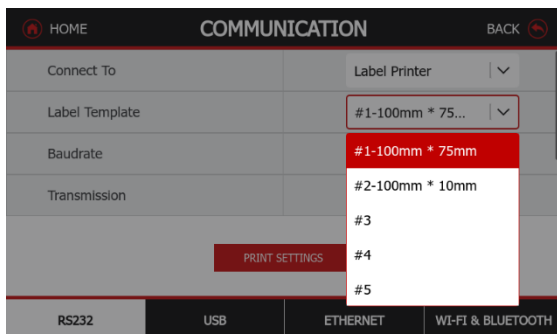

5.13.2 RS232 (DB9) Pin Connections


- Pin 2: Balance transmit line (TxD)
- Pin 3: Balance receive line (RxD)
- Pin 5: Ground signal (GND)
- Pin 7: Clear to send (hardware handshake) (CTS)
- Pin 8: Request to send (hardware handshake) (RTS)

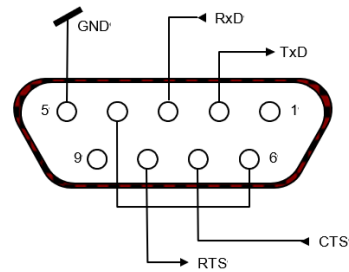

5.13.3 USB

The USB setting can individually setup the USB port to different setup appropriate peripheral device before transferring data.


- Option 1: USB Type B connection with the balance to PC or Transfer Data Direct to Microsoft Excel.
- Option 1: USB Type C connection with the balance to PC or Transfer Data Direct to Microsoft Excel.
- The format can be set in column or cell, see the detail print format in section of Print Setting xxx

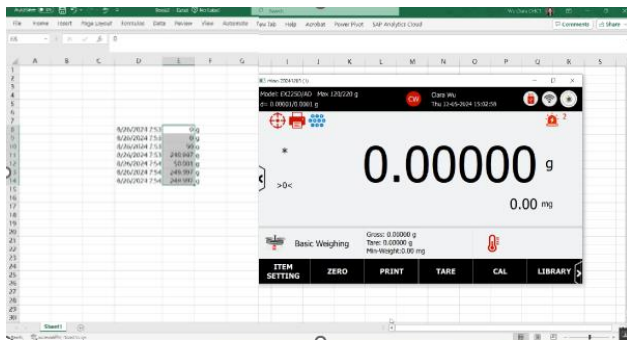

5.13.4 Ethernet

Using the Ethernet Port to connect to a Local Area Network. After successful connect ethernet cable with balance, user can setup the Ethernet options on the Communication Menu.

- **DHCP:** When connecting to a local area network (LAN) or a wireless network and the IP address is unknown, user can use DHCP to automatically obtain an IP address. Enable DHCP: the balance will automatically obtain IP, other ethernet setting will be locked.
- If you prefer to use a static IP address for the connection, you should disable DHCP and manually enter the IP address.

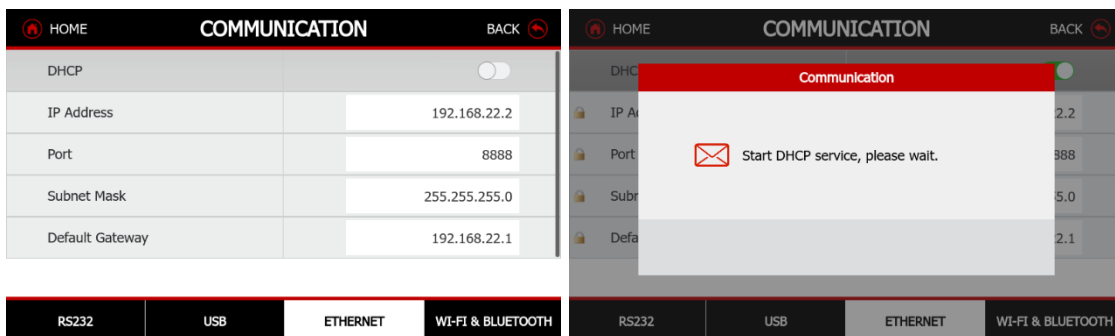


5.13.5 WI-FI & Bluetooth

The Explorer Plus balance use LMB842 USB dongle to connect BT and Wi-Fi function.

- The LM842 USB Dongle complies with Bluetooth® 5.0 (Dual Mode) and IEEE 802.11ac WiFi Standards of operation.
- See detail compliance and specification of the LM842 USB Dongle in section of Accessory Specification xxx


Bluetooth Connection

- The user needs to enter the Bluetooth name to pair with another PC's Bluetooth.
- Once the balance connects to a Bluetooth device, an icon will appear in the upper right corner of the main screen.

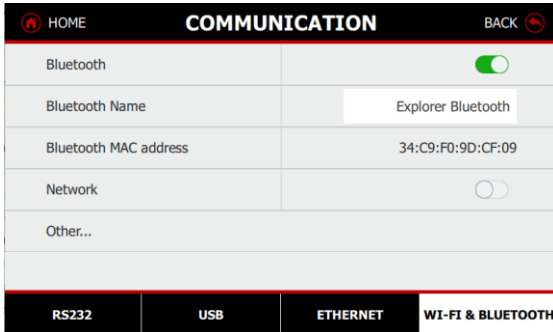

Wi-Fi Connection

- First, the user needs to enable the Network function. The balance will then search for available networks.
- Enter the password to pair and connect to the Wi-Fi.
- If the network is not visible, the user can press the "Other..." button to manually add the network.
- Once the balance connects to a Wi-Fi, an icon will appear in the upper right corner of the main screen.

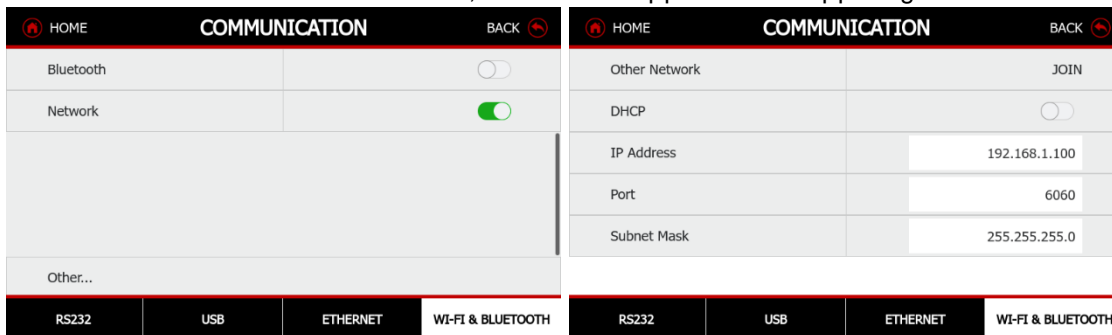

6 Print Setting

Users can configure this function individually based on their specific application requirements. The function allows users to select from the following settings:

- **Print Content:** Customize the content to be printed, including all desired results and titles along with the weighing value.
- **Connect to Printer:** Define the data to be sent to the printer.
- **Connect to PC:** Specify the data to be transmitted to the PC.
- **Data to Excel:** Format the layout of the data that is sent directly to Microsoft Excel.
- **Save to USB:** Determine how data is saved to a USB flash device, such as the data format and whether the process is manual or automatic.

Note: When the balance is turned on in the Legal For Trade mode, the print settings will immediately be changed to comply with relevant Weights and Measures Regulation, such as those of the International Organization of Legal Metrology (OIML), the National Type Evaluation Program (NTEP) and other local approves.


See detail in the section of Approved Mode in xxxx.

6.1 Print Content

Access this sub-menu to specify the content of the printed data. Users can toggle the content on or off. The Print Review function will directly provide a layout template based on the selected options.

The available print contents:

- Select All
- Date and Time, Balance ID and Balance name
- User ID, First Name/ Last Name
- Project Name
- Application Mode
- Sample Name, Sample ID
- Batch ID, Lot ID
- Customized IDs
- Result
- Gross, Net, and Tare
- Signature Line
- Verified Line
- Feeds Lines (1, 4, 10)

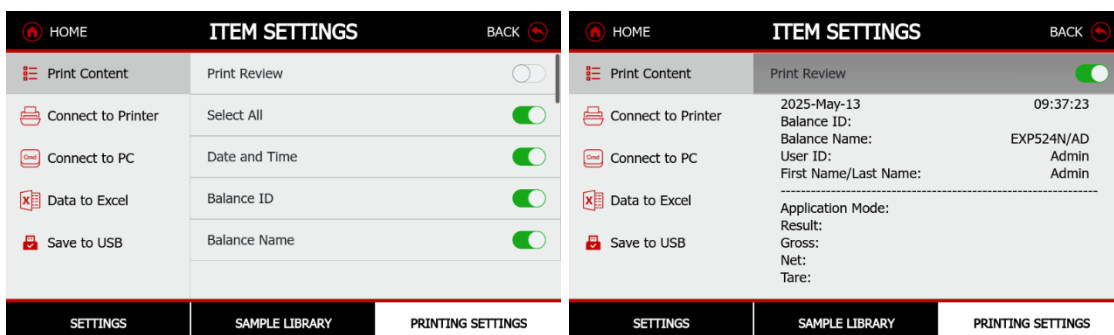

6.2 Connect to Printer

Numeric Only

On = Print only the numeric weight value.

Off = Print all the weighing value that enabled in the Print Content.

Single Header Only

When user turn it On, Header line will be printed only once each day. The header line includes the following elements:

- Date and Time, Balance ID, and Balance Name
- Project Name
- Application Mode
- Sample Name, Sample ID
- Batch ID, Lot ID
- Customized IDs

Manual Print

Stable only = Print only the stable value

All values = Print all values

Auto Print

Off = disabled

On Stability = printing occurs each time the stability criteria are met.

Interval = printing occurs at the defined time interval (1 to 3600 seconds)

Continuous = printing occurs continuously.

When ON STABLE is selected, set the conditions for printing.

- LOAD = Prints when the displayed load is stable.
- LOAD ZERO = Prints when the displayed load or zero reading is stable.
- When Approved Mode is On, users can only print stable only values

6.3 Connect to PC

Output Format

Users can choose the output format based on various output strings.

- OHAUS
- SICS = MT-SICS
- ST =ST-SICS (10)

Manual Print

Stable only = Print only the stable value

All values = Print all values

Auto Print

Off = disabled

On Stability = printing occurs each time the stability criteria are met.

Interval = printing occurs at the defined time interval (1 to 3600 seconds)

Continuous = printing occurs continuously.

When ON STABLE is selected, set the conditions for printing.

- LOAD = Prints when the displayed load is stable.
- LOAD ZERO = Prints when the displayed load or zero reading is stable.
- When Approved Mode is On, users can only print stable only values

6.4 Data to Excel

The balance supports HID (Human Interface Device) connection to a computer without the need for drivers.

Manual Print

Stable only = Print only the stable value

All values = Print all values

Auto Print

Off = disabled

On Stability = printing occurs each time the stability criteria are met.

Interval = printing occurs at the defined time interval (1 to 3600 seconds)

Continuous = printing occurs continuously.

When ON STABLE is selected, set the conditions for printing.

- LOAD = Prints when the displayed load is stable.
- LOAD ZERO = Prints when the displayed load or zero reading is stable.
- When Approved Mode is On, users can only print stable only values


Text to

Direct to Excel Function:

- Column: Print all data into a single column in Microsoft Excel.
- Cell: Print all data into a single cell in Microsoft Excel.

6.5 Save to USB

- Off = Disable save to USB
- Manual = Save to USB manually
- Auto = Auto save to USB
 - Once the user has set the auto-print interval time, the data will be sent to the USB flash drive at the specified intervals.
 - The CSV format is not allow to use in Auto save mode
- Format
 - PDF
 - CSV

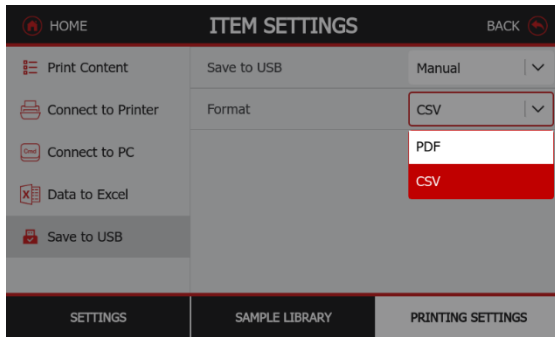

6.5.1 Application Print Out Template

| Basic Weighing | | Parts Counting | | Check Counting | |
|-----------------------|----------------|-----------------------|----------------|-----------------------|----------------|
| 12-24-2024 | 09:31:46 | 12-24-2024 | 10:13:34 | 12-24-2024 | 10:17:42 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN058 | Project Name: | PN070 | Project Name: | PN081 |
| ----- | | ----- | | ----- | |
| Application Name: | Basic Weighing | Application Name: | Parts Counting | Application Name: | Check Counting |
| Sample Name: | SN058 | Sample Name: | SN070 | Sample Name: | SN081 |
| Sample ID: | S058 | Sample ID: | S070 | Sample ID: | S081 |
| Batch ID: | B058 | Batch ID: | B070 | Batch ID: | B081 |
| Lot ID: | L058 | Lot ID: | L070 | Lot ID: | L081 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | Customer ID2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 3.5275 | Result: | 25 PCS | Result: | 74 PCS |
| oz N | | Gross: | 78.48 g | Status: | Accept |
| Gross: | 6.2960 | Net: | 78.48 g N | Gross: | 178.49 g |
| oz | | Tare: | 0.00 g T | Net: | 178.49 g N |
| Net: | 3.5275 oz N | Library: | Library 070 | Tare: | 0.00 g T |
| Tare: | 2.7690 oz T | APW: | 3.124 g | Library: | Library 081 |
| Library: | Library 058 | Samples: | 10 PCS | APW: | 2.412 g |
| Signature: _____ | | Signature: _____ | | Samples: | 81 PCS |
| Verify By: _____ | | Verify By: _____ | | Over: | 254 PCS |
| | | | | Under: | 51 PCS |
| | | | | Signature: _____ | |
| | | | | Verify By: _____ | |

| Percent Weighing | | Check Weighing | | Dynamic | |
|-----------------------|------------------|-----------------------|----------------|-----------------------|------------------|
| 12-26-2024 | 13:30:04 | 12-26-2024 | 13:33:25 | 12-26-2024 | 13:36:21 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID : | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN010 | Project Name: | PN028 | Project Name: | PN055 |
| ----- | | ----- | | ----- | |
| Application Name: | Percent Weighing | Application Name: | Check Weighing | Application Name: | Dynamic Weighing |
| Sample Name: | SN010 | Sample Name: | SN028 | Sample Name: | SN055 |
| Sample ID: | S010 | Sample ID: | S028 | Sample ID: | S055 |
| Batch ID: | B010 | Batch ID: | B028 | Batch ID: | B055 |
| Lot ID: | L010 | Lot ID: | L028 | Lot ID: | L055 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | C2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 361.69 % | Result: | Accept | Result: | 367.00 g |
| Gross: | 198.93 g | Gross: | 198.92 g | Gross: | 198.93 g |
| Net: | 198.93 g N | Net: | 198.92 g N | Net: | 198.93 g N |
| Tare: | 0.00 g T | Tare: | 0.00 g T | Tare: | 0.00 g T |
| Library: | Library 010 | Library: | Library 028 | Library: | Library 055 |
| Reference Weight: | 550.00 g | Over Limit: | 214.25 g | Averaging Time: | 3 seconds |
| Reference Factor: | 10.0 % | Under Limit: | 12.54 g | Signature: | _____ |
| Difference Weight: | 143.93 g | Signature: | _____ | Verify By: | _____ |
| Difference Factor: | 261.69 % | Verify By: | _____ | | |
| Signature: | _____ | | | | |
| Verify By: | _____ | | | | |

| Totalization | | Formulation | | Differential | |
|---------------------------|---------------------|-----------------------|---------------------|-----------------------|----------------------|
| 12-26-2024 | 15:11:42 | 12-26-2024 | 16:30:04 | 2025-Apr-08 | 19:58:46 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN072 | Project Name: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Name: | Totalization | Application Name: | Formulation | Application Mode: | Differential |
| Sample Name: | SN072 | Result: | 199.09 g N | Result: | 3.89[3] g N |
| Sample ID: | S072 | Gross: | 299.09 g | Gross: | 18.42[0] g |
| Batch ID: | B072 | Net: | 199.09 g N | Net: | 3.89[3] g N |
| Lot ID: | L072 | Tare: | 100.00 g T | Tare: | 14.52[7] g T |
| Customer ID1: | C1 | Recipe: | Recipe 001 | Start Time: | 2025-Apr-08 19:57:07 |
| Customer ID2: | C2 | Item Name | Item Weight | End Time: | 2025-Apr-08 19:58:41 |
| Customer ID3: | C3 | Item 1 | 100.00 g | Reciprocal Proportion | On |
| Customer ID4: | C4 | Item 2 | 199.08 g | Absolute Value | On |
| Customer ID5: | C5 | Start Time: | 12-26-2024 16:29:44 | Item Name: | ITEM1 |
| Customer ID6: | C6 | End Time: | 12-26-2024 16:30:01 | Initial Weight: | 0.766 g |
| Customer ID7: | C7 | -----Sample Data----- | | Final Weight: | 0.760 g |
| Customer ID8: | C8 | Item Name: | Item 1 | Difference Weight: | 0.006 g |
| Customer ID9: | C9 | Target Weight: | 100.00 g | Different Percent: | 0.8 % |
| Customer ID10: | C10 | Actual Weight: | 100.00 g | Item Name: | ITEM2 |
| Result: | 278.33 g | Difference: | 0.00 % | Initial Weight: | 1.528 g |
| Gross: | 0.00 g | Item Name: | Item 2 | Final Weight: | 1.534 g |
| Net: | 0.00 g N | Target Weight: | 199.08 g | Difference Weight: | 0.006 g |
| Tare: | 0.00 g T | Actual Weight: | 199.08 g | Different Percent: | 0.4 % |
| Library: | Library 072 | Difference: | 0.00 % | Item Name: | ITEM3 |
| Start Time: | 12-26-2024 15:08:29 | Total Weight: | 299.08 g | Initial Weight: | 2.292 g |
| End Time: | 12-26-2024 15:11:42 | Signature: _____ | | Final Weight: | 2.325 g |
| -----Sample Data (g)----- | | Verify By: _____ | | Difference Weight: | 0.033 g |
| '1: | 198.93 | | | Different Percent: | 1.4 % |
| '2: | 19.85 | | | Item Name: | ITEM4 |
| '3: | 19.86 | | | Initial Weight: | 3.094 g |
| '4: | 19.87 | | | Final Weight: | 3.110 g |
| '5: | 19.88 | | | Difference Weight: | 0.016 g |
| Samples: | 5 | | | Different Percent: | 0.5 % |
| Average: | 55.67 g | | | Item Name: | ITEM5 |
| Maximum: | 198.93 g | | | Initial Weight: | 3.871 g |
| Minimum: | 19.85 g | | | Final Weight: | 3.893 g |
| Range: | 179.08 g | | | Difference Weight: | 0.022 g |
| Standard Deviation: | 71.63 g | | | Different Percent: | 0.6 % |
| Relative Deviation(std%): | 128.68 % | | | Signature: _____ | |
| Signature: _____ | | | | Verified By: _____ | |
| Verify By: _____ | | | | | |

| Density Determination | | Peak Hold | | Pipette Adjustment | |
|-----------------------|---------------------------|-----------------------|-----------|---------------------------------------|---------------------------|
| 2025-Apr-08 | 20:01:25 | 2025-Apr-08 | 20:25:41 | 025-Apr-08 | 20:11:58 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| ----- | | ----- | | ----- | |
| Application Mode: | Density Determination | Application Mode: | Peak Hold | Application Mode: | Pipette Adjustment |
| Result: | 7.435 g/cm ³ | Result: | 355.512 g | Result: | Pass |
| Gross: | 19.44[1] g | Gross: | 0.000 g | Gross: | 24.76[4] g ? |
| Net: | 4.91[4] g N | Net: | 0.000 g N | Net: | 1.01[1] g ? N |
| Tare: | 14.52[7] g T | Tare: | 0.000 g T | Tare: | 23.75[3] g T |
| Weight in air.: | 5.67[4] g | On Stability: | No | Start Time: | 2025-Apr-08 20:09:03 |
| Weight in liquid.: | 4.91[3] g | Signature: _____ | | End Time: | 2025-Apr-08 20:11:52 |
| Auxiliary liquid.: | Water | Verified By: _____ | | Liquid Density: | 0.99823 g/cm ³ |
| Liquid Density: | 0.99823 g/cm ³ | | | Pressure: | 1.0 ATM |
| Water Temp.: | 20.0 °C | | | Pipette Test Method: | Method 001 |
| Porous Material: | Off | | | Nominal Inaccuracy Imprecision Number | 1000 ul 5.00 % 5.00 % 3 |
| Signature: _____ | | | | Pipette Test Results | |
| Verified By: _____ | | | | ----- | |
| | | | | Pipette Name: | pip001 |
| | | | | Pipette Number: | |
| | | | | Nominal Volume: | 1000 ul Pass |
| | | | | -----Sample Data----- | |
| | | | | 1 | 1023 ul 1.021 g |
| | | | | 2 | 1032 ul 1.030 g |
| | | | | 3 | 1016 ul 1.014 g |
| | | | | Inaccuracy | |
| | | | | E%: | 2.64 % |
| | | | | E% Limit: | 5.00 % |
| | | | | Average: | 1026 ul |
| | | | | Imprecision | |
| | | | | CV%: | 0.97 % |
| | | | | CV% Limit: | 5.00 % |
| | | | | Standard Deviation: | 10 ul |
| | | | | > +2S: | 0.00 % , 0 |
| | | | | > +1S: | 0.00 % , 0 |
| | | | | +1S > Mean > -1S: | 0.00 % , 0 |
| | | | | < -1S: | 0.00 % , 0 |
| | | | | < -2S: | 100.00 % , 10 |
| | | | | Signature: _____ | |
| | | | | Verified By: _____ | |

| SQC | | Fill Weight Variation | | Flow Rate Control | |
|------------------------------------|-----------------|-----------------------|-----------------------|------------------------|-------------------|
| 2025-Apr-08 | 20:13:55 | 2025-Apr-08 | 20:17:46 | 2025-Apr-08 | 19:49:45 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | | Project ID: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Mode: | SQC | Application Mode: | Fill Weight Variation | Application Mode: | Flow Rate Control |
| Result: | Accept | Sample Name: | | Result: | |
| Average: | 100.010 g | Sample ID: | | Max Flow Rate: | 0.543 ml/min |
| Minimum: | 100.010 g | Batch ID: | | Min Flow Rate: | 0.526 ml/min |
| Maximum: | 100.011 g | Lot ID: | | Average Flow Rate: | 0.533 ml/min |
| Standard Deviation: | 0.000 g | Samples: | 5 | Pump Cycle: | 3 times |
| Test Result: | | Unit1: | 100.01[1] g | Test Result: | |
| Sample ID Weight T1/T2 | | Unit2: | 100.01[1] g | 1: | 0.529 ml/min |
| Sample 001 100.011 g >=-T1 & <=+T1 | | Unit3: | 100.01[1] g | 2: | 0.526 ml/min |
| Sample 002 100.010 g >=-T1 & <=+T1 | | Unit4: | 100.01[1] g | 3: | 0.543 ml/min |
| Sample 003 100.010 g >=-T1 & <=+T1 | | Unit5: | 100.01[1] g | Pump Speed: | 0.500 ml/min |
| Batch: | Batch 01 | Difference: | | Liquid Density: | 0.99823 g/ml |
| Number in Batch: | 100% Inspection | 1: 0.00[3] g 0.00 % | | Target Volume: | 10.00 ml |
| Inspection Quantity: | 3 | 2: 0.00[3] g 0.00 % | | Number of Pump Cycles: | 3 times |
| Tare Mode: | Off | 3: 0.00[3] g 0.00 % | | Mode: | Flow Rate Control |
| Pretare: | 0.000 g | 4: 0.00[3] g 0.00 % | | Signature: | _____ |
| Sample Type: | Solid | 5: 0.00[3] g 0.00 % | | Verified By: | _____ |
| EU Criteria: | Off | Result: | | | |
| Nominal Weight: | 100.000 g | Accept: | 5 Units | | |
| +T1: | 0.300 g | Fail: | 0 Units | | |
| -T1: | -0.300 g | Acceptance Criteria: | Off | | |
| +T2: | 0.600 g | Average Weight: | 100.00[8] g | | |
| -T2: | -0.600 g | +/- Limit %: | 7.50 % | | |
| Signature: | _____ | Signature: | _____ | | |
| Verified By: | _____ | Verified By: | _____ | | |

6.6 Printout Examples

Here are the examples for each application are displayed with all items enabled in the Print Content menu. The default values for Header lines 1-5 are also indicated.

Notes:

- When the user management and electronic signature are active, the User ID will display in the Signature Line. The electronic signature is disabled by default.
- When a library is activated, "Library Name:" will appear below "Application" in the printout.

6.6.1 Calibration Report Template

| Internal Calibration | Span Calibration | Repeatability Test | | | | | | | | | | | | | | | | | | | | | |
|--|--|---|--------|-----------|-----------|---|-----------|------------|---|------------|------------|---|------------|------------|---|------------|------------|---|-----------|------------|---|-----------|------------|
| Calibration Report -----Internal Calibration----- Date and Time: 2025-May-12 10:34:02 Balance ID: Balance Name: EXP224/AD Result: 0.0000g Difference Internal Calibration Successful Signature: User ID Verified By: _____ | Calibration Report -----Span Calibration----- Date and Time: 2025-May-12 10:34:02 Balance ID: Balance Name: EXP224/AD Result: 0.0000g Difference Span Calibration Successful Signature: User ID Verified By: _____ | Routine Test: Repeatability Test Date: 2025-Apr-11 Start Time: 19:13:45 End Time: 19:15:52 Test Weights ID: Test Weights Value: 25.00000 g Test Weights Class: Test Result: <table border="1"> <thead> <tr> <th>Number</th> <th>Zero Load</th> <th>Full Load</th> </tr> </thead> <tbody> <tr> <td>1</td> <td>0.00000 g</td> <td>25.00496 g</td> </tr> <tr> <td>2</td> <td>-0.00002 g</td> <td>25.00262 g</td> </tr> <tr> <td>3</td> <td>-0.00013 g</td> <td>25.00063 g</td> </tr> <tr> <td>4</td> <td>-0.00006 g</td> <td>25.01753 g</td> </tr> <tr> <td>5</td> <td>0.00001 g</td> <td>25.00375 g</td> </tr> <tr> <td>6</td> <td>0.00002 g</td> <td>25.00014 g</td> </tr> </tbody> </table> SD (Span): 0.006441 g Signature: _____ Verified By: _____ | Number | Zero Load | Full Load | 1 | 0.00000 g | 25.00496 g | 2 | -0.00002 g | 25.00262 g | 3 | -0.00013 g | 25.00063 g | 4 | -0.00006 g | 25.01753 g | 5 | 0.00001 g | 25.00375 g | 6 | 0.00002 g | 25.00014 g |
| Number | Zero Load | Full Load | | | | | | | | | | | | | | | | | | | | | |
| 1 | 0.00000 g | 25.00496 g | | | | | | | | | | | | | | | | | | | | | |
| 2 | -0.00002 g | 25.00262 g | | | | | | | | | | | | | | | | | | | | | |
| 3 | -0.00013 g | 25.00063 g | | | | | | | | | | | | | | | | | | | | | |
| 4 | -0.00006 g | 25.01753 g | | | | | | | | | | | | | | | | | | | | | |
| 5 | 0.00001 g | 25.00375 g | | | | | | | | | | | | | | | | | | | | | |
| 6 | 0.00002 g | 25.00014 g | | | | | | | | | | | | | | | | | | | | | |

6.6.2 Application Print Out Template

| Basic Weighing | | Parts Counting | | Check Counting | |
|-----------------------|----------------|-----------------------|----------------|-----------------------|----------------|
| 12-24-2024 | 09:31:46 | 12-24-2024 | 10:13:34 | 12-24-2024 | 10:17:42 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN058 | Project Name: | PN070 | Project Name: | PN081 |
| ----- | | ----- | | ----- | |
| Application Name: | Basic Weighing | Application Name: | Parts Counting | Application Name: | Check Counting |
| Sample Name: | SN058 | Sample Name: | SN070 | Sample Name: | SN081 |
| Sample ID: | S058 | Sample ID: | S070 | Sample ID: | S081 |
| Batch ID: | B058 | Batch ID: | B070 | Batch ID: | B081 |
| Lot ID: | L058 | Lot ID: | L070 | Lot ID: | L081 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | Customer ID2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 3.5275 | Result: | 25 PCS | Result: | 74 PCS |
| oz N | | Gross: | 78.48 g | Status: | Accept |
| Gross: | 6.2960 | Net: | 78.48 g N | Gross: | 178.49 g |
| oz | | Tare: | 0.00 g T | Net: | 178.49 g N |
| Net: | 3.5275 oz N | | | Tare: | 0.00 g T |
| Tare: | 2.7690 oz T | Library: | Library 070 | | |
| | | APW: | 3.124 g | Library: | Library 081 |
| Library: | Library 058 | Samples: | 10 PCS | APW: | 2.412 g |
| | | | | Samples: | 81 PCS |
| Signature: _____ | | Signature: _____ | | Over: | 254 PCS |
| | | | | Under: | 51 PCS |
| Verify By: _____ | | Verify By: _____ | | Signature: _____ | |
| | | | | Verify By: _____ | |

| Percent Weighing | | Check Weighing | | Dynamic | |
|-----------------------|------------------|-----------------------|----------------|-----------------------|------------------|
| 12-26-2024 | 13:30:04 | 12-26-2024 | 13:33:25 | 12-26-2024 | 13:36:21 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID : | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN010 | Project Name: | PN028 | Project Name: | PN055 |
| ----- | | ----- | | ----- | |
| Application Name: | Percent Weighing | Application Name: | Check Weighing | Application Name: | Dynamic Weighing |
| Sample Name: | SN010 | Sample Name: | SN028 | Sample Name: | SN055 |
| Sample ID: | S010 | Sample ID: | S028 | Sample ID: | S055 |
| Batch ID: | B010 | Batch ID: | B028 | Batch ID: | B055 |
| Lot ID: | L010 | Lot ID: | L028 | Lot ID: | L055 |
| Customer ID1: | C1 | Customer ID1: | C1 | Customer ID1: | C1 |
| Customer ID2: | C2 | Customer ID2: | C2 | C2: | C2 |
| Customer ID3: | C3 | Customer ID3: | C3 | Customer ID3: | C3 |
| Customer ID4: | C4 | Customer ID4: | C4 | Customer ID4: | C4 |
| Customer ID5: | C5 | Customer ID5: | C5 | Customer ID5: | C5 |
| Customer ID6: | C6 | Customer ID6: | C6 | Customer ID6: | C6 |
| Customer ID7: | C7 | Customer ID7: | C7 | Customer ID7: | C7 |
| Customer ID8: | C8 | Customer ID8: | C8 | Customer ID8: | C8 |
| Customer ID9: | C9 | Customer ID9: | C9 | Customer ID9: | C9 |
| Customer ID10: | C10 | Customer ID10: | C10 | Customer ID10: | C10 |
| Result: | 361.69 % | Result: | Accept | Result: | 367.00 g |
| Gross: | 198.93 g | Gross: | 198.92 g | Gross: | 198.93 g |
| Net: | 198.93 g N | Net: | 198.92 g N | Net: | 198.93 g N |
| Tare: | 0.00 g T | Tare: | 0.00 g T | Tare: | 0.00 g T |
| Library: | Library 010 | Library: | Library 028 | Library: | Library 055 |
| Reference Weight: | 550.00 g | Over Limit: | 214.25 g | Averaging Time: | 3 seconds |
| Reference Factor: | 10.0 % | Under Limit: | 12.54 g | Signature: | _____ |
| Difference Weight: | 143.93 g | Signature: | _____ | Verify By: | _____ |
| Difference Factor: | 261.69 % | Verify By: | _____ | | |
| Signature: | _____ | | | | |
| Verify By: | _____ | | | | |

| Totalization | | Formulation | | Differential | |
|---------------------------|---------------------|-----------------------|---------------------|-----------------------|----------------------|
| 12-26-2024 | 15:11:42 | 12-26-2024 | 16:30:04 | 2025-Apr-08 | 19:58:46 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | PN072 | Project Name: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Name: | Totalization | Application Name: | Formulation | Application Mode: | Differential |
| Sample Name: | SN072 | Result: | 199.09 g N | Result: | 3.89[3] g N |
| Sample ID: | S072 | Gross: | 299.09 g | Gross: | 18.42[0] g |
| Batch ID: | B072 | Net: | 199.09 g N | Net: | 3.89[3] g N |
| Lot ID: | L072 | Tare: | 100.00 g T | Tare: | 14.52[7] g T |
| Customer ID1: | C1 | Recipe: | Recipe 001 | Start Time: | 2025-Apr-08 19:57:07 |
| Customer ID2: | C2 | Item Name | Item Weight | End Time: | 2025-Apr-08 19:58:41 |
| Customer ID3: | C3 | Item 1 | 100.00 g | Reciprocal Proportion | On |
| Customer ID4: | C4 | Item 2 | 199.08 g | Absolute Value | On |
| Customer ID5: | C5 | Start Time: | 12-26-2024 16:29:44 | Item Name: | ITEM1 |
| Customer ID6: | C6 | End Time: | 12-26-2024 16:30:01 | Initial Weight: | 0.766 g |
| Customer ID7: | C7 | -----Sample Data----- | | Final Weight: | 0.760 g |
| Customer ID8: | C8 | Item Name: | Item 1 | Difference Weight: | 0.006 g |
| Customer ID9: | C9 | Target Weight: | 100.00 g | Different Percent: | 0.8 % |
| Customer ID10: | C10 | Actual Weight: | 100.00 g | Item Name: | ITEM2 |
| Result: | 278.33 g | Difference: | 0.00 % | Initial Weight: | 1.528 g |
| Gross: | 0.00 g | Item Name: | Item 2 | Final Weight: | 1.534 g |
| Net: | 0.00 g N | Target Weight: | 199.08 g | Difference Weight: | 0.006 g |
| Tare: | 0.00 g T | Actual Weight: | 199.08 g | Different Percent: | 0.4 % |
| Library: | Library 072 | Difference: | 0.00 % | Item Name: | ITEM3 |
| Start Time: | 12-26-2024 15:08:29 | Total Weight: | 299.08 g | Initial Weight: | 2.292 g |
| End Time: | 12-26-2024 15:11:42 | Signature: _____ | | Final Weight: | 2.325 g |
| -----Sample Data (g)----- | | Verify By: _____ | | Difference Weight: | 0.033 g |
| '1: | 198.93 | | | Different Percent: | 1.4 % |
| '2: | 19.85 | | | Item Name: | ITEM4 |
| '3: | 19.86 | | | Initial Weight: | 3.094 g |
| '4: | 19.87 | | | Final Weight: | 3.110 g |
| '5: | 19.88 | | | Difference Weight: | 0.016 g |
| Samples: | 5 | | | Different Percent: | 0.5 % |
| Average: | 55.67 g | | | Item Name: | ITEM5 |
| Maximum: | 198.93 g | | | Initial Weight: | 3.871 g |
| Minimum: | 19.85 g | | | Final Weight: | 3.893 g |
| Range: | 179.08 g | | | Difference Weight: | 0.022 g |
| Standard Deviation: | 71.63 g | | | Different Percent: | 0.6 % |
| Relative Deviation(std%): | 128.68 % | | | Signature: _____ | |
| Signature: _____ | | | | Verified By: _____ | |
| Verify By: _____ | | | | | |

| Density Determination | | Peak Hold | | Pipette Adjustment | |
|-----------------------|---------------------------|-----------------------|-----------|---------------------------------------|---------------------------|
| 2025-Apr-08 | 20:01:25 | 2025-Apr-08 | 20:25:41 | 025-Apr-08 | 20:11:58 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| ----- | | ----- | | ----- | |
| Application Mode: | Density Determination | Application Mode: | Peak Hold | Application Mode: | Pipette Adjustment |
| Result: | 7.435 g/cm ³ | Result: | 355.512 g | Result: | Pass |
| Gross: | 19.44[1] g | Gross: | 0.000 g | Gross: | 24.76[4] g ? |
| Net: | 4.91[4] g N | Net: | 0.000 g N | Net: | 1.01[1] g ? N |
| Tare: | 14.52[7] g T | Tare: | 0.000 g T | Tare: | 23.75[3] g T |
| Weight in air.: | 5.67[4] g | On Stability: | No | Start Time: | 2025-Apr-08 20:09:03 |
| Weight in liquid.: | 4.91[3] g | Signature: _____ | | End Time: | 2025-Apr-08 20:11:52 |
| Auxiliary liquid.: | Water | Verified By: _____ | | Liquid Density: | 0.99823 g/cm ³ |
| Liquid Density: | 0.99823 g/cm ³ | | | Pressure: | 1.0 ATM |
| Water Temp.: | 20.0 °C | | | Pipette Test Method: | Method 001 |
| Porous Material: | Off | | | Nominal Inaccuracy Imprecision Number | |
| Signature: _____ | | | | 1000 ul 5.00 % 5.00 % 3 | |
| Verified By: _____ | | | | Pipette Test Results | |
| | | | | ----- | |
| | | | | Pipette Name: | pip001 |
| | | | | Pipette Number: | |
| | | | | Nominal Volume: | 1000 ul Pass |
| | | | | -----Sample Data----- | |
| | | | | 1 1023 ul 1.021 g | |
| | | | | 2 1032 ul 1.030 g | |
| | | | | 3 1016 ul 1.014 g | |
| | | | | Inaccuracy | |
| | | | | E%: | 2.64 % |
| | | | | E% Limit: | 5.00 % |
| | | | | Average: | 1026 ul |
| | | | | Imprecision | |
| | | | | CV%: | 0.97 % |
| | | | | CV% Limit: | 5.00 % |
| | | | | Standard Deviation: | 10 ul |
| | | | | > +2S: | 0.00 % , 0 |
| | | | | > +1S: | 0.00 % , 0 |
| | | | | +1S > Mean > -1S: | 0.00 % , 0 |
| | | | | < -1S: | 0.00 % , 0 |
| | | | | < -2S: | 100.00 % , 10 |
| | | | | Signature: _____ | |
| | | | | Verified By: _____ | |

| SQC | | Fill Weight Variation | | Flow Rate Control | |
|------------------------------------|-----------------|-----------------------|-----------------------|------------------------|-------------------|
| 2025-Apr-08 | 20:13:55 | 2025-Apr-08 | 20:17:46 | 2025-Apr-08 | 19:49:45 |
| Balance ID: | | Balance ID: | | Balance ID: | |
| Balance Name: | | Balance Name: | | Balance Name: | |
| User ID: | Admin | User ID: | Admin | User ID: | Admin |
| First Name/Last Name: | Admin | First Name/Last Name: | Admin | First Name/Last Name: | Admin |
| Project Name: | | Project ID: | | Project Name: | |
| ----- | | ----- | | ----- | |
| Application Mode: | SQC | Application Mode: | Fill Weight Variation | Application Mode: | Flow Rate Control |
| Result: | Accept | Sample Name: | | Result: | |
| Average: | 100.010 g | Sample ID: | | Max Flow Rate: | 0.543 ml/min |
| Minimum: | 100.010 g | Batch ID: | | Min Flow Rate: | 0.526 ml/min |
| Maximum: | 100.011 g | Lot ID: | | Average Flow Rate: | 0.533 ml/min |
| Standard Deviation: | 0.000 g | Samples: | 5 | Pump Cycle: | 3 times |
| Test Result: | | Unit1: | 100.01[1] g | Test Result: | |
| Sample ID Weight T1/T2 | | Unit2: | 100.01[1] g | 1: | 0.529 ml/min |
| Sample 001 100.011 g >=-T1 & <=+T1 | | Unit3: | 100.01[1] g | 2: | 0.526 ml/min |
| Sample 002 100.010 g >=-T1 & <=+T1 | | Unit4: | 100.01[1] g | 3: | 0.543 ml/min |
| Sample 003 100.010 g >=-T1 & <=+T1 | | Unit5: | 100.01[1] g | Pump Speed: | 0.500 ml/min |
| Batch: | Batch 01 | Difference: | | Liquid Density: | 0.99823 g/ml |
| Number in Batch: | 100% Inspection | 1: 0.00[3] g 0.00 % | | Target Volume: | 10.00 ml |
| Inspection Quantity: | 3 | 2: 0.00[3] g 0.00 % | | Number of Pump Cycles: | 3 times |
| Tare Mode: | Off | 3: 0.00[3] g 0.00 % | | Mode: | Flow Rate Control |
| Pretare: | 0.000 g | 4: 0.00[3] g 0.00 % | | Signature: _____ | |
| Sample Type: | Solid | 5: 0.00[3] g 0.00 % | | Verified By: _____ | |
| EU Criteria: | Off | Result: | | | |
| Nominal Weight: | 100.000 g | Accept: | 5 Units | | |
| +T1: | 0.300 g | Fail: | 0 Units | | |
| -T1: | -0.300 g | Acceptance Criteria: | Off | | |
| +T2: | 0.600 g | Average Weight: | 100.00[8] g | | |
| -T2: | -0.600 g | +/- Limit %: | 7.50 % | | |
| Signature: _____ | | Signature: _____ | | | |
| Verified By: _____ | | Verified By: _____ | | | |

6.6.3 Batch Printing Output

| Printout | Export to PDF |
|---------------------------------|---------------|
| User ID: Admin | |
| First Name/Last Name: Admin | |
| Sample Name: | |
| Sample ID: | |
| Customized ID: | |
| -----Sample Data----- | |
| 1 25.00120 g 23:19:39 | |
| 2 24.99835 g 23:19:46 | |
| 3 25.00070 g 23:19:47 | |
| 4 25.00087 g 23:19:48 | |
| 5 25.00106 g 23:19:49 | |
| Signature: | |
| _____ | |
| Verified | |
| By: _____ | |

7 Library

Each Explorer Plus balance is equipped with an advanced library function that easy to create, edit, and activate libraries within seconds. The substantial memory capacity accommodates up to 15 weighing application modes and can store around 200 libraries.

In the Library menu, users can review the list of libraries they have generated for various application modes. These libraries can be exported to PC software and imported whenever necessary.

In the upper right corner, the total library memory is displayed. For instance, 9/1% indicates that 9 libraries have been created and are occupying approximately 1% of the total capacity. When the memory usage reaches 80%, the balance will display a message prompting the user to delete or back up library records.

Note:

In some application, the library name will change to Method, Recipe an

| HOME | | LIBRARY | | | 9 / 1% |
|-------------|---------|----------------|-------------|----------------------|--------|
| LIBRARY ID | USER ID | APPLICATION | SAMPLE NAME | DATE AND TIME | |
| Library 009 | Admin | Check Weighing | | 2025-Apr-29 13:39:19 | |
| Library 008 | Admin | Check Weighing | | 2025-Apr-29 13:39:18 | |
| Library 007 | Admin | Check Weighing | | 2025-Apr-29 13:39:16 | |
| Library 006 | Admin | Basic Weighing | | 2025-Apr-29 13:39:00 | |
| Library 005 | Admin | Basic Weighing | | 2025-Apr-29 13:38:59 | |

< 1 / 2 >

| | | |
|------------|--------|------------|
| DELETE ALL | IMPORT | EXPORT ALL |
|------------|--------|------------|

7.1 Library Data

The following data is stored for the Application used:

| Application Mode | Library Function | Data Specific to Application Modes | Sample Data |
|-----------------------|------------------|---|--|
| Basic Weighing | Yes | Minimum Weight | Sample Name Sample ID Batch ID Lot ID Project Name 10 x Customer ID |
| Parts Counting | Yes | APW, Sample Size | |
| Check Counting | Yes | APW, Sample Size, Over Limit, Under Limit | |
| Percent Weighing | Yes | Reference weight, Reference factor | |
| Check Weighing | Yes | Over Limit, Under Limit, Nominal Weight, +Weight Tolerance, -Weight Tolerance, +Percent Tolerance, -Percent Tolerance | |
| Dynamic Weighing | Yes | Average Time | |
| Totalization | Yes | | |
| Density Determination | Yes | Sinker Volume, Water Temp., Liquid Density, Oil Density | |
| Peak Hold | Yes | | |
| Formulation | Recipe | Recipe Name, Item Name, Item Weight | |
| Differential Weighing | No Library | | |

| | | | |
|------------------------------|-------------------|--|--|
| Pipette Adjustment | Method | Method Name, Nominal, Pipette Name, Pipette Number, Inaccuracy, Impression, Number of Samples | N/A |
| SQC | Batch | Number in Batch, Inspection Quantity, Tare Mode, Sample Type, EU Criteria, Nominal Weight +T1, -T1, +T2, -T2 | Sample Name Sample ID Batch ID Lot ID Project Name 10 x Customer ID |
| Fill Weight Variation | No Library | | |
| Flow Rate Control | Yes | Pump Speed, Pump Time, Liquid Density, Target Volume | Sample Name Sample ID Batch ID Lot ID Project Name 10 x Customer ID |

7.2 Import and Export Library

Users can import a "library.db" file into the balance.

- Save the "library.db" file to a USB flash drive. Connect the USB drive to the balance. Once connected, the IMPORT button will become active.
- Press the IMPORT button to initiate the import process.

Users can export a "library.db" file to a USB flash device.

- Connect the USB drive to the balance. Once connected, the EXPORT button will become active.
- Press the EXPORT button to download the "library.db" file to the USB flash device.

8 Maintenance

This function serves to is user to diagnostic hardware functions, update software, review the service log file, and provide service technicians to use service menu.


Notes:

- Service Menu is locked to prevent authority change make the incorrect weighing performance.
- If you encounter any issues while using the balance, please contact Ohaus or your local Ohaus dealer for assistance.


8.1 Maintenance Menu

User can diagnostic hardware functions, such as Autodoor, IR sensor, Digital Leveling Bubble, Terminal Status Lights, Draftshield Lights, and Fingerprint Accessory if connected.


Auto Doors: Press **Auto Doors** to start testing. If this function is working properly, the doors will open and then close automatically.

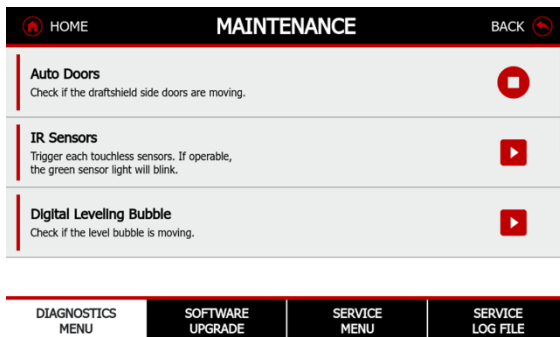

IR Sensors: Press **IR Sensors** to start testing. Waving your hand over it will trigger that function and cause the sensor light to turn green. When the function is working properly, the status icon on the right side will turn green.

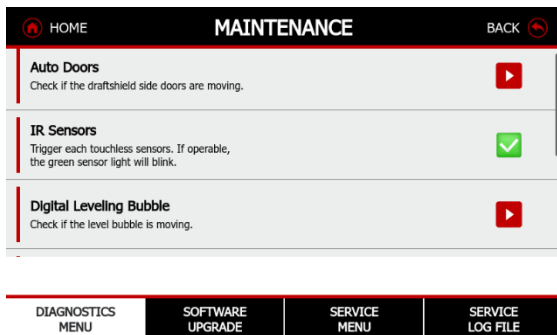


Digital Leveling Bubble: To check the bubble movement, the user needs to lift the balance. When the function is working properly, the status icon on the right side will turn green. If the leveling bubble is not moving, the status icon will display a warning sign.

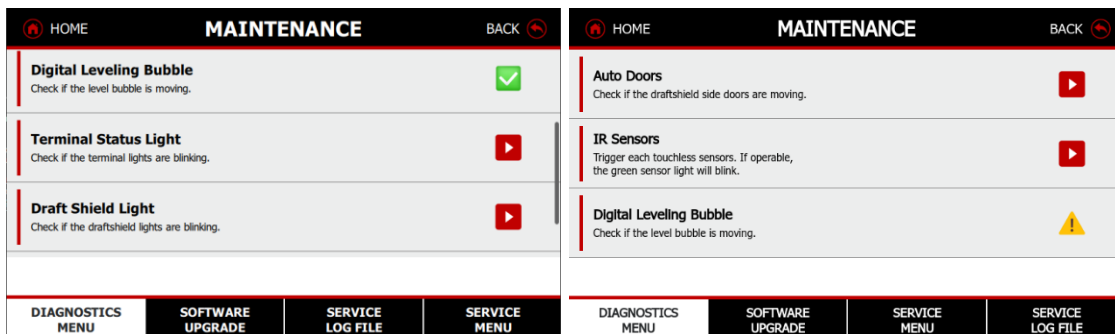

Terminal Lights: Press **Terminal Lights** button to check if the lights are functioning. The balance will cycle through all the light colors on the terminal.

Draft Shied Lights: Press **Draft Shied Lights** button to check if the lights are functioning. The balance will cycle through all the light colors in the draftshield.

Fingerprint: Press **Fingerprint** button to check if the accessory is functioning. When the function is working properly, the balance will identify the serial number of the fingerprint module.

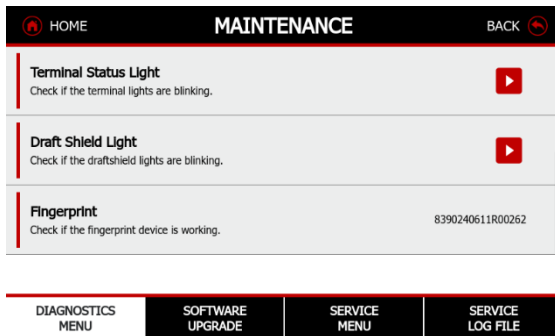


8.2 Software Upgrade

In this function, users need to obtain the correct software to upgrade the balance for feature enhancements or bug fixes. If you encounter any issues while using the balance, please contact Ohaus or your local Ohaus dealer for assistance.

8.2.1 Software Upgrade Process

- Save the upgrade file into a USB flash drive.
- Connect the USB flash driver to the Balance.
- Press START button to upgrade the system.
- When completed, the balance would reboot automatically.

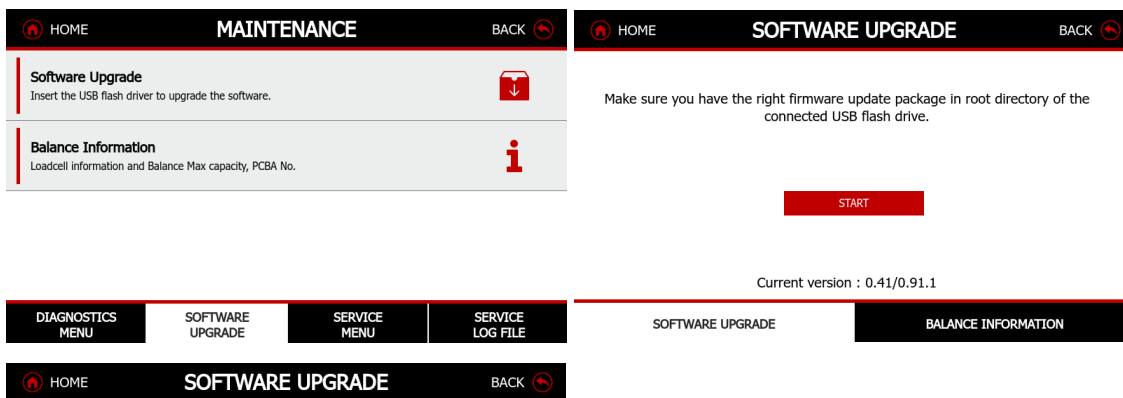

The upgrade has finished successfully. The balance will reboot.



8.2.2 Balance Information

This screen would provide essential information of the balance.

8.3 Service Menu

Service Menu is locked to prevent authority change make the incorrect weighing performance. If you encounter any issues while using or service the balance, please contact Ohaus or your local Ohaus dealer for assistance.

8.4 Service Log File

The Service Log file is record the balance Failure log, System Log, Calibration Log, Printed Data Log

- Failure Log: 100 of electronic record of record the number of times an overload /underload occurs.
- System Log: 100,000 entries which contain electronic records of the menu change on the balance, create/delete user, log in/off, date and time change and etc. Only Administrator and supervisor can be export to PDF as the un-editable format.
- Calibration Log: 3000 of electronic record of the user performed calibration reports and service performed calibration reports. Only Administrator and supervisor can be export to PDF as the un-editable format.
- Printed Data Log- 100,00 of electronic record of the routine weighing data that is sent either by manually pressing the print key or through a command from a PC.

8.5 Factory Reset

Use this sub-menu to reset the menus to their Factory default settings.

- Reset All
- Quick Setup
- Calibration
- Balance Setup
- Application Modes
- Weighing Units
- Communication
- Library

8.6 Log Off

Press this button to log out of the current user account.

8.7 Power Off

Press this button to turn off the balance.

9 Maintenance

When the balance is used in trade or a legally controlled application it must be set up, verified and sealed in accordance with local weights and measures regulations. It is the responsibility of the purchaser to ensure that all pertinent legal requirements are met.





*For Explorer Plus...N...models, please refer to Legal for Trade Settings.

9.1 Legal for Trade Setting

Before verification and sealing, perform the following steps in Order:

- Verify that the menu settings meet the local weights and measures regulations.
- Weighing Unit menu should be reviewed. Verify the units turned On meet the local weights and measures regulations.
- Perform a calibration as explained in 5.4 Calibration.
- Set the position of the Legal for Trade Switch to the locked position.

■ **Legal for Trade Switch:**

| | Explorer Plus Analytical and Precision Balance | Explorer Plus High-Capacity Balance |
|-------------|---|--|
| Position |  |  |
| Lock/Unlock |  |  |

- Set Legal for Trade to ON in the Balance Setup menu (refer to Approved Mode in section 5.3.1).

9.2 Settings Change When Approved Mode is ON

When Legal for Trade is set to ON, the menu settings are affected as follows:

| Menu | Changes |
|------------------------|---|
| Calibration Menu | <ul style="list-style-type: none"> • Internal Calibration will be locked at On. For EXP...N... models: <ul style="list-style-type: none"> ○ If Internal Calibration is Off before turning on Approved Mode, the Internal Calibration field in Calibration Settings will be locked at Off. ○ If Internal Calibration is On before turning on Approved Mode, the field and tab will remain unchanged. • AutoCal™ will be locked at its current setting. • AutoCal™ Interval (Hours) will be locked at current setting. • Span Calibration will be locked at off and greyed out. |
| Balance Setup Menu | <ul style="list-style-type: none"> • Auto Zero Tracking is limited to 0.5d or Off. • If AZT is Off before turning on Approved Mode, then the field will be locked at "Off". <ul style="list-style-type: none"> ○ For 0.01mg models, if AZT is on before turning on Approved Mode, the field will be forced to "3d" and locked. ○ For other models, if AZT is on before turning on Approved Mode, the field will be forced to "0.5d" and locked. • Gross Indicator will be locked at the current setting. • Graduations will be forced to 1 Division and the field will be hidden. • For EXP...N...models, graduations will be locked at its current setting. • In System Log menu, Auto Login will be locked at Off. |
| Maintenance Menu | <ul style="list-style-type: none"> • Software Upgrade will be locked at Off. • Service Menu will be locked at Off. |
| Weighing Units Menu | <ul style="list-style-type: none"> • Units will be restricted to metric units. • For EXP...N...models, Units will be restricted to metric and imperial units. |
| Printing Settings Menu | <ul style="list-style-type: none"> • In Connect to Printer, Numeric Only will be locked at off, and Manual Print will be locked at Stable only. • In Connect to PC, Numeric Only will be locked at off, and Manual Print will be locked at Stable only. • In Data to Excel, Manual Print will be locked at Stable only |
| Terminal Setting: | <ul style="list-style-type: none"> • After turning on Approved Mode, it will not be allowed to switch terminal. |
| Applications | <ul style="list-style-type: none"> • In Item Settings menu, Auto Tare will be locked at Off. |

9.3 Verification

A weights and measures official or authorized service agent must perform the verification procedure.

9.4 Sealing

After the Balance has been verified, it must be sealed to prevent undetected access to the legally controlled settings. Before sealing the device, ensure that the security switch is in the Locked position and the Legal for Trade setting in the Balance Setup menu has been set to ON.

If using a wire seal, pass the sealing wire through the holes in the security switch and Bottom Housing as shown.

If using a paper seal, place the seal over the security switch and Bottom Housing as shown.

- Explore Plus Semi-Micro, Analytical and Precision Balance



- Explore Plus High Capacity Balance



9.5 Output Format

Strings Definition

| Field: | Label ¹ | Space ² | Weight ³ | Space ² | Unit ⁴ | Space | Stability ⁵ | Space | G/N ⁶ | Space | Term. Characters ⁷ |
|---------|--------------------|--------------------|---------------------|--------------------|-------------------|-------|------------------------|-------|------------------|-------|-------------------------------|
| Length: | | 1 | 11 | 1 | | 1 | ≤ 1 | ≤ 1 | ≤ 3 | 0 | ≤ 8 |

- In certain cases, a Label field of up to 11 characters is included.
- Each field is followed by a single delimiting space (ASCII 32).
- The Weight field is 9 right justified characters. If the value is negative, the “-” character is located at the immediate left of the most significant digit.
- The Unit field contains the unit of measure abbreviation up to 5 characters.
- The Stability field contains the “?” character if the weight reading is not stable. The Stability field and the following Space field are omitted if the weight reading is stable.
- The G/N field contains the net or gross indication. For net weights, the field contains “NET”. For gross weights, the field contains nothing, “G” or “B”, depending on the GROSS INDICATOR menu setting.
- The Termination Characters field contains CRLF, Four CRLF or Form Feed (ASCII 12), depending on the LINE FEED menu setting.

10 MAINTENANCE

10.1 Calibration

Periodically verify calibration by placing an accurate weight on the balance and viewing the result. If calibration is required, perform a Balance internal calibration.

10.2 Cleaning



WARNING: Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.



Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The housing may be cleaned with a cloth dampened with a mild detergent if necessary.

10.3 Battery Power



CAUTION: Battery is to be replaced only by an authorized Ohaus service dealer. Risk of explosion can occur if the rechargeable battery is replaced with the wrong type or if it is not properly connected. Dispose of the rechargeable battery according to local laws and regulations.

10.4 Troubleshooting

| Symptom / Display | Possible Cause | Remedy |
|-----------------------------|--|---|
| Balance will not turn on | <ul style="list-style-type: none"> No power supply to the Balance | <ul style="list-style-type: none"> Verify connection and voltage |
| Inaccuracy weighing results | <ul style="list-style-type: none"> Improper calibration Unstable environment | <ul style="list-style-type: none"> Perform calibration Move balance to suitable location |
| Cannot calibrate | <ul style="list-style-type: none"> Calibration Menu locked Approved Mode set to on Instable environment Incorrect calibration masses | <ul style="list-style-type: none"> Turn Calibration menu lock off Turn Approved Mode off Move balance to suitable location Use correct calibration masses |
| Cannot change menu settings | <ul style="list-style-type: none"> Sub-menu locked Approved Mode set to on | <ul style="list-style-type: none"> Unlock sub-menu Turn Approved Mode off |
| Low Reference Weight | <ul style="list-style-type: none"> Reference weight too small The weight on the pan is too small to define a valid reference weight. | <ul style="list-style-type: none"> Increase sample size |
| Invalid Piece Weight | <ul style="list-style-type: none"> Average piece weight is too small | <ul style="list-style-type: none"> Increase average piece weight |
| Operation Timeout | <ul style="list-style-type: none"> Weight reading is not stable | <ul style="list-style-type: none"> Move balance to suitable location |
| ----- | <ul style="list-style-type: none"> Busy (tare, zero, printing) | <ul style="list-style-type: none"> Wait until completion |

10.5 End of Life Instruction

OHAUS electronic balances are precision instruments consisting of a metal housing, coated aluminum weighing cells, stainless steel parts, ABS/PC plastic parts, cardboard and foam packaging, and other materials. Kindly comply with the sustainable instructions for handling the balance when it is not in use or reaches the end of its operational lifespan.

Data Management: Before disposing of the electronic balance, ensure that all sensitive data or user information is erased from the device. Follow the balance instruction manual for data removal or seek OHAUS service engineer provider assistance.

Reuse or Donate: Consider donating your electronic balance to a school, community center, or charitable organization that might have a use for it. Reusing the equipment is the most sustainable option.

Recycling: If the electronic balance is no longer functional or cannot be reused, consider recycling it. Look for electronic recycling centers in your area that accept electronic equipment. Make sure to choose a reputable recycling center that follows proper e-waste disposal regulations. In both the US and the EU, aluminum and stainless steel are considered readily recyclable while ABS/PC plastic can be recycled, but not as easily. In the UK, aluminum, stainless steel, and ABS/PC plastics are considered readily recyclable.

Disposal: If recycling is not an option, dispose of the electronic balance responsibly. Do not throw it in the regular trash, as it can potentially harm the environment. Check with your local waste management authorities for guidance on properly disposing of electronic equipment, such as EU directive 2002/96/EC (WEEE), refer to www.ohaus.com/weee.

Packaging: When transporting the electronic balance for recycling or disposal, use minimal packaging material and consider using eco-friendly packaging options. Avoid single-use plastics and opt for recyclable or biodegradable materials.

Sustainable Alternatives: When purchasing a new electronic balance, choose a product from a manufacturer that prioritizes sustainability.

10.5.1 Material Composition of Draftshield Models

- Material Composition for Explorer Draftshield Models

| | Material Composition | Main Usage | Percentage of Row Materials | DS Model Weight (kg) | % |
|--------------------------|----------------------|-----------------------------------|-----------------------------|----------------------|----------------|
| Products | Metal Parts | Housing, Loadcell | 100% Aluminum | 4.54 | 41% |
| | Plastic Parts | Leveling feet, wing rings, covers | 50% ABS | 1.34 | 12% |
| | Glass | Draftshield | | 1.58 | 14% |
| | Electronic Parts | PCB | | 0.74 | 7% |
| | Fastener | Screws | 70% stainless steel | 0.08 | 1% |
| | Cables | Connections | copper+rubber skin | 0.12 | 1% |
| Packaging | Paper Parts | Instruction Manuals, carton box | 100% recycle paper | 1.88 | 17% |
| | Foams | Shipping Foams | 100% EPP | 0.92 | 8% |
| Total Weight (kg) | | | | 11.19 | 100.00% |

10.6 Service Information

OHAUS prioritizes energy-efficient design, eco-friendly packaging, and other sustainable objectives, reflected in all its laboratory weighing products that incorporate sustainable product goals. By following these sustainable end-of-life instructions, you can ensure that your electronic balance is disposed of in an environmentally friendly and sustainable manner. Thank you for choosing the OHAUS instrument and taking the necessary steps to protect the environment.

If the troubleshooting section does not resolve your problem, contact an Authorized Ohaus Service Agent. Please visit our website www.ohaus.com to locate the Ohaus office nearest you. An Ohaus Product Service Specialist will be available to assist you.

11 TECHNICAL DATA

11.1 Service Information

Ambient Conditions

- Indoor use only
- Altitude: 2000 m
- Specified Temperature range: 10°C to 30°C. Operability is assured at ambient temperatures between 5°C and 40°C
- Humidity: Maximum relative humidity 80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40°C.
- Electrical supply:
 - 12VDC, 1.5A. (For use with certified or approved power supply, which must have a SELV and limited energy output.). (For the models powered by external power adapter.)
 - 100 - 240V~, 0.5A, 50/60Hz. (For Explorer Plus High-Capacity models)
- Mains supply voltage fluctuations: up to ±10% of the nominal voltage
- Installation category II
- Pollution degree: 2

Materials

- Bottom Housing;
 - Die-cast Aluminum, Painted
 - IP54 waterproof protected base (EXP24001, EXP35001, and EXP65001)
- Top Housing: Die-cast Aluminum, Painted
- Terminal: Glass, Die-cast Aluminum
- Weighing Platforms:
 - Plastic (PC)
 - Die-cast Aluminum (EXP24001, EXP35001, and EXP65001)
- Weighing Pan:
 - Zink alloy (0.01mg models)
 - 316 SST (0.1mg, 1mg, 0.01g, 0.1g models)
 - 304 SST (EXP24001, EXP35001, and EXP65001)
- In-use Cover: plastic (PET)
- Draft Shield; Glass, Aluminum, Plastic

11.2 Model Specifications

| MODEL | EXP125D/AD | EXP125/AD | EXP225D/AD | EXP225/AD |
|---|--|------------|---------------|------------|
| | EXP125DM/AD | EXP125M/AD | EXP225DM/AD | EXP225M/AD |
| Capacity (g) | 82 g / 120 g | 120 g | 120 g / 220 g | 220 g |
| Readability d, fine range (g) | 0.01 mg | 0.01 mg | 0.01 mg | 0.01 mg |
| Readability d, full range (g) | 0.1 mg | 0.01 mg | 0.1 mg | 0.01 mg |
| Repeatability (sd.), ≤5% of full load | 0.01 mg | 0.01 mg | 0.01 mg | 0.01 mg |
| Repeatability (sd.), 5% of full load to fine range max | 0.02 mg | 0.02 mg | 0.02 mg | 0.02 mg |
| Repeatability (sd.), fine range max to full range | 0.1 mg | 0.02 mg | 0.1 mg | 0.02 mg |
| Linearity deviation typical | ± 0.06 mg | ± 0.06 mg | ± 0.06 mg | ± 0.06 mg |
| Linearity deviation | ± 0.1 mg | ± 0.1 mg | ± 0.1 mg | ± 0.1 mg |
| Verification interval, e (EXP...M..models only) | 1 mg | 1 mg | 1 mg | 1 mg |
| Approval Class (EXP...M.. models only) | Class I | Class I | Class I | Class I |
| Span Calibration Points (g) | 25g, 50g, 75g, 100g | | | |
| Weighing units* | 15 weighing units: g, mg, ct, N, oz, ozt, GN, dwt, mo, msg, tcl, tola, baht, Customer 1, Customer 2 | | | |
| Weighing units* (EXP...M.. models only) | g, mg, ct | | | |
| Applications | 15 application modes: Basic Weighing, Parts Counting, Check Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Totalization, Formulation, Differential, Density Determination, Peak Hold, Pipette Adjustment, SQC, Fill Weight Variation, Flow Rate Control | | | |
| Stabilization time (typical) | 0.1 mg: ≤2 seconds, 0.01 mg: ≤ 5 seconds | | | |
| Sensitivity Temperature Drift (ppm/K) | 0.8 | | | |
| Min-Weight (typical) (USP, K=2, U=0.10%) | 20 mg | | | |
| Min-Weight (optimal) (USP, K=2, U=0.10%, SRP≤0.41d) | 8.2 mg | | | |
| Terminal Display | 7-inch TFT 16.7M Color Press Screen with Glass Overlay, Detachable Terminal | | | |
| Number of Dots | 800 RGB x 480 DOTS | | | |
| Communication | USB Host (Type A) x 2, USB Device (Type B) x 1, USB Device (Type C) x 1 Ethernet (RJ45) x 1, RS232 x 1; Optional Wi-Fi, Bluetooth Dongle | | | |
| Leveling System | Motorized Leveling System with Digital Leveling Bubble | | | |
| Power Supply | AC Adapter Output: 12 VDC 1.5A | | | |
| Weighing Pan Size | Aero Pan 80 x 80 mm | | | |
| Build-in Ionizer | Standard Equipped | | | |
| Environmental Sensor RH, Air pressure Vibration Sensor | Temperature Sensor, Relative Humidity Sensor (RH), Air Pressure Sensor, Vibration Sensor | | | |
| IR Sensor | 2 Touchless Sensors on the terminal | | | |
| Anti-static glass panel | Standard Equipped | | | |
| Automatic Draftshield Door Function | Standard Equipped, Open Range: 1/2 open and fully open | | | |
| Draftshield & Terminal Status Lights | Standard Equipped | | | |
| Terminal Housing Dimensions (W x D x H) | 203mm x 124mm x 65mm | | | |
| Base Housing Dimensions (W x D x H) | | | | |
| Shipping Dimensions (W x D x H) | 630 x 415 x 516 mm | | | |
| Net Weight | 7825 g | | | |

| Shipping Weight | 120010 g | | | |
|---|---|---------------------------------|-------------------------|----------------------------------|
| MODEL | EXP124/AD | EXP224/AD | EXP324/AD | EXP524/AD |
| | EXP124M/AD | EXP224M/AD | EXP324M/AD | EXP524M/AD |
| | EXP124N/AD | EXP224N/AD | EXP324N/AD | EXP524N/AD |
| Capacity (g) | 120 g | 220 g | 320 g | 520 g |
| Readability d, full range (mg) | 0.1 mg | 0.1 mg | 0.1 mg | 0.1 mg |
| Repeatability (sd.), ≤5% of full load | 0.07 mg | 0.07 mg | 0.07 mg | 0.07 mg |
| Repeatability (sd.), 5% of full load to fine range max | 0.1 mg | 0.1 mg | 0.1 mg | 0.1 mg |
| Linearity deviation typical | 0.06 mg | 0.06 mg | 0.06 mg | 0.06 mg |
| Verification interval, e (EXP...M..models only) | 1 mg | 1 mg | 1 mg | 1 mg |
| Approval Class (EXP...M.. models only) | Class I | Class I | Class I | Class I |
| Span Calibration Points (g) | 25g, 50g, 75g, 100g | 50g, 100g, 150g, 200g | 100g, 200g, 300g | 200g, 300g, 400g, 500g |
| Weighing units* | 18 weighing units: g, mg, ct, N, oz, ozt, GN, dwt, mo, msg, tl H, tl S, tl T, tcl, tola, baht, Customer 1, Customer 2 | | | |
| Weighing units* (EXP...M.. models only) | g, mg, ct | | | |
| Applications | 15 application modes: Basic Weighing, Parts Counting, Check Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Totalization, Formulation, Differential, Density Determination, Peak Hold, Pipette Adjustment, SQC, Fill Weight Variation, Flow Rate Control | | | |
| Stabilization time (typical) | 2 seconds | | | |
| Sensitivity Temperature Drift (ppm/K) | 1.5 | 1.5 | 1.5 | 1.5 |
| Min-Weight (typical) (USP, K=2, U=0.10%) | 160 mg | | | |
| Min-Weight (optimal) (USP, K=2, U=0.10%, SRP≤0.41d) | 82 mg | | | |
| Terminal Display | 7-inch TFT 16.7M Color Press Screen with Glass Overlay, Detachable Terminal | | | |
| Number of Dots | 800 RGB x 480 DOTS | | | |
| Communication | USB Host (Type A) x 2, USB Device (Type B) x 1, USB Device (Type C) x 1 Ethernet (RJ45) x 1, RS232 x 1; Optional Wi-Fi, Bluetooth Dongle | | | |
| Leveling System | Motorized Leveling System with Digital Leveling Bubble | | | |
| Power Input | 12 VDC, 1.5A | | | |
| Power Supply | AC Adapter Input: 100-240 VAC 0.5A 50-60 Hz AC Adapter Output: 12 VDC 1.5A | | | |
| Weighing Pan Size | Square Pan 90 x 90 mm | | | |
| Build-in Ionizer | Standard Equipped | | | |
| Environmental Sensor RH, Air pressure Vibration Sensor | Temperature Sensor, Relative Humidity Sensor (RH), Air Pressure Sensor, Vibration Sensor | | | |
| IR Sensor | 2 Touchless Sensors on the terminal | | | |
| Anti-static glass panel | Standard Equipped | | | |
| Automatic Draftshield Door Function | Standard Equipped, Open Range: 1/2 open and fully open | | | |
| Draftshield & Terminal Status Lights | Standard Equipped | | | |
| Terminal Housing Dimensions (W x D x H) | 203mm x 124mm x 65mm | | | |
| Base Housing Dimensions (W x D x H) | | | | |
| Shipping Dimensions (W x D x H) | | | | |
| Net Weight | | | | |
| Shipping Weight | | | | |

| MODEL | EXP223/AD | EXP423/AD | EXP623/AD | EXP1203/AD | EXP3203/AD | EXP5203/AD |
|---|--|-------------------------------|--------------------------------|--------------------------------|--|-----------------------------------|
| | EXP223M/AD | EXP423M/AD | EXP623M/AD | EXP1203M/AD | EXP3203M/AD | EXP5203M/AD |
| | EXP223N/AD | EXP423N/AD | EXP623N/AD | EXP1203N/AD | EXP3203N/AD | EXP5203N/AD |
| Capacity (g) | 220 g | 420 g | 620 g | 1200 g | 3200 g | 5200 g |
| Readability d (mg) | 1mg | 1mg | 1mg | 1mg | 1mg | 1mg |
| Repeatability (sd.), ≤5% of full load (mg) | 0.7 mg | 0.7 mg | 0.7 mg | 0.7 mg | 0.7 mg | 0.7 mg |
| Repeatability (sd.), 5% of full load to fine range max (mg) | 1 mg | 1 mg | 1 mg | 1 mg | 1 mg | 1 mg |
| Linearity Deviation typical | ± 0.6 mg | ± 0.6 mg | ± 0.6 mg | ± 0.6 mg | ± 0.6 mg | ± 0.6 mg |
| Verification interval, e (EXP...M models and EXP...N models only) | 10 mg | 10 mg | 10 mg | 10 mg | 10 mg | 10 mg |
| Approval Class (EXP...M.. models and EXP..N..models only) | Class II | Class II | Class II | Class I | Class I | Class I |
| Span Calibration Points (g) | 50g, 100g, 150g, 200g | 100g, 200g, 300g, 400g | 300g, 400g, 500g , 600g | 400g, 600g, 800g, 1000g | 1000g, 1500g, 2000g, 2500g, 3000g | 2000g, 3000g, 4000g, 5000g |
| Weighing units* | 20 weighing units: g, mg, ct, N, oz, ozt, GN, dwt, mo, msg, tl H, tl S, tl T, tcl, tola, baht, Customer 1, Customer 2, lb (models with capacity ≥ 620 g), kg (models with capacity ≥ 1200 g) | | | | | |
| Weighing units* (EXP...M.. models and EXP..N..models only) | EXP...M.. models: g, mg, ct, kg (models with capacity ≥ 1200 g) EXP...N.. models: g, mg, ct, oz, ozt, GN, dwt, lb (models with capacity ≥ 620 g), kg (models with capacity ≥ 1200 g) | | | | | |
| Applications | 15 application modes: Basic Weighing, Parts Counting, Check Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Totalization, Formulation, Differential, Density Determination, Peak Hold, Pipette Adjustment, SQC, Fill Weight Variation, Flow Rate Control | | | | | |
| Stabilization time (typical) | ≤ 1.5 seconds | | | | | |
| Sensitivity Temperature Drift (ppm/K) | 3 | | | | | |
| Min-Weight (typical) (USP, K=2, U=0.10%) | 2 g | | | | | |
| Min-Weight (optimal) (USP, K=2, U=0.10%, SRP≤0.41d) | 0.82 g | | | | | |
| Terminal Display | 7-inch TFT 16.7M Color Press Screen with Glass Overlay, Detachable Terminal | | | | | |
| Number of Dots | 800 RGB x 480 DOTS | | | | | |
| Communication | USB Host (Type A) x 2, USB Device (Type B) x 1, USB Device (Type C) x 1 Ethernet (RJ45) x 1, RS232 x 1; Optional Wi-Fi, Bluetooth Dongle | | | | | |
| Leveling System | Motorized Leveling System with Digital Leveling Bubble | | | | | |
| Power Input | 12 VDC, 1.5A | | | | | |

| | |
|---|---|
| Power Supply | AC Adapter Input: 100-240 VAC 0.5A 50-60 Hz AC Adapter Output: 12 VDC 1.5A |
| Weighing Pan Size | Square Pan 130 x 130 mm |
| Build-in Ionizer | Standard Equipped |
| Environmental Sensor RH, Air pressure Vibration Sensor | Temperature Sensor, Relative Humidity Sensor (RH), Air Pressure Sensor, Vibration Sensor |
| IR Sensor | 2 Touchless Sensors on the terminal |
| Anti-static glass panel | Standard Equipped |
| Automatic Draftshield Door Function | Standard Equipped, Open Range: 1/2 open and fully open |
| Draftshield & Terminal Status Lights | Standard Equipped |
| Terminal Housing Dimensions (W x D x H) | 203mm x 124mm x 65mm |
| Base Housing Dimensions (W x D x H) | |
| Shipping Dimensions (W x D x H) | |
| Net Weight | |
| Shipping Weight | |

| MODEL | EXP2202 | EXP4202 | EXP6202 | EXP8202 | EXP12202 |
|---|--|-----------------------------------|---|-----------------------------------|-------------------------------------|
| | EXP2202M | EXP4202M | EXP6202M | EXP8202M | EXP12202M |
| | EXP2202N | EXP4202N | EXP6202N | EXP8202N | EXP12202N |
| Capacity (g) | 2200 g | 4200 g | 6200 g | 8200 g | 12200 g |
| Readability d, full range (g) | 0.01 g | 0.01 g | 0.01 g | 0.01 g | 0.01 g |
| Repeatability (sd.), ≤5% of full load (g) | 0.007 g | 0.007 g | 0.007 g | 0.007 g | 0.007 g |
| Repeatability (sd.), 5% of full load to fine range max (g) | 0.01 g | 0.01 g | 0.01 g | 0.01 g | 0.01 g |
| Linearity Deviation Typical | ± 0.006 g | ± 0.006 g | ± 0.006 g | ± 0.006 g | ± 0.006 g |
| Linearity Deviation | ± 0.02 g | ± 0.02 g | ± 0.02 g | ± 0.02 g | ± 0.02 g |
| Verification interval, e (EXP...M.. models and EXP...N..models only) | 0.1 g | 0.1 g | 0.1 g | 0.1 g | 0.1 g |
| Approval Class (EXP...M.. models and EXP...N..models only) | Class II | Class II | Class II | Class II | Class I |
| Span Calibration Points (g) | 500g, 1000g, 1500g, 2000g | 1000g, 2000g, 3000g, 4000g | 2000g, 3000g, 4000g, 5000g , 6000g | 2000g, 4000g, 6000g, 8000g | 6000g, 8000g, 10000g, 12000g |
| Weighing units* | 19 weighing units: g, ct, N, oz, ozt, GN, dwt, mo, msg, tl H, tl S, tl T, tcl, tola, baht, lb, kg Customer 1, Customer 2, | | | | |
| Weighing units* (EXP...M.. models and EXP...N..models only) | EXP...M.. models: g, mg, ct, kg EXP...N.. models: g, mg, ct, oz, ozt, GN, dwt, lb, kg | | | | |
| Applications | 15 application modes: Basic Weighing, Parts Counting, Check Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Totalization, Formulation, Differential, Density Determination, Peak Hold, Pipette Adjustment, SQC, Fill Weight Variation, Flow Rate Control | | | | |
| Stabilization time (typical) | ≤ 1 second | | | | |
| Sensitivity Temperature Drift (ppm/K) | 3 | | | | |
| Min-Weight (typical) (USP, K=2, U=0.10%) | 20 g | | | | |
| Min-Weight (optimal) (USP, K=2, U=0.10%, SRP≤0.41d) | 8.2 g | | | | |
| Terminal Display | 7-inch TFT 16.7M Color Press Screen with Glass Overlay, Detachable Terminal | | | | |
| Number of Dots | 800 RGB x 480 DOTS | | | | |
| Communication | USB Host (Type A) x 2, USB Device (Type B) x 1, USB Device (Type C) x 1 Ethernet (RJ45) x 1, RS232 x 1; Optional Wi-Fi, Bluetooth Dongle | | | | |
| Leveling System | Motorized Leveling System with Digital Leveling Bubble | | | | |
| Power Input | 12 VDC, 1.5A | | | | |
| Power Supply | AC Adapter Input: 100-240 VAC 0.5A 50-60 Hz AC Adapter Output: 12 VDC 1.5A | | | | |
| Weighing Pan Size | Trapezoid Pan 178 x 201 mm | | | | |
| Build-in Ionizer | Standard Equipped | | | | |
| Environmental Sensor RH, Air pressure Vibration Sensor | Temperature Sensor, Relative Humidity Sensor (RH), Air Pressure Sensor, Vibration Sensor | | | | |
| IR Sensor | 2 Touchless Sensors on the terminal | | | | |
| Terminal Status Lights | Standard Equipped | | | | |
| Terminal Housing Dimensions (W x D x H) | 203mm x 124mm x 65mm | | | | |
| Base Housing Dimensions (W x D x H) | | | | | |
| Shipping Dimensions (W x D x H) | | | | | |
| Net Weight | | | | | |
| Shipping Weight | | | | | |

| MODEL | EXP6201 | EXP8201 | EXP10201 |
|---|--|----------------------------|-----------------------------|
| | EXP6201M* | EXP8201M* | EXP10201M* |
| | EXP6201N* | EXP8201N* | EXP10201N* |
| Capacity (g) | 6200 g | 8200 g | 10200 g |
| Readability d, full range (g) | 0.1 g | 0.1 g | 0.1 g |
| Repeatability (sd.), ≤5% of full load | 0.07 g | 0.07 g | 0.07 g |
| Repeatability (sd.), 5% of full load to fine range max | 0.1 g | 0.1 g | 0.1 g |
| Linearity Deviation Typical | ± 0.06 g | ± 0.06 g | ± 0.06 g |
| Linearity Deviation | ± 0.2 g | ± 0.2 g | ± 0.2 g |
| Verification interval, e (EXP...M..models and EXP...N..models only) | 0.1 g | 0.1 g | 0.1 g |
| Approval Class (EXP...M.. models and EXP...N..models only) | Class II | Class II | Class I |
| Span Calibration Points (g) | 2000g, 3000g, 4000g, 5000g, 6000g | 2000g, 4000g, 6000g, 8000g | 2500g, 5000g, 7500g, 10000g |
| Weighing units* | 19 weighing units: g, ct, N, oz, ozt, GN, dwt, mo, msg, tl H, tl S, tl T, tcl, tola, baht, lb, kg, Customer 1, Customer 2 | | |
| Weighing units* (EXP...M.. models and EXP...N..models only) | EXP...M.. models: g, mg, ct, kg EXP...N.. models: g, mg, ct, oz, ozt, GN, dwt, lb, kg | | |
| Applications | 15 application modes: Basic Weighing, Parts Counting, Check Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Totalization, Formulation, Differential, Density Determination, Peak Hold, Pipette Adjustment, SQC, Fill Weight Variation, Flow Rate Control | | |
| Stabilization time (typical) | ≤ 1 second | | |
| Sensitivity Temperature Drift (ppm/K) | 5 | 3 | 3 |
| Min-Weight (typical) (USP, K=2, U=0.10%) | 160 g | | |
| Min-Weight (optimal) (USP, K=2, U=0.10%, SRP≤0.41d) | 82 g | | |
| Terminal Display | 7-inch TFT 16.7M Color Press Screen with Glass Overlay, Detachable Terminal | | |
| Number of Dots | 800 RGB x 480 DOTS | | |
| Communication | USB Host (Type A) x 2, USB Device (Type B) x 1, USB Device (Type C) x 1 Ethernet (RJ45) x 1, RS232 x 1; Optional Wi-Fi, Bluetooth Dongle | | |
| Leveling System | Motorized Leveling System with Digital Leveling Bubble | | |
| Power Input | 12 VDC, 1.5A | | |
| Power Supply | AC Adapter Input: 100-240 VAC 0.5A 50-60 Hz AC Adapter Output: 12 VDC 1.5A | | |
| Weighing Pan Size | Trapezoid Pan 178 x 201 mm | | |
| Environmental Sensor RH, Air pressure Vibration Sensor | Temperature Sensor, Relative Humidity Sensor (RH), Air Pressure Sensor, Vibration Sensor | | |
| IR Sensor | 2 Touchless Sensors on the terminal | | |
| Terminal Status Lights | Standard Equipped | | |
| Terminal Housing Dimensions (W x D x H) | 203mm x 124mm x 65mm | | |
| Base Housing Dimensions (W x D x H) | | | |
| Shipping Dimensions (W x D x H) | | | |
| Net Weight | | | |

| MODEL | EXP24001 | EXP35001 | EXP65001 |
|--|--|---------------------------------------|---------------------------------------|
| | EXP24001M* | EXP35001M* | EXP65001M* |
| | EXP24001N* | EXP35001N* | EXP65001N* |
| Capacity (g) | 24000 | 35000 | 65000 |
| Readability d, full range (g) | 0.1 g | 0.1 g | 0.1 g |
| Repeatability (sd.), ≤5% of full load | 0.08 g | 0.08 g | 0.08 g |
| Repeatability (sd.), 5% of full load to fine range max | 0.1 g | 0.1 g | 0.1 g |
| Linearity deviation | ± 0.07 g | ± 0.07 g | ± 0.07 g |
| Linearity deviation | ± 0.2 g | ± 0.2 g | ± 0.2 g |
| Verification interval, e (EXP...M models and EXP...N..models only) | 1 g | 1 g | 1 g |
| Approval Class (EXP...M.. models and EXP..N..models only) | Class II | Class II | Class II |
| Span Calibration Points (g) | 10000g, 15000g, 20000g, 24000g | 10000g, 20000g, 30000g, 35000g | 20000g, 40000g, 60000g, 65000g |
| Weighing units* | 19 weighing units: g, ct, N, oz, ozt, GN, dwt, mo, msg, tl H, tl S, tl T, tcl, tola, baht, lb, kg, Customer 1, Customer 2 | | |
| Weighing units* (EXP...M.. models and EXP..N..models only) | EXP...M.. models: g, mg, ct, kg EXP...N.. models: g, mg, ct, oz, ozt, GN, dwt, lb, kg | | |
| Applications | 14 application modes: Basic Weighing, Parts Counting, Check Counting, Percent Weighing, Check Weighing, Dynamic Weighing, Totalization, Formulation, Differential, Density Determination, Peak Hold, SQC, Fill Weight Variation, Flow Rate Control | | |
| Stabilization time (typical) | ≤ 1 second | | |
| Sensitivity Temperature Drift (ppm/K) | 5 | | |
| Min-Weight (typical) (USP, K=2, U=0.10%) | 200 g | | |
| Min-Weight (optimal) (USP, K=2, U=0.10%, SRP≤0.41d) | 82 g | | |
| Terminal Display | 7-inch TFT 16.7M Color Press Screen with Glass Overlay, Detachable Terminal | | |
| Number of Dots | 800 RGB x 480 DOTS | | |
| Communication | USB Host (Type A) x 2, USB Device (Type B) x 1, USB Device (Type C) x 1 Ethernet (RJ45) x 1, RS232 x 1; Optional Wi-Fi, Bluetooth Dongle | | |
| Leveling System | 4 leveling feet with Digital Leveling Bubble | | |
| Power Input | 12 VDC, 1.5A | | |
| Power Supply | AC Adapter Input: 100-240 VAC 0.5A 50-60 Hz AC Adapter Output: 12 VDC 1.5A | | |
| Weighing Pan Size | Square Pan 377 mm x 311 mm | | |
| Environmental Sensor RH, Air pressure Vibration Sensor | Temperature Sensor, Relative Humidity Sensor (RH), Air Pressure Sensor, Vibration Sensor | | |
| IR Sensor | 2 Touchless Sensors on the terminal | | |
| Terminal Status Lights | Standard Equipped | | |
| Terminal Housing Dimensions (W x D x H) | | | |
| Base Housing Dimensions (W x D x H) | | | |
| Shipping Dimensions (W x D x H) | | | |
| Gross Weight | | | |
| Net Weight | | | |

Note 1: M = EC Type approved

N = NTEP certified and Measurement Canada approved

C1= Customer Unit 1

C2= Customer Unit 2

Note 2: Default Calibration weights shown in Bold

Note 3: * Availability is dependent on region.

11.3 Accessory Specifications

The LM842 USB Adapter (Dongle)

TECH SPEC



| | |
|------------------------|--|
| WIRELESS STANDARD | 802.11ac |
| BACKWARD COMPATIBILITY | 802.11n, 802.11g, 802.11b, 802.11a, 802.11n, 802.11g, 802.11b, 802.11a |
| FREQUENCY | 2.4 GHz and 5 GHz |
| ADAPTER TYPE | Host Controller Interface (HCI) |
| INTERFACES | USB |
| ANTENNA | 2 x Metal Frame Antennas, SMA Connector |
| ANTENNA OPTION | 1 x Metal Frame Antenna + 1 x SMA Connector |
| COMPATIBLE ANTENNAS | LM256 2dBi, LM255 1.5dBi, LM251 2dBi |
| DIMENSIONS | 32-37mm x 17.1mm x 94mm |
| OPERATING TEMPERATURE | -20°C to +85°C |
| BLUETOOTH TECHNOLOGY | Bluetooth Classic, Bluetooth Low Energy (LE) |
| COMPATIBILITY | Android, Linux, Win 10, Win 7 |
| CERTIFICATIONS | |
| COMPLIANCE | |

11.4 Drawings and Dimensions

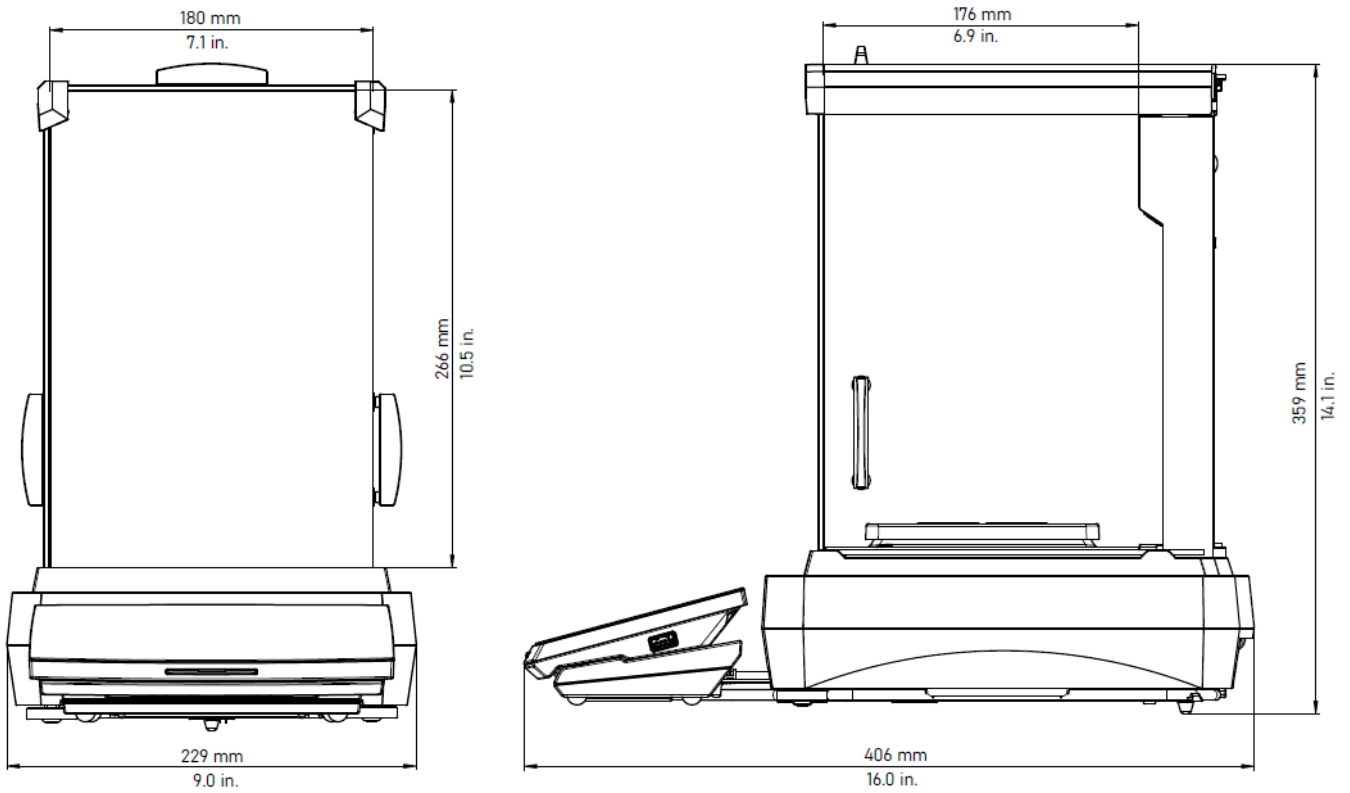


Figure 9-1. Draft Shield models

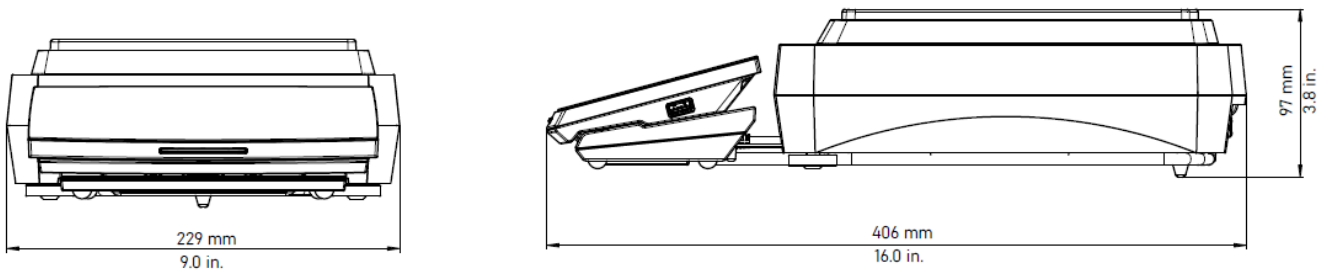


Figure 9-2. Non-Draft Shield models

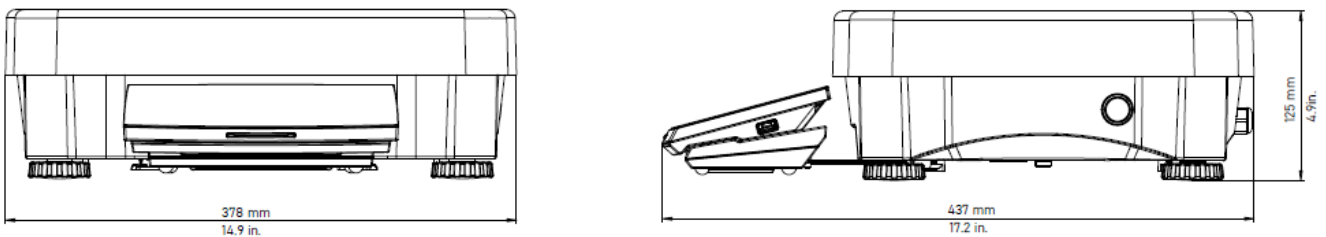


Figure 9-3. EXP24001, EXP35001 and EXP65001 models

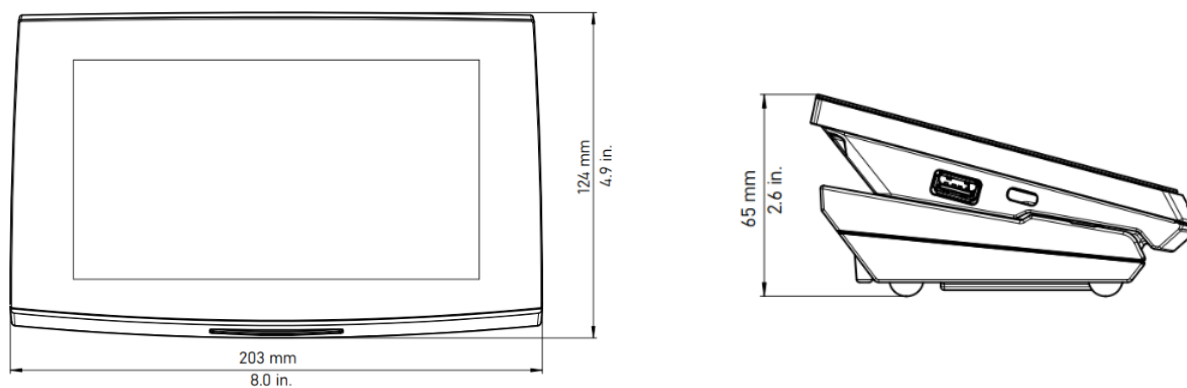











Figure 9-4. EXP Terminal

11.5 Accessories

| Image | P.N | Accessory Name |
|---|----------|--|
|  | 30095929 | ION-100A - Ionizer |
|  | 80253384 | Density Kit for Solids |
|  | 83034024 | Sinker Glass for Density Determination |
|  | 31059237 | Weighing Kit |

| | | |
|---|-----------------|--|
|  | <p>31059238</p> | <p>Fingerprint Module FIN-100</p> |
|  | <p>30252145</p> | <p>Pipette Evaporation Trap</p> |
| | <p>31059239</p> | <p>BT dongle, Wi-Fi dongle LM832</p> |
|  | <p>30064202</p> | <p>Impact Printer SF40A</p> |
|  | <p>30944359</p> | <p>Bluetooth Impact Printer SF40A/BT</p> |
| | <p>12120799</p> | <p>SF40A Paper roll (57,5mm 2pcs)</p> |
| | <p>12120798</p> | <p>SF40A Ink ribbon cassette</p> |
|  | <p>30808539</p> | <p>Tower Kit for High-Capacity Model</p> |
| | <p>30078078</p> | <p>Terminal extension cable (9 m)</p> |

11.6 Interface Commands

Commands listed in the following table will be acknowledged by the balance.

- Commands sent to the indicator must be terminated by a carriage return-line feed (CRLF).
- Data output is always terminated with a carriage return-line feed (CRLF).
- The balance will return “ES” for invalid commands.
- There is 40-second timeout control for print under stable requirement. If the unstable condition continues over 40 seconds, balance will respond “ES”.

| Command | Function |
|---------|--|
| SI | Print weighing value without weighing unit |
| IP | Immediate Print of displayed weight (stable or unstable). Attention: for EXP...N...models, when APPROVED MODE IS ON ON, IP could only print stable displayed weight. |
| P | Print displayed weight according to “Stable only” setting in the communication menu. Attention: when APPROVED MODE IS ON ON, P could only print stable displayed weight. |
| CP | Continuous Print. For EXP...N...models, when APPROVED MODE IS ON ON, CP could not work. |
| SP | Print on Stability. |
| P | Interval Print x = Print Interval (1-3600 sec) OP ends interval Print. Attention: the corresponding settings in the communication menu are modified too. |
| Z | Same as pressing Zero Key |
| T | Same as pressing Tare Key. |
| M | Print current application mode |
| U | Print Current weighing unit: g, Kg, lb, oz, etc... |
| ON | Brings out of Standby |
| OFF | Goes to Standby. |
| C3 | Begin internal Calibration, same as trigger from calibration menu. |
| WI | Left door open, right door closed. |
| PSN | Print Serial Number. |
| PV | Print terminal software version, base software version and Approved Mode is set to ON. |
| # | Set Counting APW (x) in grams. (must have APW stored) |
| % | Set Percent application reference weight (x) in grams. (must have reference weight stored) |
| CO | Set Check weighing Over Limit in grams x. |





| | |
|-----|--|
| CU | Set Check weighing Under Limit in grams x. |
| CU | Print Check weighing Under Limit. |
| TIM | Print current time. |
| DAT | Print current date. |
| TIM | Set Time, x format: hh:mm:ss. |
| DAT | Set Date, x format: mm/dd/yyyy. |

Note:

There is 40-second timeout control for print under stable requirement. If the unstable condition continues over 40 seconds, balance will respond "ES".

12 COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

| Mark | Standard |
|---|---|
|  | This product complies with the applicable harmonized standards of EU Directives 2011/65/EU (RoHS), 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at www.ohaus.com/ce . |
|  | This product complies with the EU Directive 2012/19/EU (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee . |
|  | EN 61326-1 |
|  | CAN/CSA-C22.2 No. 61010-1 UL 61010-1 |
| LM842 USB Dongle | Compliant to; IEEE 802.11ac, abgn additional standards, Compliant to; Bluetooth® 5.02 and backward compatible with Bluetooth v2.1+EDR / v3.0 / v3.0+HS / v4.0, 4.1, & 4.2 with both BR/EDR. Classic and LE can operate simultaneously. The Realtek IC, RTL8822CU uses a USB main interface. LM842 offers high throughput for both WiFi and Bluetooth® connections, connected via a (HCI) USB 2.0 TYPE A interface |

Important notice for Explorer Plus...M verified weighing instruments in the EU.

When the instrument is used in trade or a legally controlled application it must be set up, verified and sealed in accordance with local weights and measures regulations. It is the responsibility of the purchaser to ensure that all pertinent legal requirements are met.

Weighing Instruments verified at the place of manufacture must bear one of the following supplementary metrology marking on the descriptive plate.



Weighing Instruments to be verified in two stages have no supplementary metrology marking on the descriptive plate. The second stage of conformity assessment must be carried out by the applicable weights and measures authorities.

If national regulations limit the validity period of the verification, the user of the weighing instrument must strictly observe the re-verification period and inform the weights and measures authorities.

As verification requirements vary by jurisdiction, the purchaser should contact their local weights and measures office if they are not familiar with the requirements.

Industry Canada Note

CAN ICES-003(A) / NMB-003(A)

ISO 9001 Registration

The management system governing the production of this product is ISO 9001 certified.

Notes:

All the icons used for terminal design are sourced from a free platform: (<https://icons8.com/icons>).

Note: All the sound used for volume design are sourced from a free platform: (<https://pixabay.com/sound-effects>).

FCC Supplier Declaration of Conformity

Unintentional Radiator per 47CFR Part B

Trade Name: OHAUS CORPORATION

Model: Explorer Plus™ EXP...

Party issuing Supplier's Declaration of Conformity:

Ohaus Instruments (Changzhou) Co., Ltd.

Building C, No. 6 Zhengqiang Road, Xuejia Town, Xinbei District, Changzhou

Jiangsu 213022,

China

Phone: +86 519 85287270

Responsible Party – U.S. Contact Information:

Ohaus Corporation

8 Campus Drive, Suite 105

Parsippany, NJ 07054

United States

Phone: +1 973 377 9000

Web: www.ohaus.com

FCC Compliance Statement:

Note: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense. Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

13 LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or because of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.



Ohaus Corporation
8 Campus Drive, Suite 105
Parsippany, NJ 07054
United States
Phone: +1 973 377 9000
With offices worldwide.
Web: www.ohaus.com



* 3 1 0 6 2 8 6 5 *

P/N 31082865 A © 2025 Ohaus Corporation, all rights reserved.